

CATALOGUE PRODUCTS 1 1 2 0 - 1 0 0

SERIES OF REVERSE OSMOSIS APPARATUSES FOR PURE WATER PRODUCTION FROM DRINK WATER



ver. 004-21/05/2009-EN

IWA 20-100

IWA 20-100 ro, roi, roip, rot, roit, EDI

Apparatuses of raw IWA is designed for very pure water production from drink water by reverse osmosis with conducitivity below 1 mS/cm and producing 20 - 100 lph.

Apparatus contain four parts of treatment:

- mechanical filtration = removing mechanical particular parts from water
- **dechloration** = removing chlorine from waterodstranění volného chlóru
- **reverse osmotics modul** = removing anorganic ionts with efficiency 95 99%, organics parts up 99% and bacteriums and microorganisms
- **finish treatment** = by ionexchange resin value of conducitivty below 1 mS/cm (apparatus series roi, roip a roit)

or

= **electrodeionization** = A new modern technology of ultrapure water production. Under the action of electric field, ions are continuously transferred via ion exchanger grains to the corresponding ion-selective membranes. The key benefit of this technology is obtaining output water of a constant quality over a period of time (in EDI and EDI-T type apparatuses).

Priority apparatus:

- preatrement water before the reverse osmoses (mechanical filtrtion and dechloration
- reverse osmoses modul
- automatic operationin of connection with the level switch of water in tank
- function alloving of the exit of product after rearching of quality water
- periodic automatic rinsing to obstruct production of deposits and growing of microorganism
- cleaning making possible the process to clean
- alfanumeric display showing represent measuring vallues, autodiagnostic parameters and other function
- finish cleaning of product (ionexchange resin od EDI module)



SERIES OF REVERSE OSMOSIS
APPARATUSES FOR PURE
WATER PRODUCTION FROM
DRINK WATER



Principles of operation of individual device types

Demiwater production

Raw water flows in through catch pit, pressure-reducing valve (adjusted entrance pressure is cca 400 kPa) and elmag. valve into column of mechanica filtering [5a], removes with gentle filtering on filter sleeve water-insoluble impurities larger than 1µm. Flow through second section of filtration [5b] - dechlorination column - removes main part of organic compounds and active chlorine. Pump [6] forces water into RO module [7]. Osmotic modulus placed here removes cca 95% of ion type compounds from water.

Demiwater further:

- *type ro*: pressureless flow from RO module [7] to drawoff spot **B** and from here flows through tube straight for use $\underline{Product}$ = water with conductivity up to 50 μ S/cm. Apparatus capacity from 20 to 100 l/h.
- type rop: flows out from RO module [7] into pressure reservoir [21], then to drawoff spot B and from here flows through tube straight for use. <u>Product</u> = pressure outlet of water with conductivity up to 50 μS/cm. Apparatus capacity from 20 to 100 l/h.
- type roi: flows from RO module [7] into ion exchanger column [8], conductivity is further decreased here to values from 0,2 to 5 μS/cm. Demiwater flows out pressureless from ion exchanger column [8] through coarse filter, filter serves as catch pit of loose particles of ion exchange resin that could escape from column to drawoff spotB. From here flows through tube for use. Product = water with conductivity up to 5 μS/cm. Apparatus capacity from 20 to 100 l/h.
- type roip: flows from RO module [7] into ion exchanger column [8], conductivity is further decreased here to values from 0,2 to 5 μS/cm. Demiwater flows out from ion exchanger column [8] through coarse filter that serves as catch pit of ion exchange resin loose particles which could escape from column into pressure reservoir [21], further to drawoff spot B. From here flows through tube for use. <u>Product</u> = pressure outlet of water with conductivity up to 5 μS/cm. Apparatus capacity from 20 to 100 l/h.
- type rot: flows from RO module [7] into gravitational reservoir [30], further to drawoff spot B and from here flows through tube for use. Product = pressure outlet of water with conductivity up to 50 μS/cm. Apparatus capacity from 20 to 100 l/h.
- *type roit*: from RO module [7] flows into ion exchange column [8] conductivity is further decreased here to values from 0,2 to 5 μS/cm. Demiwater flows out from ion exchanger column [8] through coarse filter that serves as catch pit of ion exchange resin loose particles which could escape from column into gravitational reservoir β0], further to drawoff spot B. From here flows through tube straight for use.


<u>**Product**</u> = pressure outlet of water with conductivity of up to $5 \mu S/cm$. Apparatus capacity from 20 to 100 l/h.



• type **EDI**: is discharged from the RO module [7] in the absence of pressure through the flow sensor into the EDI electrodeionization unit for final treatment and conductivity reduction. From the EDI module demiwater is further discharged into drawoff spot **B** for immediate use.

<u>Product</u> = water of <u>constant quality</u> with conductivity of up to $5 \mu S/cm$. Apparatus capacity from 40 to 100 l/h.

Apparatus main parts (series of ro, roi, rop, roip)

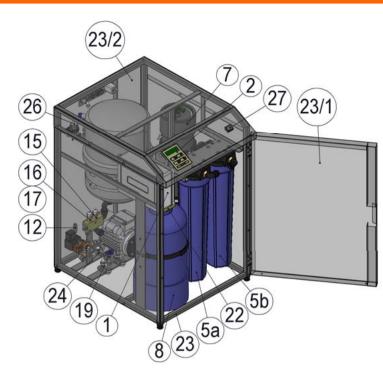

Connecting and drawoff spots of demiwater

- A inlet of pressure raw water into apparatus
- **B** outlet of demiwater
- **C** waste water


Apparatus main parts IWA ro, roi, rop, roip

- 1 control system
- 2 keyboard
- 3 catch pit
- 4 inlet elmag. valve
- 5a mechanical filtration
- 5b dechloration column
- 6 pump
- 7 RO module
- 8 ion exch. column
- 9 elmag. Valve of quick circulation
- 10 three-way plast elmag. valve
- 11 check valve
- 12 check valve
- 13 pressure sensor
- 14 pressure sensor
- 15 conductivity measuring probe of inlet water


16 conductivity measuring probe of demiwater

17 reading of demiwater temperature

- 18 pressure sensor
- 19 restrictor R
- 20 shutoff valve
- 21 pressure reservoir
- 22 ion exch.column
- 23 housing
- 23/1 doors
- 23/2 upper cover
- 24 check valve
- 25 control valve
- 26 coarse filter
- 27 main switch
- 29 pressure regulator

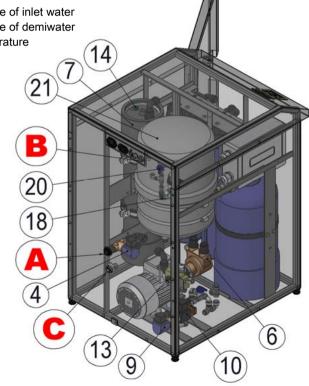

Apparatus main parts (series of rot, roit)

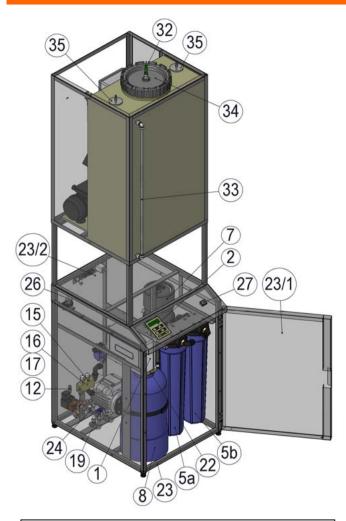


Connecting and drawoff spots of demiwater


- A inlet of pressure raw water into apparatus
- **B** outlet of demiwater
- **C** waste water

Apparatus main parts IWA rot, roit:

- 1 control system
- 2 keyboard
- 3 catch pit
- 4 inlet elmag. valve
- 5a mechanical filtration
- 5b dechloration column
- 6 pump
- 7 RO module
- 8 ion exch. column
- 9 elmag. Valve of quick circulation
- 10 three-way plast elmag. valve
- 11 check valve
- 12 check valve
- 13 pressure sensor
- 14 pressure sensor
- 15 conductivity measuring probe of inlet water
- 16 conductivity measuring probe of demiwater
- 17 reading of demiwater temperature
- 19 restrictor
- 23 housing
- 23/1 doors
- 23/2 upper cover
- 24 check valve ZK2
- 25 control valve V4
- 26 coarse filter
- 27 main switch
- 29 pressure regulator
- 30 reservoir 100 l
- 31 sensor min. level
- 32 sensor max, level
- 33 level gauge
- 34 closure reservoir
- 35 air microbial filter
- **36** pump with pressure tank
- 37 check valve
- 38 shutoff valve V9
- 39 shutoff valve V8
- 40 pressure switch of pump station
- 41 electrical instalation of tank

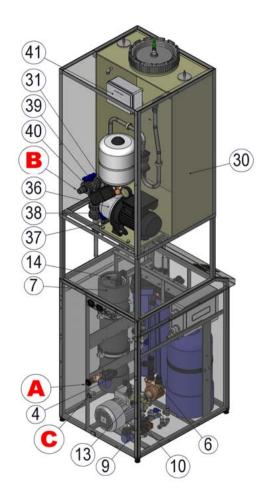

Apparatus main parts (series of EDI)

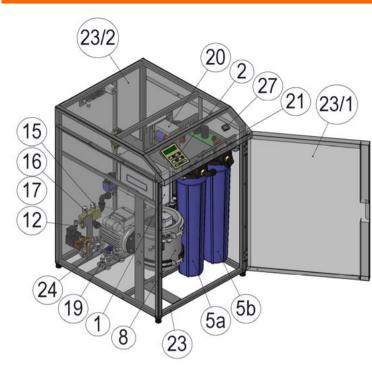

Connecting and drawoff spots of demiwater

- A inlet of pressure raw water into apparatus
- **B** outlet of demiwater
- C waste water

Apparatus main parts IWA XX EDI:

- 1 control system
- 2 keyboard
- 3 catch pit
- 4 inlet elmag. valve
- 5a mechanical filtration
- 5b dechloration column
- 6 pump
- 7 RO module
- 8 electrodeionization module EDI
- 9 elmag. valve of quick circulation
- 10 three-way plast elmag. valve
- 11 check valve
- 12 check valve
- 13 pressure sensor
- 14 pressure sensor
- 15 conductivity measuring probe of inlet water

16 conductivity measuring probe of demiwater


17 reading of demiwater temperature


- 18 restrictor
- 19 restrictor
- 20 EDI unit power supply
- 21 EDI unit control panel
- 22 flow sensor
- 23 housing
- 23/1 doors
- 23/2 upper cover
- 24 check valve ZK2
- 25 control valve V4
- 27 main switch
- 29 pressure regulator


This process utilizes exchanger resins under the action of electric field. The key benefit of this technology is the constant quality of output water over a period of time.

Compared to the conventional DI technology, EDI:


- nullifies the use of chemicals (as compared to DI resin regeneration)
- does not require shutdowns
- produces water of constant quality
- requires minimum energy
- is an economical option

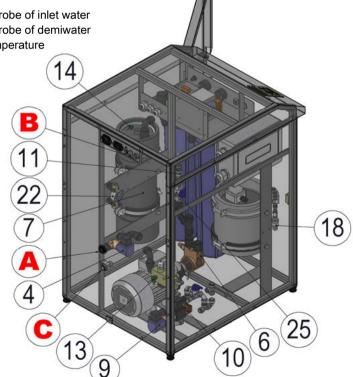

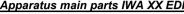

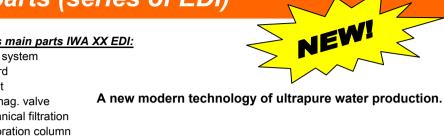

Apparatus main parts (series of EDI-T)

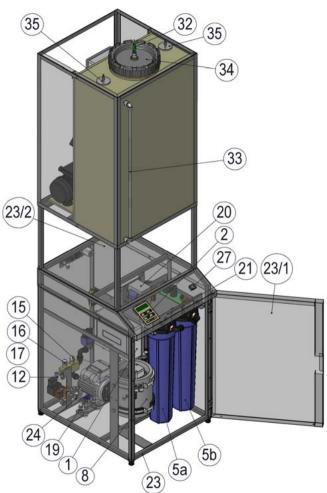

Connecting and drawoff spots of demiwater


- A inlet of pressure raw water into apparatus
- **B** outlet of demiwater
- C waste water

Apparatus main parts IWA XX EDI-T:

- 1 control system
- 2 keyboard
- 3 catch pit
- 4 inlet elmag. valve
- 5a mechanical filtration
- 5b dechloration column
- 6 pump
- 7 RO module
- 8 electrodeionization module EDI
- 9 elmag. Valve of quick circulation
- 10 three-way plast elmag, valve
- 11 check valve
- 12 check valve
- 13 pressure sensor
- 14 pressure sensor
- 15 conductivity measuring probe of inlet water
- **16** conductivity measuring probe of demiwater
- 17 reading of demiwater temperature
- 18 restrictor
- 19 restrictor
- 20 EDI unit power supply
- 21 EDI unit control panel
- 22 flow sensor
- 23 housing
- 23/1 doors
- 23/2 upper cover
- 24 check valve
- 25 control valve
- 27 main switch
- 29 pressure regulator
- 30 reservoir 100 l
- 31 sensor min. level
- 32 sensor max. level
- 33 level gauge
- **34** closure reservoir
- 35 air microbial filter
- 36 pump with pressure tank
- 37 check valve
- 38 shutoff valve V9
- 39 shutoff valve V8
- **40** pressure switch of pump station
- 41 electrical instalation of tank
- 42 valve



A new modern technology of ultrapure water production.

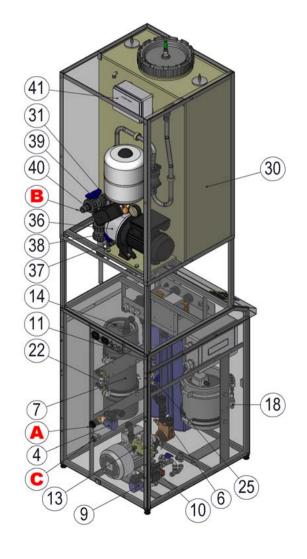


| Ordering Numbers | | | | | | | | | | | | | |
|------------------|------------|--------------|------------|--|---------------|------------|------|--|--|--|--|--|--|
| Apparatus | Order. No. | Apparatus | Order. No. | | Apparatus | Order. No. | | | | | | | |
| IWA 20 ro | D09010000 | IWA 20 roi | D09060000 | | IWA 40 EDI | D29310000 | | | | | | | |
| IWA 40 ro | D09020000 | IWA 40 roi | D09070000 | | IWA 60 EDI | D09320000 | | | | | | | |
| IWA 60 ro | D09030000 | IWA 60 roi | D09080000 | | IWA 80 EDI | D09330000 | | | | | | | |
| IWA 80 ro | D09040000 | IWA 80 roi | D09090000 | | IWA 100 EDI | D09340000 | | | | | | | |
| IWA 100 ro | D09050000 | IWA 100 roi | D09100000 | | | | EW! | | | | | | |
| | | | | | | | | | | | | | |
| IWA 20 rop | D09110000 | IWA 20 roip | D09160000 | | IWA 40 EDI-T | D09350000 | | | | | | | |
| IWA 40 rop | D09120000 | IWA 40 roip | D09170000 | | IWA 60 EDI-T | D09360000 | | | | | | | |
| IWA 60 rop | D09130000 | IWA 60 roip | D09180000 | | IWA 80 EDI-T | D09370000 | | | | | | | |
| IWA 80 rop | D09140000 | IWA 80 roip | D09190000 | | IWA 100 EDI-T | D09380000 | | | | | | | |
| IWA 100 rop | D09150000 | IWA 100 roip | D09200000 | | | ₹N | EWIS | | | | | | |
| | | | | | | | | | | | | | |
| IWA 20 rot | D09210000 | IWA 20 roit | D09260000 | | | | | | | | | | |
| IWA 40 rot | D09220000 | IWA 40 roit | D09270000 | | | | | | | | | | |
| IWA 60 rot | D09230000 | IWA 60 roit | D09280000 | | | | | | | | | | |
| IWA 80 rot | D09240000 | IWA 80 roit | D09290000 | | | | | | | | | | |
| IWA 100 rot | D09250000 | IWA 100 roit | D09300000 | | | | | | | | | | |


On special request can apparatuses be equipped with different demiwater reservoir with possibility to guard maximal level, if need be also with water house that is fed from apparatus and guarded minimal level in reservoir.

On special request can apparatuses be equipped with UV lamp and microbial filter at outlet. Service life of UV lamp is 8 000 hour run and its replacement is done by manufacturer or delegated establishment serviceman. Microbial filter is replaced when clogged, latest after one year of service.

Dimension sketch

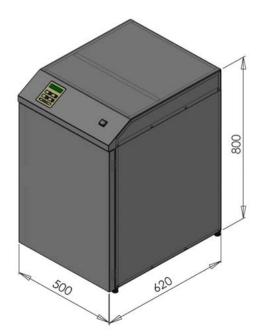

Apparatus series of type ro, roi, rop, roip, EDI

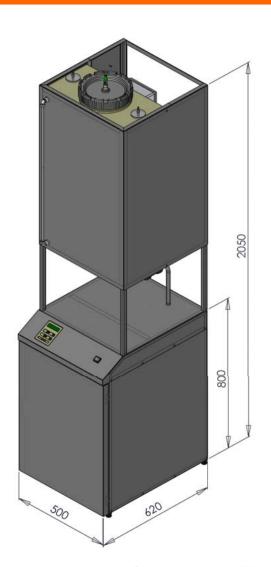

Apparatus series of type rot, roit, EDI-T

Spare Parts

| | | IWA | | | | | | | | | | | | | | | | | | | | | | | | | | | | |
|--------------------------------------|---------------|------|------|------|------|------|------|--------|------|---------|------|------|------------------|------|----|--------|-------|--------|-------|-------|-----|------|------|-----|-----|---------|------|---------|--------|------|
| Parts | Order. No. | ro | | | roi | | | rop | | | | roip | | | | rot | | | | roit | | | | EDI | | | | | | |
| | NO. | 20 | 40 | 60 | 80 | 100 | 20 | 40 | 60 E | 30 100 | 20 | 40 | 60 | 30 1 | 00 | 20 40 | 60 | 80 1 | 00 20 | 40 | 60 | 80 | 100 | 20 | 40 | 60 80 | 100 | 40 60 | 80 | 100 |
| Microprocesor system DEMI 2 | N00005017 | 1рс | 1рс | 1рс | 1рс | 1рс | 1рс | 1pc 1 | pc 1 | рс 1рс | 1pc | 1pc | 1pc 1 | pc 1 | рс | 1pc 1p | с 1ро | 1pc 1 | ос 1р | с 1рс | 1pc | 1pc | 1рс | 1рс | 1рс | 1pc 1pc | 1pc | 1pc 1pc | 1pc | 1рс |
| DEMI 2 - PIO | N00005017-PIC | 1рс | 1рс | 1рс | 1pc | 1pc | 1рс | 1pc 1 | pc 1 | рс 1рс | 1pc | 1рс | 1pc | pc 1 | рс | 1pc 1p | с 1ро | 1pc 1 | ос 1р | с 1рс | 1pc | 1pc | 1рс | 1рс | 1pc | 1pc 1pc | 1pc | 1pc 1pc | 1рс | 1pc |
| DEMI 2 - CU | N00005017-CU | 1рс | 1pc | 1рс | 1рс | 1pc | 1рс | 1pc 1 | рс 1 | рс 1рс | 1pc | 1рс | 1pc | pc 1 | рс | 1pc 1p | с 1ро | 1pc 1 | ос 1р | с 1рс | 1pc | 1рс | 1рс | 1рс | 1pc | 1pc 1pc | 1pc | 1pc 1pc | 1рс | 1pc |
| DEMI 2 - IAN | N00005017-IAN | 1pc | 1pc | 1рс | 1рс | 1pc | 1рс | 1pc 1 | рс 1 | рс 1рс | 1pc | 1рс | 1pc | pc 1 | рс | 1pc 1p | с 1ро | 1pc 1 | ос 1р | с 1рс | 1pc | 1рс | 1рс | 1рс | 1pc | 1pc 1pc | 1pc | 1pc 1pc | 1рс | 1pc |
| Mechanical filtration FCPS1-L | | 1рс | 1pc | 1рс | 1рс | 1pc | 1рс | 1pc 1 | рс 1 | рс 1рс | 1pc | 1рс | 1pc | pc 1 | рс | 1pc 1p | с 1ро | 1pc 1 | ос 1р | с 1рс | 1pc | 1рс | 1рс | 1рс | 1pc | 1pc 1pc | 1pc | 1pc 1pc | 1рс | 1pc |
| Dechloration FCCBL | D99440000 | 1рс | 1pc | 1рс | 1рс | 1pc | 1рс | 1pc 1 | рс 1 | рс 1рс | 1pc | 1рс | 1pc | pc 1 | рс | 1pc 1p | с 1ро | 1pc 1 | ос 1р | с 1рс | 1pc | 1рс | 1рс | 1рс | 1pc | 1pc 1pc | 1pc | 1pc 1pc | 1рс | 1pc |
| RO module FILMTEC XLE 2521 | NOXLE2521 | 1рс | - | - | - | - | 1рс | - | - | | 1pc | - | - | - | - | 1рс - | - | - | - 1p | с - | - | - | - | 1рс | - | | - | | - | - |
| Ro module HYDR. ESPA 1 4014 | | | 1рс | - | - | - | | 1рс | - | | | 1pc | - | - | - | 1р | с - | - | - | 1рс | - | - | - | | 1pc | | - | 1рс - | - | - |
| Ro module HYDR. ESPA 2 4021 | | | - | 1рс | - | - | | - 1 | рс | | | - | 1pc | - | - | - | 1pc | - | - | - | 1рс | - | - | | - | 1pc - | - | - 1pc | - | - |
| Ro module HYDR. ESPA 1 4021 | | | - | - | 1 pc | 1 pc | | - | - 1 | рс 1 рс | | - | - 1 | pc 1 | рс | - | - | 1 pc 1 | рс | - | - | 1 pc | 1 pc | | - | - 1 pc | 1 pc | - | 1 pc | 1 pc |
| Pressure switch KTS 618 | N00001588 | 2 pc 2 | pc 2 | pc 2pc | 3 pc | 3рс | 3pc 3 | рс 3 | рс | 3рс 3р | с Зро | 3pc 3 | oc 2p | с 2рс | 2рс | 2рс | 2рс | 2рс | 2рс | 2рс 2рс | 2рс | 2pc 2pc | 2pc | 2рс |
| Sensor SV1 | D99100000 | 1рс | 1рс | 1рс | 1рс | 1рс | 1рс | 1pc 1 | рс 1 | рс 1рс | 1pc | 1рс | 1pc | pc 1 | рс | 1pc 1p | с 1ро | 1pc 1 | ос 1р | с 1рс | 1рс | 1рс | 1рс | 1рс | 1pc | 1pc 1pc | 1рс | 1pc 1pc | 1рс | 1рс |
| Sensor SVT4 | D99130000 | 1рс | 1рс | 1рс | 1рс | 1рс | 1рс | 1pc 1 | рс 1 | рс 1рс | 1pc | 1рс | 1pc | pc 1 | рс | 1pc 1p | с 1ро | 1pc 1 | ос 1р | с 1рс | 1рс | 1рс | 1рс | 1рс | 1pc | 1pc 1pc | 1рс | 1pc 1pc | 1pc | 1рс |
| Solenoid valve CEME G3/8", typ 8323 | N00015220 | 2рс | 2рс | 2рс | 2рс | 2рс | 2рс | 2pc 2 | pc 2 | pc 2pc | 2pc | 2рс | 2pc 2 | pc 2 | рс | 2рс 2р | с 2рс | 2pc 2 | oc 2p | с 2рс | 2рс | 2рс | 2рс | 2рс | 2рс | 2рс 2рс | 2рс | 2pc 2pc | 2pc | 2рс |
| Three way valve SIRAI | NOD344F57 | 1рс | 1рс | 1рс | 1рс | 1рс | 1рс | 1pc 1 | рс 1 | рс 1рс | 1pc | 1рс | 1pc | pc 1 | рс | 1pc 1p | с 1ро | 1pc 1 | ос 1р | с 1рс | 1рс | 1рс | 1рс | 1рс | 1pc | 1pc 1pc | 1рс | 1pc 1pc | 1pc | 1рс |
| Pump PO404X | N00PO404X | 1рс | 1рс | - | - | - | 1рс | 1рс | - | | 1pc | 1рс | - | - | - | 1pc 1p | с - | - | - 1p | с 1рс | - | - | - | 1рс | 1pc | | - | 1рс - | - | - |
| Pump PO901 | | - | - | 1рс | 1рс | 1рс | - | - 1 | рс 1 | рс 1рс | - | - | 1pc ⁻ | pc 1 | рс | | 1pc | 1pc 1 | oc - | - | 1рс | 1рс | 1рс | - | - | 1pc 1pc | 1рс | - 1pc | 1рс | 1рс |
| MOTOR 1LF7073-4AB17-ZA31 | N00000989 | 1рс | 1рс | 1рс | 1рс | 1рс | 1рс | 1pc 1 | рс 1 | рс | 1pc | 1рс | 1pc | pc 1 | рс | 1pc 1p | с 1ро | 1pc | 1р | с 1рс | 1рс | 1рс | 1рс | 1рс | 1pc | 1pc 1pc | | 1pc 1pc | 1рс | 1рс |
| lonex module 0717 (8I) | | - | - | - | - | - | 1рс | 1pc 1 | рс 1 | рс | - | - | - | - | - | 1pc 1p | с 1ро | 1pc | - | - | - | - | - | 1рс | 1рс | 1pc 1pc | | 1pc 1pc | 1pc | 1рс |
| EDI module | | - | - | - | - | - | - | - | - | | - | - | - | - | - | | - | - | | - | - | - | - | - | - | | - | 1pc 1pc | 1pc | 1рс |
| Pressure tank | N00011412 | - | - | - | - | - | - | - | - | - | 1pc | 1рс | 1pc | pc 1 | рс | 1pc 1p | с 1ро | 1pc 1 | рс - | - | - | - | - | - | - | | - | - | - | - |
| Supplier water pump NP MCX 80/48 M | N00010029 | - | - | - | - | - | - | - | - | - | - | - | - | - | - | | - | - | 1р | с 1рс | 1pc | 1рс | 1рс | 1рс | 1рс | 1pc 1pc | 1pc | - | - | - |
| Air filter MILEX FG 50 | SLFG55010 | - | - | - | - | - | - | - | - | - | - | - | - | - | - | | - | - | 2р | с 2рс | 2рс | 2рс | 2рс | 2рс | 2рс | 2рс 2рс | 2ls | - | - | - |
| Sensor of level RC5-30200S-M0M5C | N01900116 | - | - | - | - | - | - | - | - | - | - | - | - | - | - | | - | - | 1р | с 1рс | 1pc | 1рс | 1рс | 1рс | 1рс | 1pc 1pc | 1pc | | - | - |
| Analog. sensor UB800-18GM40-I-V1 | | - | - | - | - | - | - | - | - | - | - | - | - | - | - | | - | - | 1р | с 1рс | 1pc | 1рс | 1рс | 1рс | 1рс | 1pc 1pc | 1pc | - | - | - |
| Check valve ZK5 | 0ACHV0404 | 1рс | 1рс | 1рс | 1рс | 1рс | 1рс | 1pc 1 | pc 1 | рс 1рс | 1pc | 1pc | 1pc ² | pc 1 | рс | 1pc 1p | с 1ро | 1pc 1 | ос 1р | с 1рс | 1рс | 1рс | 1рс | 1рс | 1рс | 1pc 1pc | 1pc | | - | - |
| Keyboard | 000030111 | 1рс | 1рс | 1рс | 1рс | 1рс | 1рс | 1pc 1 | pc 1 | рс 1рс | 1pc | 1pc | 1pc ² | pc 1 | рс | 1pc 1p | с 1ро | 1pc 1 | ос 1р | с 1рс | 1рс | 1рс | 1рс | 1рс | 1рс | 1pc 1pc | 1pc | 1pc 1pc | pc 1pc | 1pc |
| Microbiological filter SARTOBRAN 300 | 523130745 | | | | | | | | | | | | | | | | | | | | | | | | | | | | | |
| UV lamp GCL307T5VH | NTL205068 | | | | | | | | | | | | | | | | | | | | | | | | | | | | | |
| Power UV lampy | 000030099 | | | | | | | | | | | | | | | | | | | | | | | | | | | | | |



Ledeč nad Sázavou **ZAŘÍZENÍ PRO ÚPRAVU VOD**

Jiřího Wolkera 852 584 01 LEDEČ nad Sázavou

tel. +420569/720533, +420569/720534

<u>fax.</u> +420569/721099

info: www.watek.cz