# **MOMOLINE**

Technical furniture for laboratories

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**AK** is an automatic chemical fume cupboard designed and tested to ensure full protection to lab operators in every situation, even the most challenging and advanced.

Thanks to its design, **AK** finds a place in every modern and innovative laboratory environment adding value and efficiency.

**AK** is characterized by a light appearance, thanks to a careful location of the utilities below the worktop, allowing a full employment of the available space, optimizing the efficiency and the overall usability of the fume cupboard.

Both the lateral jambs of **AK** are realized using anti-acid HPL sheets, providing at the same time stiffness and lightness to its structure, making it different from the conventional profiles available on market for chemical fume cupboards.

In case of fume hood for aggressive acids, its structure is entirely realized in PVC or PP.

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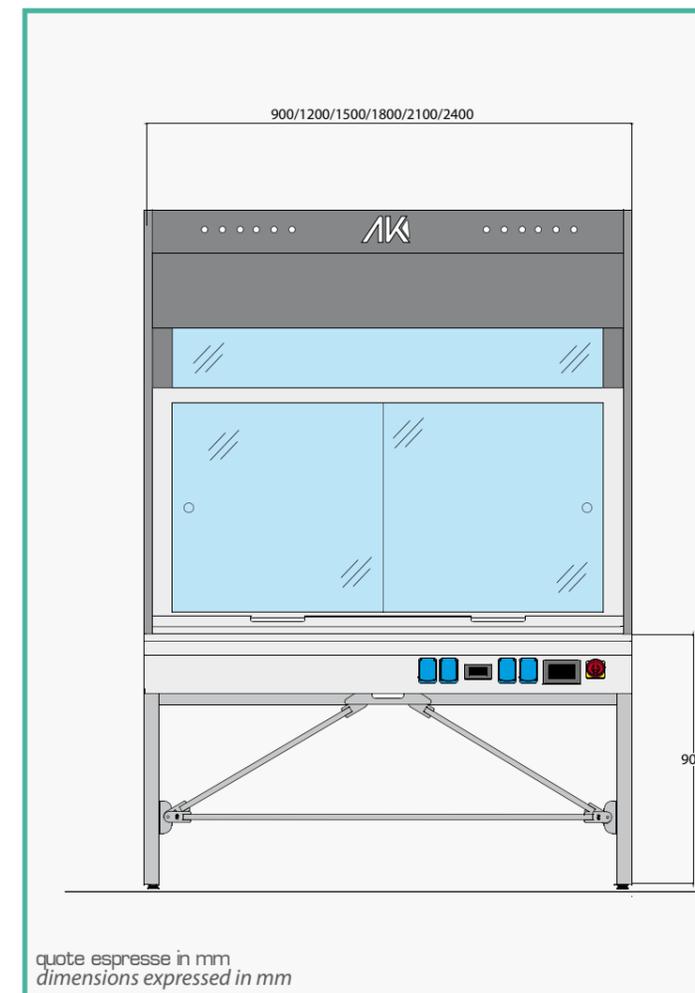
quote espresse in mm  
dimensions expressed in mm

The lateral windows, standard fitting for all models, are realized in safety glass and contribute to make the structure of **AK** very light and sustainable, providing a smooth and futuristic design, together with an optimal illumination of the internal working area.

**AK** fume cupboard has been certified by a third party for its compliance to the Norm EN 14175:2003, and the certificate has been issued by Internationally recognized third authorities, so to ensure the efficiency and safety Standards are always met.

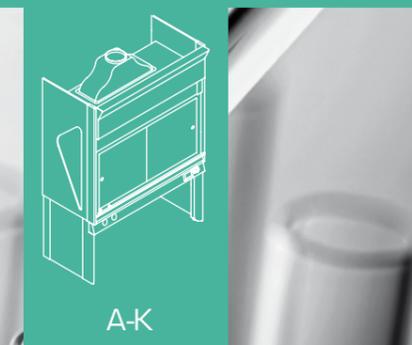
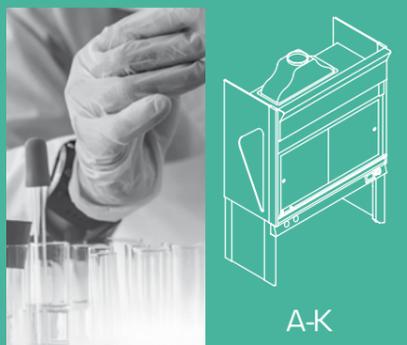
In order to obtain the EN 14175 Certification, **AK** has been tested in the most adverse and challenging conditions to grant full safety of the operators in all applications, proving superior features of containment and robustness of containment.

Together with the wide choice of materials for worktop, many different accessories and optional items are available: motorized sash, many automation optionals or employment of explosion proof components are only some examples.

**AK** fume cupboard is equipped as a standard with a highly technologic electronic control software, specifically designed for its management, and usable by the operator trough a very intuitive touch screen interface. The control software embeds a wide range of alarms and detecting probes, aimed to continuously monitor the working conditions of the fume cupboard, and ready to promptly intervene in case of necessity, so that operators are always safeguarded.

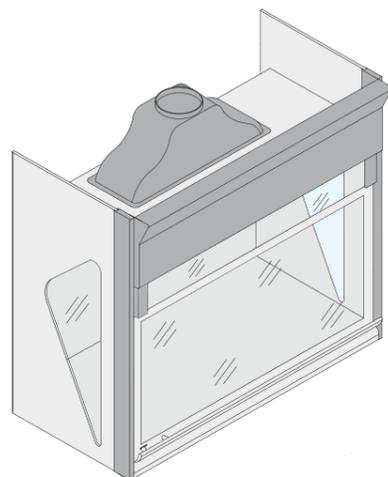
In addition to the standard software, fume cupboard can also embed some advanced airflow control features, designed to implement active safety and energy saving logics, such as the VAV, CLIMA and the TWIN SET systems.

Detailed specifications regarding all available accessories can be found in this Paragraph.

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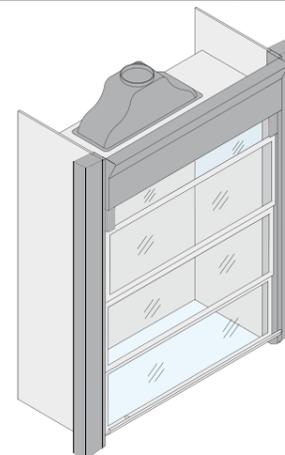
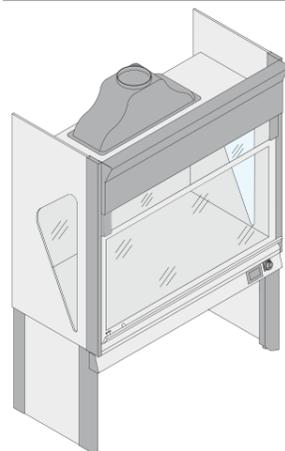
Equipements of **AK** walk-in can be positioned on side columns. Details at pag.162

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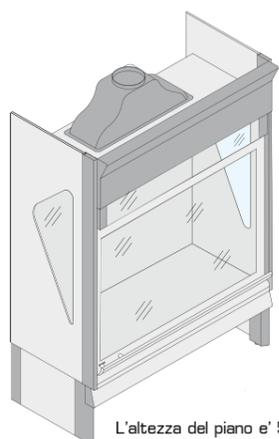
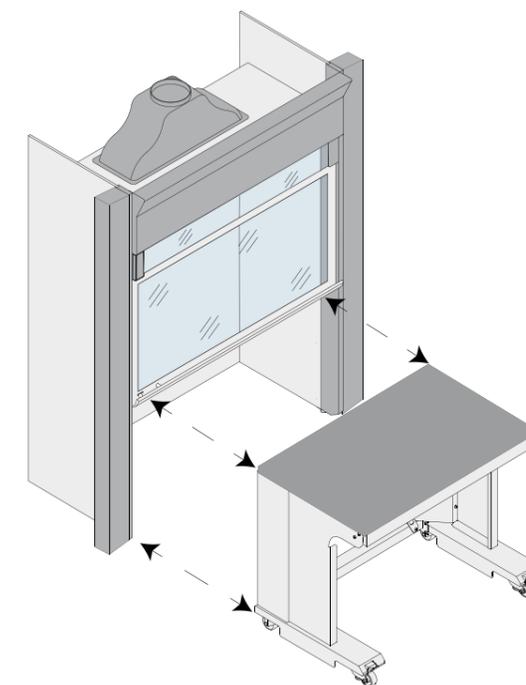
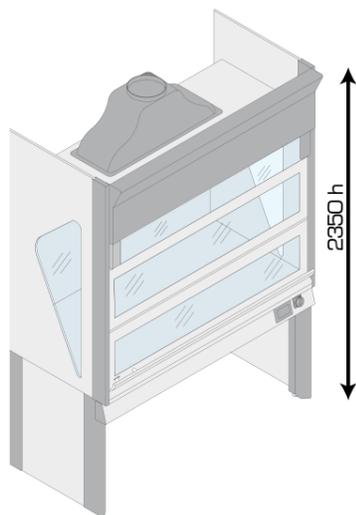
Walk-in version is not equipped with the conventional worktop to allow the insertion under the fume cupboard of tall equipment and trolley benches. Walk-in fume cupboards are often coupled to **Skylab** line trolley benches (code: BCAR), described in lab benches technical catalogue.

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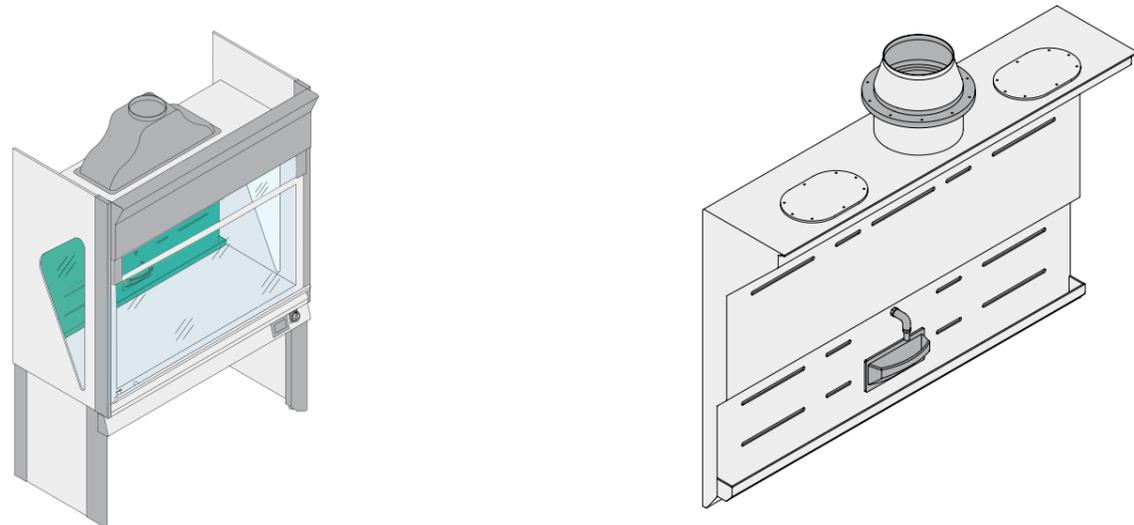
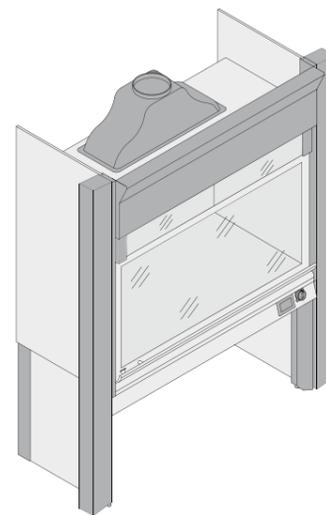
*Upon request it is possible to equip fume cupboards with a fume washing system embedded in the plenum, in the back of the hood.*

*This system consist in a device spraying water in counter-current to the exhaust air, which is sent through some swirlers to increase the interface surface. Water can be collected into a tank and recirculated or directly properly disposed.*

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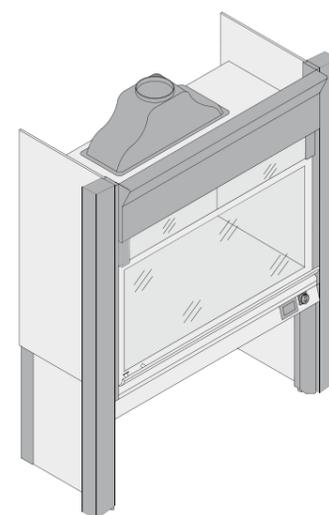
*The upper part of this hood is chemical-resistant and entirely realized in PVC. It is suitable for application where highly acidic chemicals are needed. Sash sliding mechanism and his anti-fall safety device are positioned outside the working chamber in order to be protected from acid attacks.*

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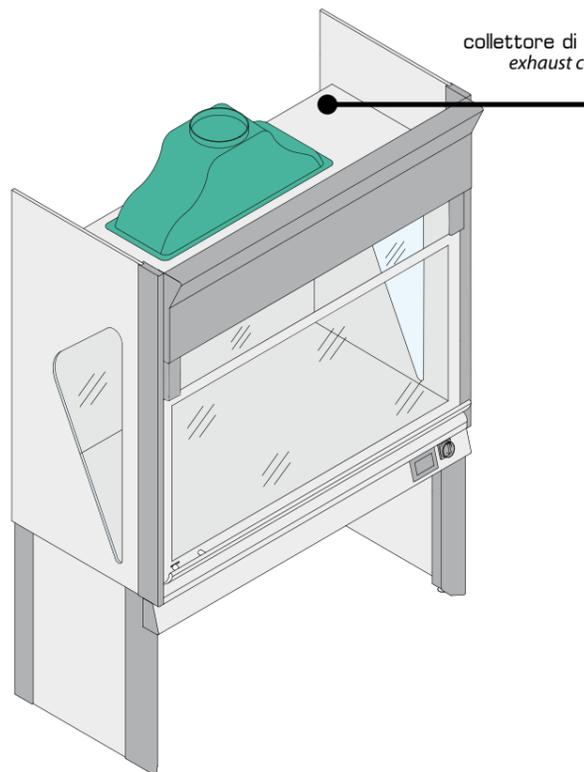
*The structure for use of highly acidic and thermal load differs from the structure for use of highly acidic for the presence of a ceramic lining inside the chamber.*

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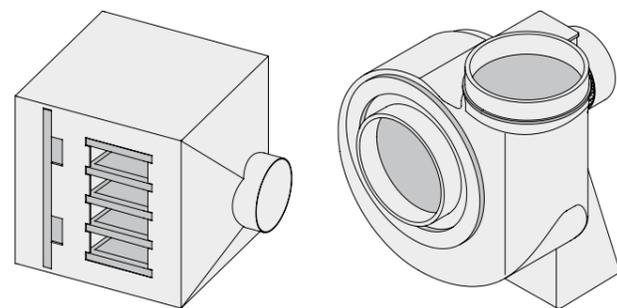
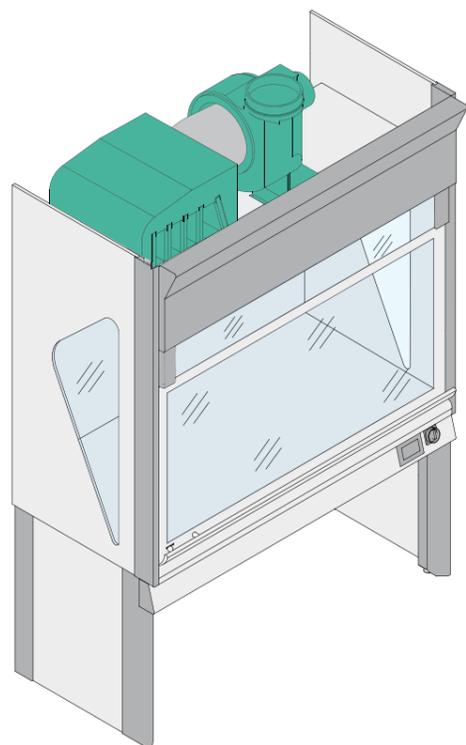
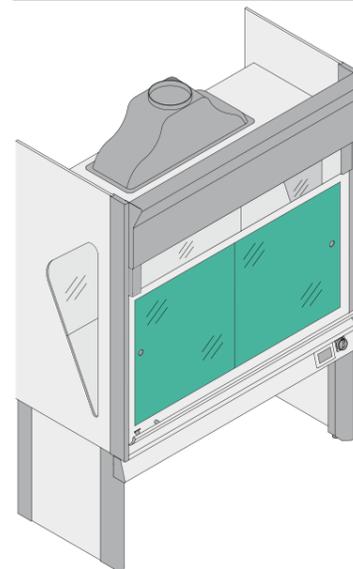
**AK** is equipped, as a standard, with an exhaust collector maximising its containment performances.  
If a filter box or an exhaust fan are provided, these are installed either on ceiling or on the roof.

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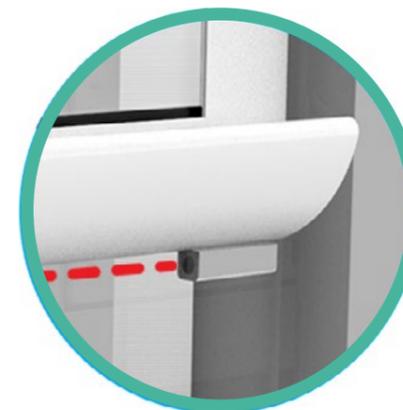
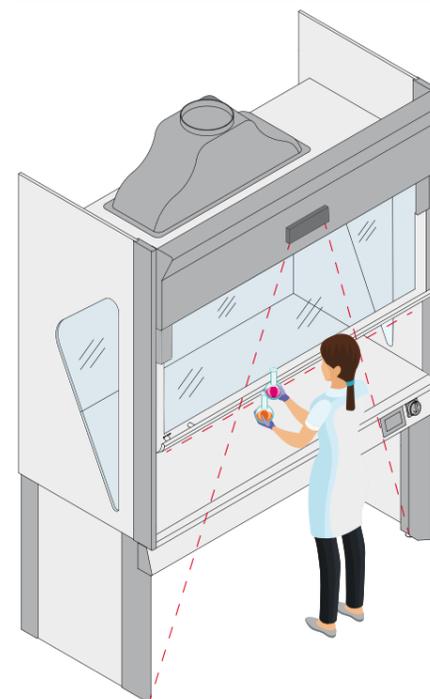
**AK** is always equipped with a vertically sliding frontal safety glass, framed into a counter-balanced sash.

Upon request, it is possible to provide a vertically sliding frame incorporating up to 4 glass panels, horizontally sliding.

Frontal sash is always equipped with a safety device granting its lock in case of failure of one or both its suspending chords, so that it is prevented from falling.

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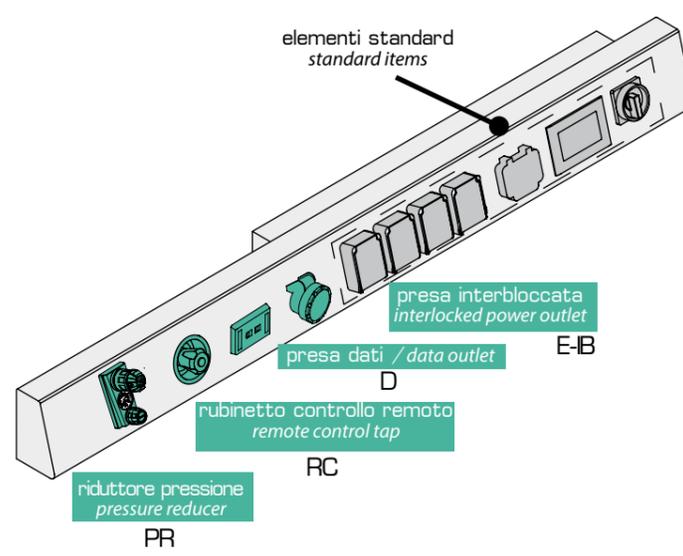
An electric engine can be provided to operate the movement of the frontal sash. The sash can be either operated manually (through dedicated buttons on the touch-screen panel) or automatically, through the operator's presence detection system. In case the operator's hands or any object is placed under the sash, the sensor automatically blocks the sash and raises it again. Even in case of motorized sash, the operator can always operate it manually.

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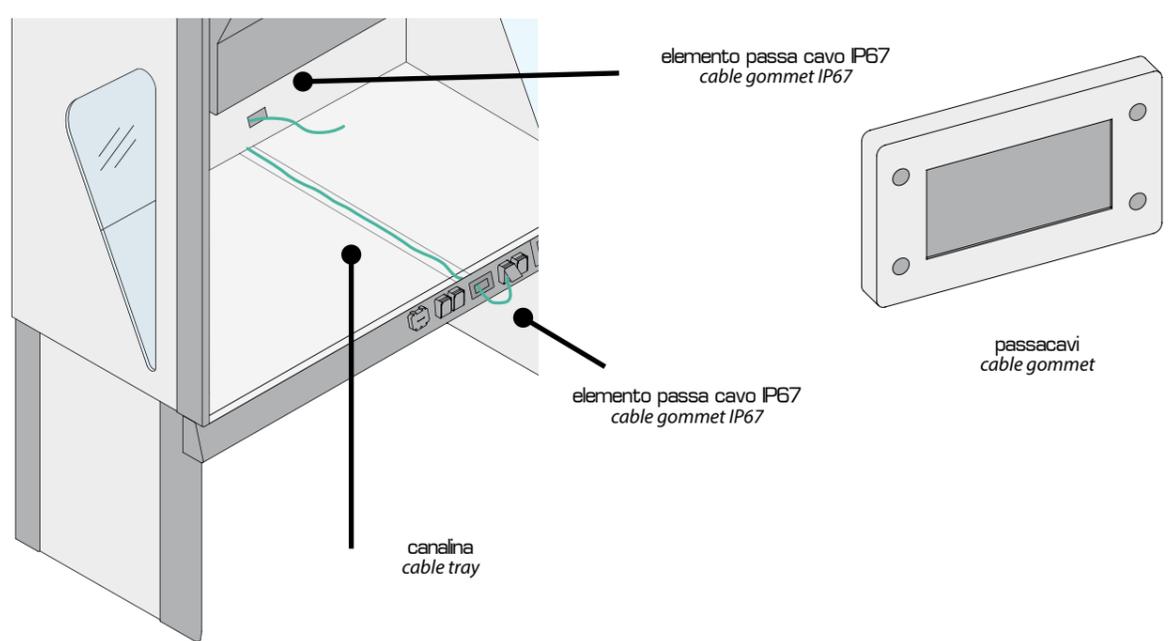
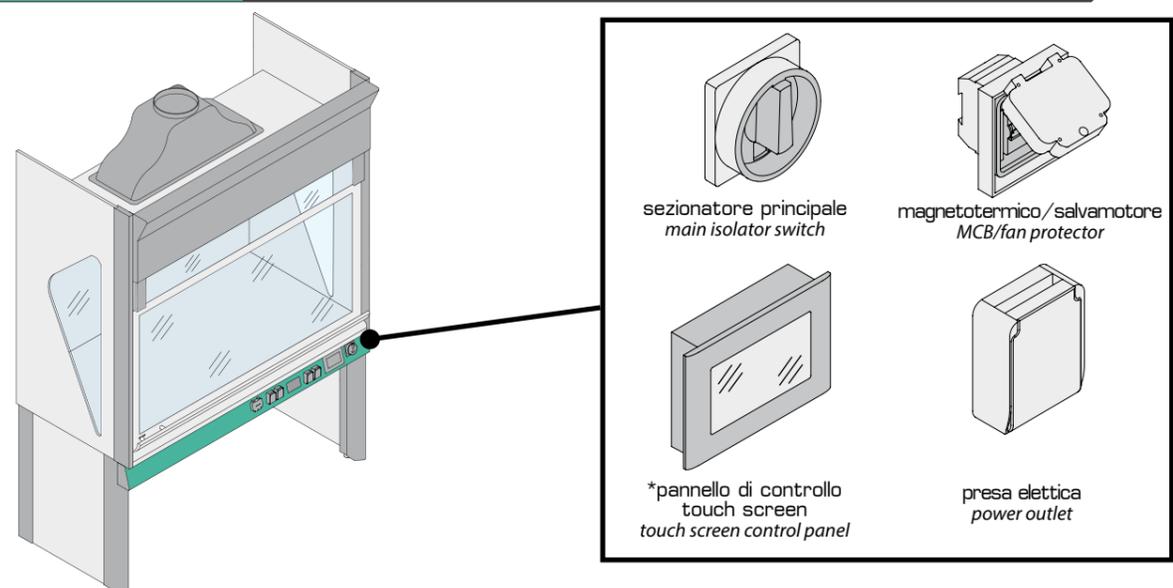
The frontal panel of **AK** fume cupboard can be equipped with many accessories, as shown above. In standard configuration, basic provisions are: master control switch for turning ON and OFF the fume cupboard, MCB, 4 power outlets and a cable grommet, to be used for supplying power to all instruments placed on the worktop, with no risk of interfering with the frontal sash movement or with the extraction performances. Right side element also incorporates a technical compartment under the worktop area.

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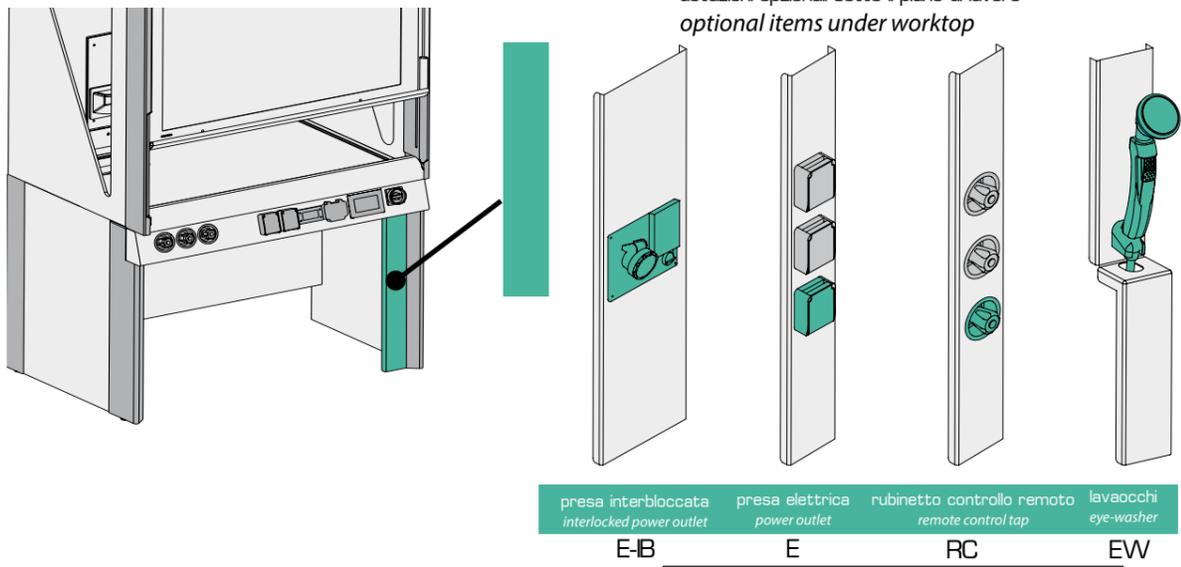
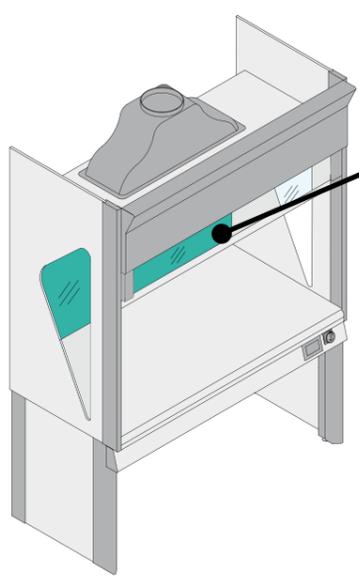
presa interbloccata  
interlocked power outlet  
E-B

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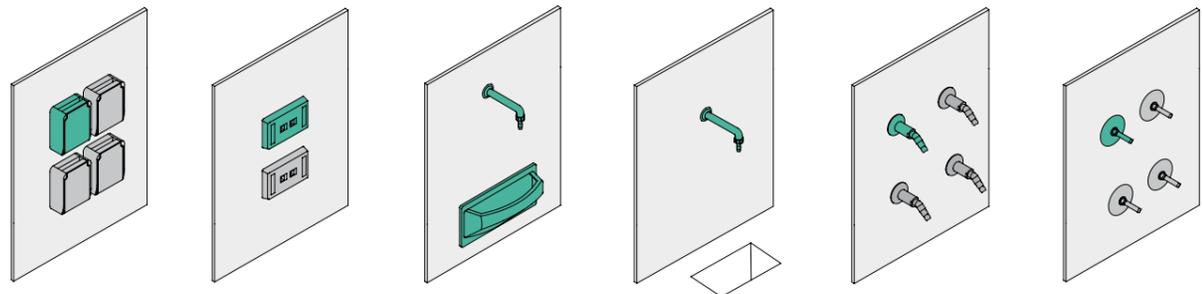
rubinetto controllo remoto  
remote control tap  
RC

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*The rear panel of **AK** is made up of modules that can be easy de-assembled. These panels can be equipped with various utilities (as below detailed) including water drip cup. If necessary, the rear panel can fit a rack, as detailed at pag. 198.*

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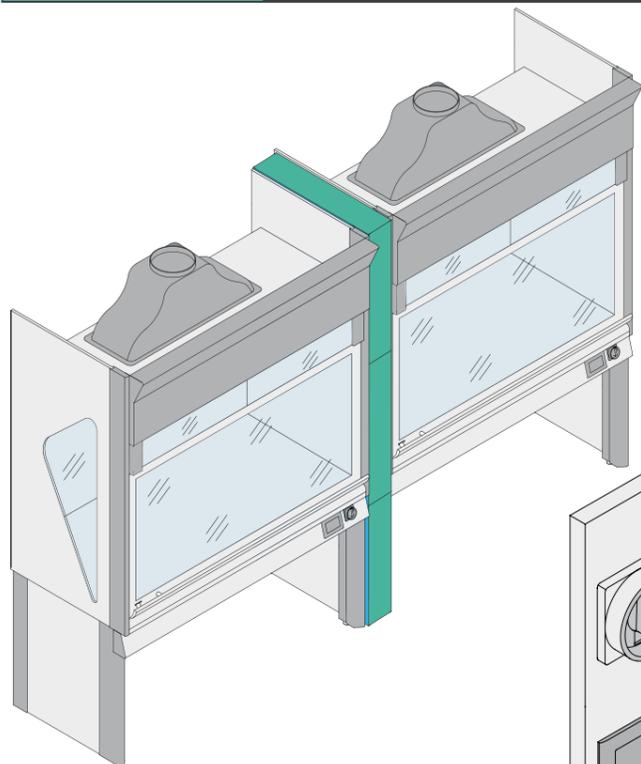
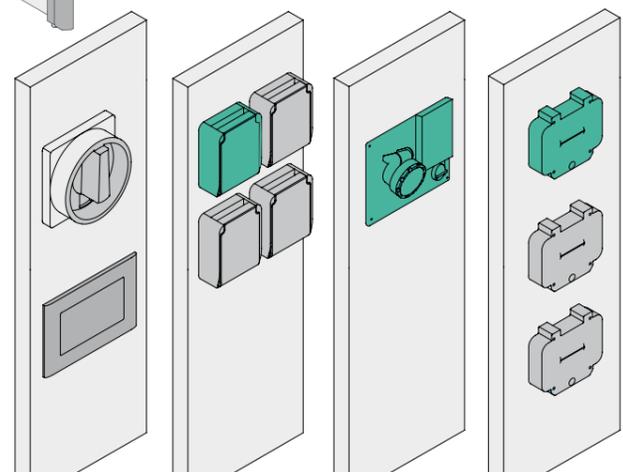
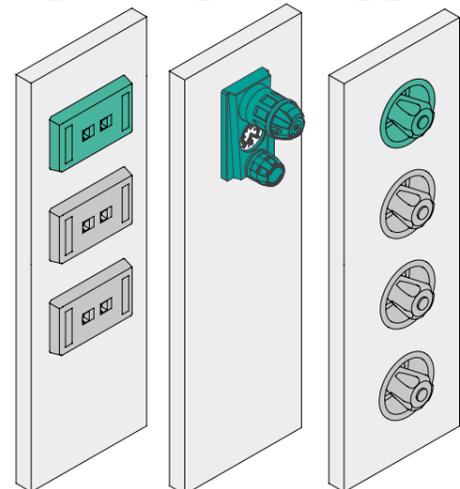
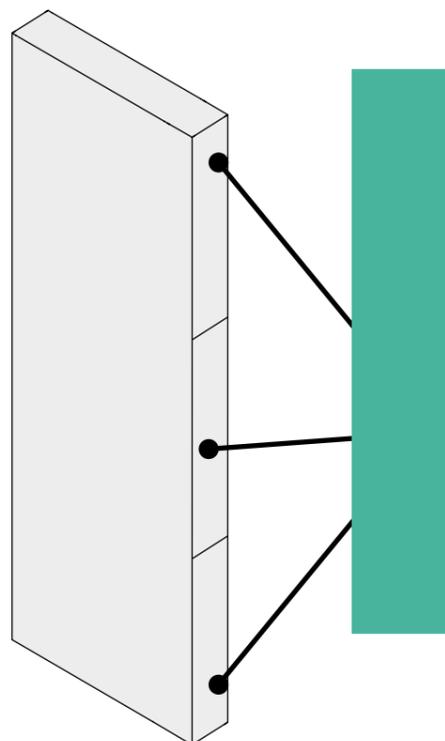
*In case two different fume cupboard, with inter-communicating working area are needed, a central supporting element is provided. Middle element can incorporate up to 3 customizable panels.*

pannello controllo control panel CP  
presa elettrica power outlet E  
presa interbloccata interlocked power outlet E-IB  
magnetotermico MCB M

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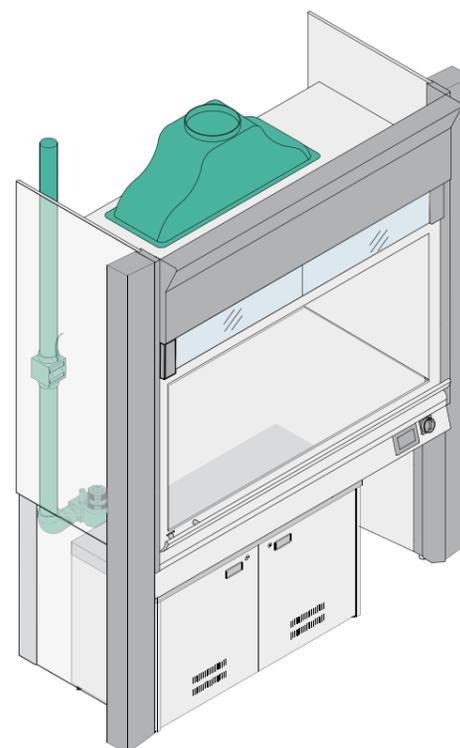
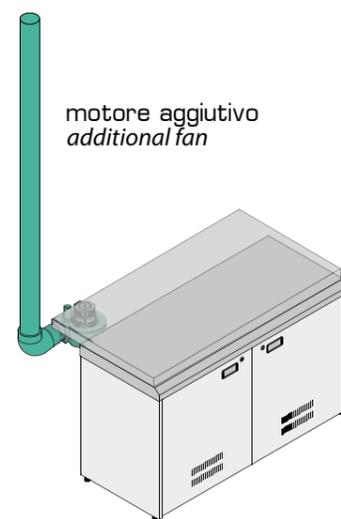
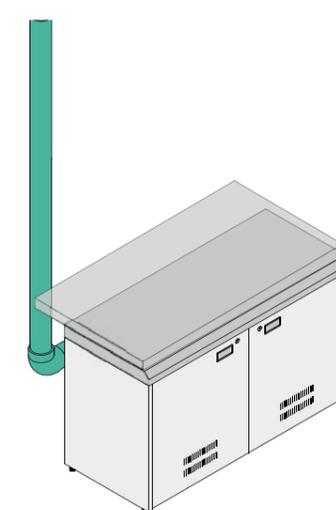
*All safety cabinets installed under the worksurface can be connected to the exhaust system. Exhaust can be either common to the fume cupboard or independent. Upon specific request, also a dedicated filter embedded in the structure of the fume cupboard can be provided for safety cabinets.*

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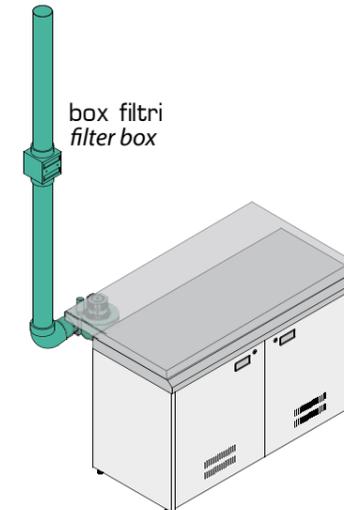
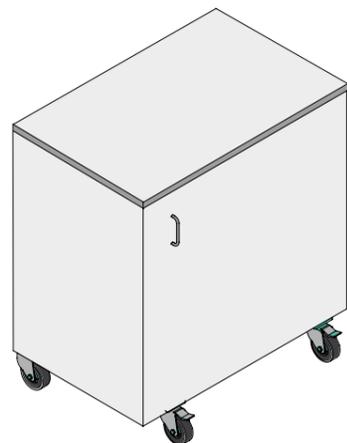
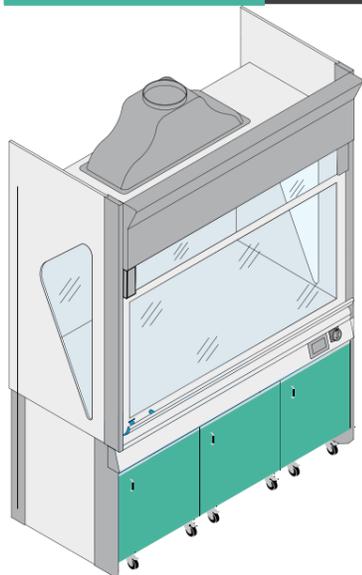
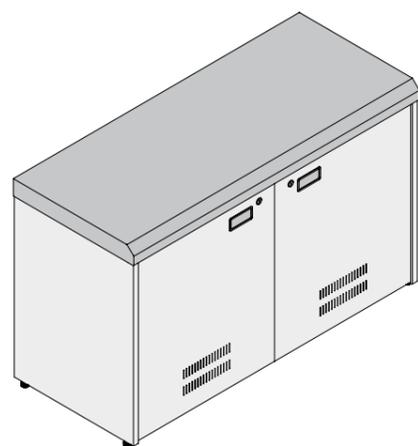
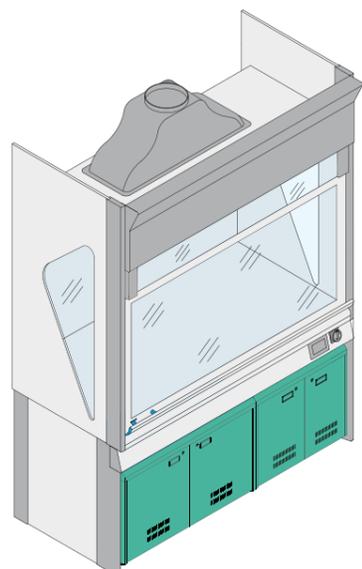
A wide range of storage units can be installed under the fume cupboard, for the user's convenience.

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In case of necessity a safety cabinet for chemicals storage can be installed under the fume cupboard. The safety cabinet can also be equipped with extraction system upon request. For details and specifications of safety cabinets please refer to chapter 4 "safety cabinets".

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If equipped with VAV (Variable Airflow Volume) fume cupboards embed a controller monitoring their face velocity in real time: once the face velocity has been set, VAV system keeps it constant with the variation of sash position. VAV system can be obtained either using an inverter (cod. VAV-I-AK) or using an air dumper (cod. VAV-V-AK). More details at pag. 123

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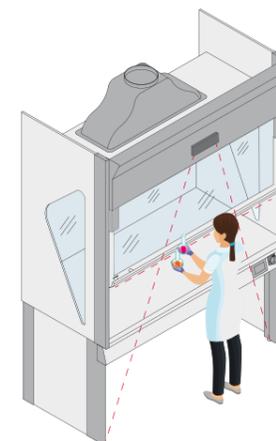
Upon specific request, the fume cupboard can be equipped with an automatic system for fire detection and suppression. The development of a fire inside the working area of the fume cupboard determines the sudden activation of the system (through a pressure rise in the probe duct) and the subsequent release of extinguishing agent through specific nozzles. The extinguishing agent is stored in a dedicated cylinder located outside of the fumehood and can be chosen according to the specific need (standard CO<sub>2</sub>)

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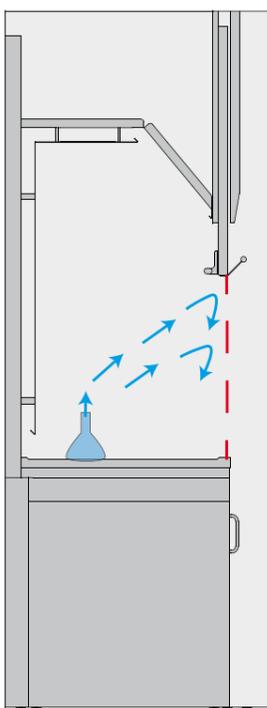
An electric engine can be provided to operate the movement of the frontal sash. The sash can be either operating manually (through dedicated buttons on the touch-screen panel) or automatically, through the operator presence detection system. In case the operator's hands or any object is placed under the sash, the sensor automatically blocks the sash and raises it again. Even in case of motorized sash, the operator can always operate it manually.

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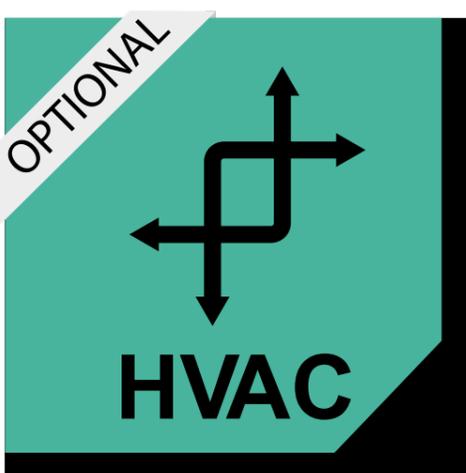
*TWIN SET is an exclusive device able to integrate energy savings to operator's safeguard.  
 If TWIN SET system is installed, in case of necessity the operator can decrease the face velocity, in order to save energy and to reduce the total airflow extracted from the room.  
 In this case, to keep the containment performances unvaried, one or more additional blowers automatically activate, so that an air barrier is created.  
 This barrier allows to completely shield the working environment under the fume cupboard, so that airborne contaminants are retained inside.*

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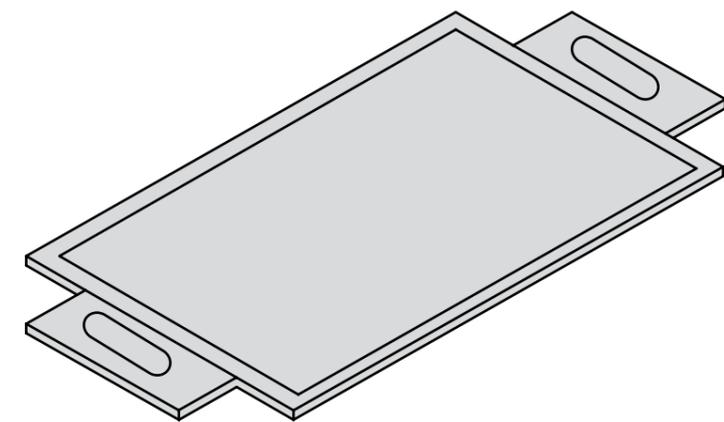
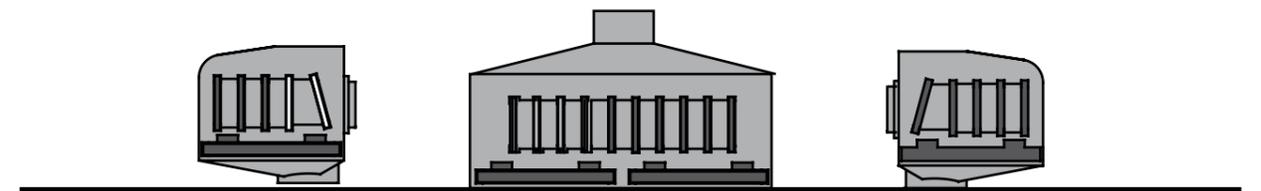
*The high flexibility and modularity of **MOMOLINE** production line allow to equip all fume cupboards with many different accessories and items, in order to meet all customer's wishes.*

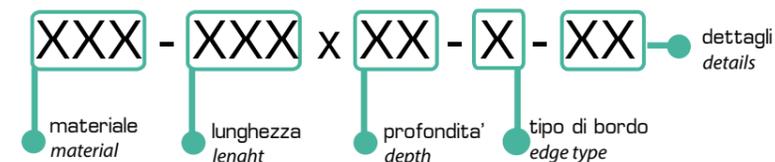
*Accessories reported here are the standard ones, but custom-made and specially designed solutions can be provided upon specific request, so that every customer can fulfill its own desires and expectations from its new fume cupboard.*

**MOMOLINE** is always pleased to receive detailed and specific enquiries from customers, in order to provide the best solution for every occasion.

*The considerable experience of **MOMOLINE** in setting up laboratories for every field and application allows to support customers in the choice of the most suitable items, both from the aesthetic and technical point of view.*

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es: COR120x75-X4

		materiale / worktop						
		COR	PP	HPL	EP	SS	GR	SP
applicazione / application	Utilizzo reagenti non aggressivi <i>Employment of non-aggressive chemicals</i>	●	●	●	●	● <small>AISI 316</small>	●	●
	Utilizzo reagenti molto aggressivi <i>Employment of very aggressive chemicals</i>	●	●	●	●	●	●	●
	Utilizzo coloranti <i>Employment of dyes</i>	●	●	●	●	● <small>AISI 316</small>	●	●
	Settore petrolchimico / Utilizzo solventi <i>Oil&amp;gas / Employment of solvents</i>	●	●	●	●	● <small>AISI 316</small>	●	●
	Settore alimentare <i>Food industry</i>	●	●	●	●	● <small>AISI 316</small>	●	●
	Settore farmaceutico <i>Pharmaceutical</i>	●	●	●	●	● <small>AISI 316</small>	●	●
	Microbiologia <i>Microbiology</i>	●	●	●	●	● <small>AISI 316</small>	●	●
	Elevata igiene <i>High hygiene standard</i>	●	●	●	●	● <small>AISI 316</small>	●	●
	Appoggio strumentazione / Scrivania <i>Equipment support / Desk</i>	●	●	●	●	●	●	●

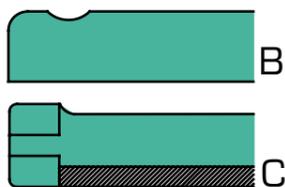
● idoneo suitable    ● poco idoneo not recommended    ● non idoneo unsuitable

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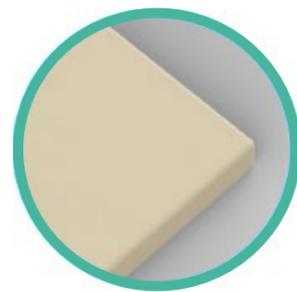
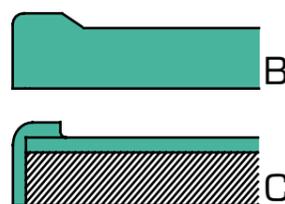
*Corian is a synthetic material, composed by 1/3 of acrylic resin and by 2/3 of natural minerals. Corian has an incredibly uniform and homogeneous structure, does not delaminate, is not porous and is completely hygienic and stain-resistant. Corian also resists high temperature (up to 100°C) and most common chemicals. This material has excellent decontamination performances against viruses and bacteria, and its particular junction method allows to obtain virtually infinite surfaces, with no evident junction grooves. For this reason, it is particularly suitable for microbiology and high hygiene applications. MOMOLINE is part of Corian Quality Network, meaning all the products are officially recognized and guaranteed by DuPont, sole authorized manufacturer of Corian.*

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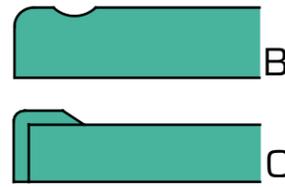
*Polypropylene worktops are injection moulded and stand out for their extreme lightness. PP worktops provide very good resistance against acids, bases and alkalis.*

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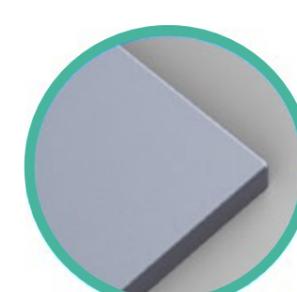
*High Pressure Laminate worktops (such as Trespa) are composed by cellulose fiber layers, soaked in phenolic resin and then heavily compressed through mechanical presses. The surface of the panel is covered with cellulose fiber soaked in melaminic resin with wide choice of color. For aesthetic reasons the external surface can be specifically treated in order to obtain high resistance to chemicals, heat and scratches. Upon specific request, the phenolic resin can be additioned of special substances able to provide the worktop with flame retardant properties.*

EP

piano in RESINA EPOSSIDICA/*EPOXY RESIN worktop*

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*Epoxy resin worktops (such as Durcon) are obtained from molding and then properly shaped using CNC (Computer Numerical Control) machines. These worktops are very easy to clean and decontaminate, are perfectly smooth and non-porous, very resistant to stains and to laboratory reagents. Together with ceramic worktops, they represent one of the best choices when it comes to employment with very aggressive acids.*

accessori  
accessories

accessori  
accessories

SS

piano in acciaio INOX / stainless steel worktop

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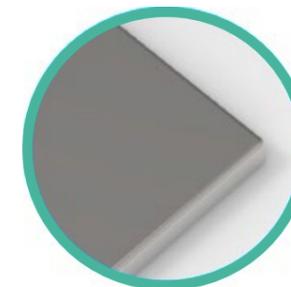
*This worktop can be realized using different SS alloys (AISI 304 or AISI 316), according to the required application. SS worktops are realized using first choice steel sheets, cold bended, and can be provided with raised edge all around. The surface is completely seamless, perfectly shining and with no joints.*

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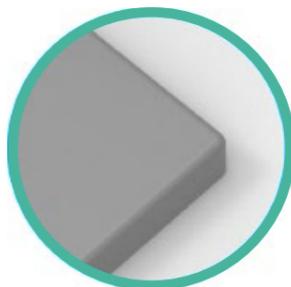
*Enamelled steel is obtained by applying special ceramic enamel on a steel substratum. These worktops provide good resistance to most of common chemicals.*

GR

piano in GRES MONOLITICO / CERAMIC worktop

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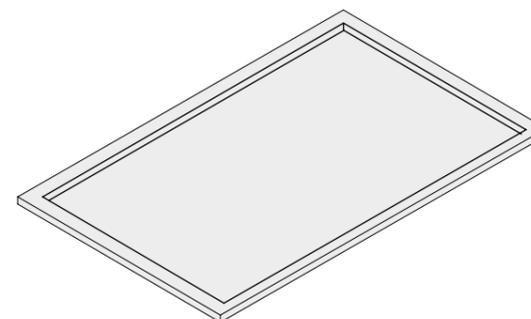
*Ceramic worktops are manufactured through sintering of clay, feldspar and quartz, at a temperature higher than 1200°C. The worktops are perfectly seamless, so that an unique flat surface is obtained, with good resistance against mechanical shocks. Ceramic worktops are recommended for all applications where a superior resistance to chemicals and heat is required.*

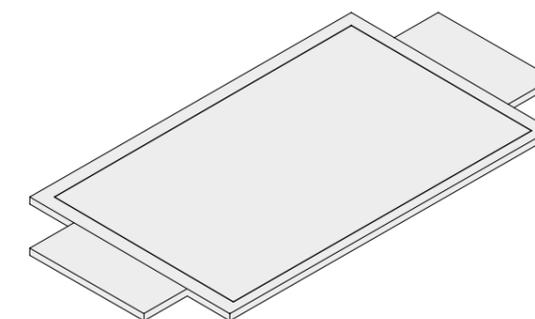
X4

bordo antidebordante 4 lati  
4 sides raised edge

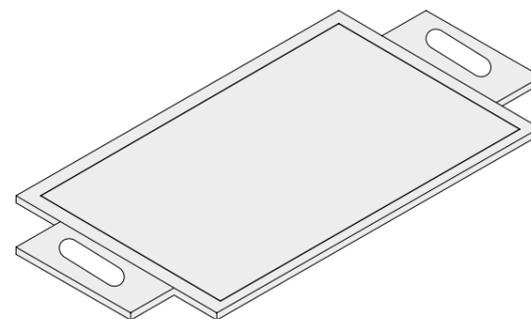
XW

ali laterali  
lateral wings

XWV

ali laterali con vasca  
lateral wings with drip cup

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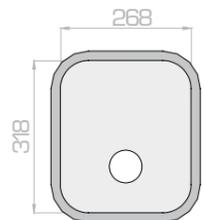
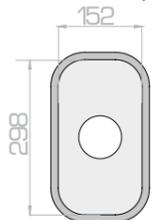
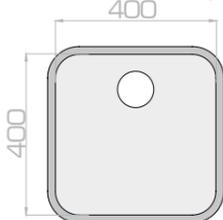
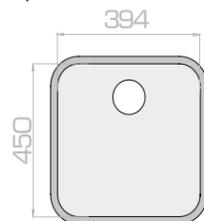
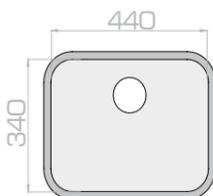
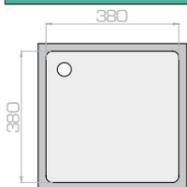
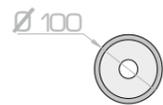
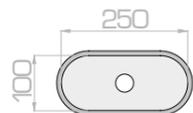
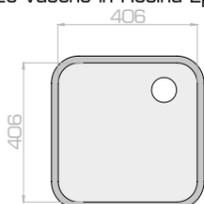
profondita'/  
depth 120 mm

VSC-DC-\_\_\_

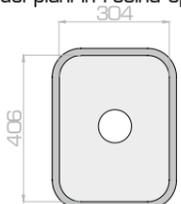
vasche in Resina Epossidica/Epoxy Resin sinks

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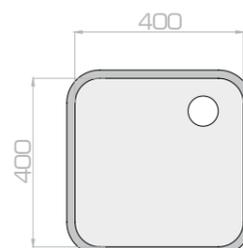
profondita'/  
depth 222 mm

VSC-SS-\_\_\_

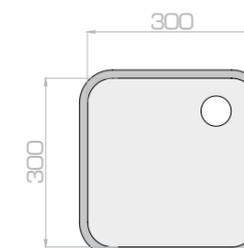
vasche in Acciaio inox/Stainless steel sinks

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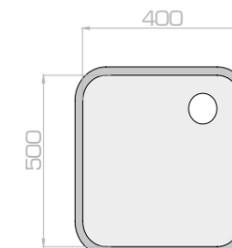
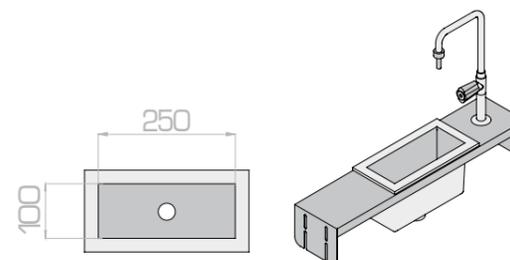
profondita'/  
depth 300 mm

vasche in Polipropilene/Polypropylene sinks

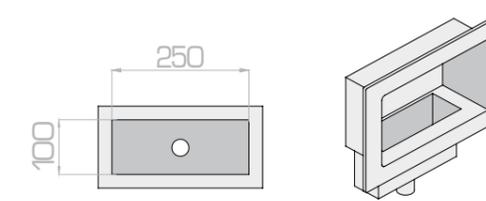
Le vasche in Polipropilene sono accessori dei piani in Polipropilene, HPL, PVC e laminato. Polypropylene sinks are suitable for Polypropylene, HPL, PVC and laminate worktops.

codice/code  
PS

profondita'/  
depth 150 mm

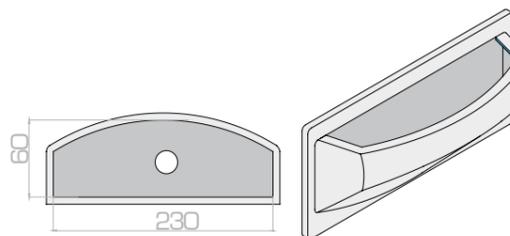
Vaschetta per fascia laterale cappa Belair56  
Drip cup for lateral bracket of Belair56 fume cupboard

codice/code  
TW

profondita'/  
depth 150 mm

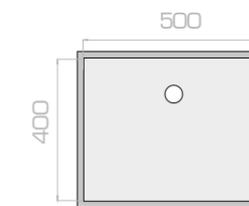
Vaschetta per spalla laterale  
Drip cup for side element

codice/code  
TX

profondita'/  
depth 150 mm

Vaschetta per parete laterale cappa Belair56  
Drip cup for lateral panel of Belair56 fume cupboard

codice/code  
VSC-PP-543L

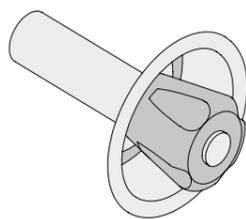
profondita'/  
depth 300 mm

accessori  
accessories

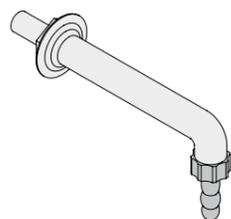
Q-\_\_\_

rubinetti e miscelatori acqua / water taps and mixers

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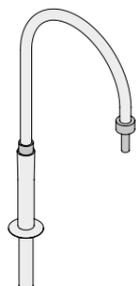
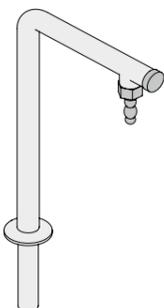
codice/code  
Q-PP-2

beccuccio da piano in PP  
PP worktop nozzle

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All taps are manufactured according to DIN 12918 norm.

The threaded demountable nozzle is manufactured according to DIN 12898 norm.

All handles are made of anti-acid polypropylene, and the colors comply to EN 13792 norm.

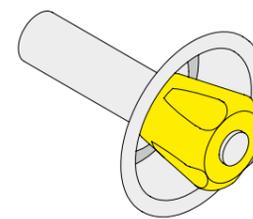
Maximun operating pressure is 7 bar. Temperature operating range 30°C to 150°C.

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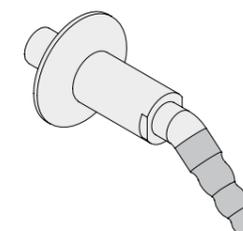
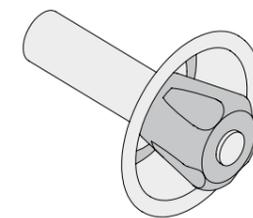
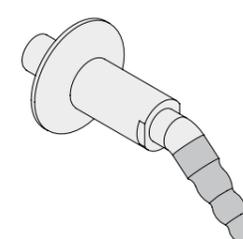
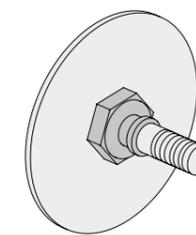
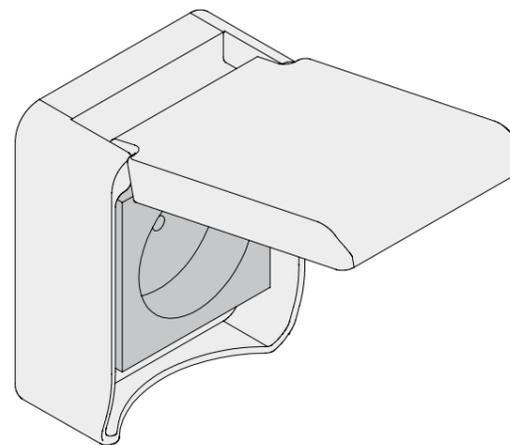
*Single stage pressure reducer for laboratory. Flow pressure gauge compliant to EN 837-1 norm, embedded needle type valves on outlet NY66. Temperature operating range : -20/+60°C.*

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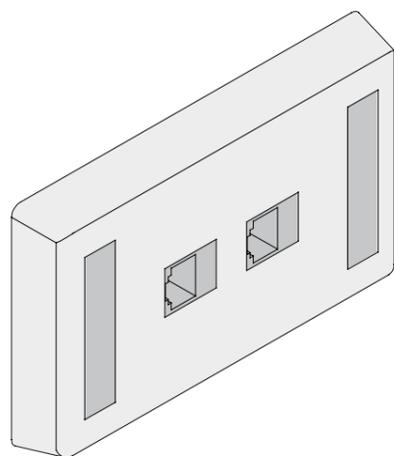
*IP65 Power outlet protection grade, according to IEC 60529 norm, IK 09 protection grade against mechanical shocks, according to IEC 62262 norm; heat resistance 850°C, according to IEC 60695-2-1 norm. 10/16A, 250V. Different socket standards available.*

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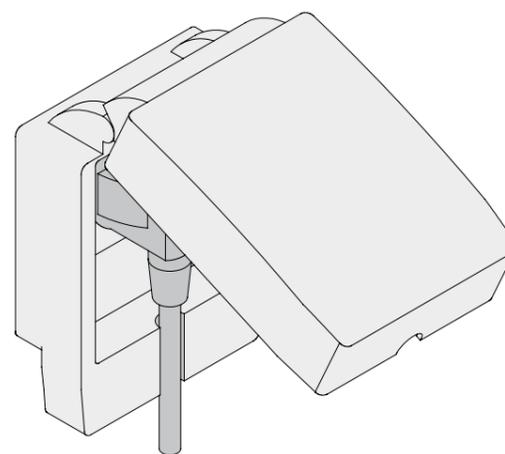
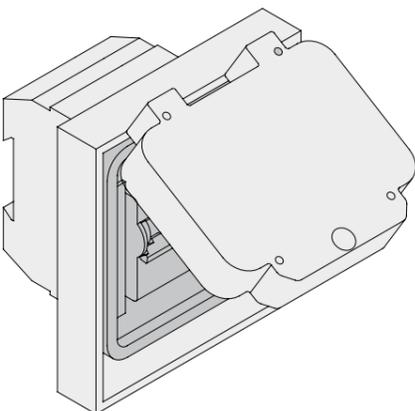
*IP66 Power outlet with fitted plug protection grade, according to IEC 60529 norm, IK 08 protection grade against mechanical shocks, according to IEC 62262 norm; heat resistance 850°C, according to IEC 60695-2-1 norm. 10/16A, 250V. Different socket standards available.*

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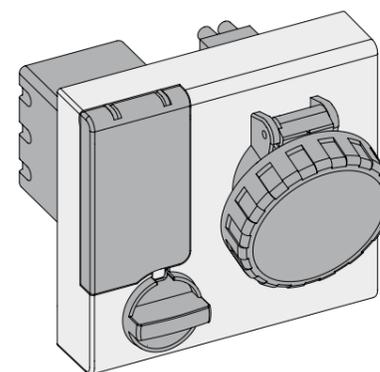
*The provided MCB is certified by CE and IMQ for compliance to CEI 23-3(EN60898) and EN 60947-2 norms. Intervention curve C. 16A, 230V. Interrupting capacity: 4,5kA a 230Vca (EN 60898).*

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*IP67 interlocked power outlet, with embedded protection fuse and switch. 16 A, 50/60 Hz. Sockets are marked IMQ and CE, and complies to EN 60309-1, EN 60309-2 norms.*

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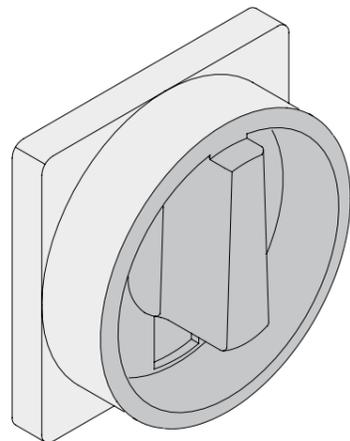
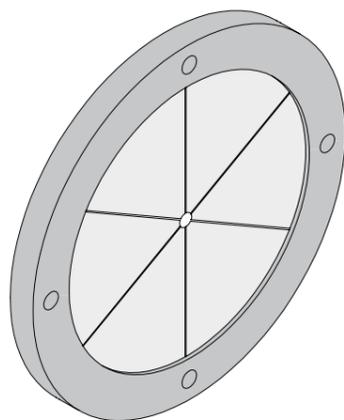
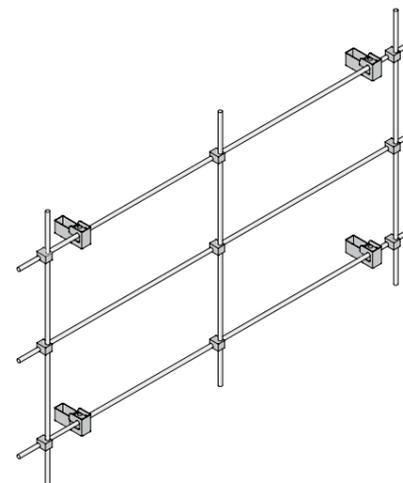
*The cable grommet has a 100mm diameter, and can be used to connect all the equipments/instruments under the fume cupboard to a power supply, not interfering with the sash movement and with the airflow patterns. The cable grommet is entirely realized in soft anti-acid PVC, in order to grant a comfortable and safe employment, as well as to keep the containment performances unvaried.*

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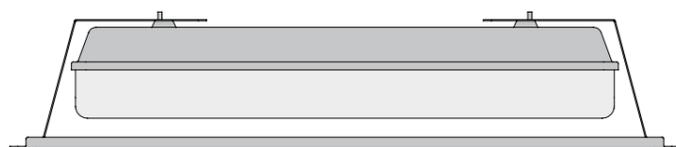
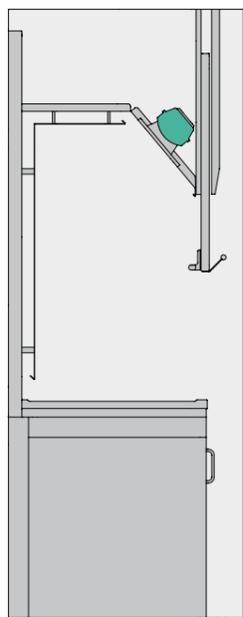
*Stainless steel rack made up of properly shaped bars and fastened to the rear of the fume cupboard. The rack can be used for hanging/fastening objects or instruments, such as distillation equipment.*

codice / code*	cappa da / cupboard length
RST-90	900
RST-120	1200
RST-150	1500
RST-180	1800
RST-220	2200
RST-240	2400

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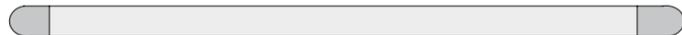
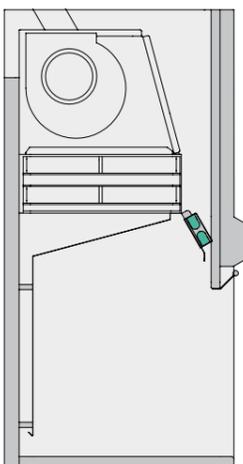
*Lamps are positioned outside the working area, so that the risk of triggering potentially explosive vapors is removed. The lamps are indeed separated from the working area by a dedicated wired safety glass. Standard illumination makes use of a IP65 water-proof lamp with 2x18 W neon tubes having suitable power for granting approved enlightenment. Upon specific request, other items can integrate or substitute the standard items:*

- IP65 water-proof lamp with LED lamps;
- further adjustable light;
- UV lamp with timer.
- ex-proof lamp.

codice/code  
**K-LED**  
lampada a led  
led lamp

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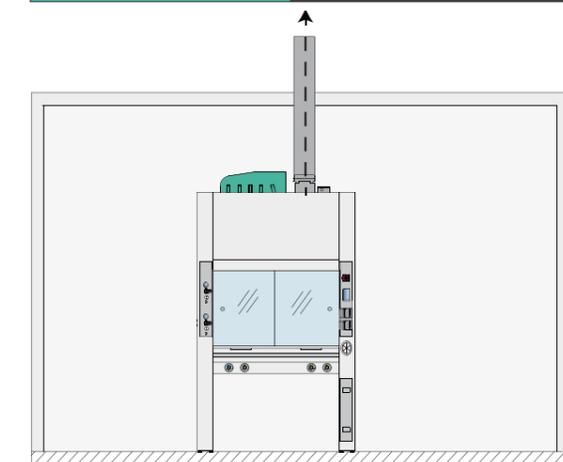
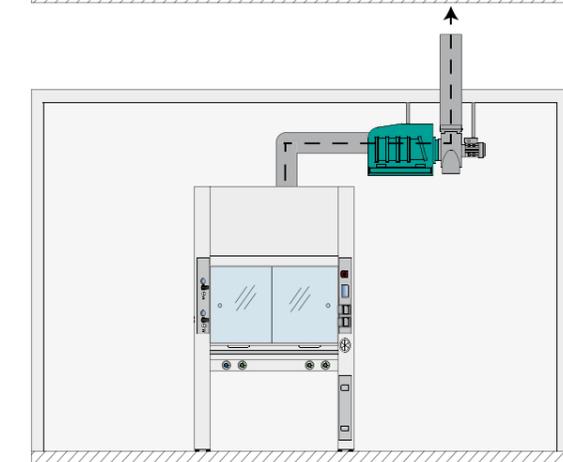
*In CARBOCHEM model the lamps are positioned outside the working area, so that the risk of triggering potentially explosive vapors is removed. The lamps are indeed separated from the working area by a dedicated wired safety glass. Standard illumination makes use of two led lamps 220V, 50Hz, IP65, 4500 K. VDE omologation, CE compliant.*

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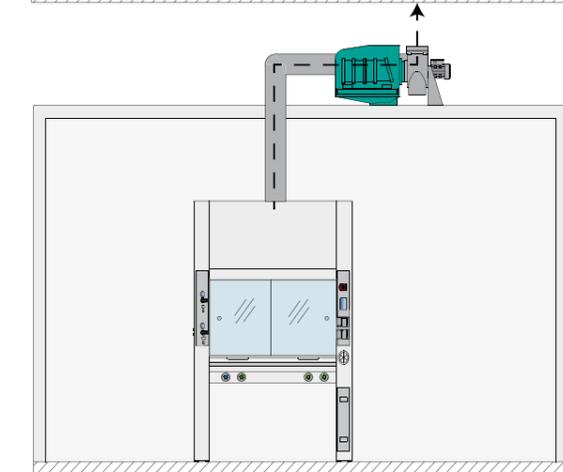
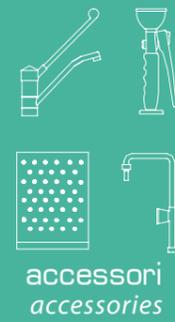
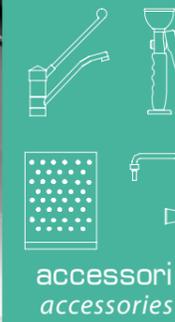
*All hood models can be equipped, upon request, with filter boxes on board, wall/ceiling, or in a remote location outside the building. The filter box is selected according to customer requirements and installation characteristics.*

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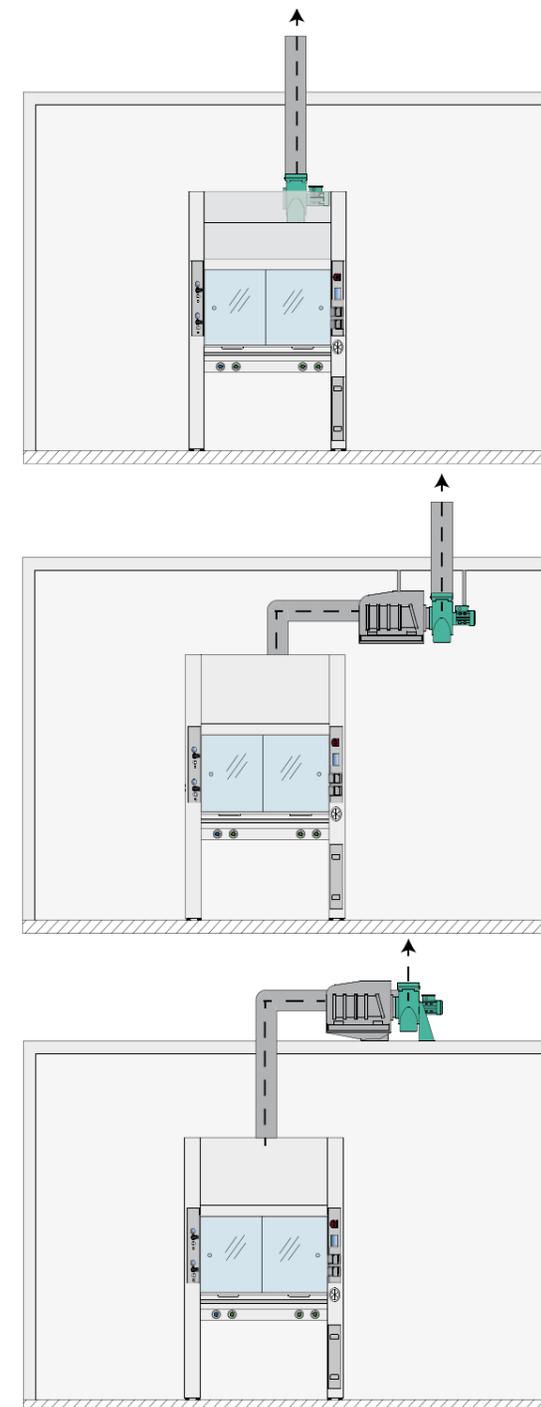
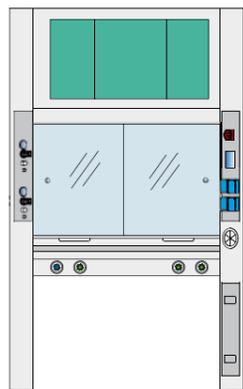
*Exhaust fan is chosen according to the technical feature of the specific fume cupboard installed as well as on the base of the exhaust duct. The aspiration and containment parameters recommended by international norms are always guaranteed. ATEX (ex-proof) extraction system is available on request. In case of blower installation outside the building, is available a kit inclusive of anchoring base, anti-weather covering, anti-birds net, IP65 electric breaker for maintenance operations (code ASP-EST).*

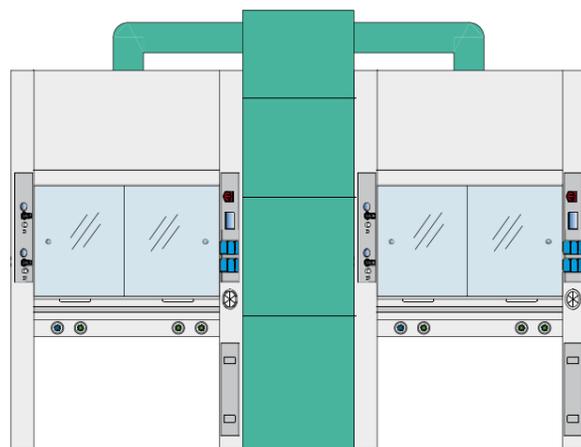
SCR

scrubber / scrubber

scrubber a bordo cappa  
scrubber on the hood

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*The installation of a gas scrubber is recommended for use in laboratories where toxic, corrosive and odorous gases are released.*

*The laboratory gas scrubbers are used mainly for the absorption of strong acids. Inside the scrubber the contaminated exhaust air is brought into contact with an aqueous absorption solution.*

*To achieve a high efficiency it is important to ensure a large interfacial area between the air and the liquid. This is achieved by distributing the liquid over a packed bed.*

*All scrubbers are equipped with a level control. If the liquid falls below a certain level, fresh water is added automatically. The electrical conductivity of the washing solution is measured. If the adjustable set value is exceeded, a portion of the washing solution is discharged and replaced by fresh water.*

*Is also available with a pH measuring and control facility. By controlled addition of caustic soda or acid, the pH-value of the washing solution is kept at a constant value. This optimizes the performance of the scrubber.*

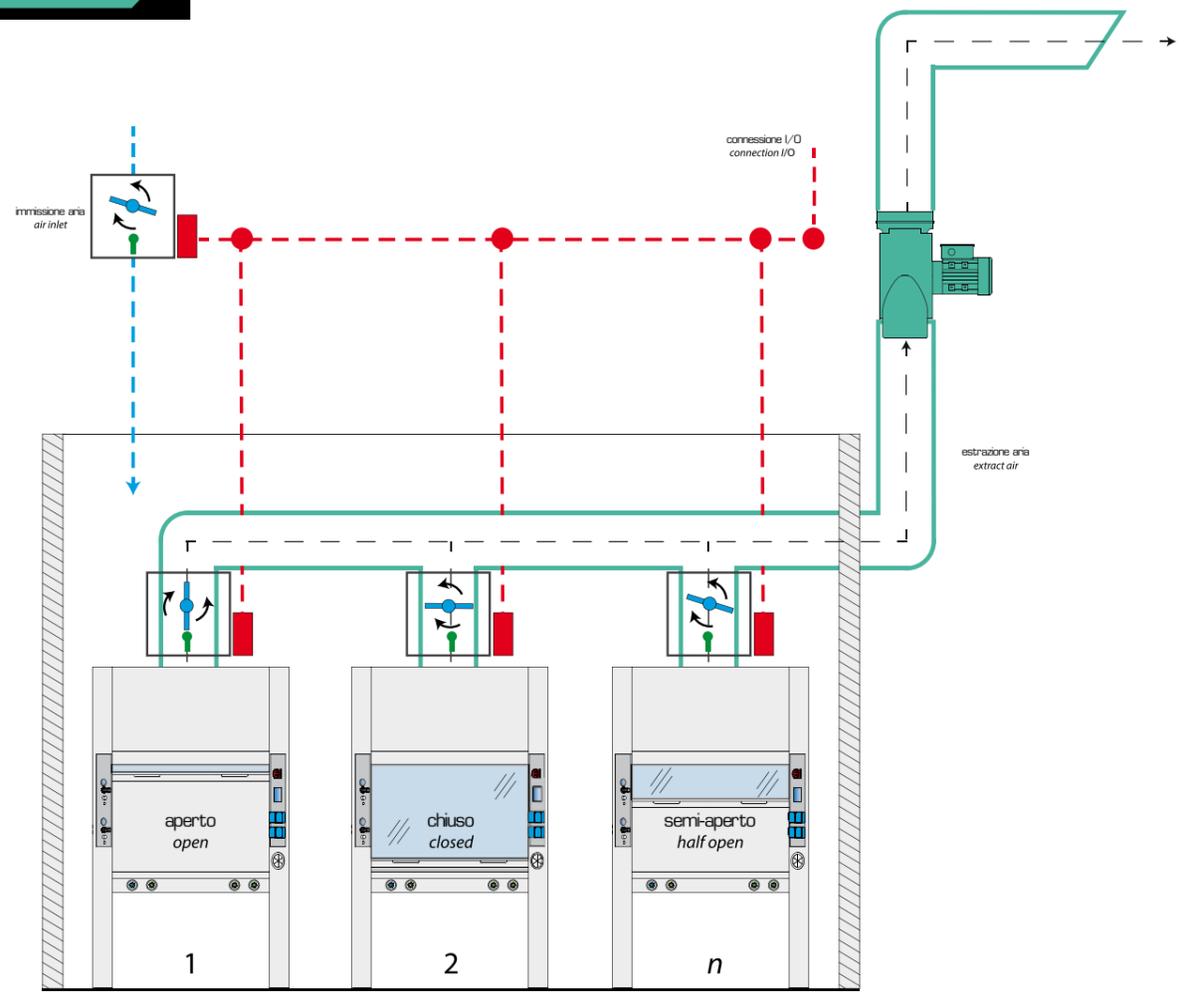
*Gas scrubber can be installed on the top of the fume cupboard, for dedicated employment, or on the floor, besides the fume cupboard.*

*Floor installation allows to increase the filtered airflow and is often employed for connecting more fume hoods to the same scrubber, optimizing the installation.*

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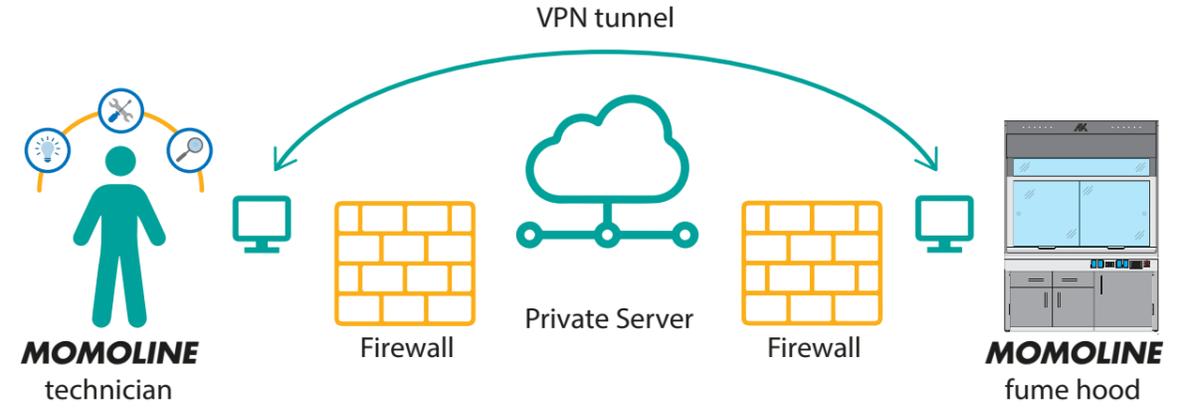
The control system of fume cupboard **BELAIR56** can be connected to the air conditioning system (HVAC) of the building. The interface grants an optimal management of inlet and outlet airflows, ensuring at the same time the efficiency of the fume cupboards and an optimization of conditioned air volumes. This solution can determine substantial energy savings.

**UBIQUITY** assistenza tecnica remota / remote technical assistance

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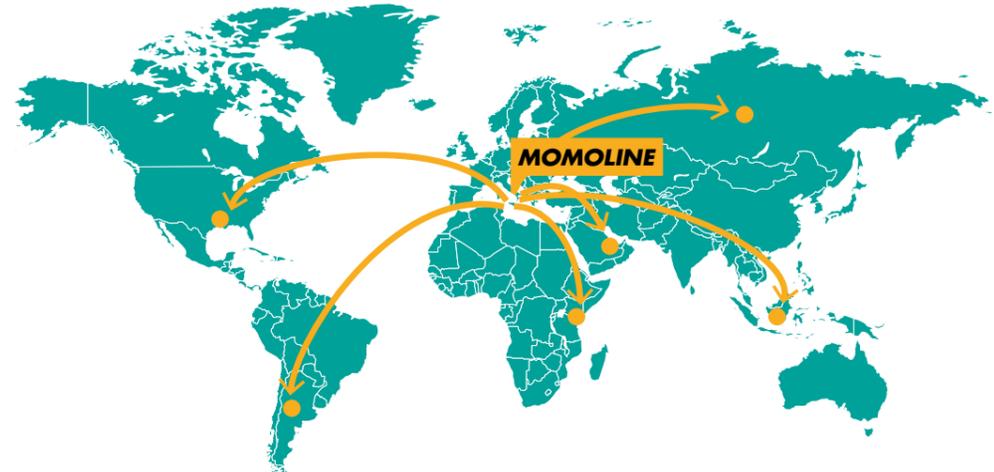
Fume hoods can be equipped with a remote assistance system to manage post-sales services such as technical assistance, software updating, maintenance, debug.

This solution enables **MOMOLINE** technicians to access the remote device (the single fume hood) connected to the network through a VPN based on proprietary technology having procedures with end-to-end sessions without any intermediate and optimized for industrial communication. Reliability and service continuity are granted thanks to a redundant and distributed server infrastructure. The server infrastructure is globally distributed (Europe, United States, South America, Asia). The system does not require additional hardware and allows technicians to operate on remote devices (after the authorization of the customer) eliminating the distance, reducing the need for on-site assistance, dramatically reducing postsale service costs.

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*Within its facilities **MOMOLINE** has a modern test laboratory for fume cupboards realized according to EN14175-3:2019 and fully equipped to carry out all tests from standard EN and ASHRAE.*

*All models of fume cupboards sold by **MOMOLINE** have been tested (type test), and test results have been validated by Bureau Veritas, who has therefore issued third-party certification.*

*Upon request, specific tests on the single fume cupboard can be executed before the supply and test report accordingly to EN14175 can be provided.*

**MOMOLINE** is also certified by a third party according to EN14175-4 for the execution of in situ tests with the issuance of test reports in accordance with EN 14175-3:2019.