General catalogue

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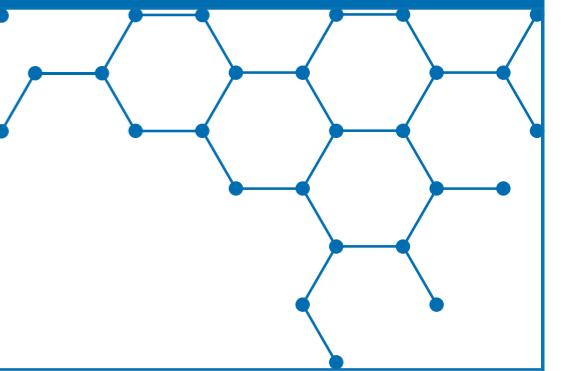
Momoline Company was founded in 1978, being at that time mainly focused on the sale of analytical instruments and consumables, but standing since the beginning as a complete and polyvalent partner for the laboratory world.

Nowadays, Momoline stands as an undoubted leader in the field of lab designing, as a reliable and experienced supplier for the set-up of the most advanced analytical facilities, with significant references in all the main fields where the presence of a laboratory is determining.

Along with its multifaceted attitude, from the beginning the strength of the company has been represented by the centralized care and management of the entire industrial process; the experienced staff of Momoline takes care of customers in all steps of the laboratory set-up: since the designing stage, the Company provides assistance through the activities of manufacture, supply, installation and after-sales assistance.

The manufacturing process of laboratory furniture is entirely controlled and managed by Momoline: as a corporate strategy, the Company does not outsource any of its activities, taking direct involvement from design to installation, also in case of customized requests from customers.

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Momoline is structured to provide its customers the best service, and to stand as the ideal partner for the set-up of your laboratory!

The architecture of the Company has evolved along the years, shaping its structure on the goal of assisting customers in all aspects: experienced technical staff provides the most specific and advanced solutions, which are shown to the customers through realistic representations.

The chosen solutions are then brought to reality in factory through the different production departments, and finally installed at customer's site by Momoline's technicians. Manufacture is articulated in the Departments of Metal Carpentry, Wooden Carpentry, Polymers Carpentry, Painting, Mechanical & electric Assembling.

Sales Office can assist the customer in all phases, from preliminary design up to after-sale services. In Sales Office there is a dedicated team caring about public tenders, with highly specialized staff, learned about the most updated regulations. All Momoline's departments work in symbiosis with the aim of always ensuring best quality and customer satisfaction.

It is Momoline's primary and steady goal achieving high quality standards by daily researching for the most modern and innovative materials, and by adopting manufacturing approaches in the van of technology. Besides, the care for environment safeguard and safety laws fulfillment, in the light of reaching a sustainable development production line, is always a must. These principles have been validated by the achievement of ISO 9001 - ISO 14001 - ISO 45001-ISO50001 certifications.

Within its facilities, Momoline has a modern test laboratory for fume cupboards realized accordingly to EN14175-3:2019 and fully equipped to carry out all tests from standard EN and ASHRAE.

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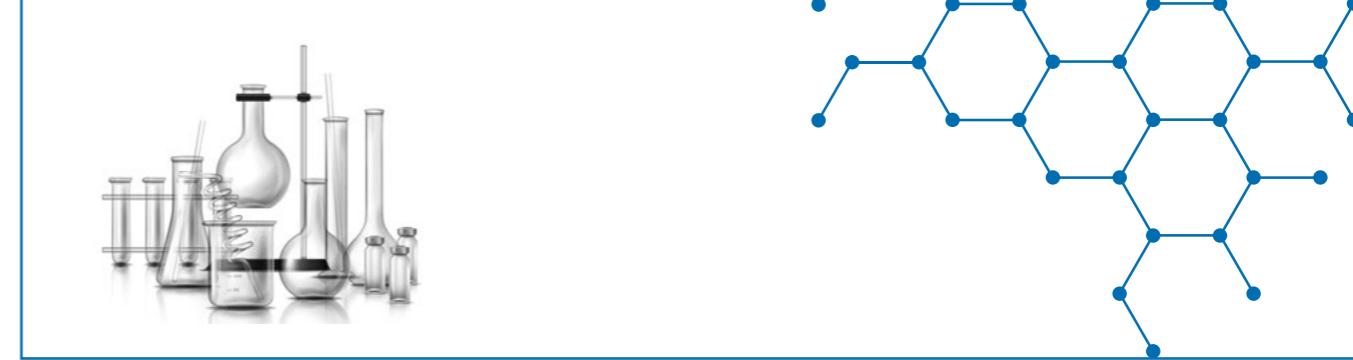
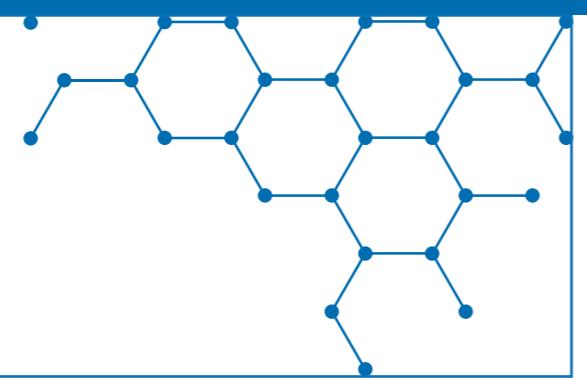
In the last decade the Company has embarked in a new thrilling adventure, putting in place an ambitious plan of Internationalization, which has allowed to enter many new stimulating markets and to get satisfying results by exporting the quality and the style of Made in Italy products towards new and broader horizons.

The company can boast a wide portfolio of references, both in Italy and abroad, in the most diversified fields: many satisfied customers can be found, for example, in the fields of pharmaceuticals, industry, chemistry, hospitals, Oil & Gas, research, didactics.

The range of products offered by Momoline to its customer is constantly updated and improved, so to keep the pace of an ever-changing market, offering the possibility of choosing among many different product, options and configurations.

Momoline is also able to execute, upon request, very special jobs and installations such as mobile laboratories and innovative solutions which can be hardly found in the Standard catalogues of main laboratory furniture manufacturers.

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Momoline is highly specialized in the design and installation of all technical systems and facilities which are essential to the laboratory, always in compliance to all applicable Standards and according to the most updated design techniques.

In the following the essential facilities to the good working of the laboratory are listed:

Facilities for the handling of exhaust air

These facilities are essential for a correct and efficient management of the laboratory, as well as to grant the safety of all its operators. The airflow extracted from the laboratory through fume cupboards, safety cabinets, extraction arms and other items should be carefully estimated and properly determined together with the number of hourly air-exchanges in the room, so to get an efficient and safe place of work. To obtain an exhaustive ventilation into the lab, also extraction fans and exhaust ducts should be suitably dimensioned by skilled and experienced personnel.

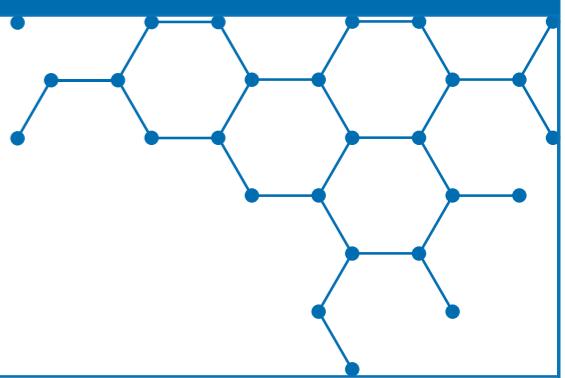
Specialty Gas Distribution Systems

Momoline can take care of the realization of the entire distribution network, starting from pressurized cylinders up to the hook-up points installed on furniture, also caring about the implementation of manifolds, valves, pressure regulators, together with safety devices and electronic controllers for the monitoring of the cylinders level.

Gas detection system

In very advanced work environments (such as Pharmaceutical or Oil fields) the installation of sensors for detecting possible leaks of toxic/flammable gases as well as smoke and flames might be necessary.

The installation of these sensors, along with its supporting network, monitoring units and alarm systems, can be carried out by Momoline personnel and carefully planned and evaluated according to the specific needs.

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MOMOWORK laboratory benches and desks represent the perfect synthesis of simplicity, strength and design, being the best solution for setting-up a modern and technologic laboratory, with a powerful visual impact.

The structure is entirely composed by epoxy powder coated steel profiles, granting the 'Class O' fire reaction, as well as high resistance against corrosion.

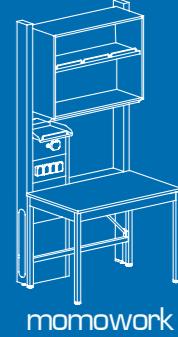
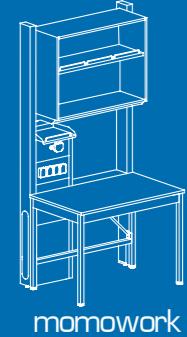
The side frames of these benches, shaped as a "reversed U", are connected by three horizontal elements ensuring high solidity and reliability.

MOMOWORK laboratory benches and desks hold Certificate stating the compliance to the European Norm EN 13150:2001. The compliance certificate has been issued by Bureau Veritas, an authorized third body officially recognized at international level.

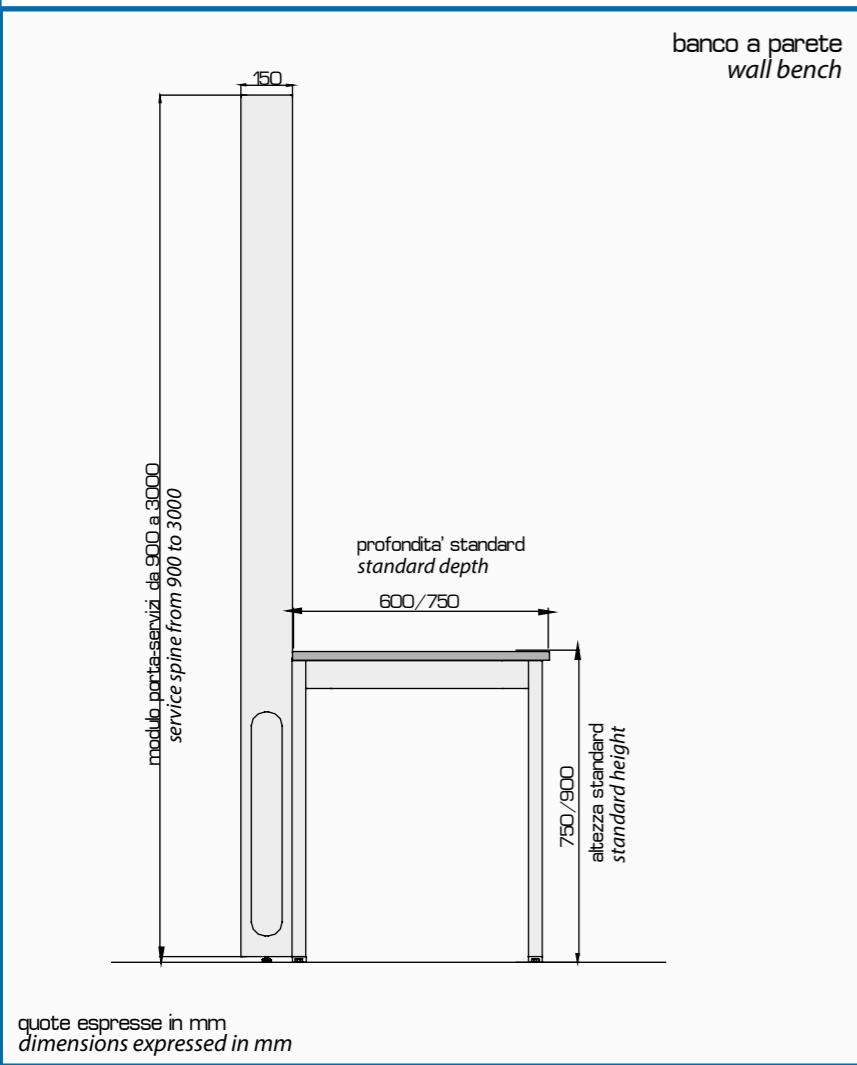
What is more, the structures have been subjected to strict and severe load tests to demonstrate and verify their effective reliability: both experimental tests and software aided simulations (Finite Elements Method - F.E.M.) have been performed, and both have led to excellent results.

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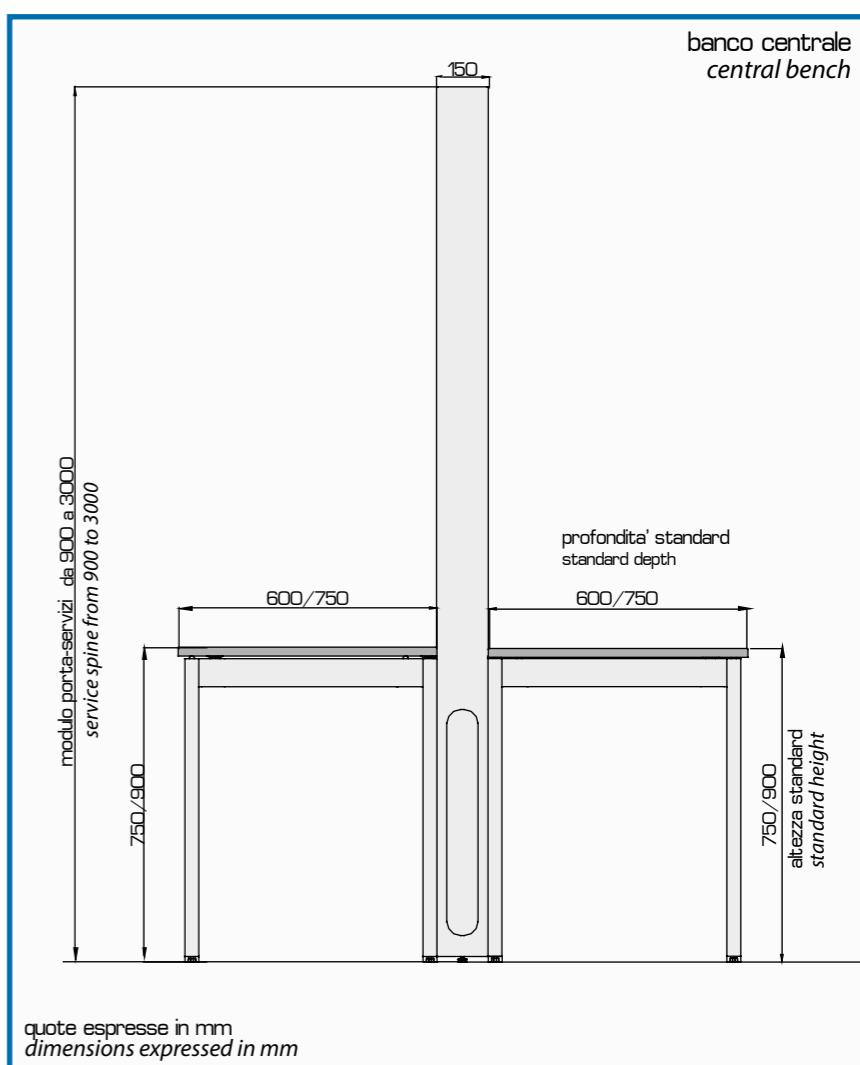
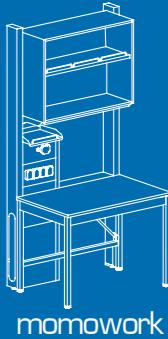
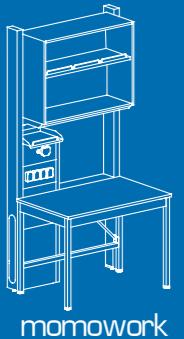
MOMOWORK structure has dedicated compartments to allow the routing of utilities and their connection to the delivery points mounted on furniture.

Benches and the desks can be equipped with special technical panels, housing all the necessary services and pick-up points for technical gases, power, data etc. The pick-up points are installed on interchangeable steel panels, fitted to the main structure through quick fitting devices, allowing to perform inspection and maintenance operations with no need of tools, as well as making future changes easier.

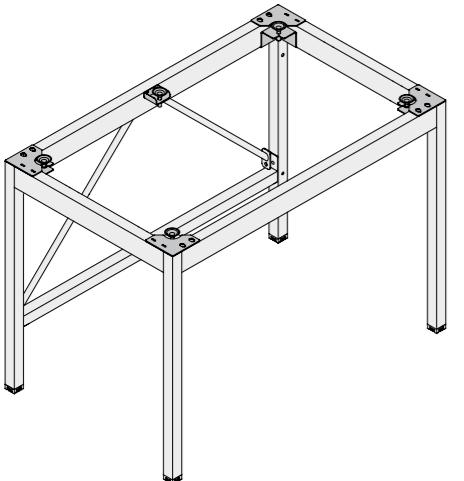
The basic structure can be then completed by a wide range of items such as upper cabinets, shelves, reagent racks and underbench cabinets.

Although **MOMOWORK** line offers a range of standard dimensions, even the most demanding needs can be fulfilled, realizing products completely customized in terms of dimensions, structure and installed equipments and accessories.

The standard height of the worktop is 900 mm in case of conventional laboratory benches, to be used in standing position, and 750 mm in case of desk application, to be used in seated position. Upon request it is possible to produce height adjustable benches, for employment with wheelchair.

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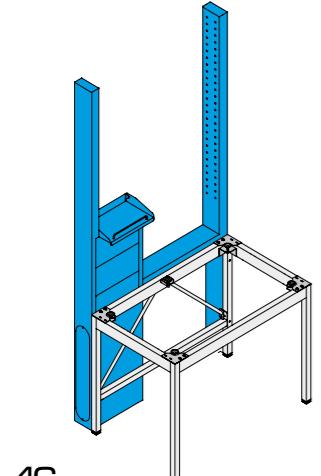
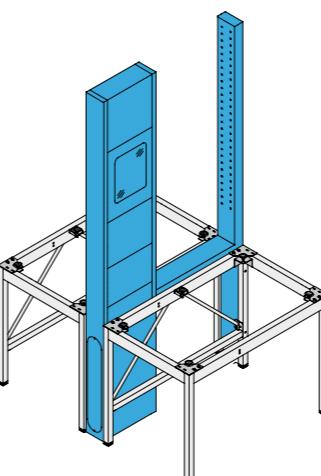
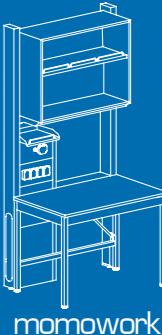
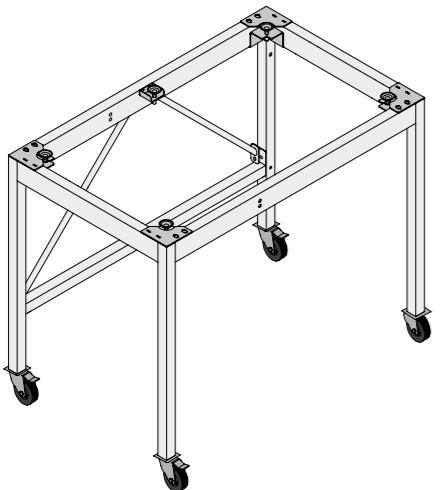
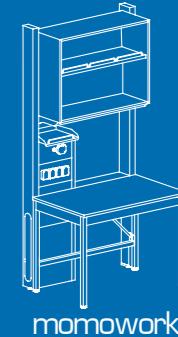
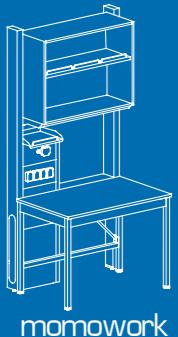
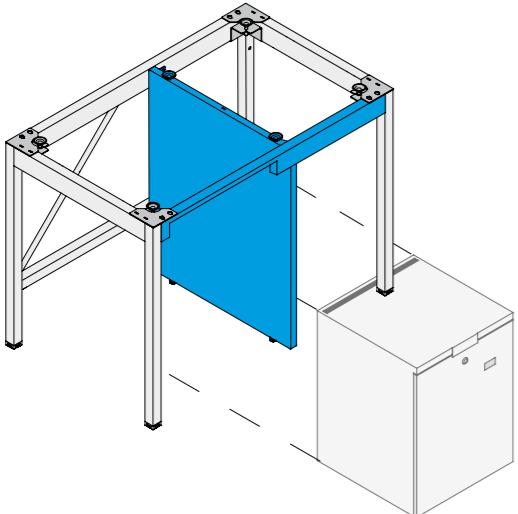
The antivibrating bench is equipped with an insulating column laying on the floor. The insulating column, which can be provided with further ballast upon request, sustains a 40 mm thick HPL pad through four adjustable rubber vibration dumpers (also different materials for the pad available upon request). The column is installed under a standard bench.

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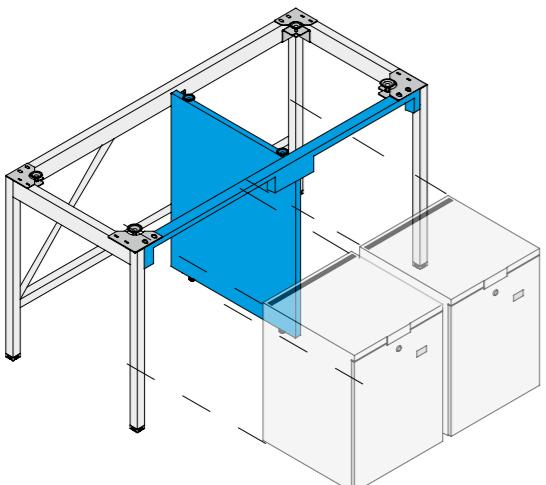
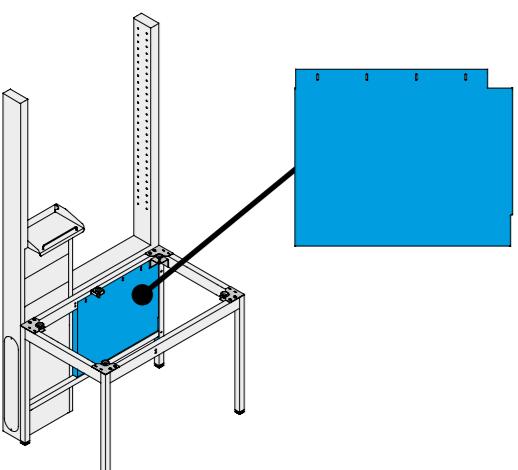
Epoxy-powder painted steel panel for concealing utility lines and services.

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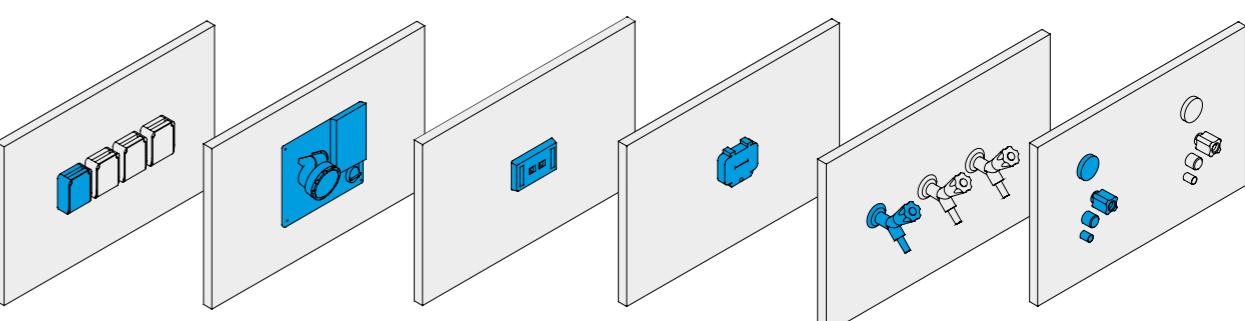
The installation of panels in the rear side of the bench allows an easy access to the utilities in case of instrumental applications (e.g. gas chromatography).

codice/code	per banco da/for bench length
PP-60-MW	600
PP-80-MW	800
PP-90-MW	900
PP-100-MW	1000
PP-120-MW	1200
PP-150-MW	1500
PP-180-MW	1800

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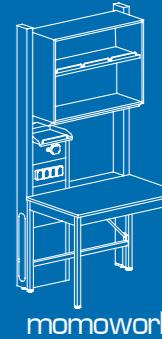
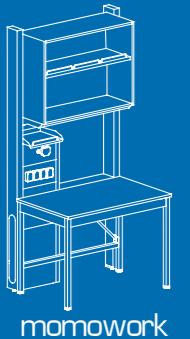
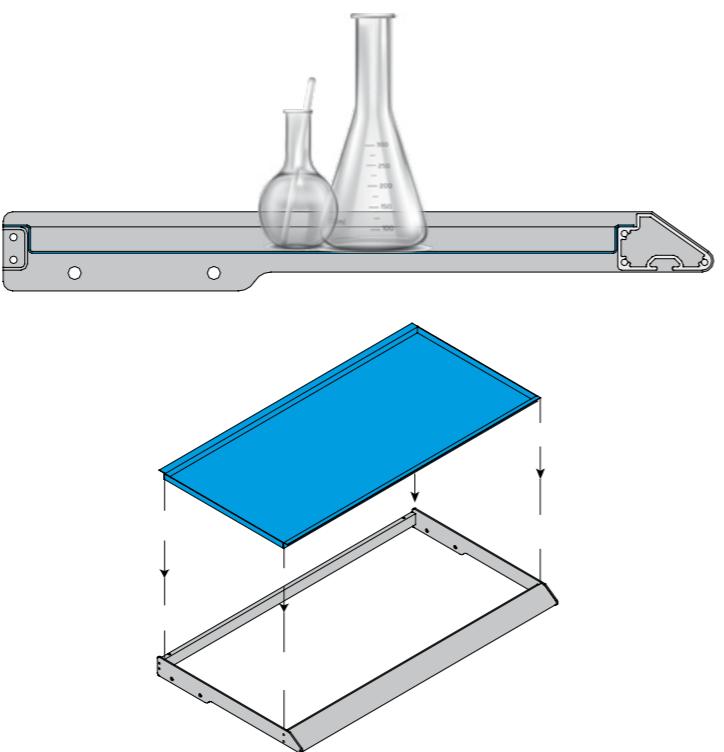
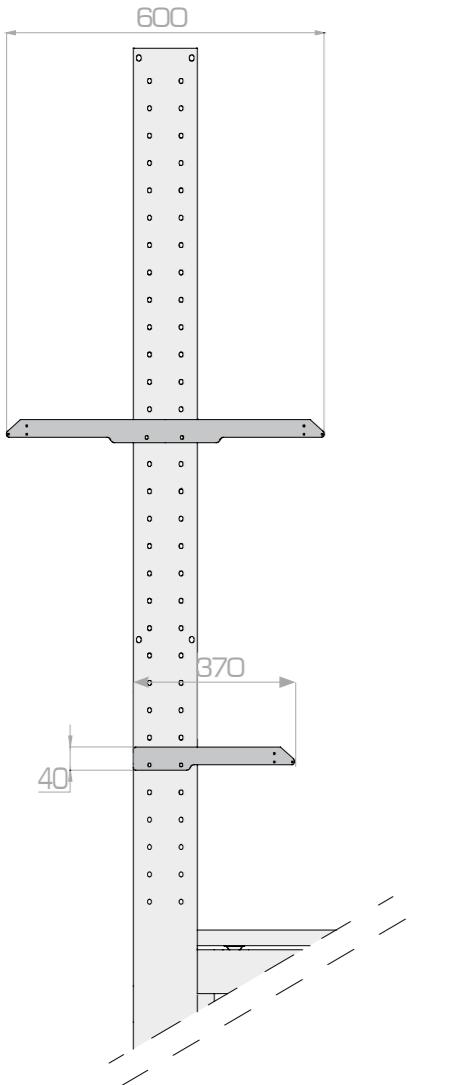
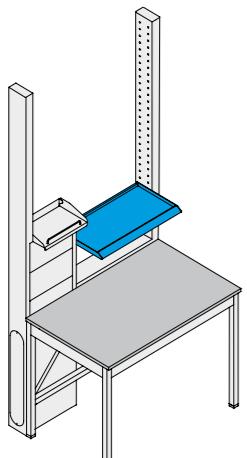
*This shelf is realized in AISI 304 stainless steel, and is equipped with a tray preventing any possible liquid spillage.
The tray is supported by an epoxy powder coated aluminium frame, which can be fitted to the service spine at different heights.*

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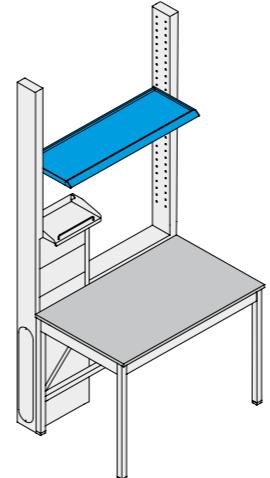
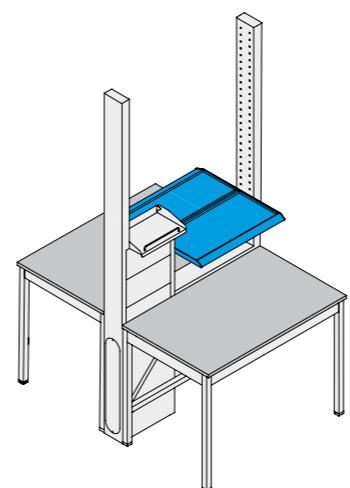
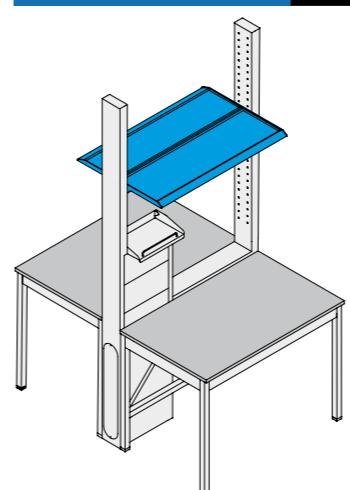
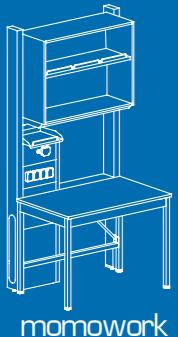
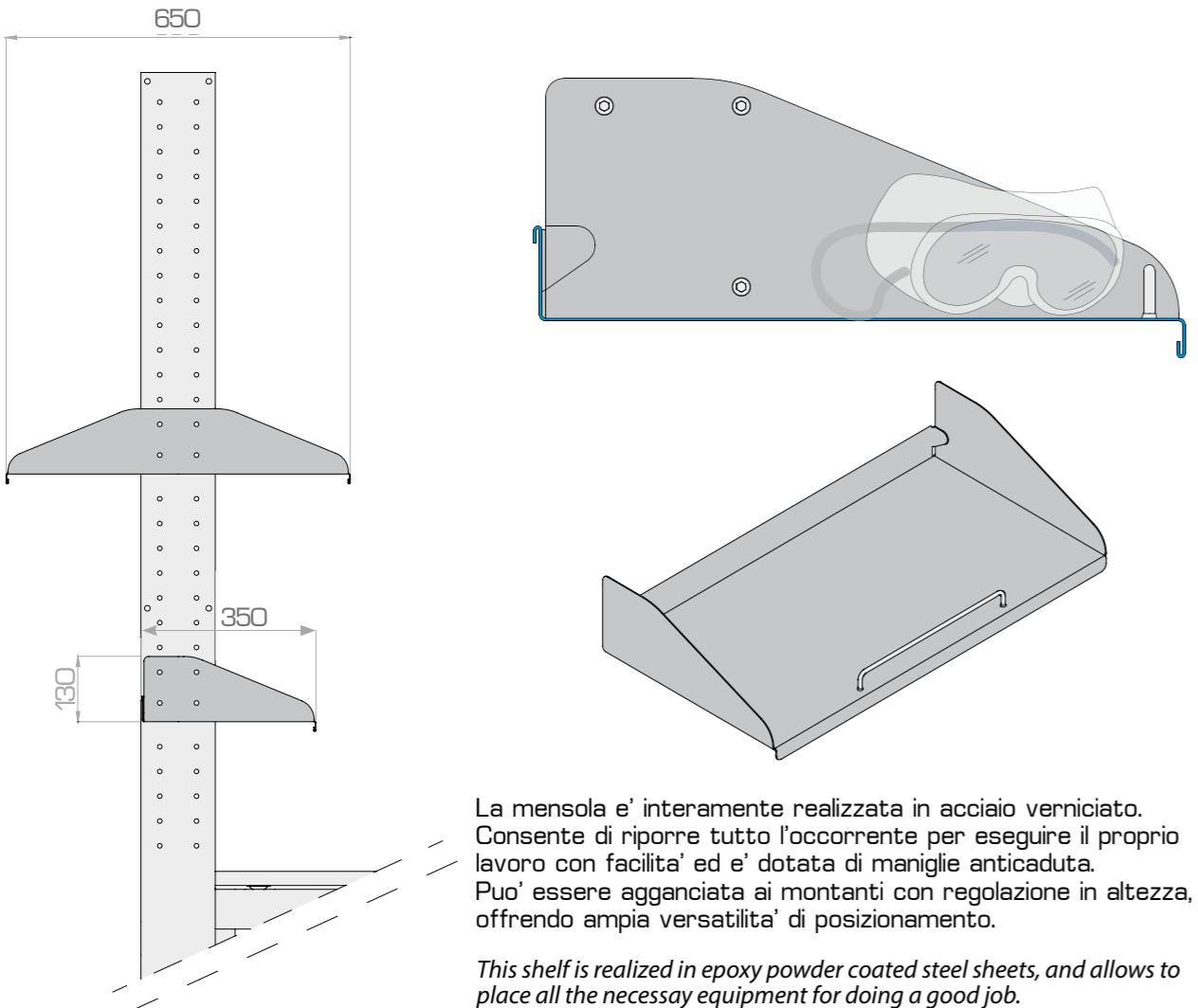
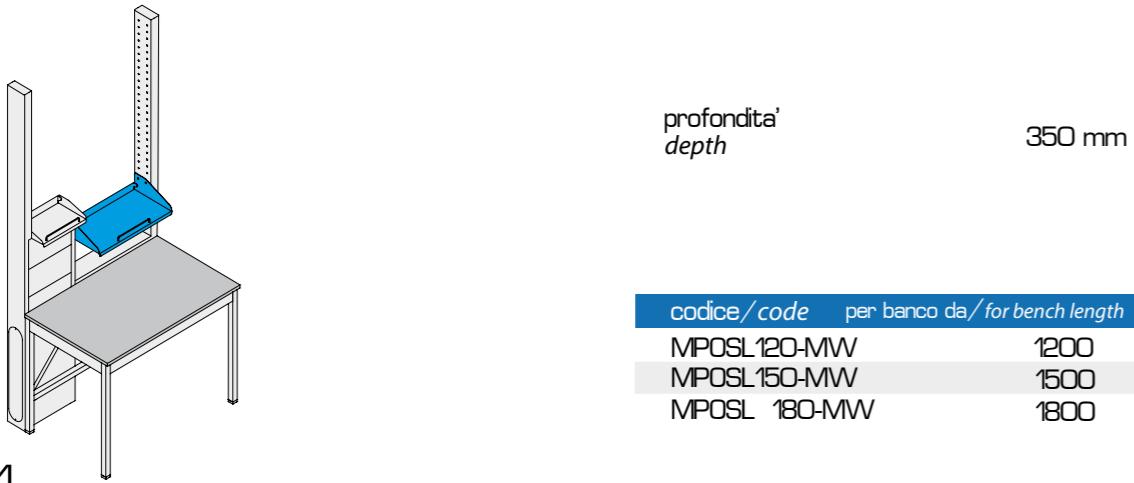
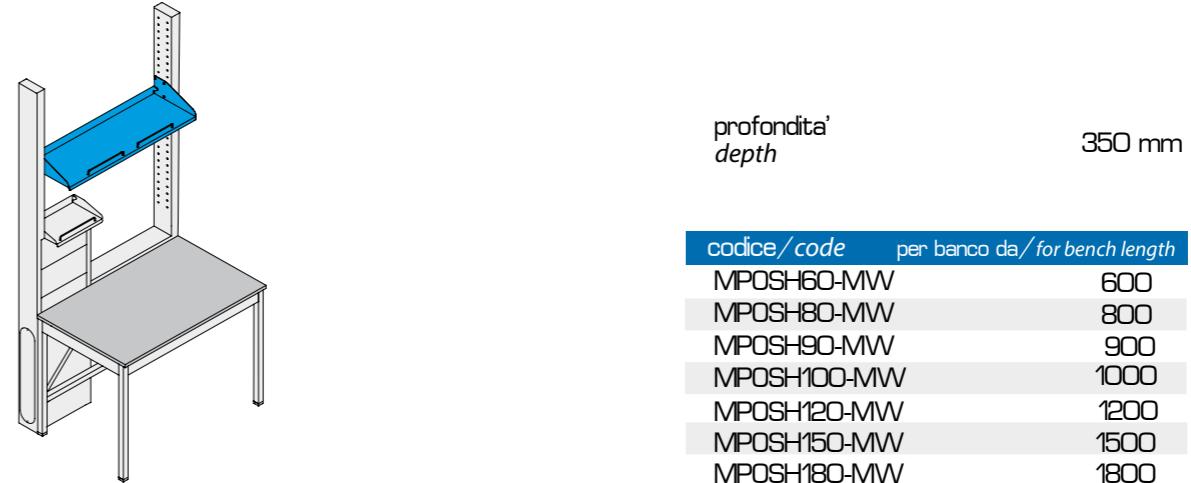
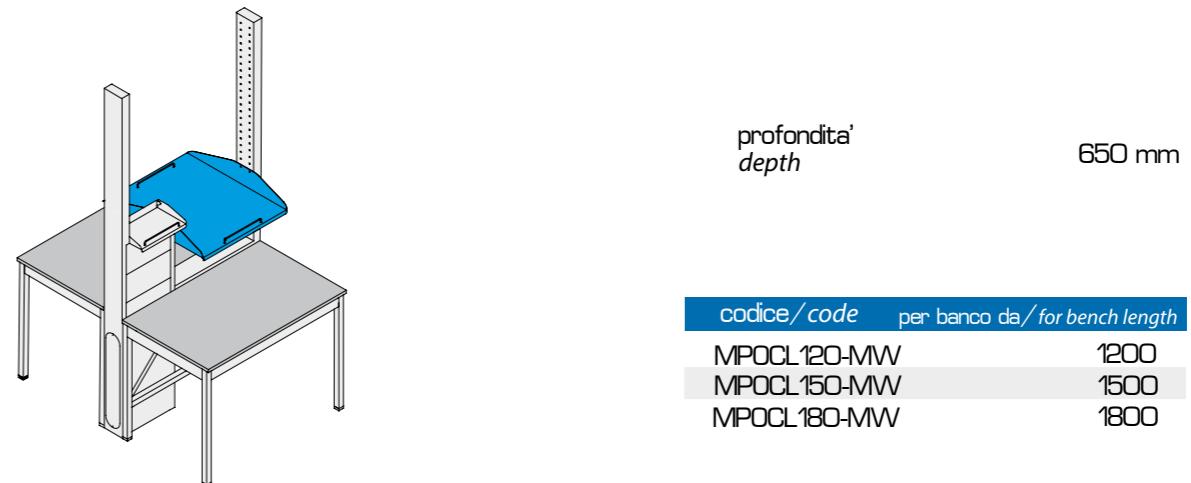
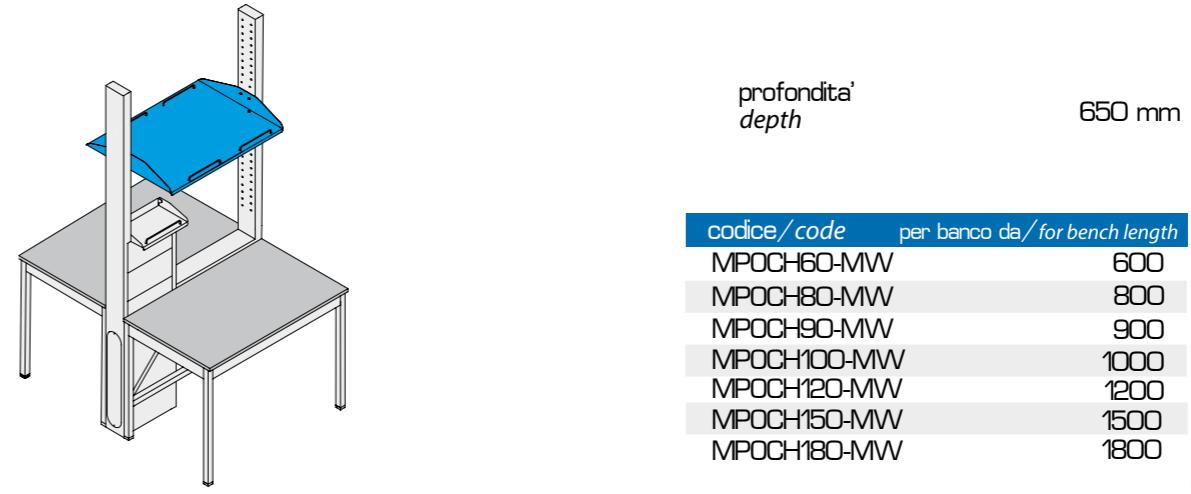
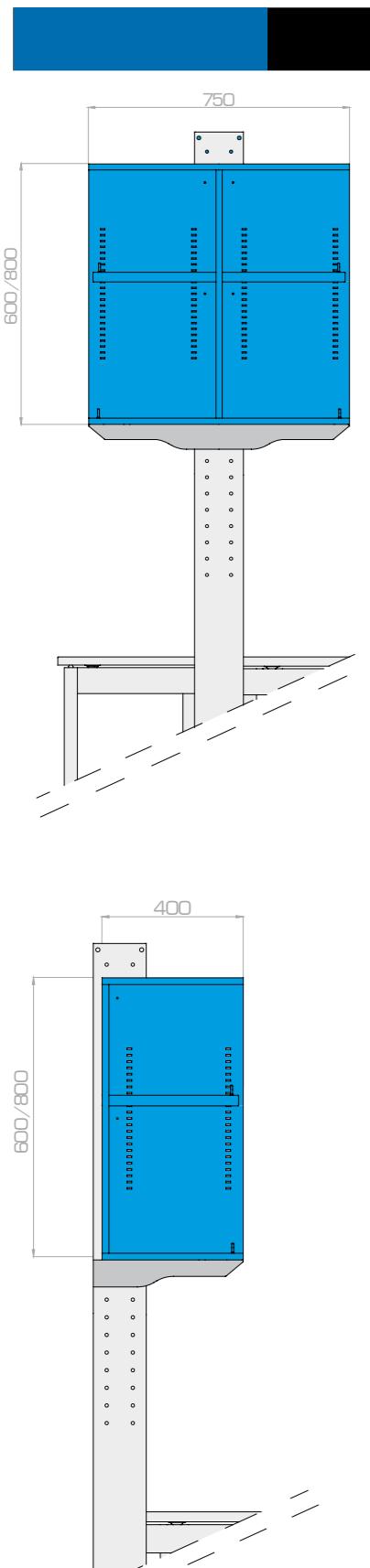
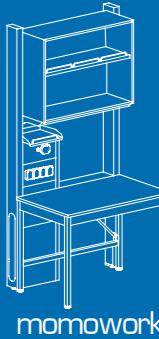
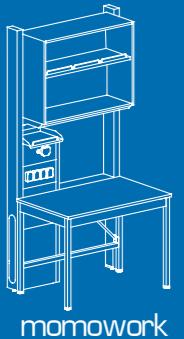
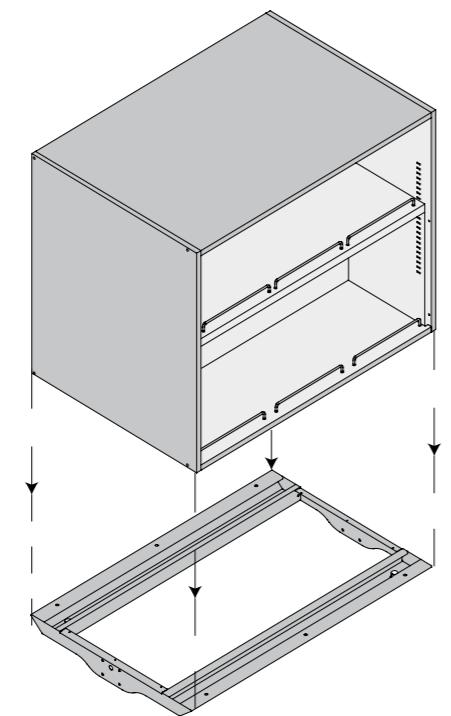
Overhead cabinets can be made either of wood or steel.

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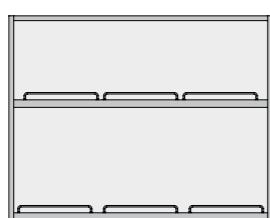
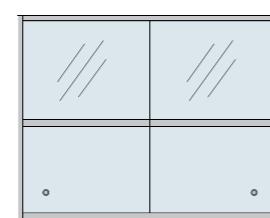
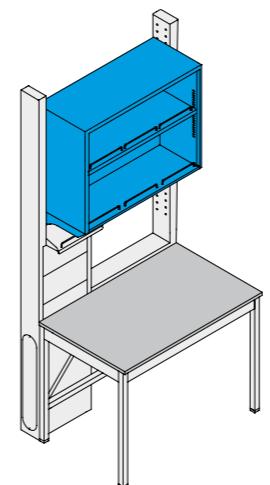
Standard heights for overhead cabinets are 600 and 800 mm. Different dimensions are available upon specific request. Cabinets have an internal height-adjustable shelf. Starting from the overhead cabinet for bench length 1500, cabinets have an internal partition.

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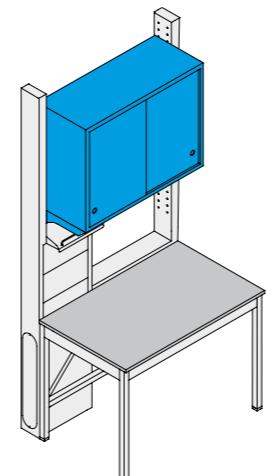
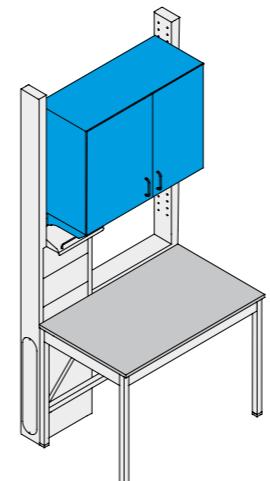
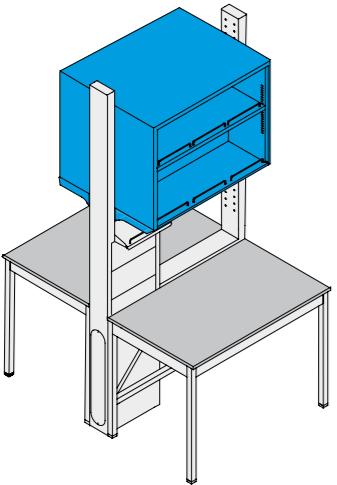
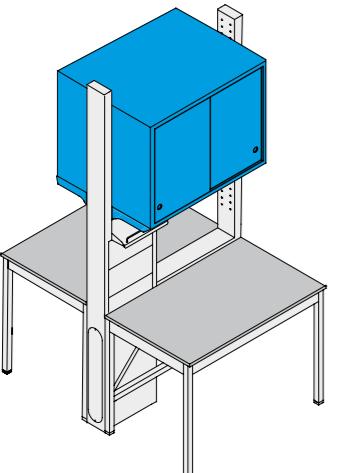
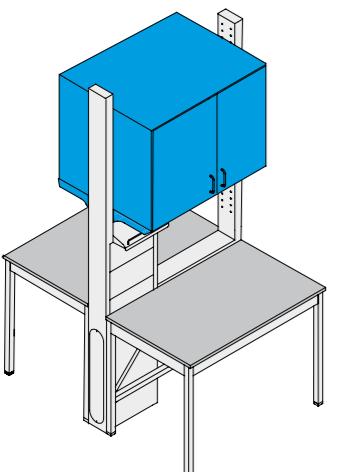
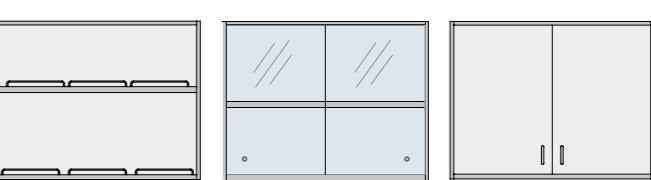
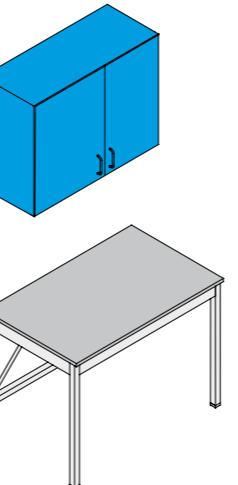
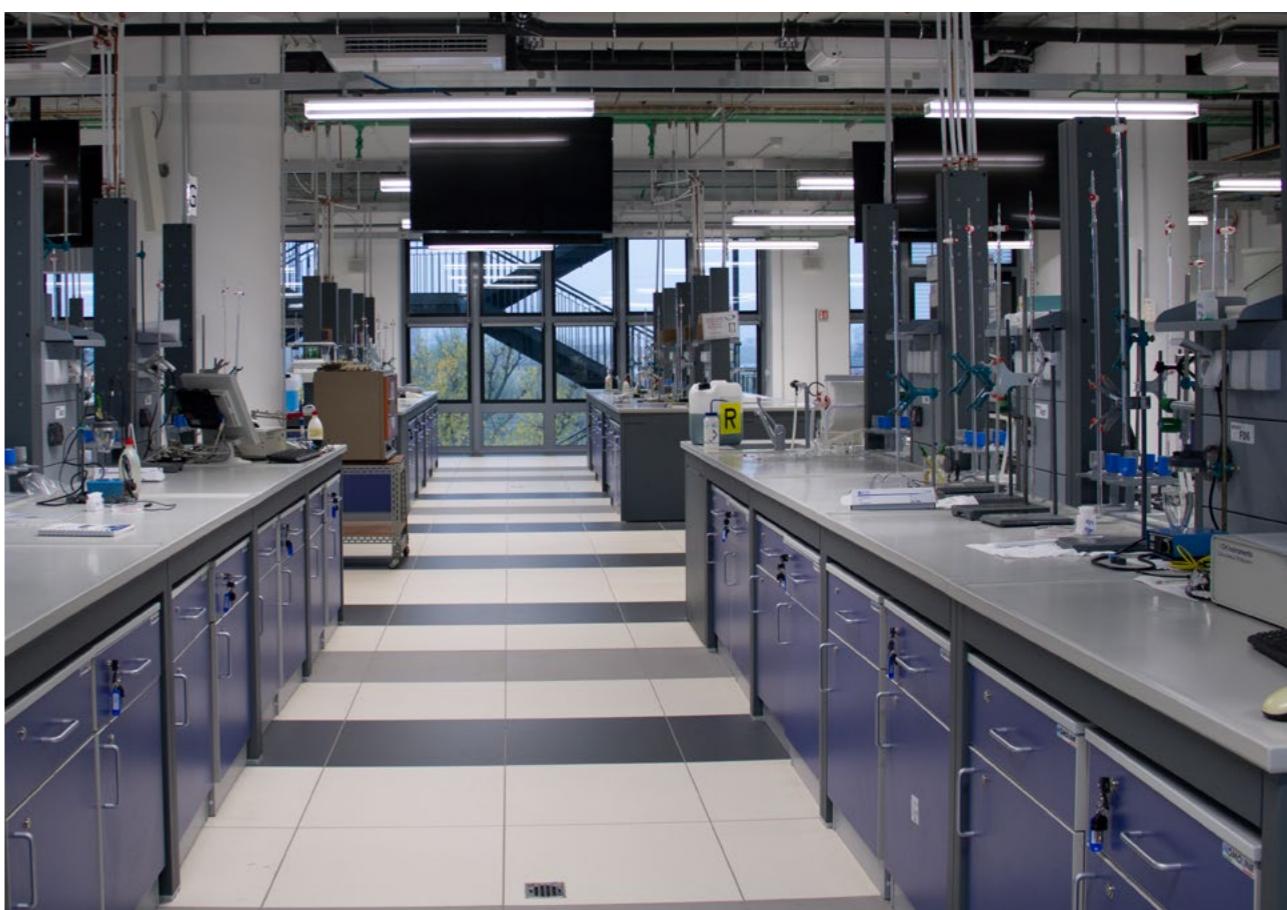
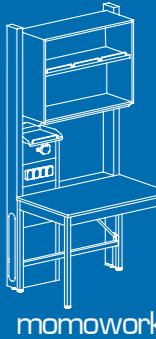
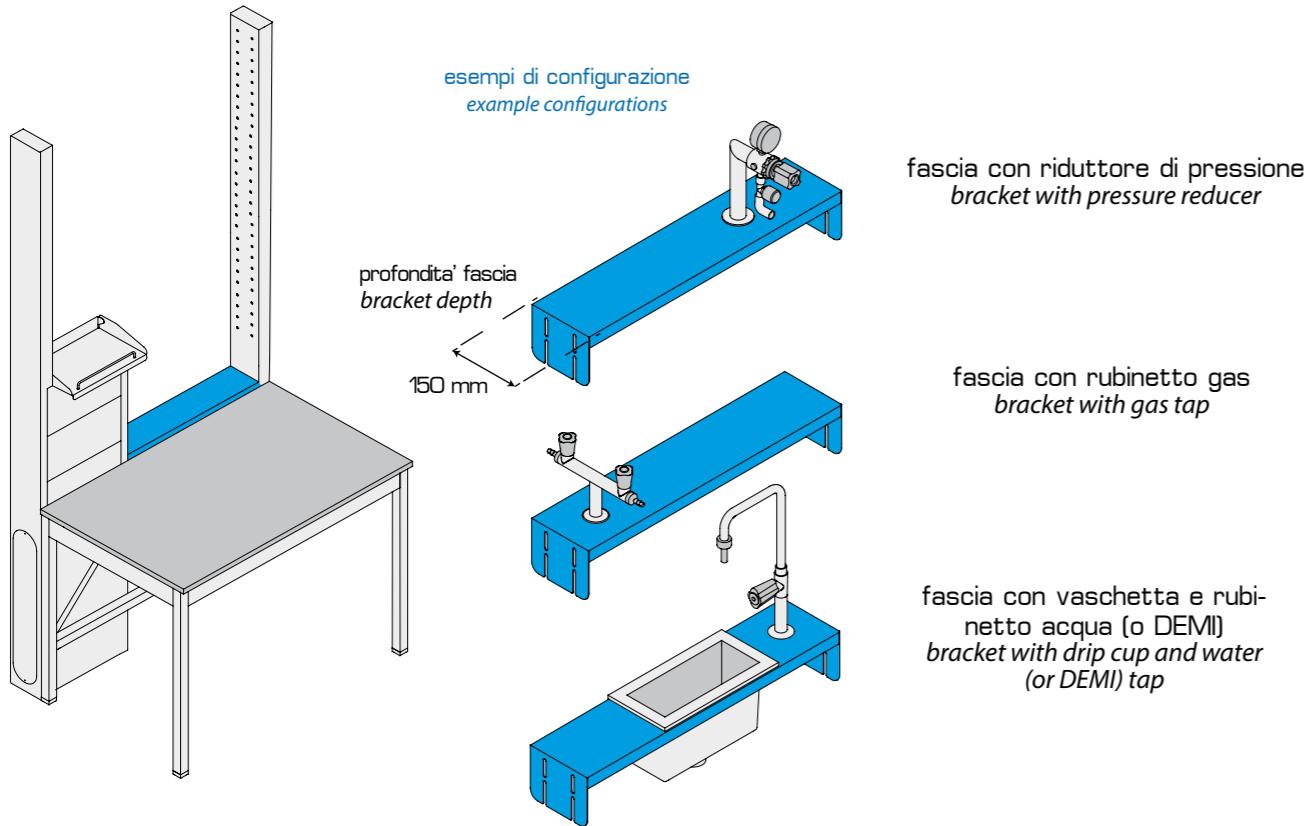
P___-WLL-

pensile a muro / overhead cabinet for wall fastening

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Shown configurations for brackets are only suggestive. Customer can choose the desired configuration and all accessories can be placed to the right or to the left of the technical panels.

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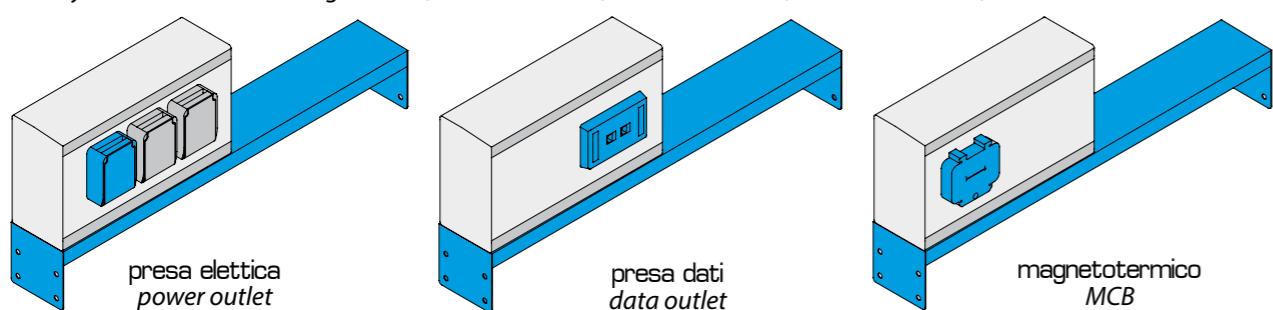
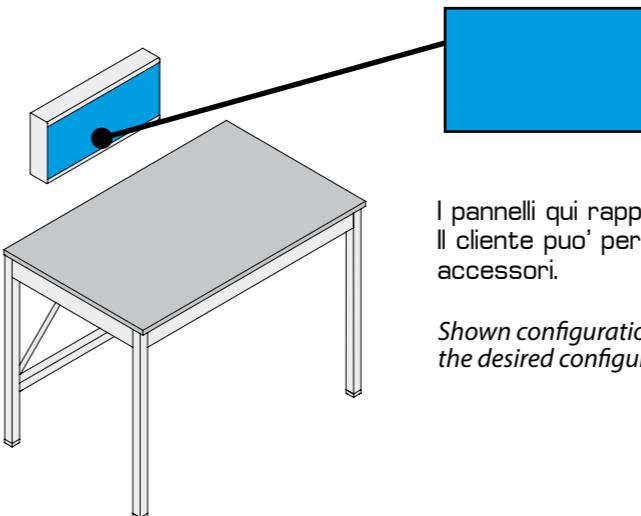
Utility boxes can be either single-sided(for wall benches) or double sided (for island benches).

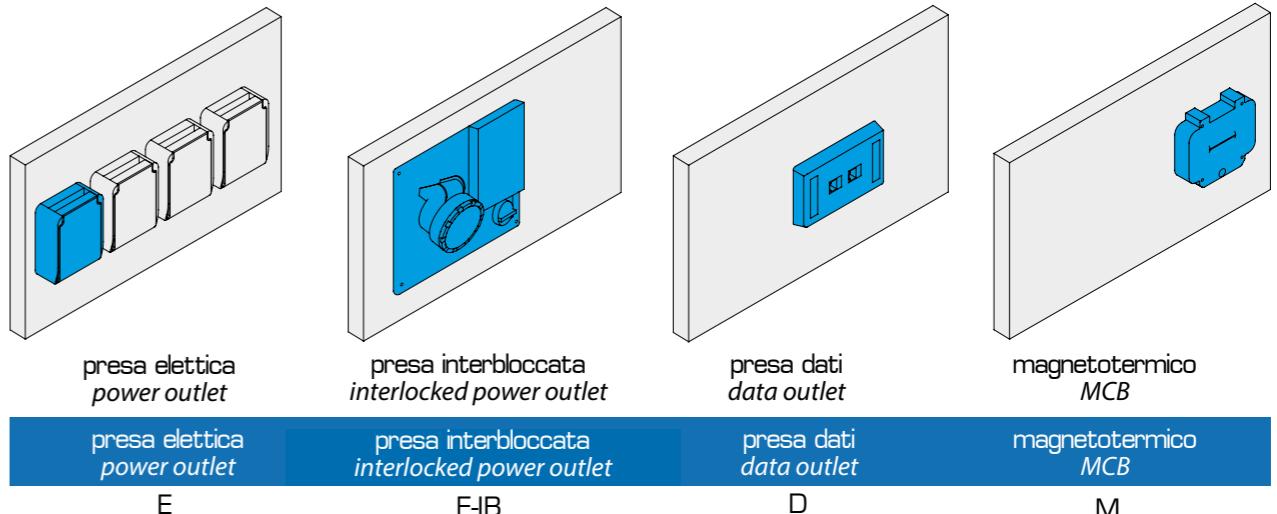
torretta utenze a muro /utility box installed on wall

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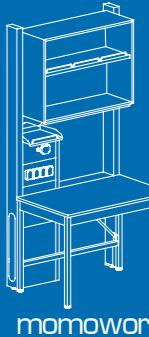
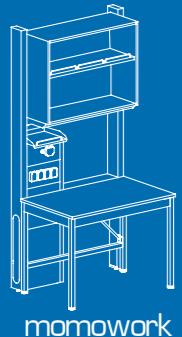
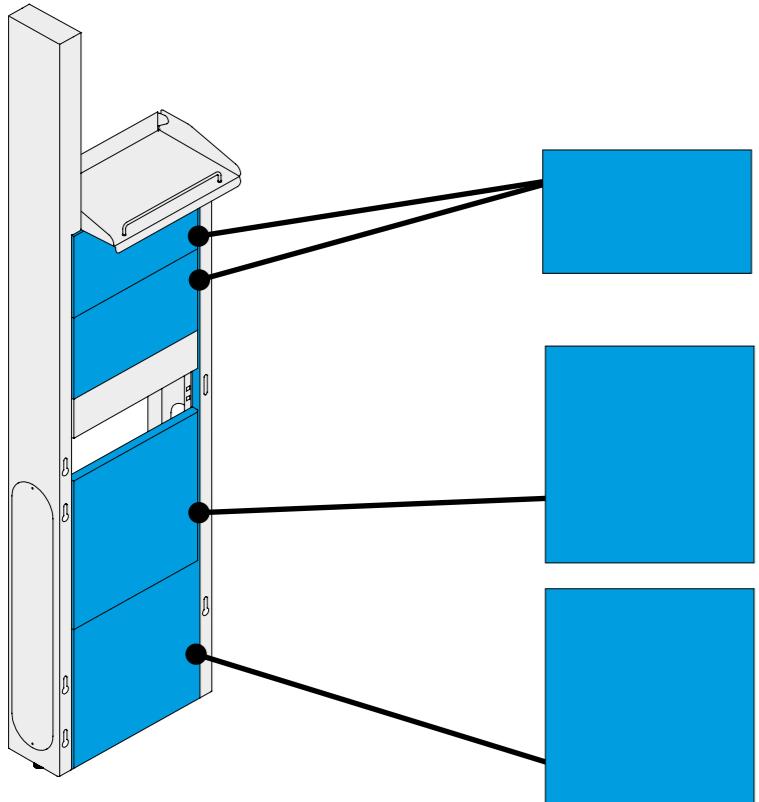
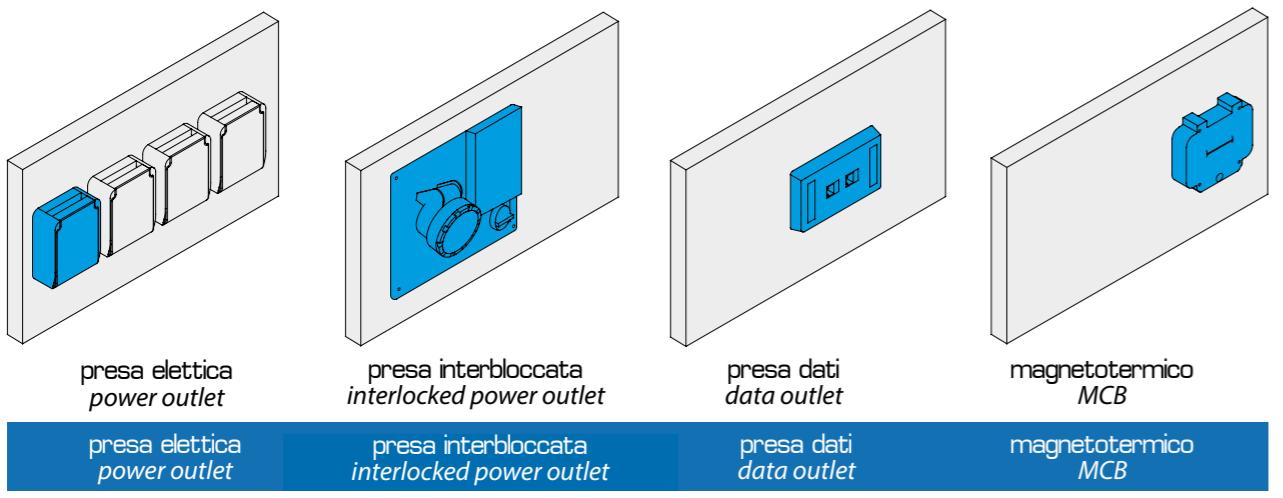
benches con parete tecnica / benches with service panels

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*Service panels come with 380x250 mm shelf, and can be equipped with all services and outlets shown in this page.
All panels are interchangeable (modular) and are fastened without the need of any tool.*

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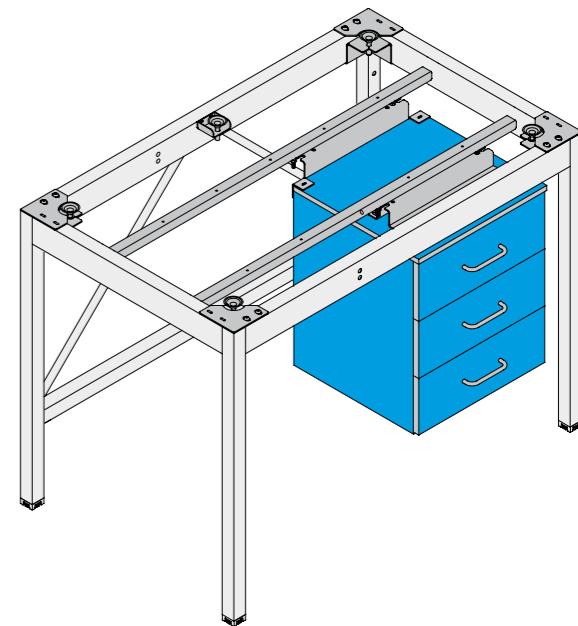
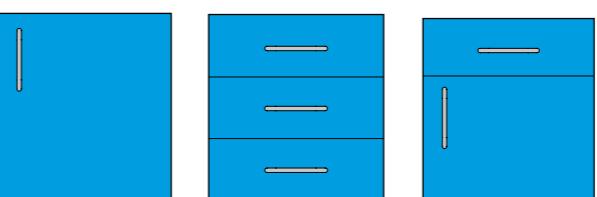
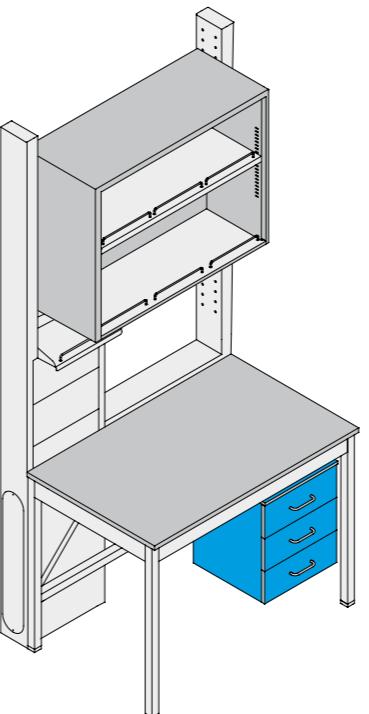
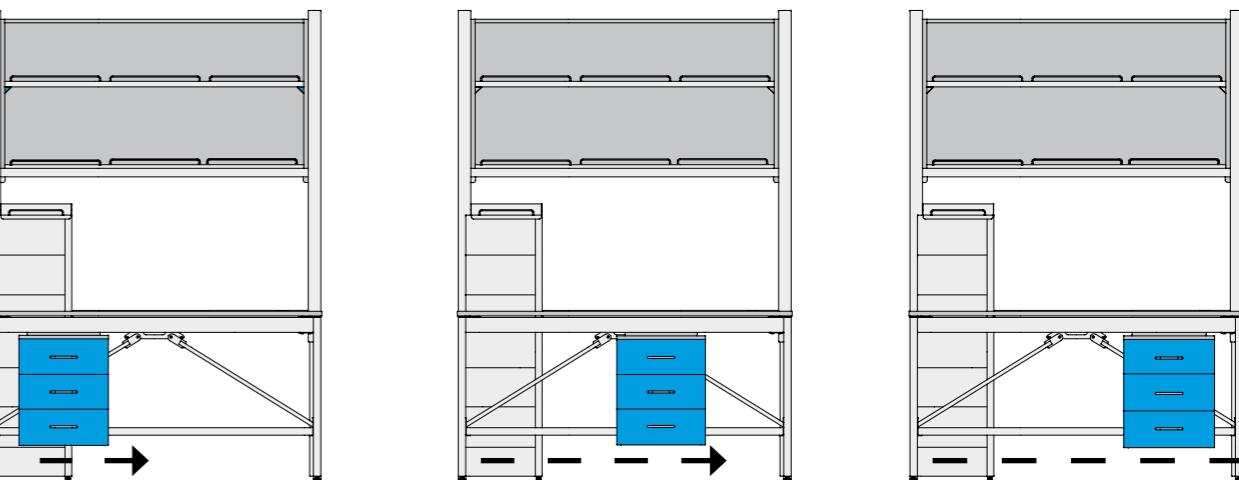
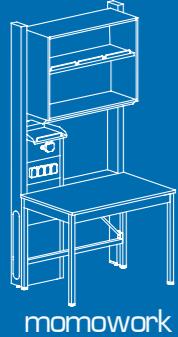
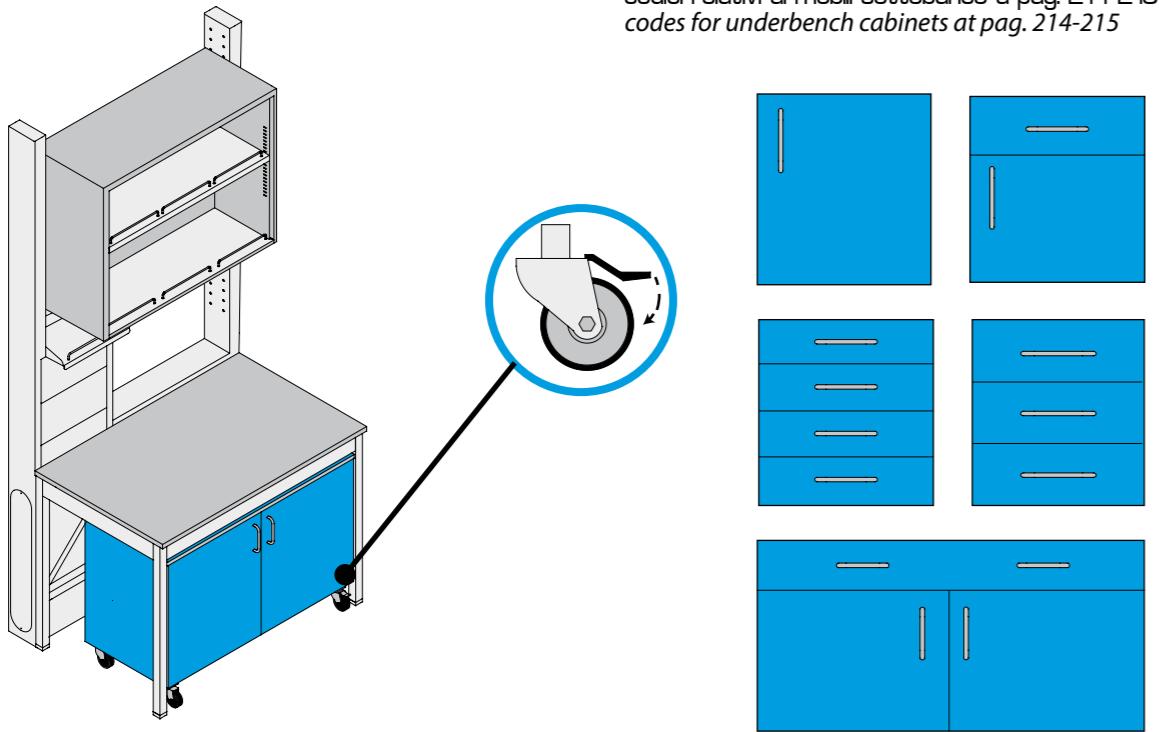
*Suspended underbench cabinet can be installed on the
MOMOWORK bench line.
Bench structure is equipped with two rails to allow the cabinet's
movement below the working surface.*

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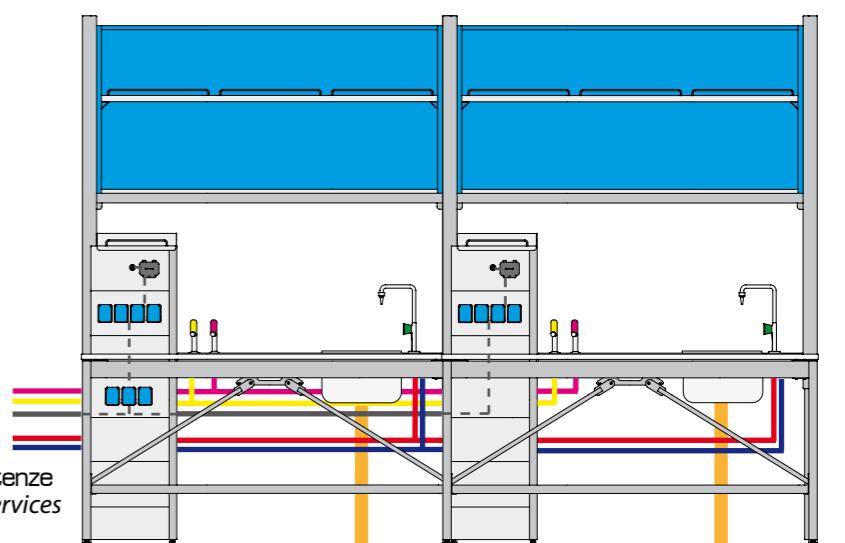
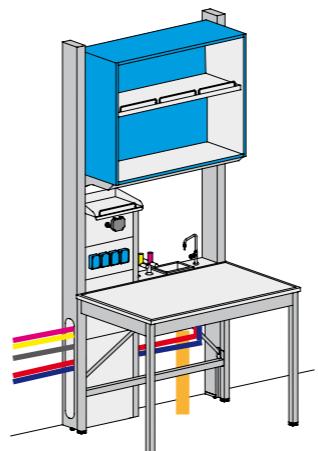
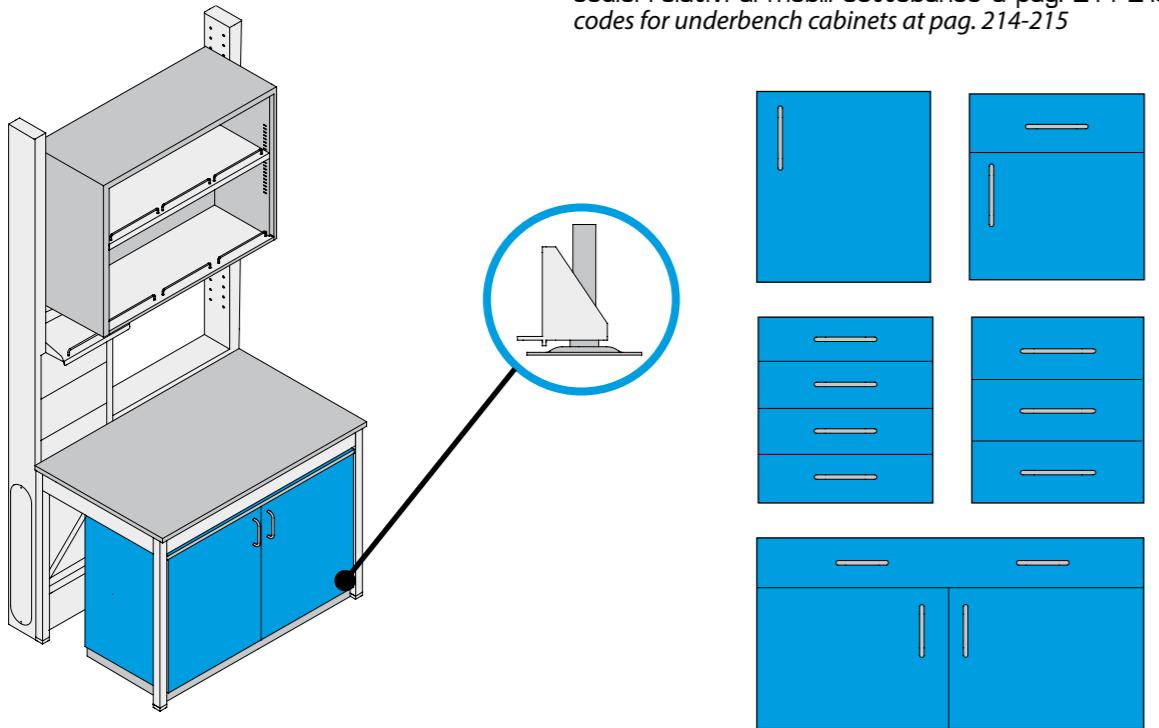
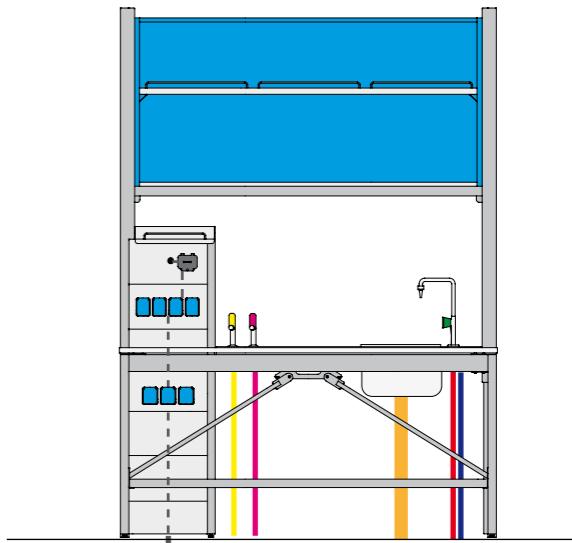
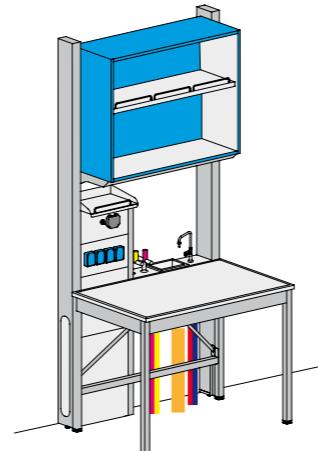
*Example of furniture installation with inlet of services from side.
In this case a steel back panel is provided to conceal the lines (see pag.21).*

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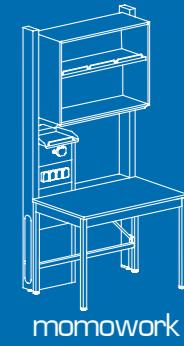
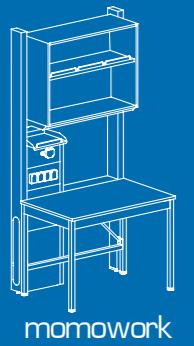
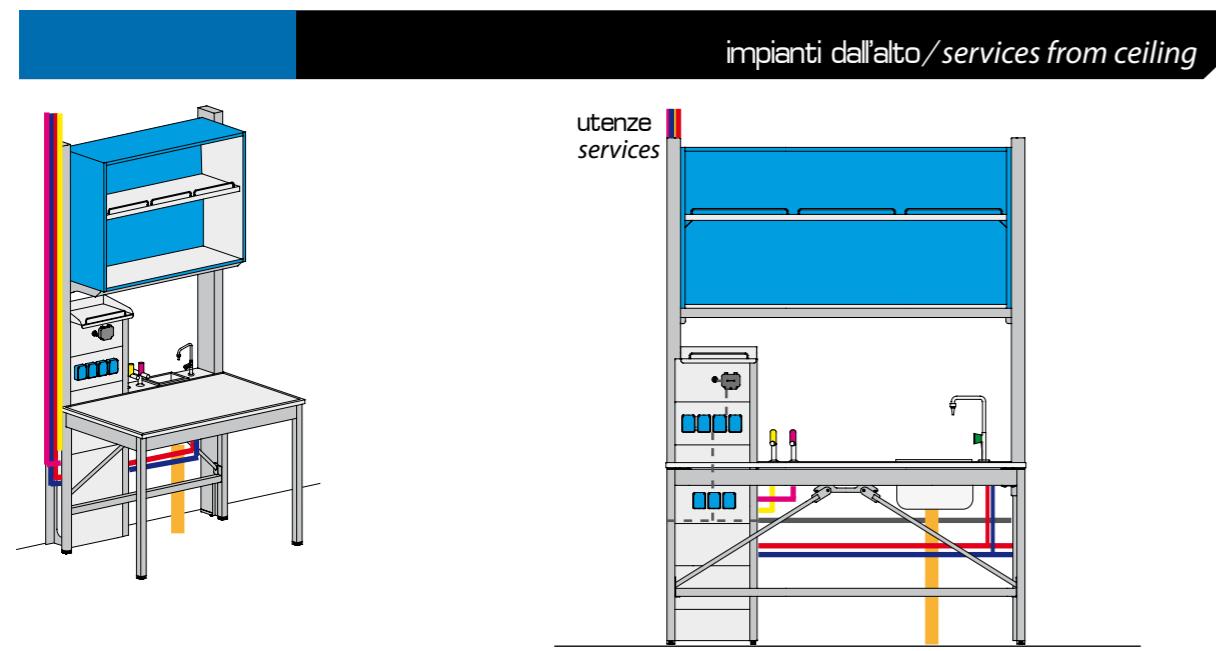
*Example of furniture installation with inlet of services from floor.
In this case a steel back panel is provided to conceal the lines (see pag.21).*

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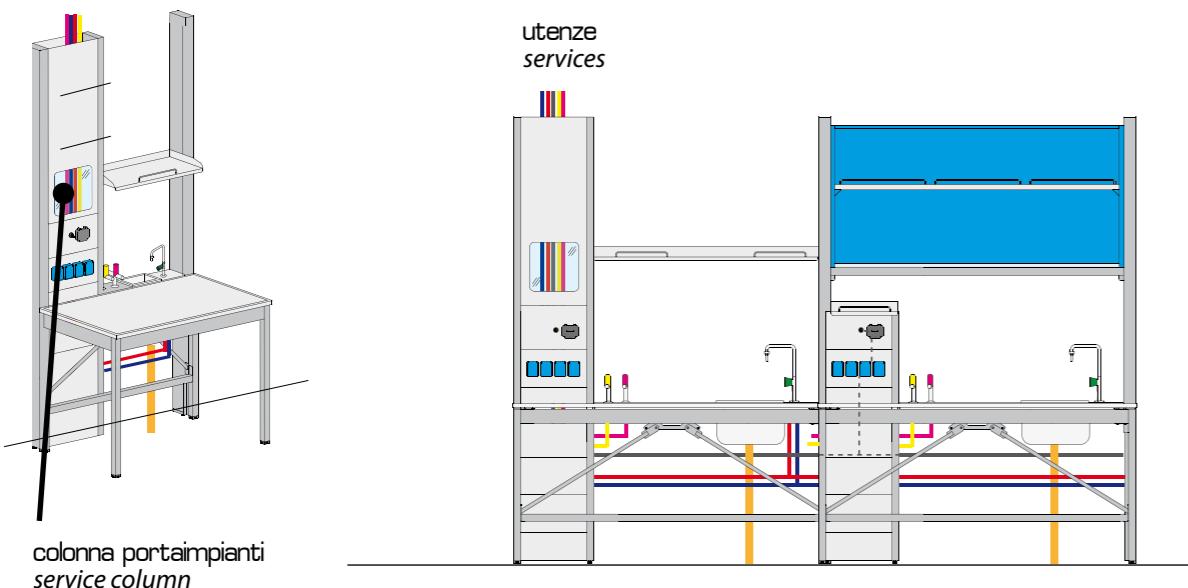
Example of furniture installation with inlet of services from ceiling. In this case a dedicated carter is fitted to the pillar to conceal the lines.

A steel back panel (see pag. 16) can be provided to conceal the lines.

Alternatively, a service column up to the ceiling with inspection window can be provided.

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MOMOWORK Plus laboratory benches and desks represent the best solution for the most demanding customers, aiming to set-up an elegant and ergonomic laboratory, at the cutting edge of technology, being at the same time safe, reliable and efficient.

MOMOWORK Plus benches have "C" shaped side frames, entirely realized using laser-cut steel sheets, bended through cold processes and then coupled together. The side frames are then connected through three horizontal elements, and the structure is finally completed and strengthened by three connecting rods, forming a triangle, in order to absorb the stress components in all directions. The steel structure is then epoxy powder coated: this grants the 'Class O' fire reaction, as well as high resistance against corrosion.

The resulting structure is absolutely solid and reliable, making it possible to accommodate even the heaviest instruments and machines: **MOMOWORK Plus** laboratory benches and desks hold Certificates stating the compliance to the European Norm EN 13150:2001. The compliance certificates have been issued by Bureau Veritas, authorized third organization officially recognized at international level.

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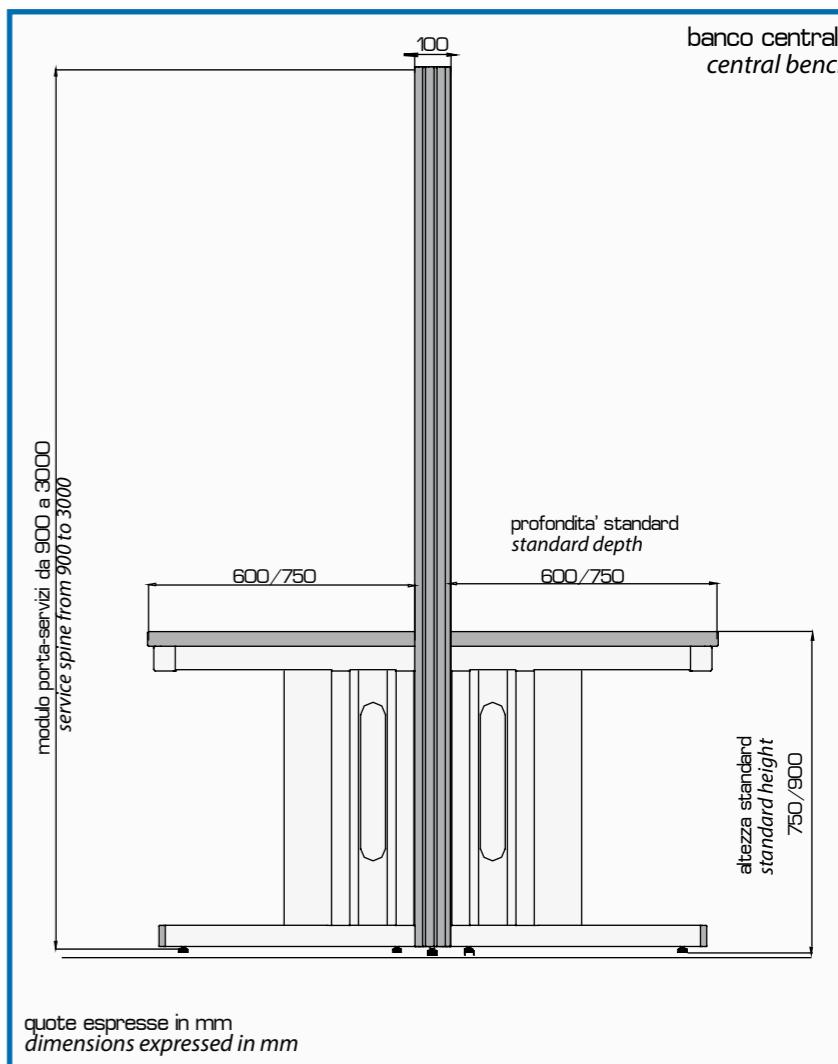
The compliance certificate of **MOMOWORK Plus** laboratory benches states the resistance to a 800 kg load placed on the working surface, considerably exceeding the minimum prescriptions of the EN 14175:2001 Norm, recommending a 200 kg maximum load.

MOMOWORK Plus laboratory benches and desks have a fully modular structure, very easy to assemble and de-assemble, so that even future improvements or moving are possible.

The structure has dedicated compartments to allow the routing of utilities and their connection to the delivery point mounted on furniture.

The basic structure can be equipped with special technical panels, housing all the necessary services and pick-up points for technical gases, power, data etc. The pick-up points are installed on interchangeable steel panels, fitted to the main structure through quick fitting devices, allowing to perform of inspection and maintenance operations with no need of tools, as well as making future changes easier.

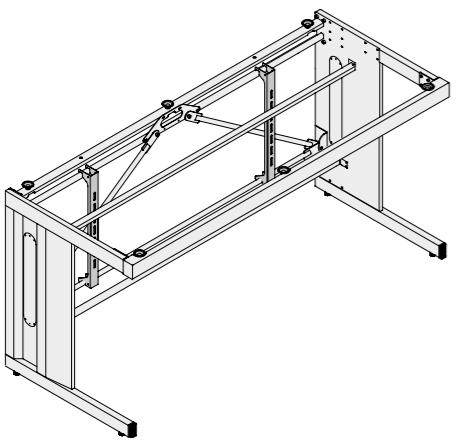
Benches and desks can be then completed by a wide range of items such as upper cabinets, shelves, reagent racks and underbench cabinets.

MOMOWORK Plus laboratory benches are also available in the anti-vibrating version, ensuring best active and passive insulation performances against vibrations: this is the optimal solution in case highly precise instruments have to be employed, or even for the most demanding microscopy applications.

The standard height of the worktop is 900 mm in case of conventional laboratory benches, to be used in standing position, and 750 mm in case of desk application, to be used in seated position.

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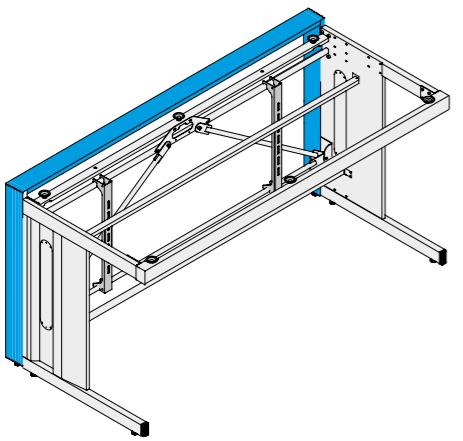
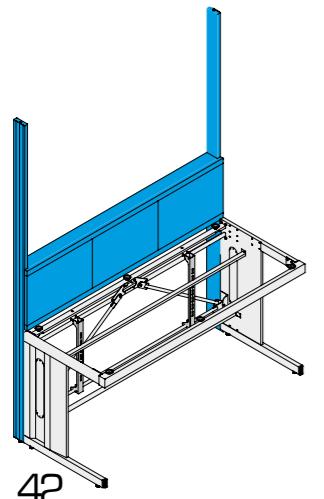
codice/ <i>code</i>	lunghezza/ <i>length</i>
BP60-S-MWP	600
BP80-S-MWP	800
BP90-S-MWP	900
BP100-S-MWP	1000
BP120-S-MWP	1200
BP150-S-MWP	1500
BP180-S-MWP	1800

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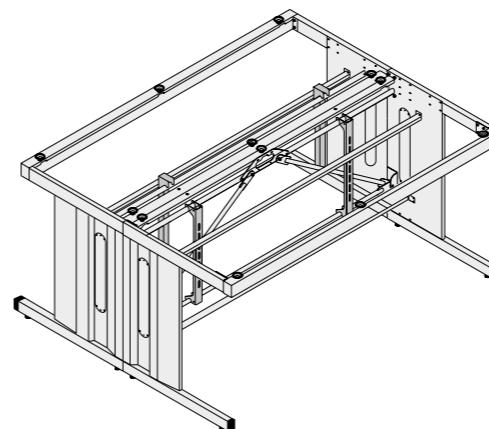
codice/ <i>code</i>	lunghezza/ <i>length</i>
BP60-F-MWP	600
BP80-F-MWP	800
BP90-F-MWP	900
BP100-F-MWP	1000
BP120-F-MWP	1200
BP150-F-MWP	1500
BP180-F-MWP	1800

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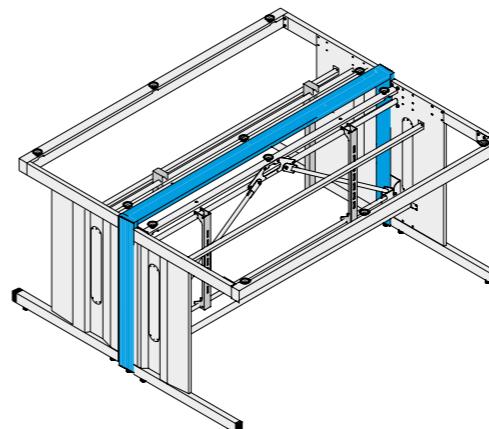
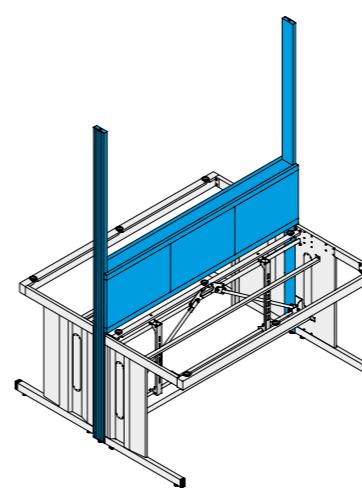
codice/ <i>code</i>	lunghezza/ <i>length</i>
BC60-S-MWP	600
BC80-S-MWP	800
BC90-S-MWP	900
BC100-S-MWP	1000
BC120-S-MWP	1200
BC150-S-MWP	1500
BC180-S-MWP	1800

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codice/ <i>code</i>	lunghezza/ <i>length</i>
BC60-F-MWP	600
BC80-F-MWP	800
BC90-F-MWP	900
BC100-F-MWP	1000
BC120-F-MWP	1200
BC150-F-MWP	1500
BC180-F-MWP	1800

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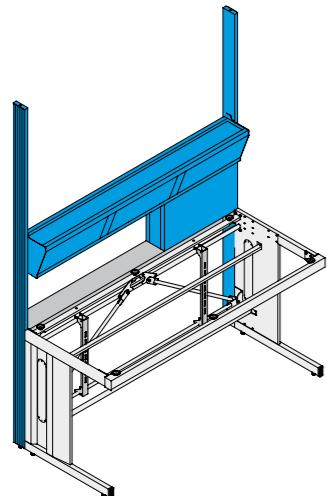
BP___-BR-MWP

struttura banco bridge/bridge bench structure

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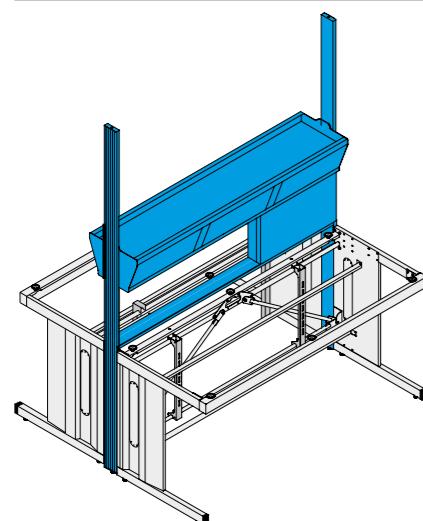
BC___-BR-MWP

struttura banco centrale con bridge/central bench structure with bridge

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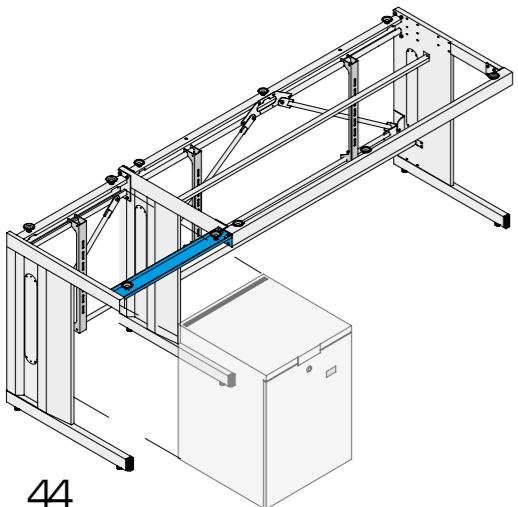
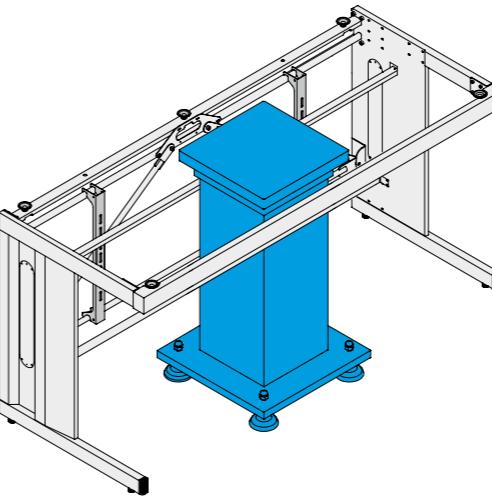
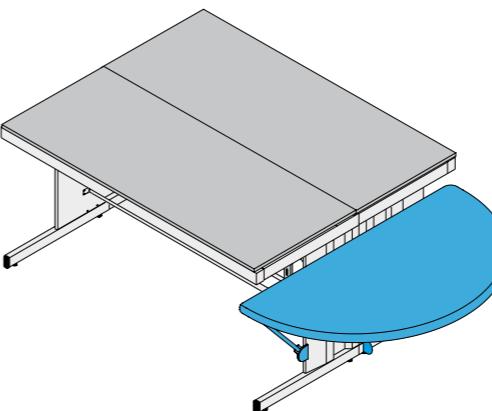
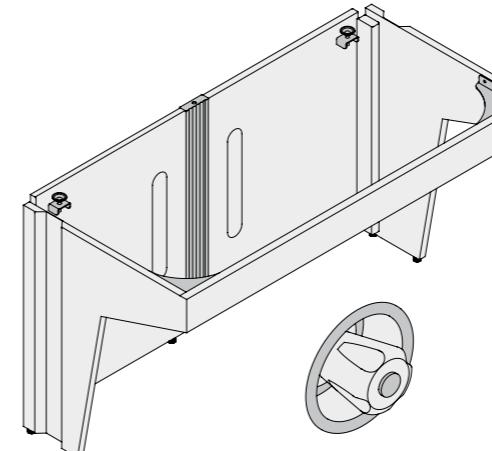
Table for desk or storage employment, for installation besides central bench, series Momowork Plus.

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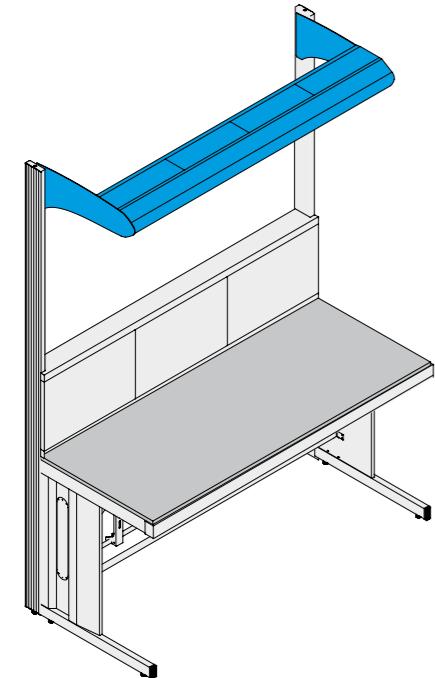
This module is for installation besides island benches. If a sink is requested, it is possible to install a remote control for the tap (code Q-RC).

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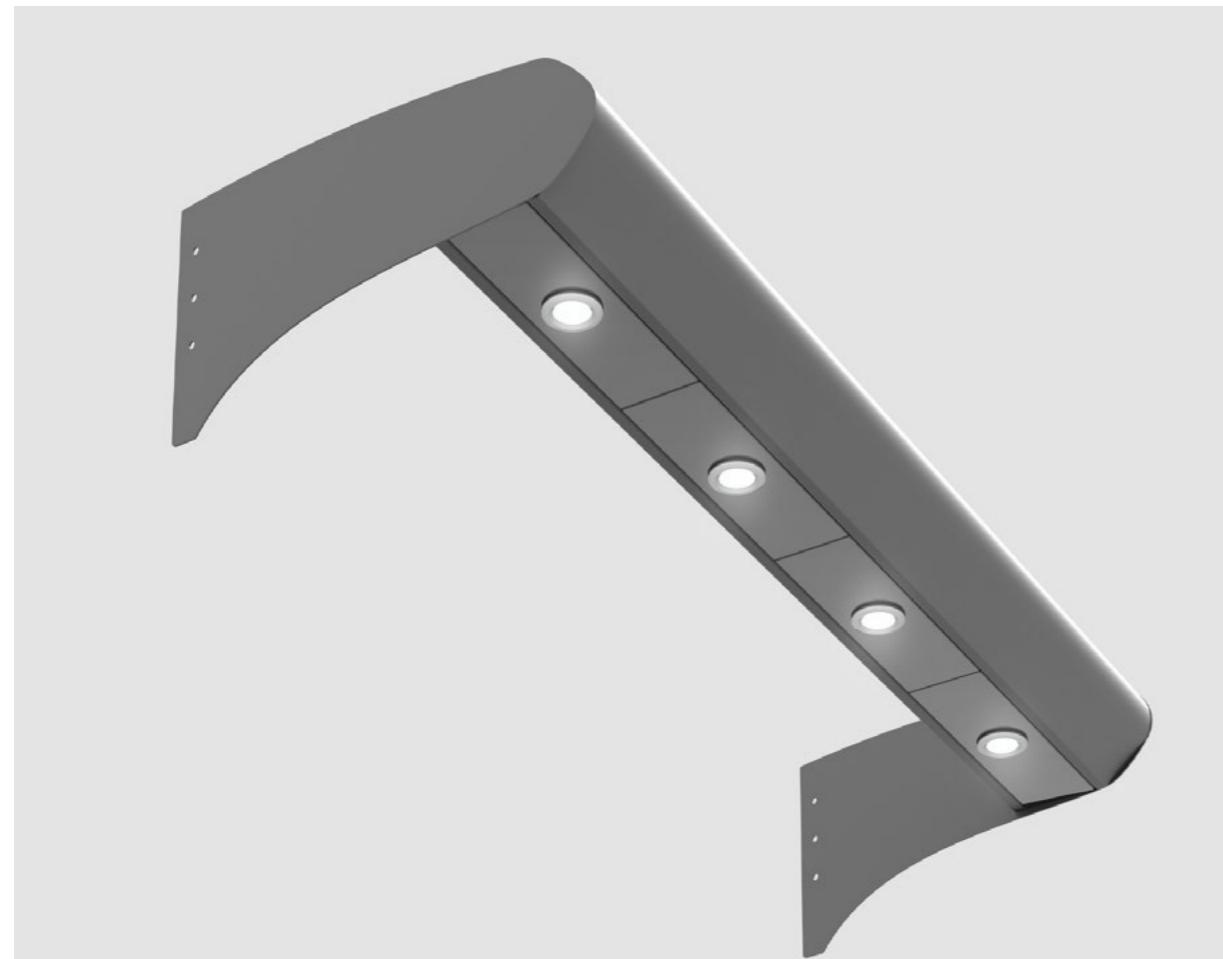
Skylight is a module fixed to the pillars of the bench embedding some lights for the working area of the bench.

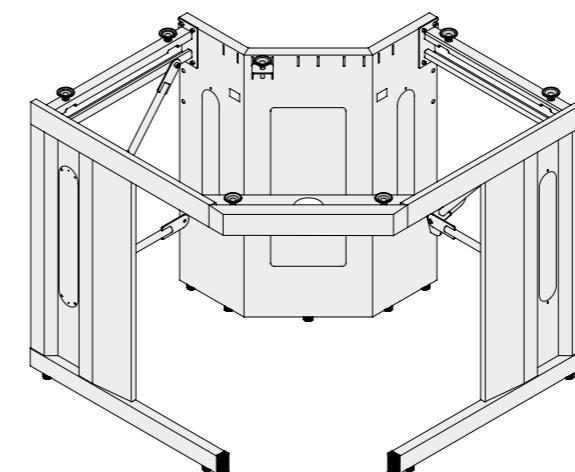
Skylight is realized in epoxy powder coated steel and is equipped with modular panels with LED lights.

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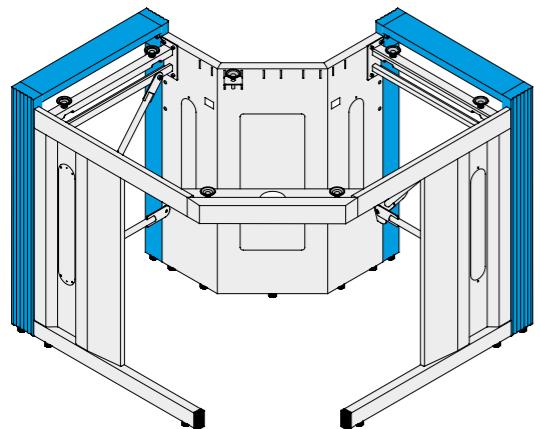
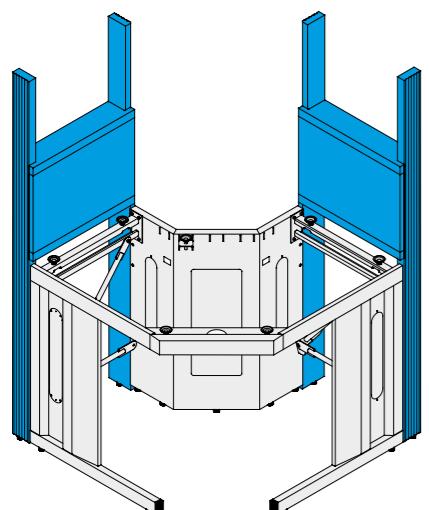
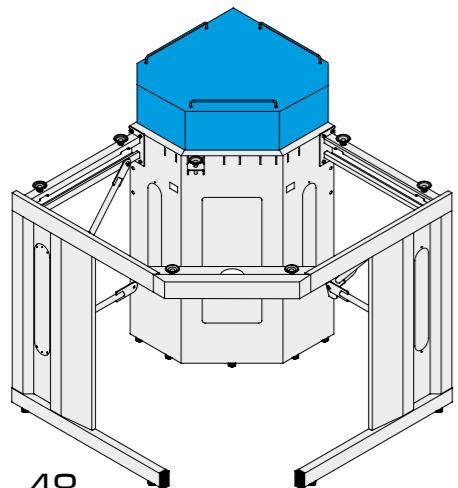
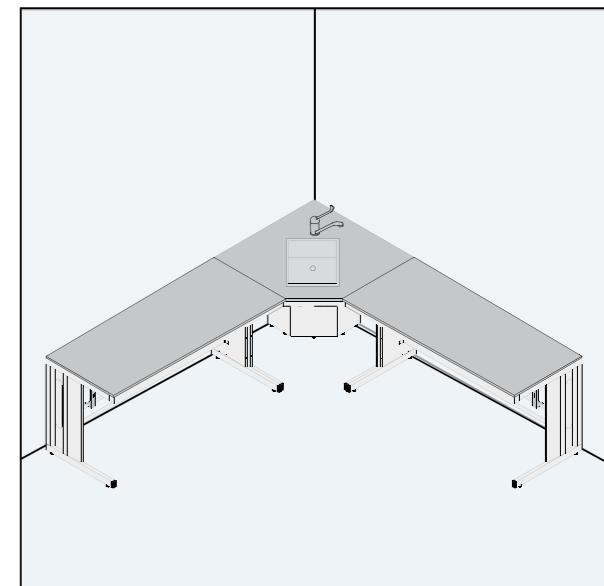
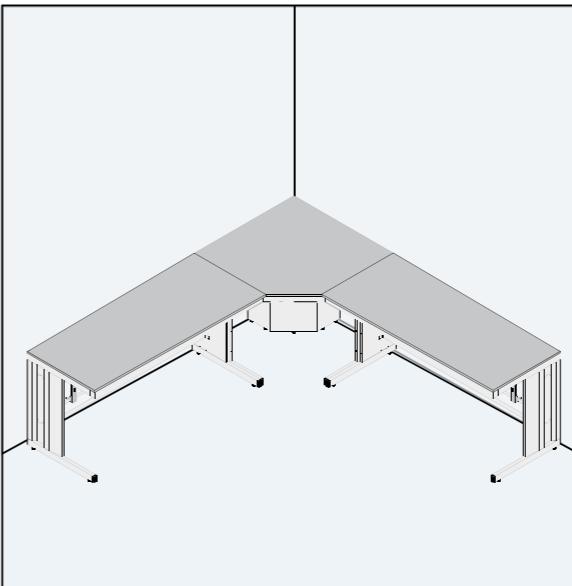
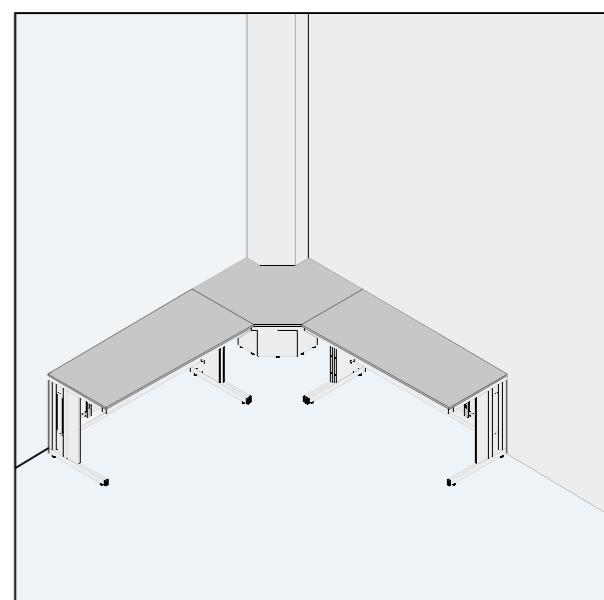
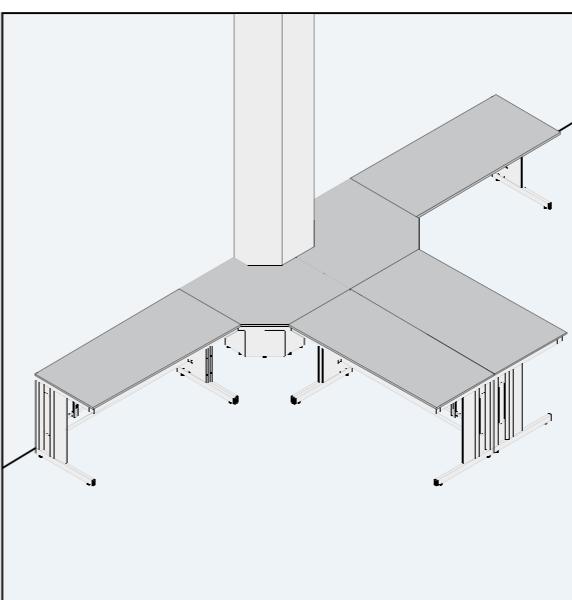
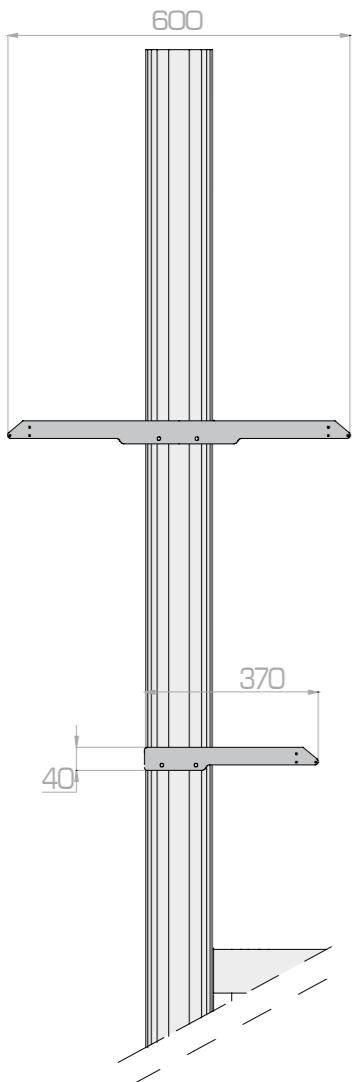
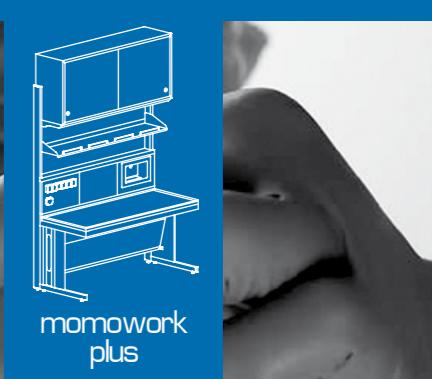
The special design of the corner bench allows to maximize the employment of worktop space on the benches, avoiding any "dead area" with difficult access. An example of optimized employment for this bench is the installation of sinks.

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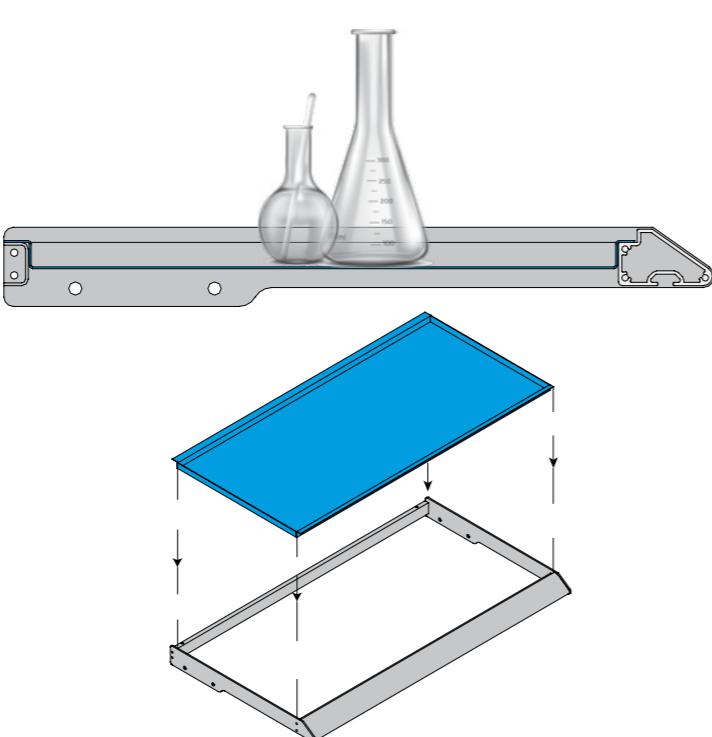
In case of necessity, the corner bench structure can be used to circumvent an architectural constraint imposed by the building (e.g: a pillar).

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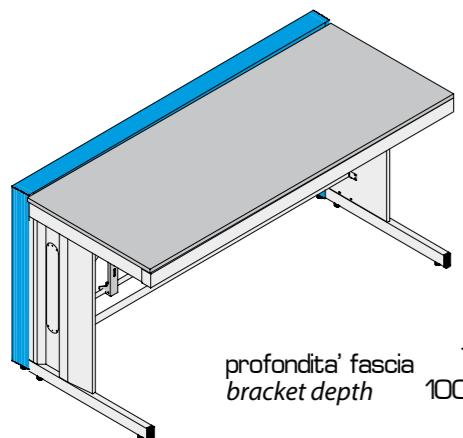
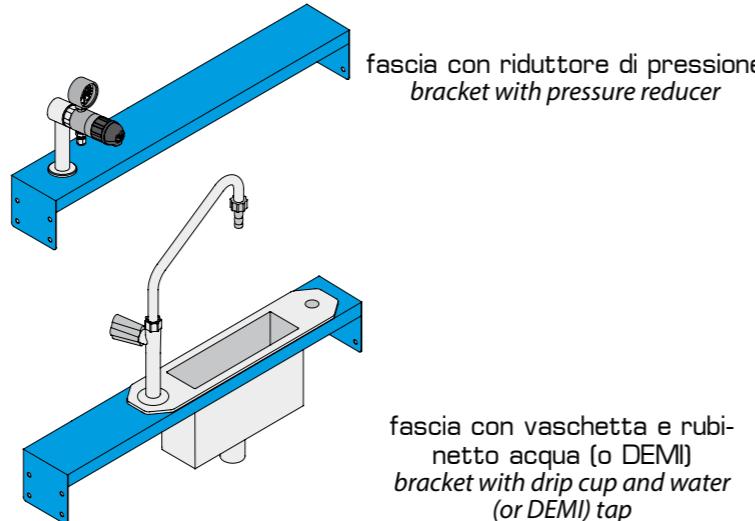
benches 1

benches con fascia / benches with bracket

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The shown configurations for brackets are only suggestive. Customer can choose the desired configuration and all accessories can be placed to the right or to the left of the spine.

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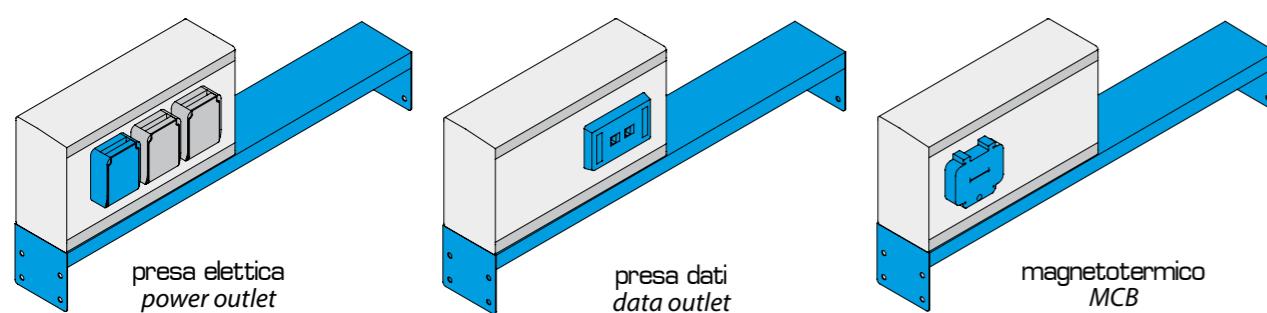
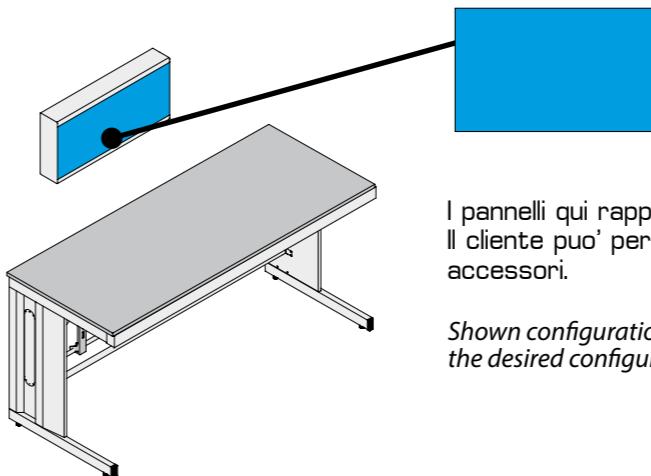
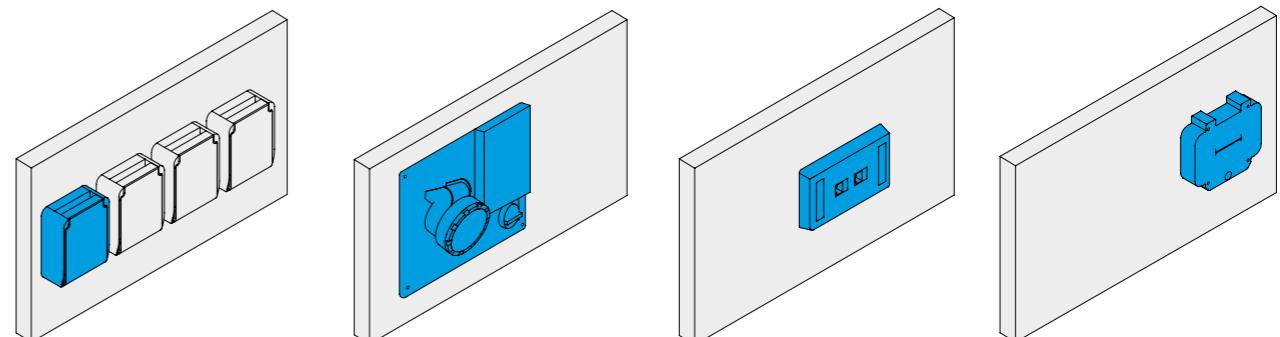
Utility boxes can be either single-sided (for wall benches) or double-sided (for island benches).

torretta utenze a muro / utility box installed on wall

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Shown configurations for panels are only suggestive. Customer can choose the desired configuration.

esempi di allestimento dei pannelli
example configurations for panelspresa elettrica
power outletpresa interbloccata
interlocked power outletpresa dati
data outletmagnetotermico
MCBpresa elettrica
power outletpresa interbloccata
interlocked power outletpresa dati
data outletmagnetotermico
MCB

E

E-IB

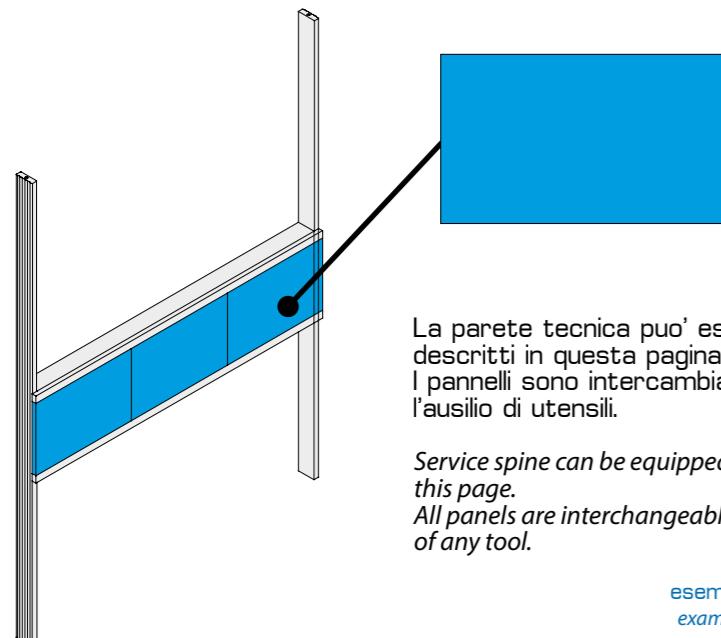
D

M

benches 1

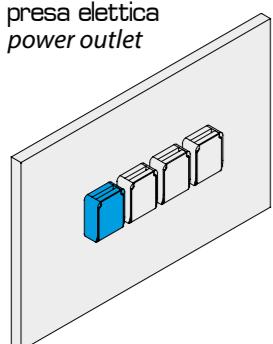
banchi con parete tecnica / benches with service panels

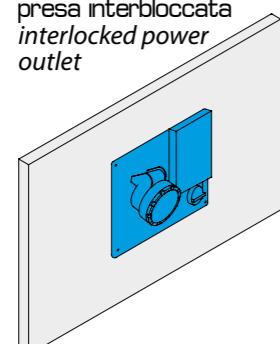
*Service spine can be equipped with all services and outlets reported in this page.
All panels are interchangeable (modular) and fastened without the need of any tool.*

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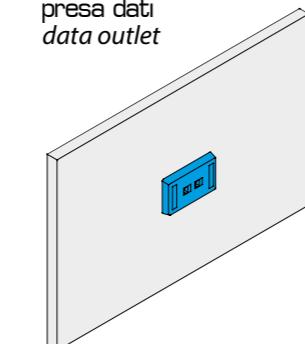
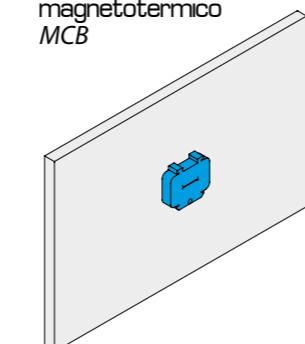
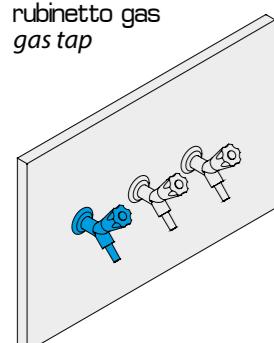
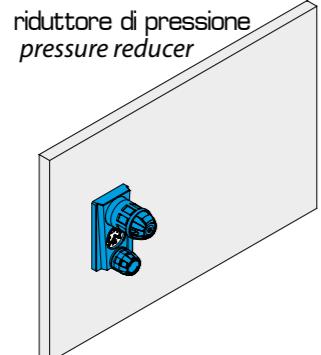
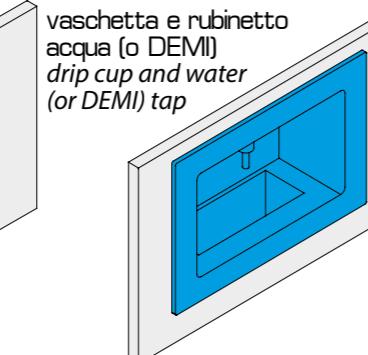
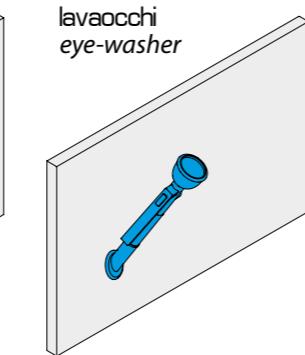
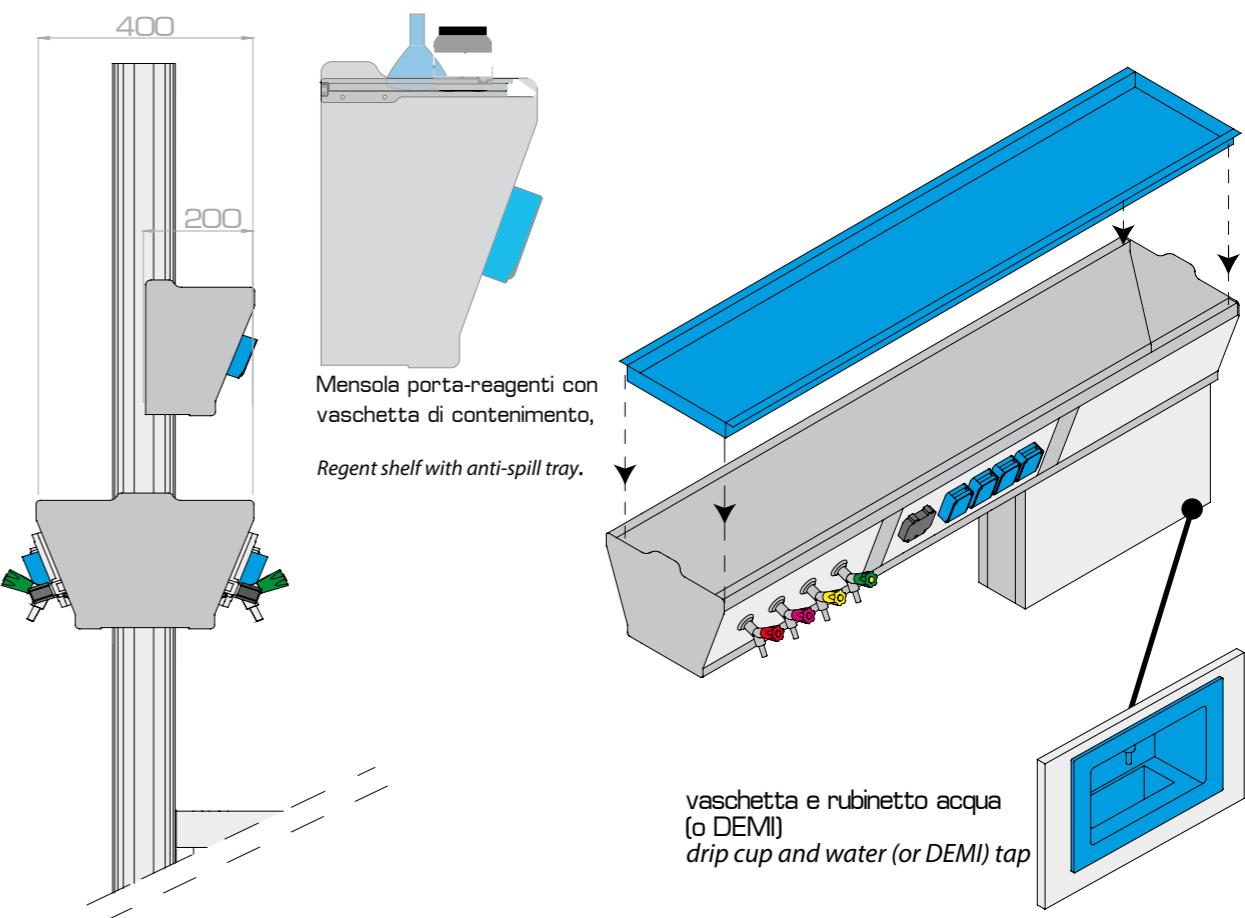
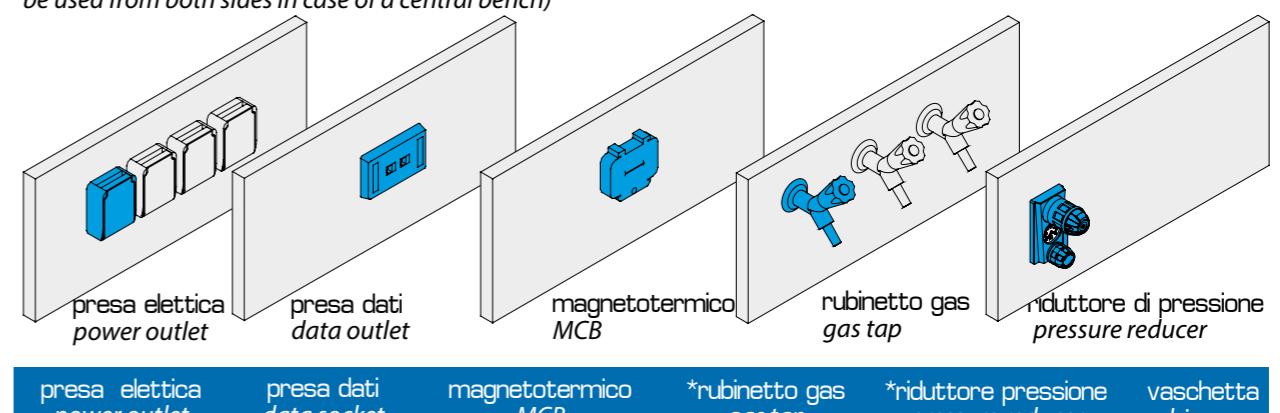
Benches with Bridge module are equipped with a dedicated technical compartment allowing the vertical passage of utilities. The standard compartment is plain, but it can be upon request fitted with drip cup and tap (the drip cup can be used from both sides in case of a central bench)

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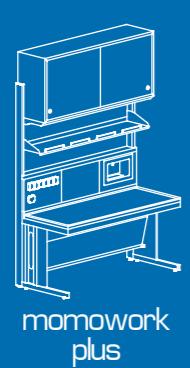
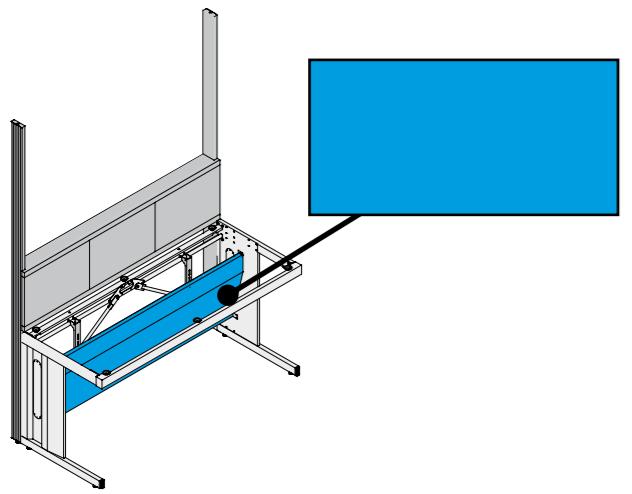
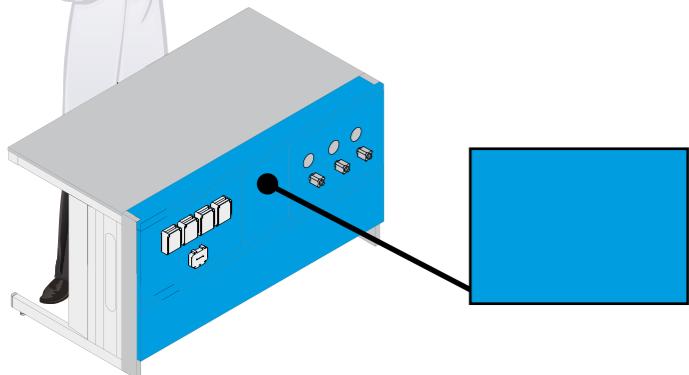
Epoxy-powder painted steel panel for concealing lines and services.

codice/code	per banco da/for bench length
FD60-MWP	600
FD80-MWP	800
FD90-MWP	900
FD100-MWP	1000
FD120-MWP	1200
FD150-MWP	1500
FD180-MWP	1800

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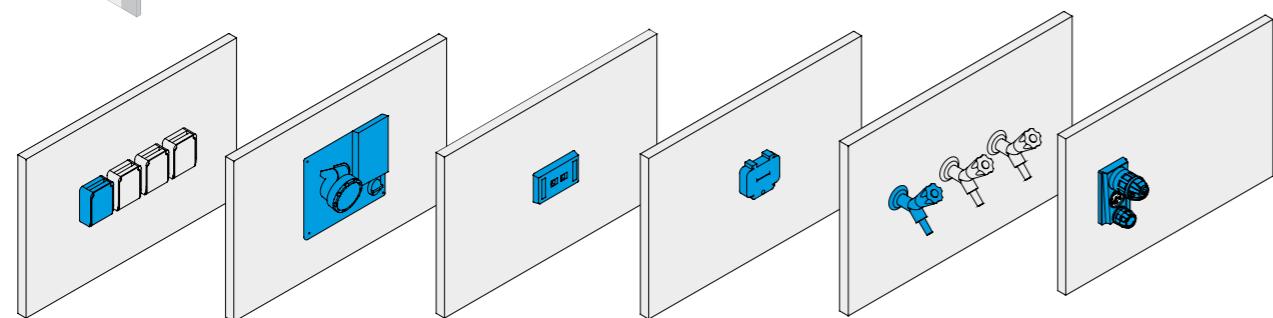
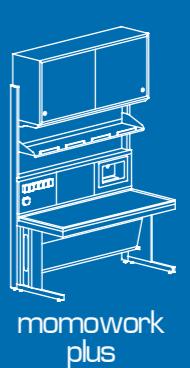
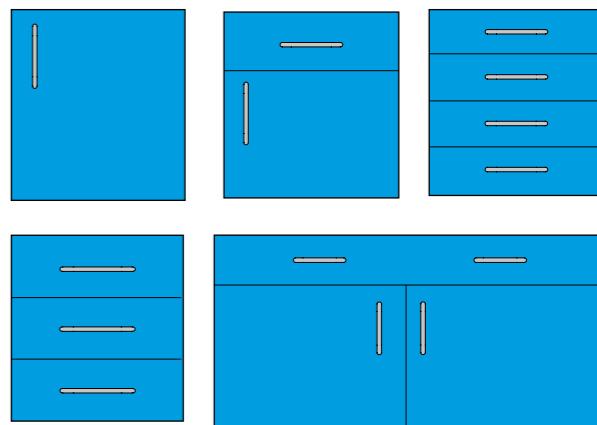
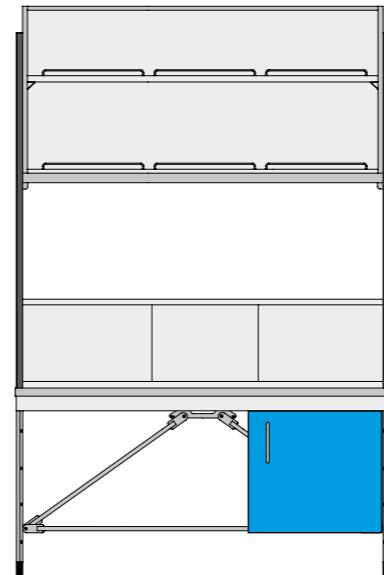
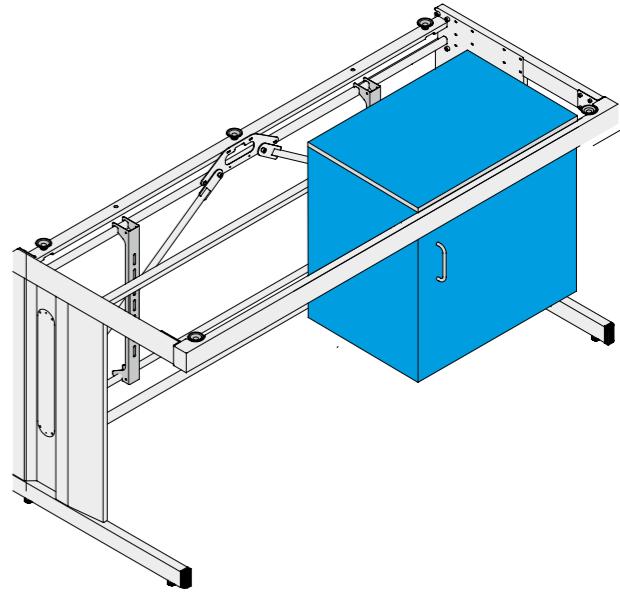
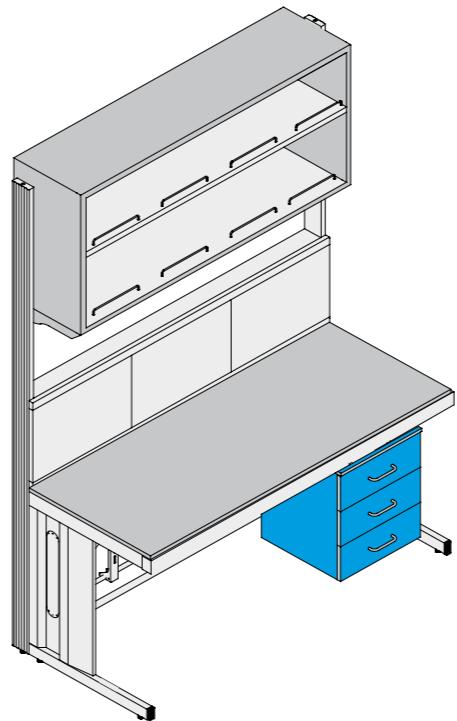
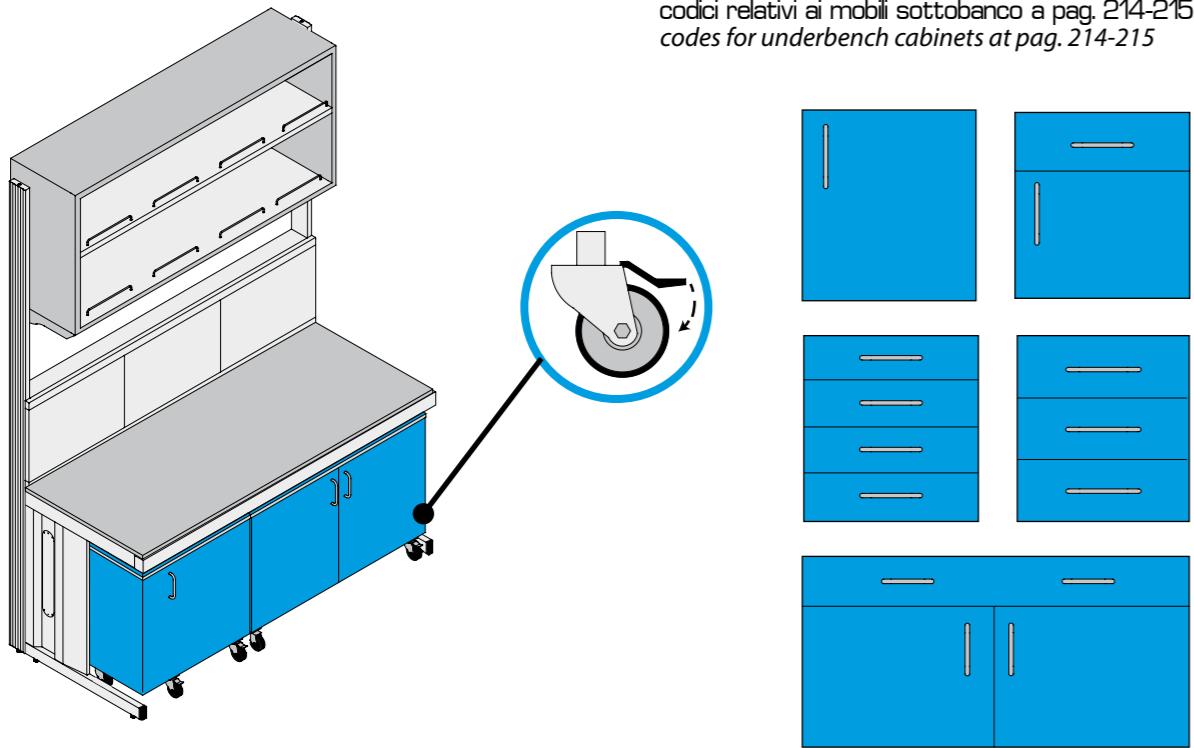
Suspended underbench cabinet can be installed on the **MOMOWORKPlus** bench line.
Bench structure of this line is equipped with two rails to allow the cabinet's movement below the working surface.

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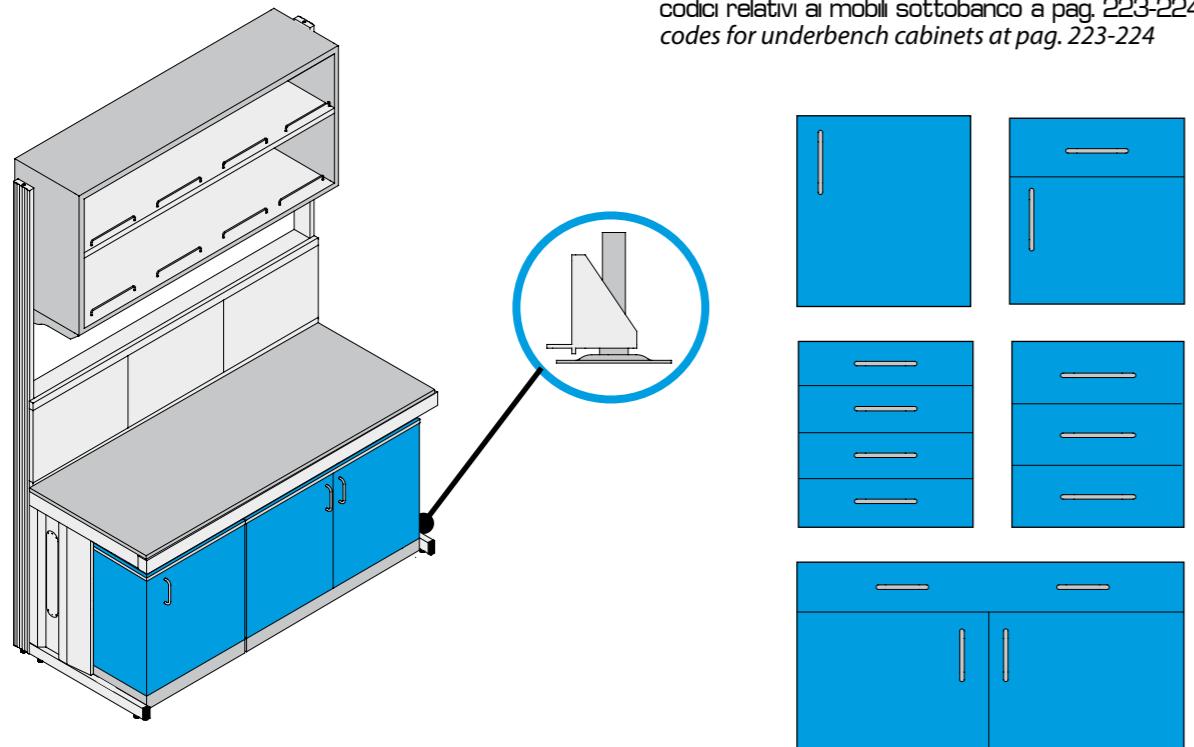
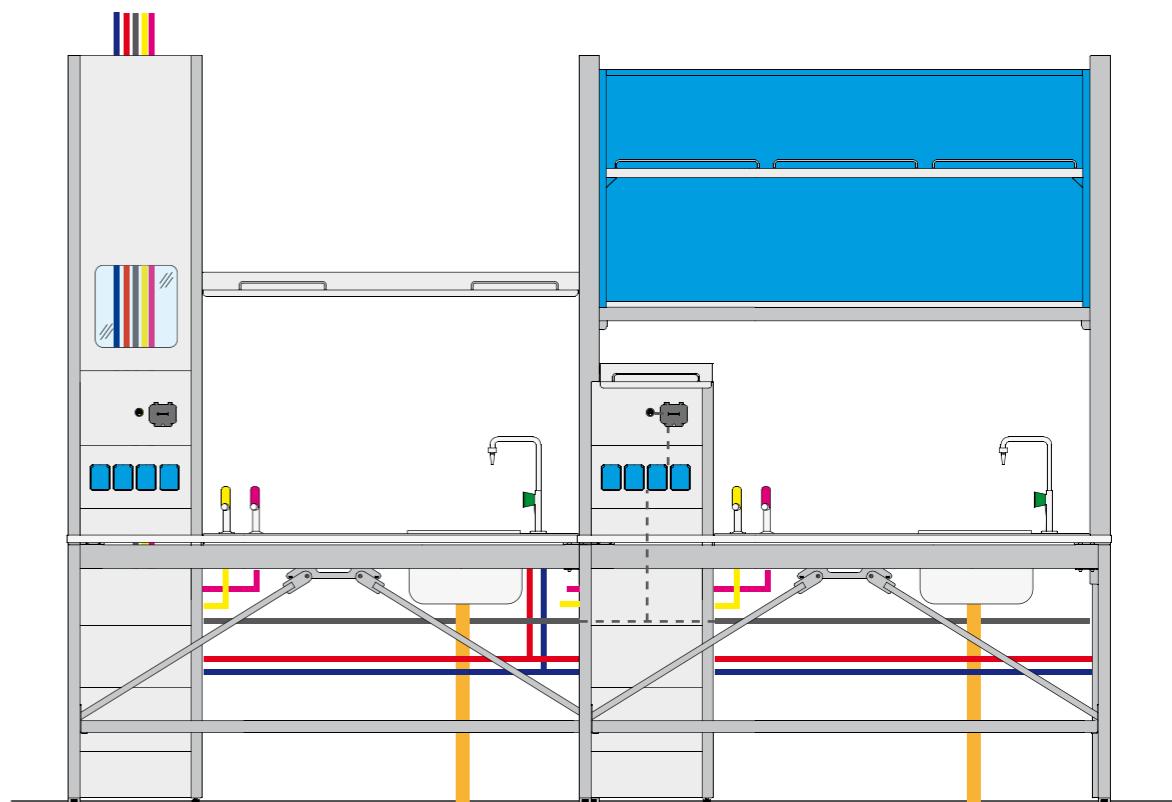
In case of services coming from ceiling the connection to the utilities on furniture is achieved through a special service column equipped with inspection window.
Alternatively, a side carter (trunk) for connection of utilities can be added (see pag. 67).

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Service columns allow the tie-in of the utilities installed on the benches to the services into ceiling/false ceiling.
Service columns are fitted, as a standard, with a compartment for utilities located at operator's height. This compartment, equipped with a window in safety glass, allows easy access for any maintenance operation, as well as for control/regulation on the utilities, since main insulation valves and pressure regulators are installed behind the glass for easy access and employment.
If necessary, the service column can be equipped with an internal partition used to separate electric and fluid utilities.

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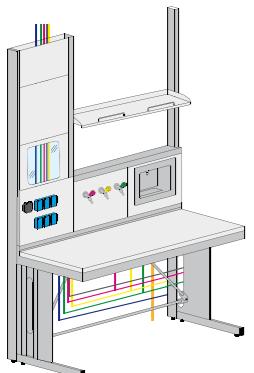
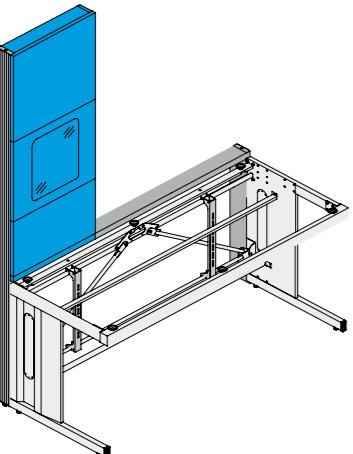
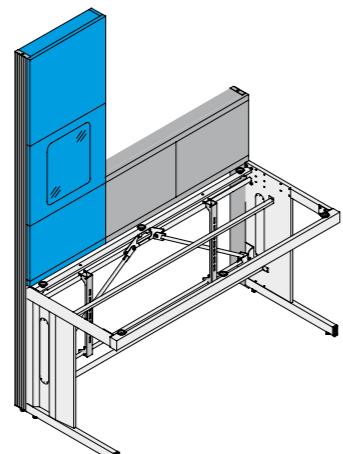
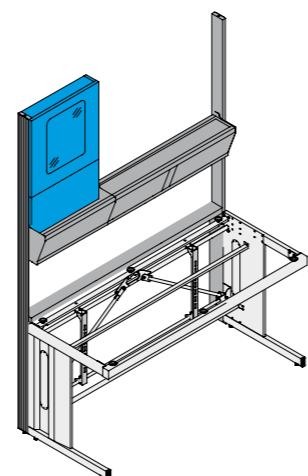
*Service column can be installed on all moduls of **MOMOWORK Plus** benches, fitted with bracket, technical panels or bridge module.
In case os island benches, service columns are equipped with a rear panel to cover the connections.*

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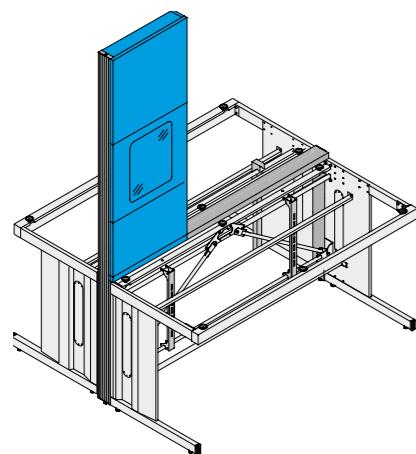
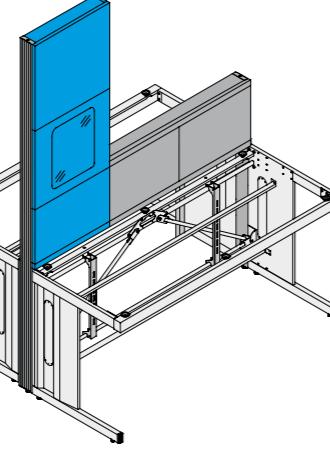
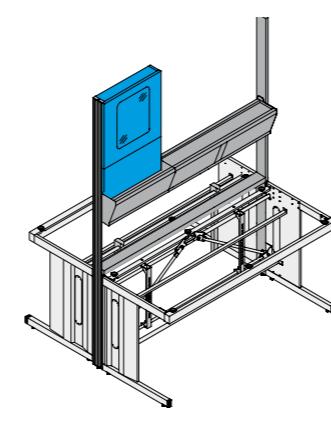
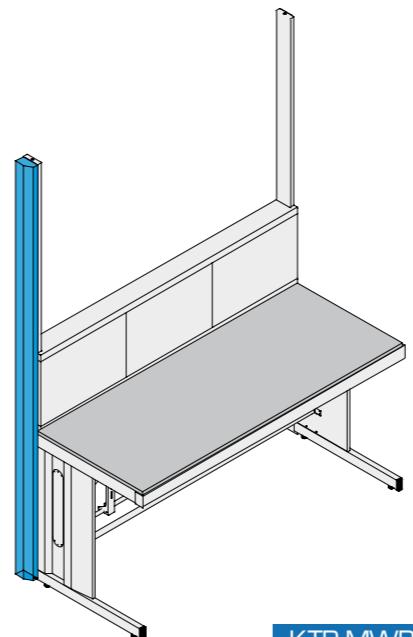
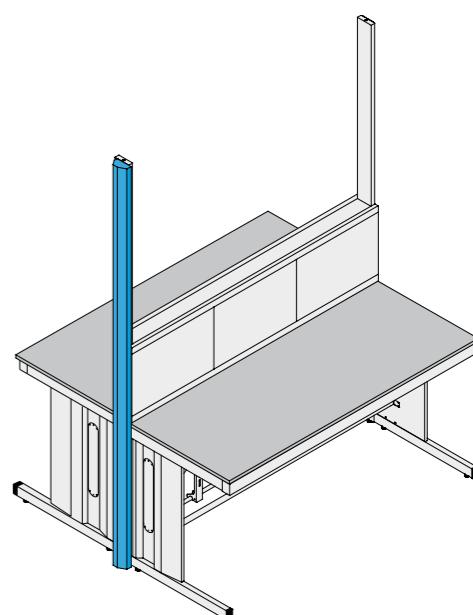
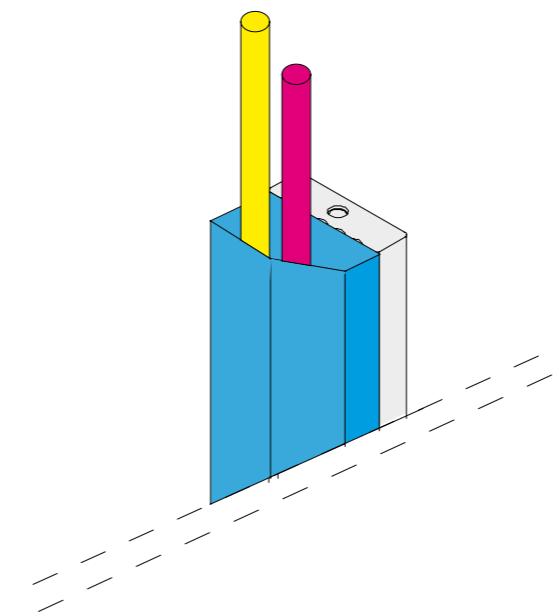
As an alternative to the service column, in case of simpler connections with no need of installing valves and regulators, a side carter in epoxy powder coated steel can be installed on the bench, needing the utility.

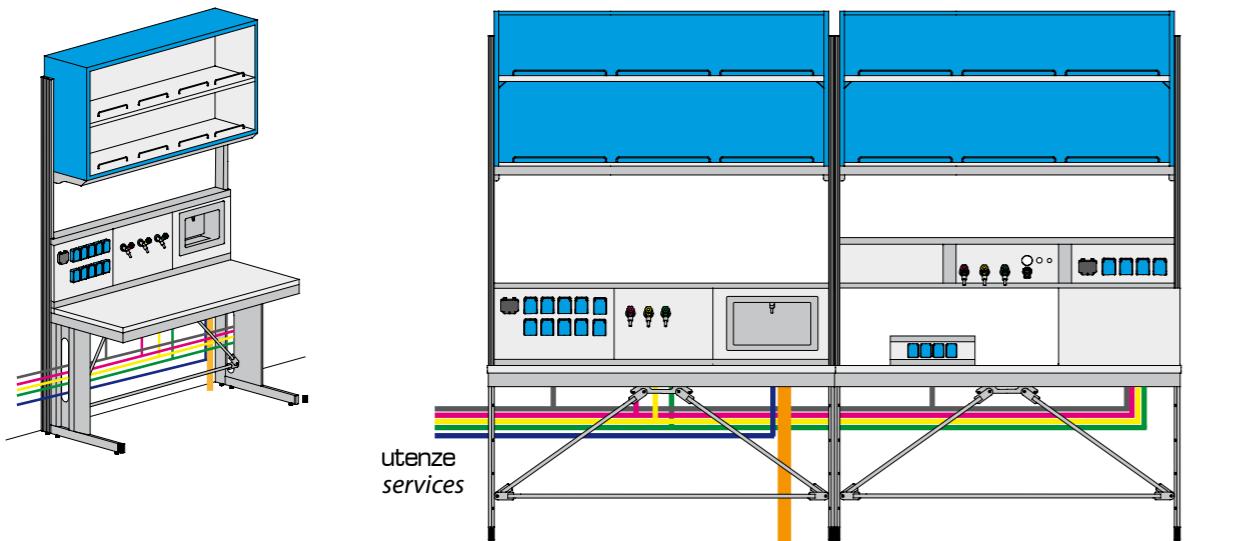
banchi
benches 1

MOMOLINE

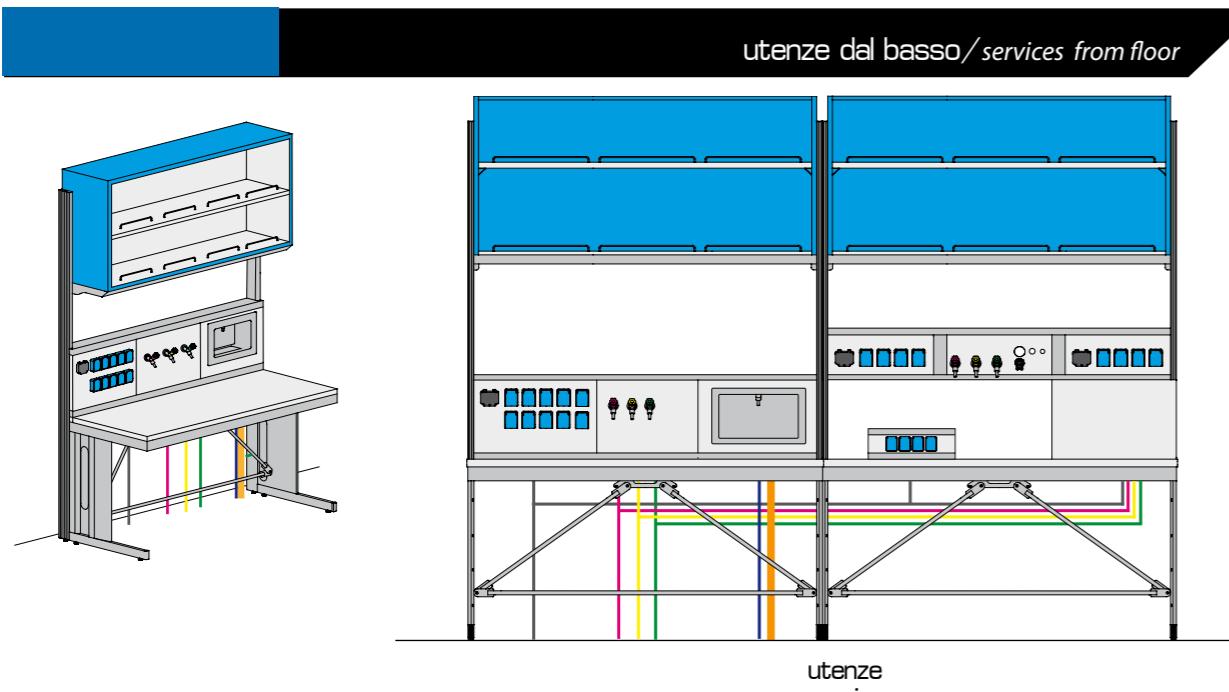
banchi
benches 1

utenze laterali / services from side

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*Example of furniture installation with services from side.
In this case a steel back panel is provided to conceal the lines (see pag.62).*

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*Example of furniture installation with services from floor.
In this case a steel back panel is provided to conceal the lines (see pag.62).*

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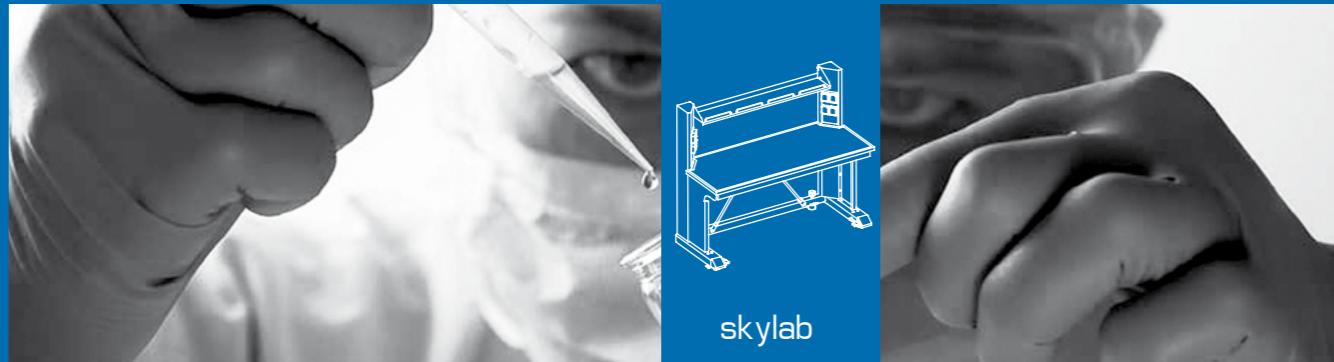
SkyLab line stands for the quintessence of flexibility and handiness, applied to laboratories.

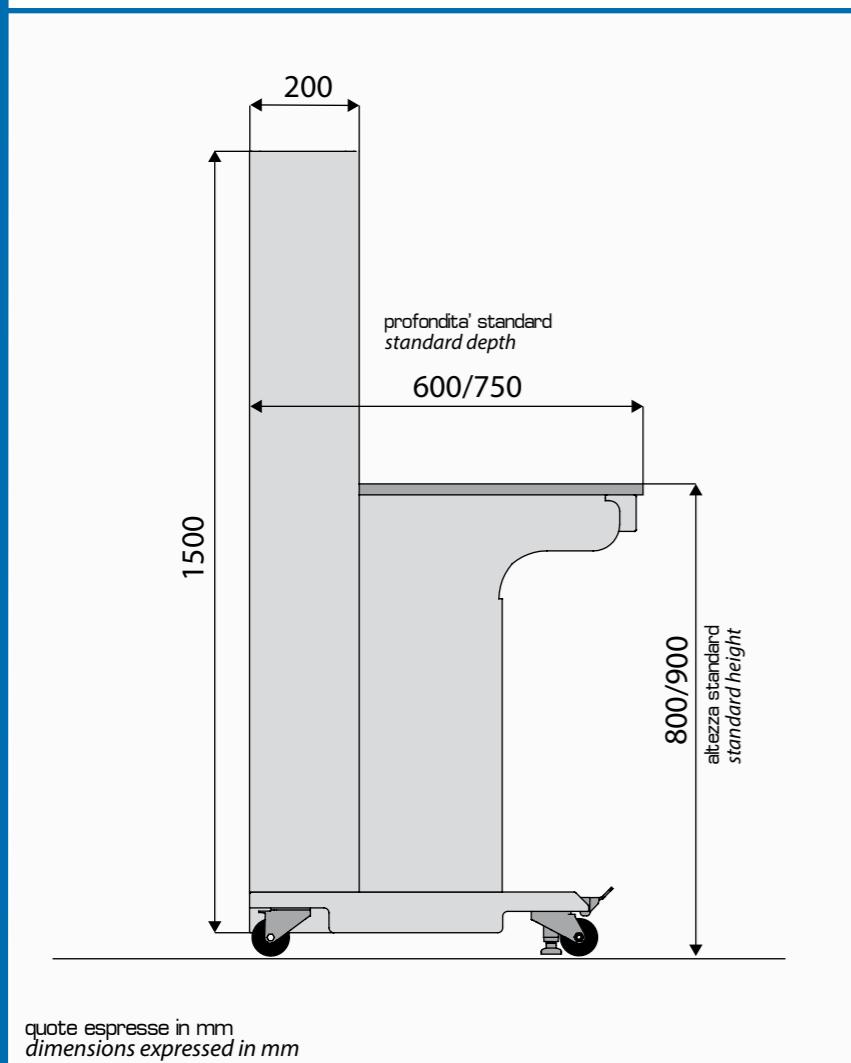
The main feature of the **SkyLab** line is the possibility of obtaining an extremely dynamic laboratory, which can be easily modified according to changing needs.

The core element of **SkyLab** line is represented by the movable trolley bench: together with the same solidity and reliability typical of all **MOMOLINE** products, the trolley benches also have the advantage of moving, so to be relocated with ease.

Thanks to special devices allowing the wheels to be lifted up once the bench has been located in the desired position, **SkyLab** trolleys act exactly the same way as traditional benches: also **SkyLab** laboratory benches hold Certificates stating the compliance to the European Norm EN 13150:2001 issued by Bureau Veritas, authorized third organization officially recognized at international level.

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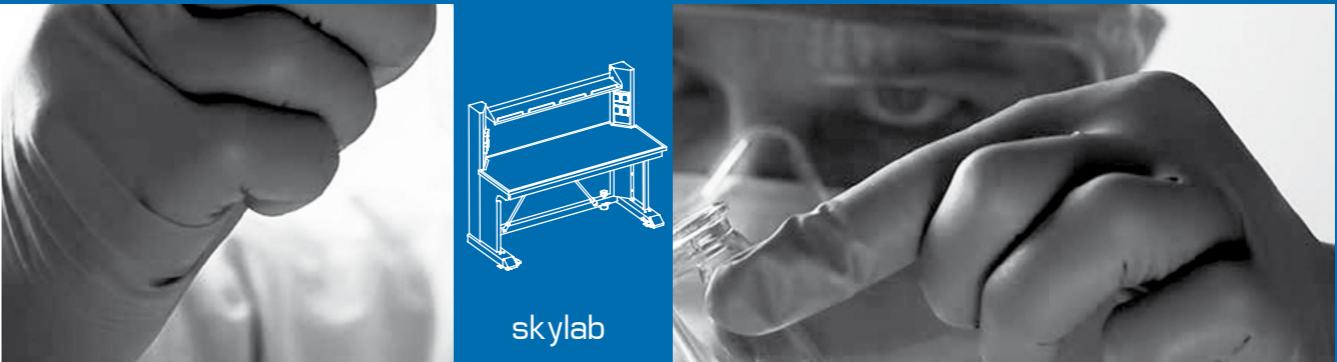
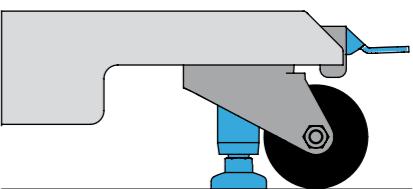
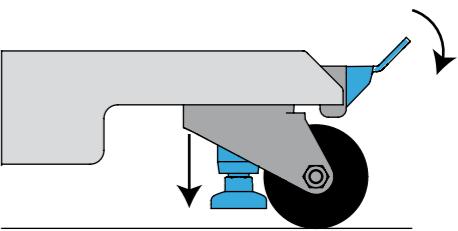
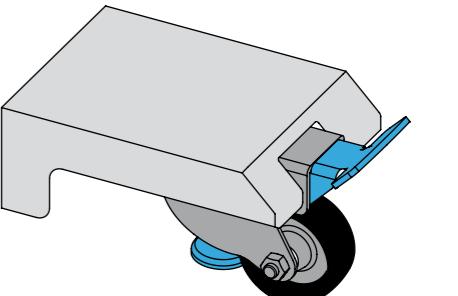
The compliance certificate states the resistance of **SkyLab** trolley benches to a 800 kg load placed on the working surface, considerably exceeding the minimum prescriptions of the EN 14175:2001 Norm, which recommends a 200 kg resistance.

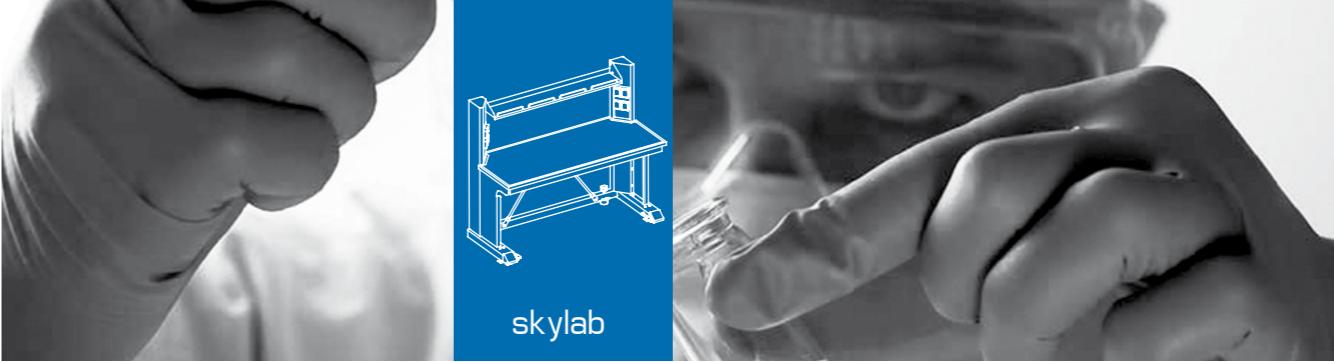
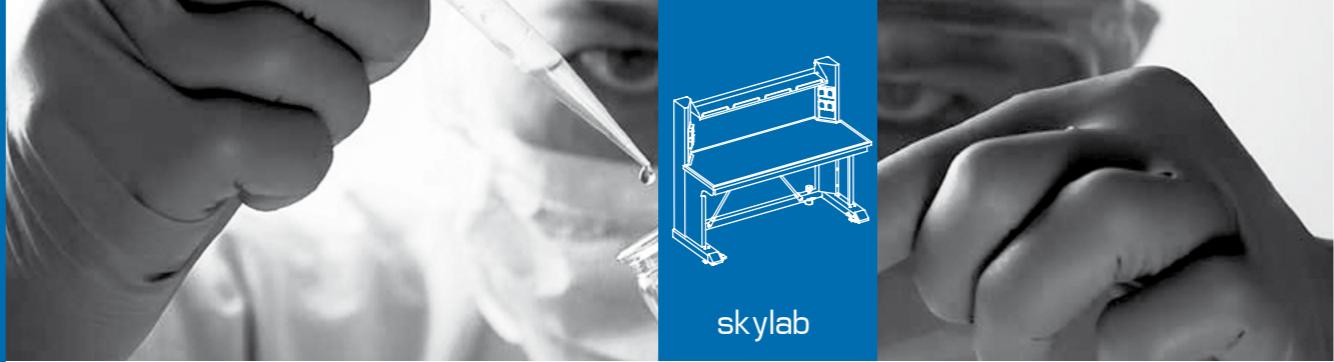
The trolley bench can be either directly connected to a power outlet (through its dedicated cable) or free of services: **SkyLab** line is completed by the 'Totem' 'Sky' and 'Service' modules, allowing to obtain an extremely efficient and functional distribution of the services to all the benches.

All the services come from the top or from the sides of the benches, and are completely independent, so that the trolley benches can be easily moved and transferred. The great flexibility coming from this particular system allows to fully customize the configuration of the laboratory, so to successfully meet any specific demand.

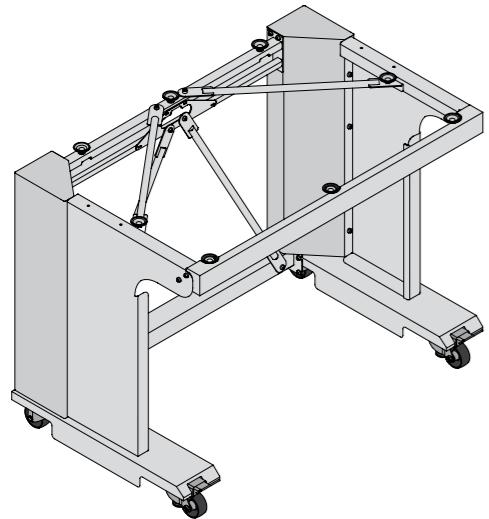
The 'Satellite' module allows to considerably optimize and enhance the performances of the gas distribution system, so that considerable savings can be achieved on the number of pressure reducers.

SkyLab line has been specifically designed for all applications where a frequent movement of the benches is required, and particularly when heavy and bulky instruments, difficult to handle in case of traditional benches, have to be employed.

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750 mm

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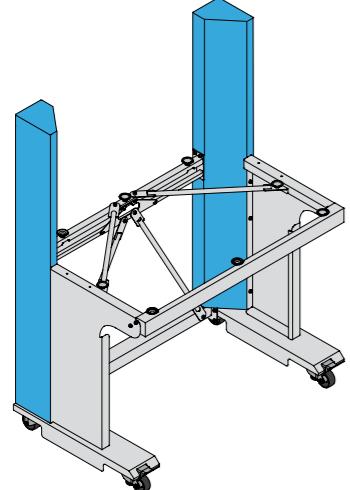
750/900 mm

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BCAR90-S-SKY	900
BCAR100-S-SKY	1000
BCAR120-S-SKY	1200
BCAR150-S-SKY	1500
BCAR180-S-SKY	1800

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750/900 mm

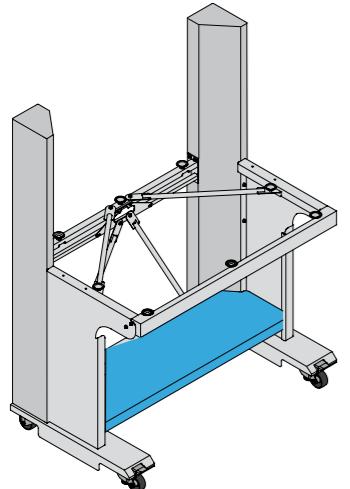
altezza service spine/
service spine's height

1500 mm

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BCAR90-T-SKY	900
BCAR100-T-SKY	1000
BCAR120-T-SKY	1200
BCAR150-T-SKY	1500
BCAR180-T-SKY	1800

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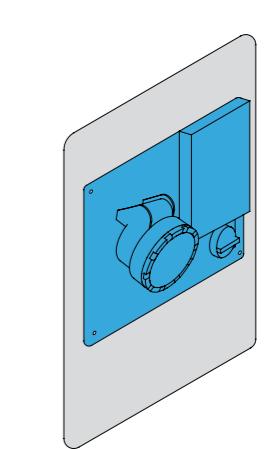
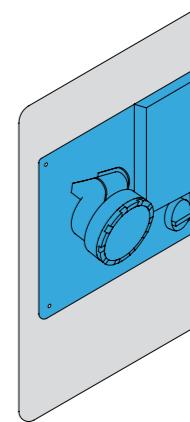
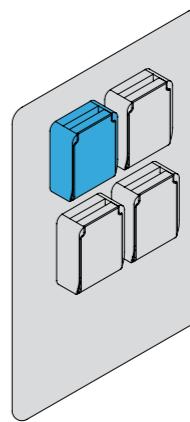
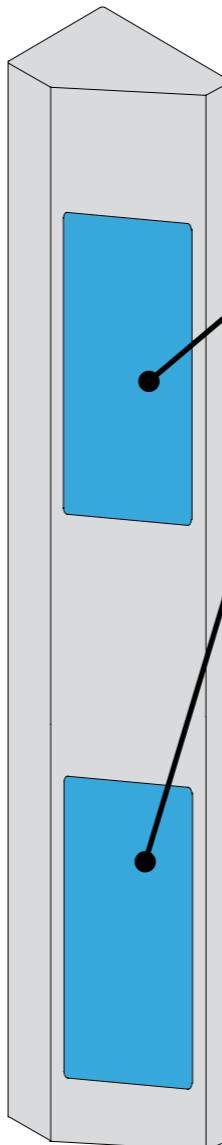
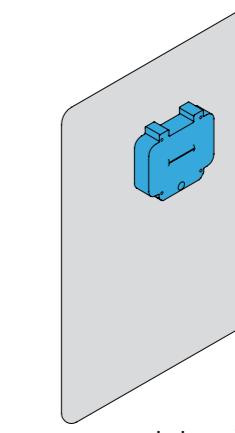
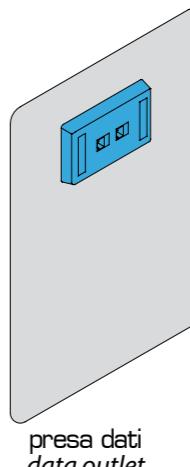
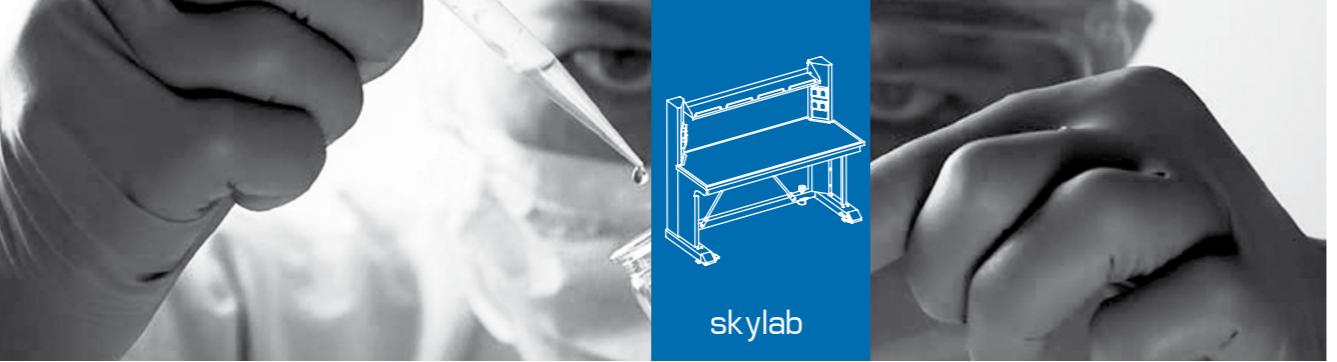
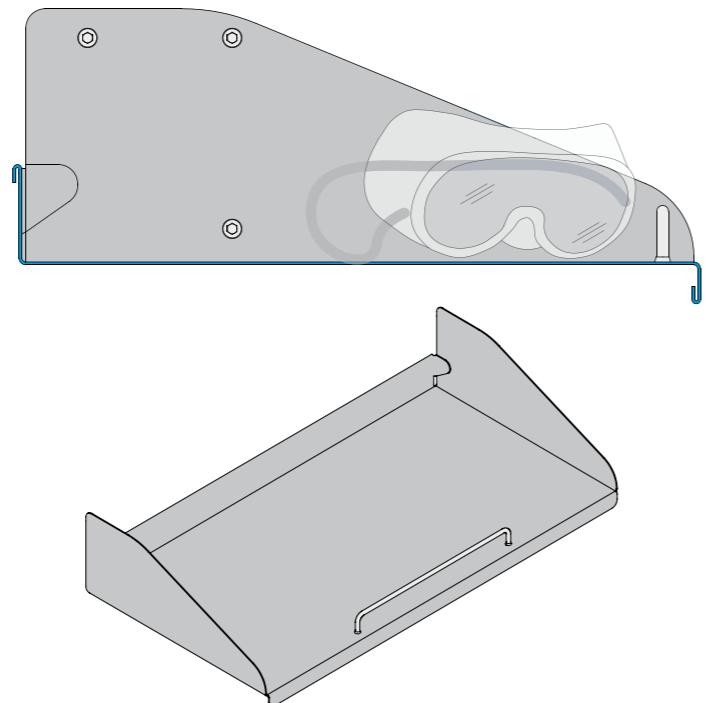
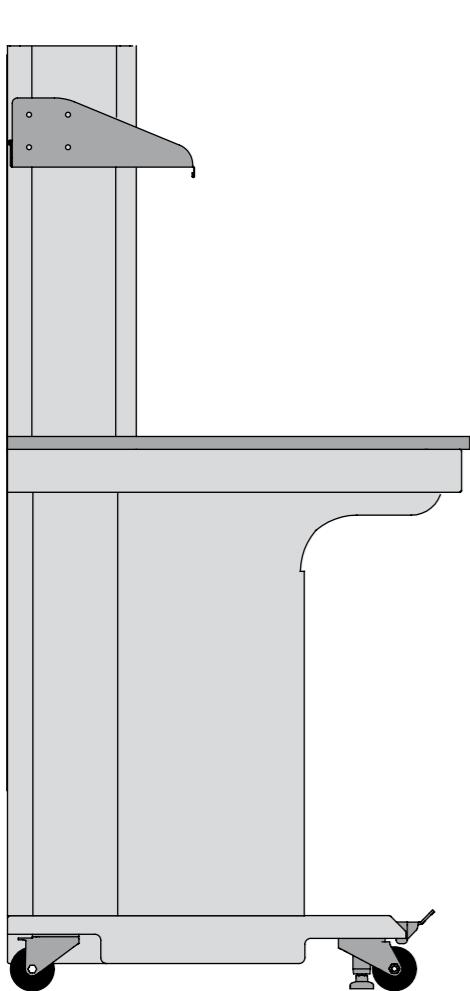
Service spine for **SkyLab** trolley bench is made up of two vertical elements located at the sides of the bench.

Electric services are fitted on interchangeable panels, and can be placed both above and under the worktop level.

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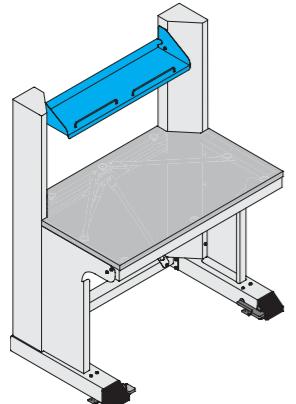
This shelf is realized in epoxy power coated steel sheets, and allows to place all the necessary equipment for doing a good job. The shelf is equipped with fall-protection handles, and can be fitted to the service spine at different heights.

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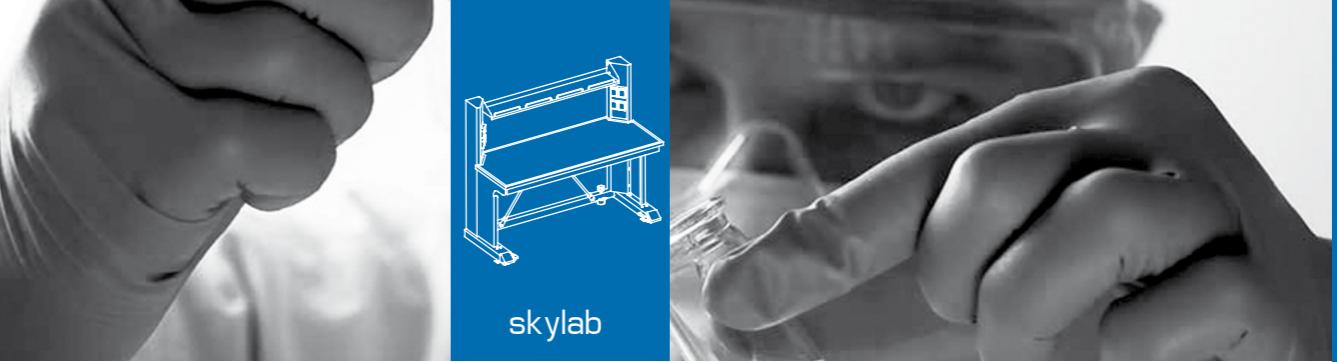
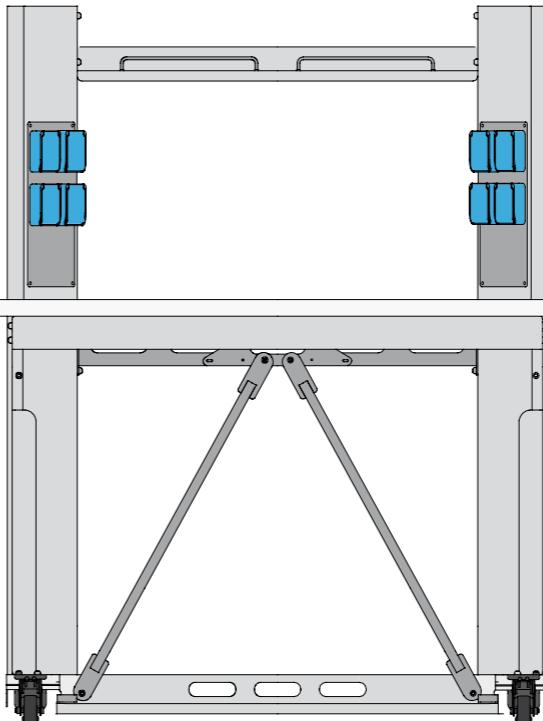
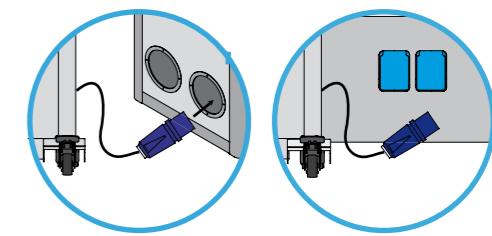
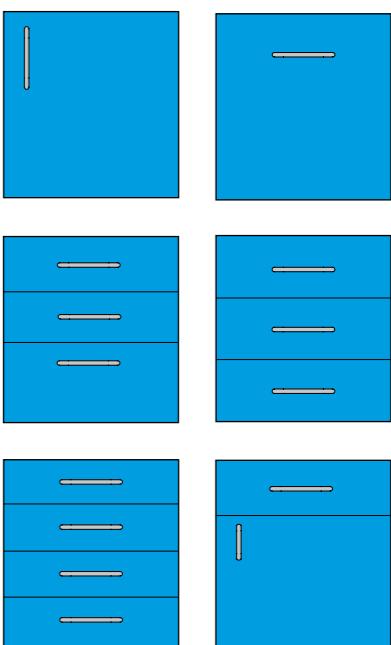
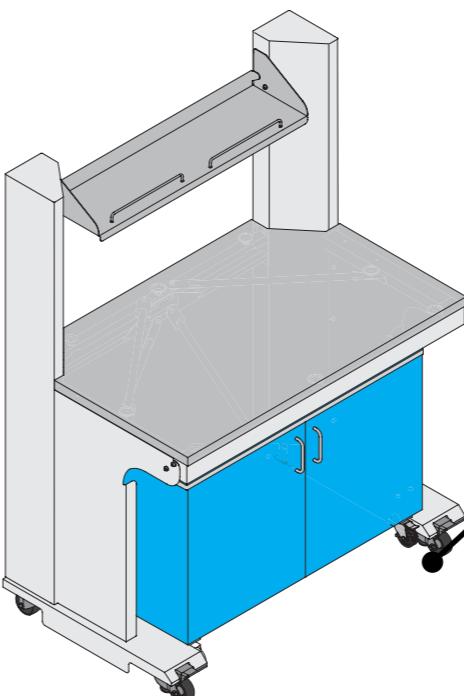
SkyLab trolley bench, designed to ensure great flexibility, can be connected to a power outlet through a dedicated cable, already fitted to the bench.

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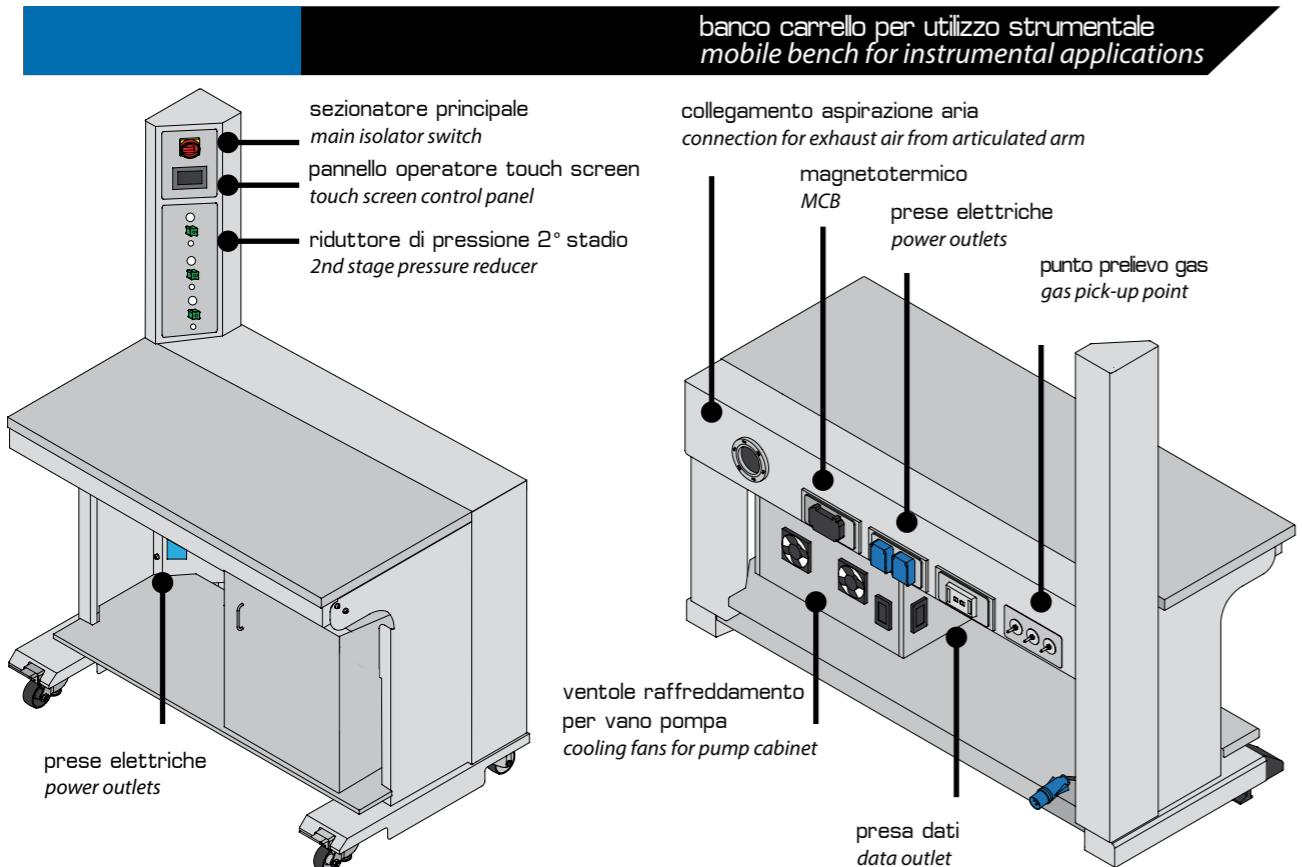
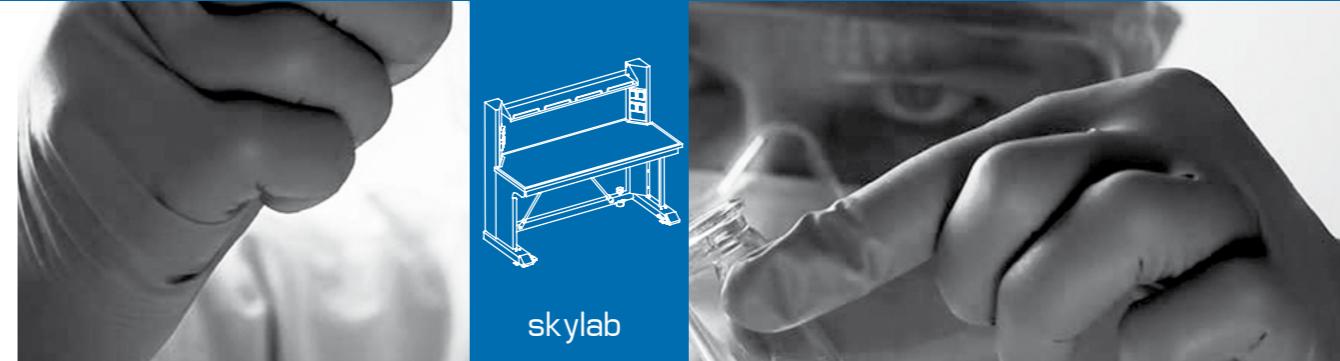
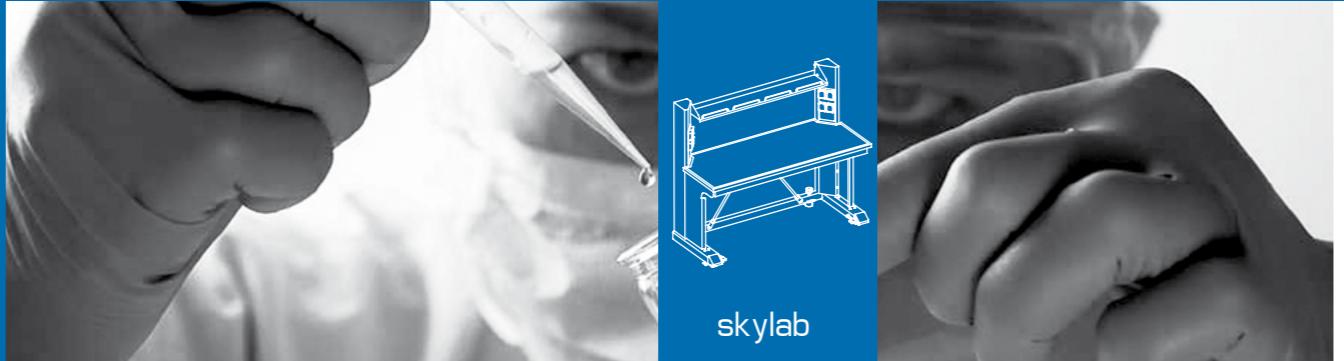
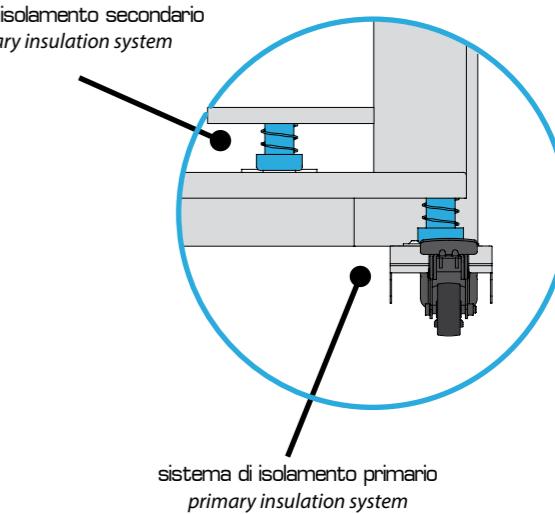
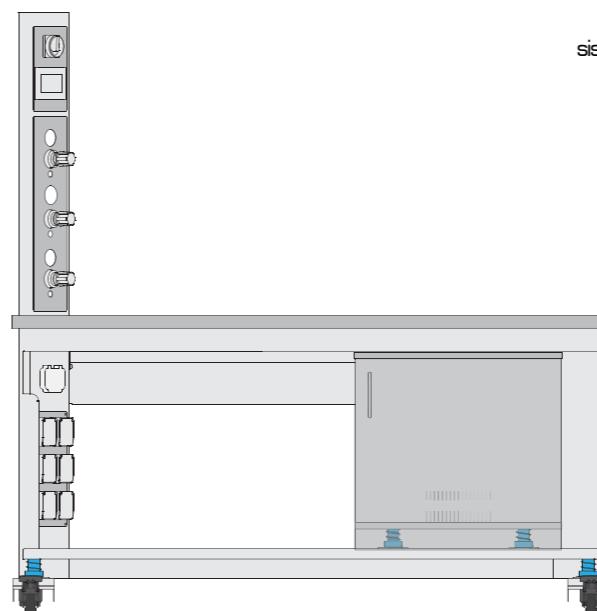
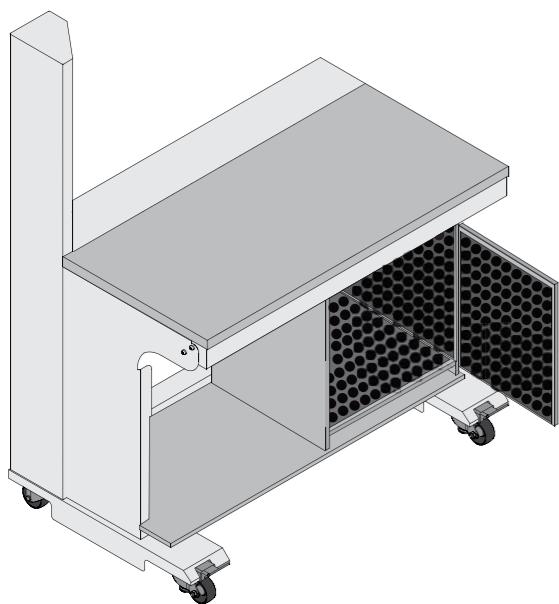
Trolley benches from **SkyLab** line can be required in a special version for analytical applications. In this case, the bench is designed for the installation of analytical instruments requiring the connection to vacuum pumps and to a PC with all its accessories. The bench is also equipped with exhaust system for the waste vapors produced by the instruments. This special bench is equipped with a sound-proofed technical compartment (approx. dimensions: 600x600x600 mm), designed for the installation of a vacuum pump to be connected to the instruments on the bench. The compartment is located under the worktop and is fitted with a ventilation cooling system. The technical compartment is also fitted with a pull-out tray used for collecting possible oil spilled by the pump during its operation or during maintenance.

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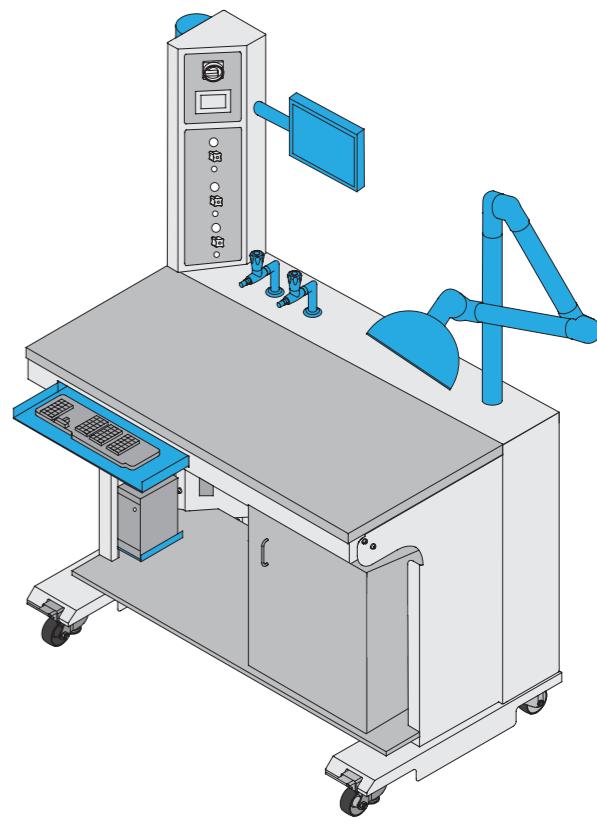
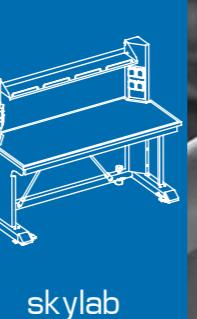
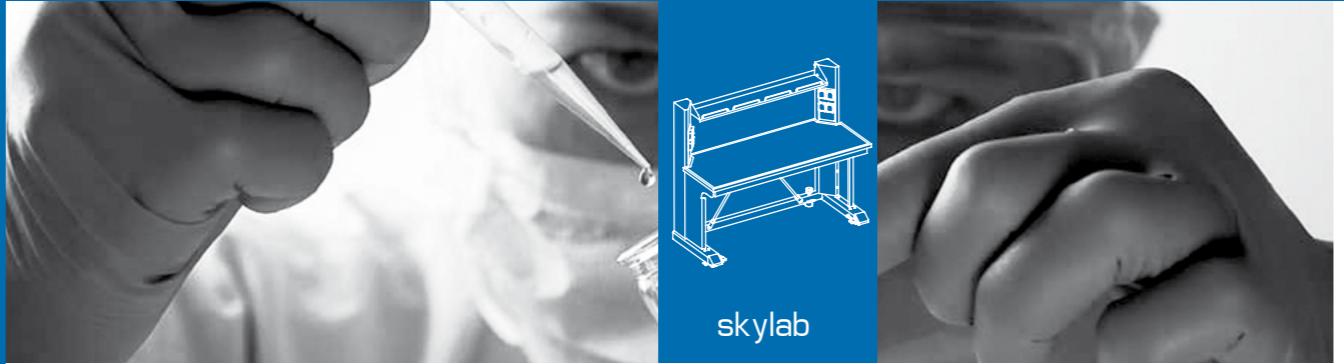
In order to grant perfect insulation from vibrations, the technical compartment for vacuum pumps is fitted with a double insulation system. Primary system is made-up of anti-vibrating elements installed between the compartment and the bench structure, while additional springs installed between the cabinet and the pump's support grant the total absence of vibrations. The bench can be completed by a wide range of accessories, such as extraction arms, gas taps, PC holders.

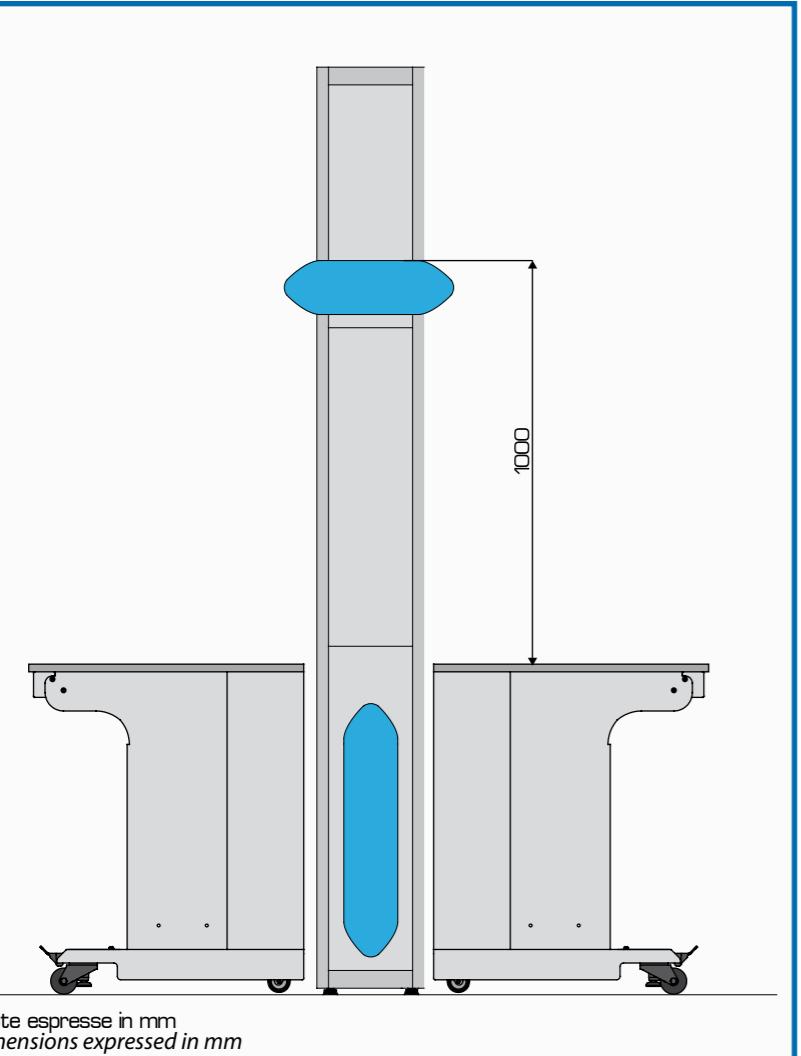
porta-tastiera keyboard holder	porta-PC PC holder	cappa a braccio articulated arm local extractor
TRAY-KB	PCH	AF-C

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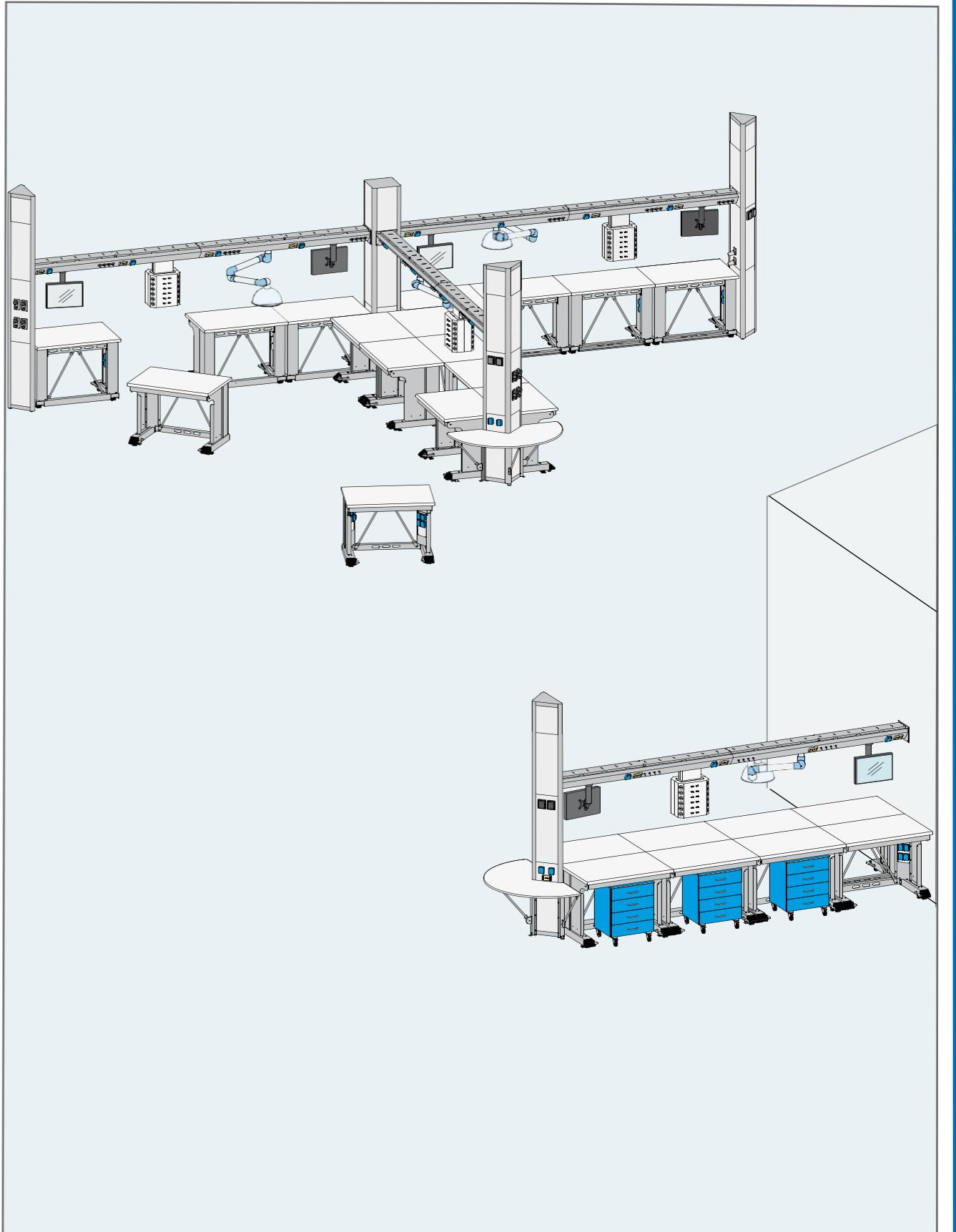
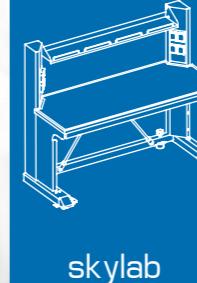
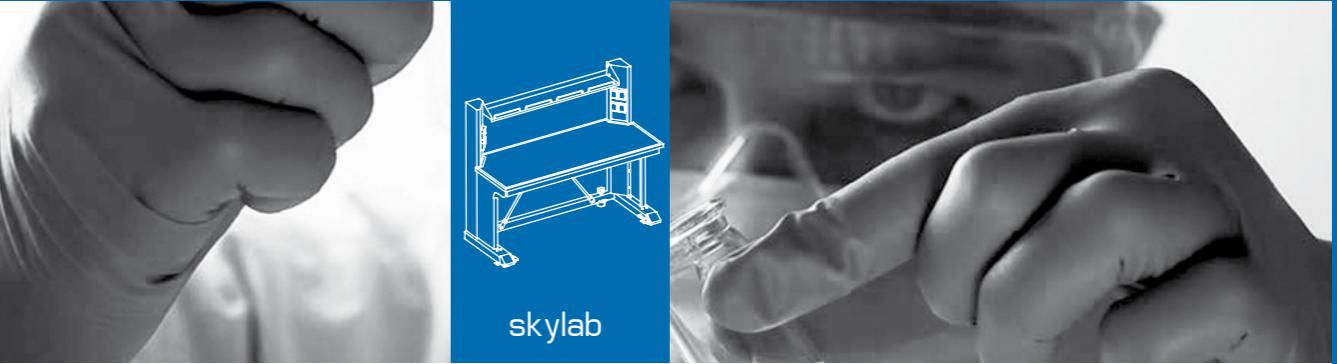
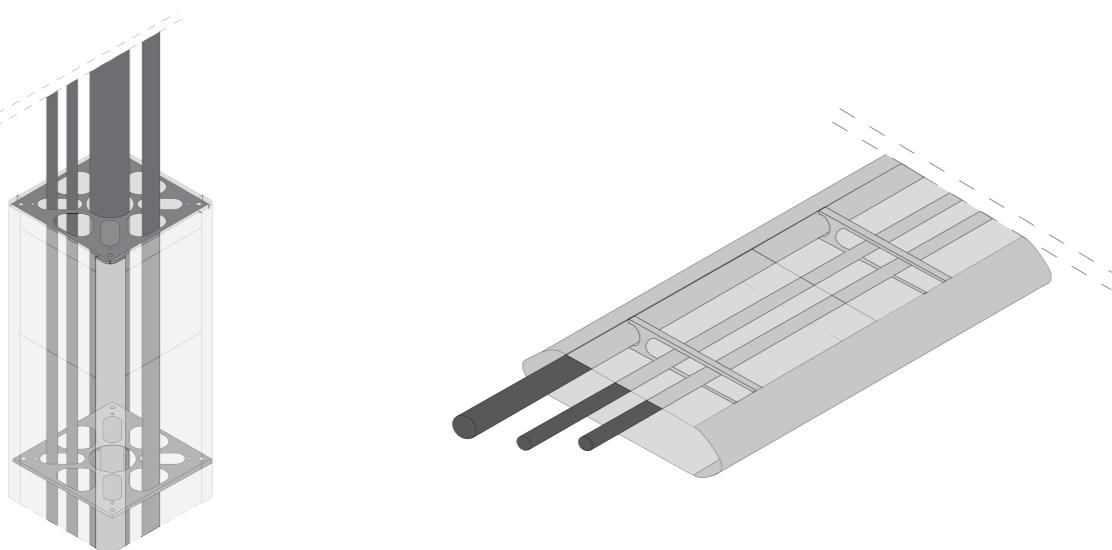
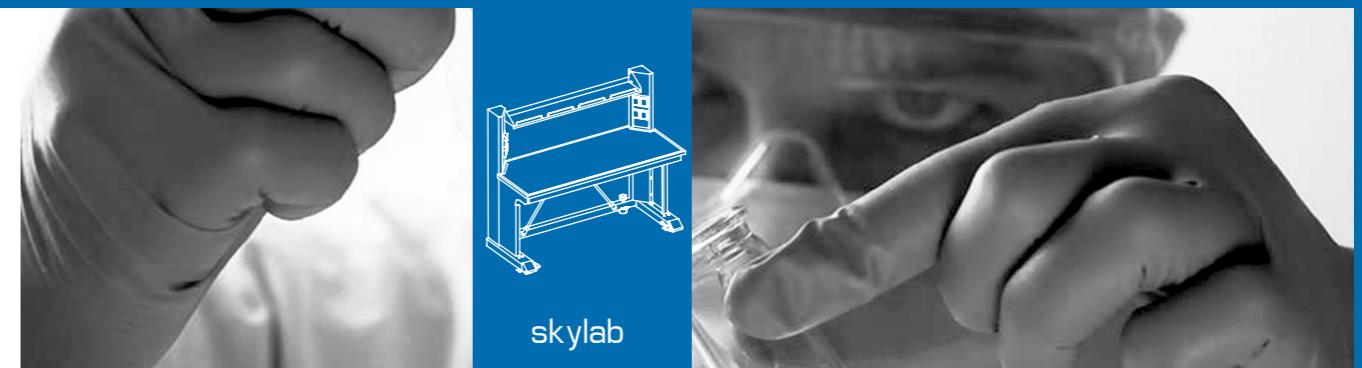
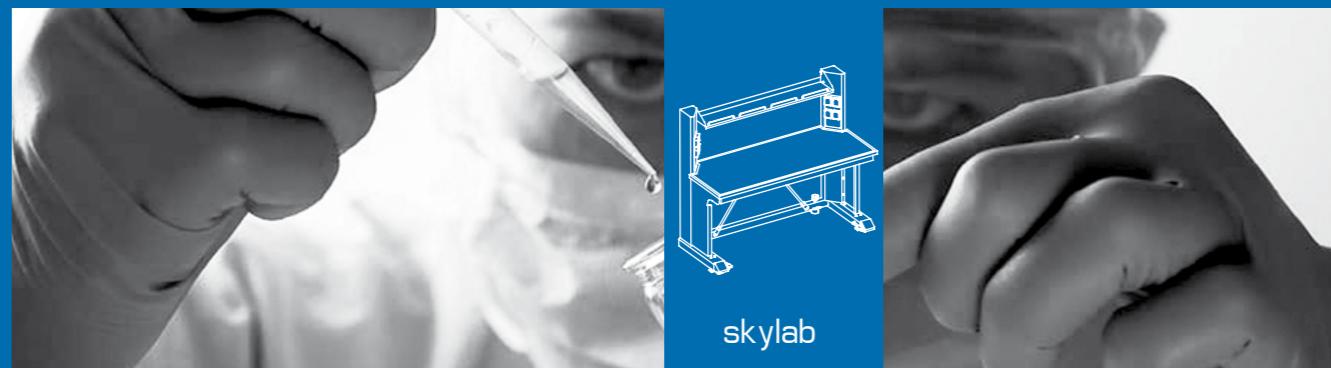
SkyLab line is completed by the 'Totem' 'Sky' and 'Service' modules, allowing to obtain an extremely efficient and functional distribution of the services to all the benches.

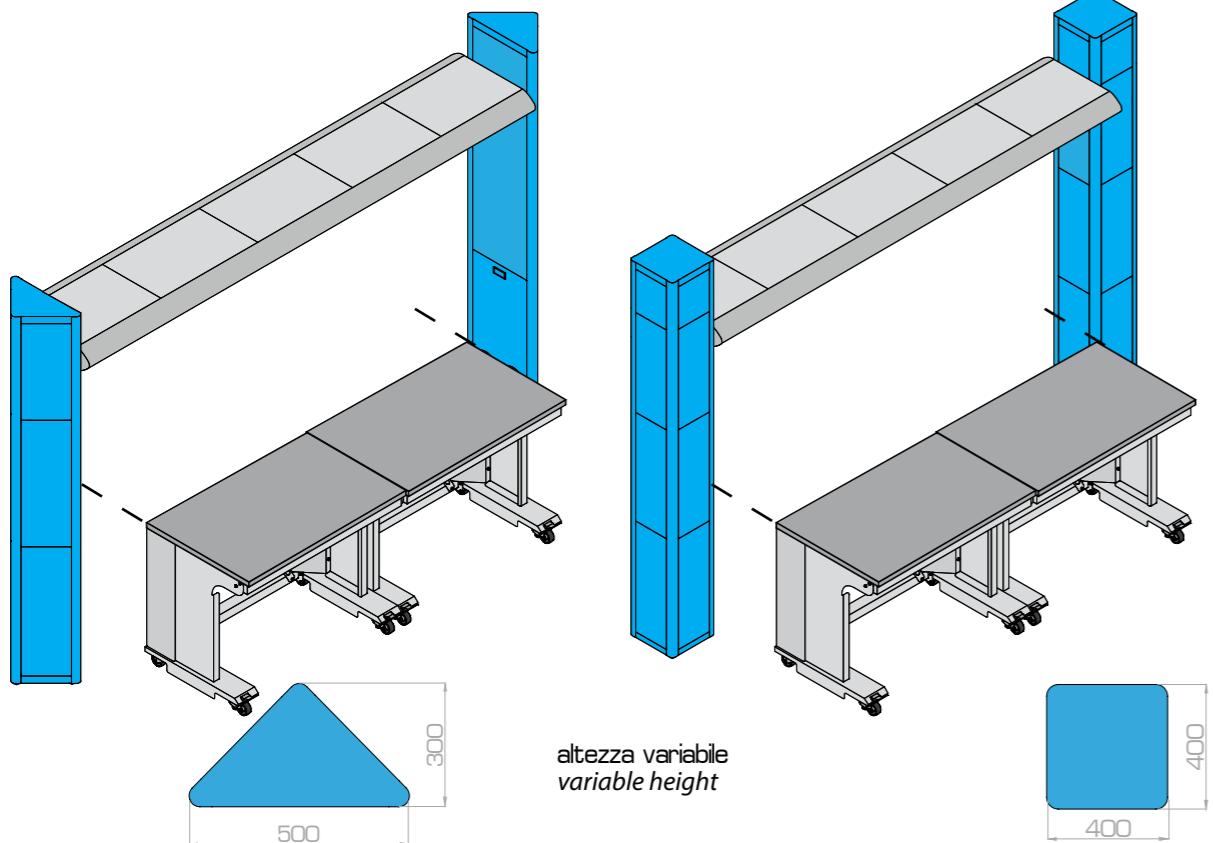
All the services come from the top or from the sides of the benches, and are completely independent, so that the trolley benches can be easily moved and transferred.

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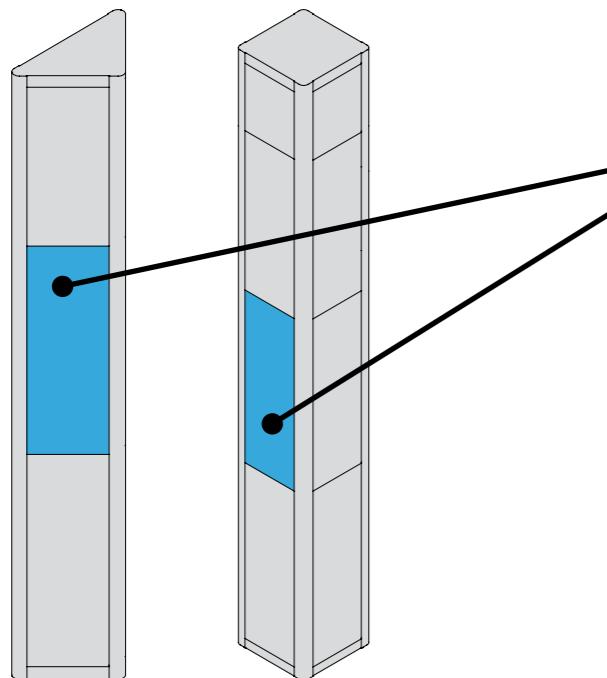
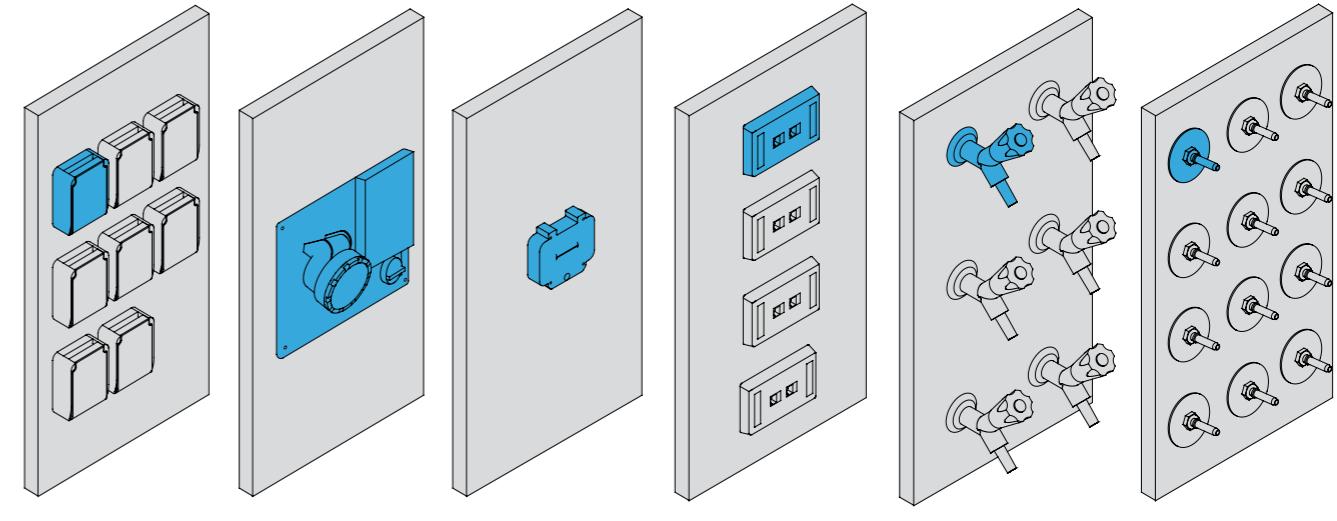
Totem module is a vertical column positioned on the sides of benches, with squared or triangular cross section. Totem stands as a supporting element for the Sky module, and can be also equipped with many accessories, fitted on interchangeable panels.

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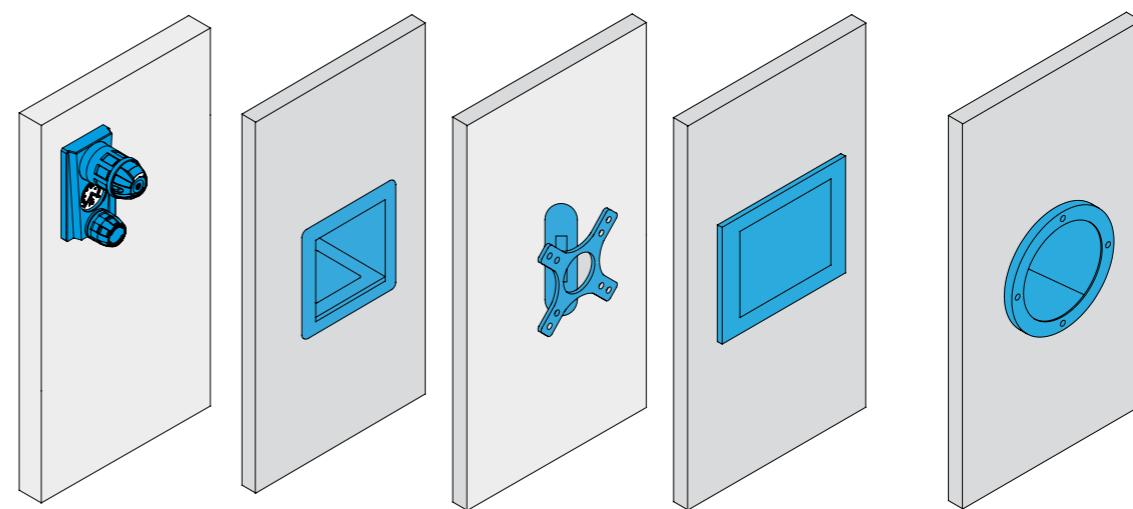
vaschetta
drip cup

braccio monitor
monitor supporting arm

monitor GDS
GDS monitor

collegamento asp.aria
connection for exhaust

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vaschetta
drip cup

braccio monitor
monitor supporting arm

monitor GDS
GDS monitor

collegamento asp. aria
connection for exhaust control panel

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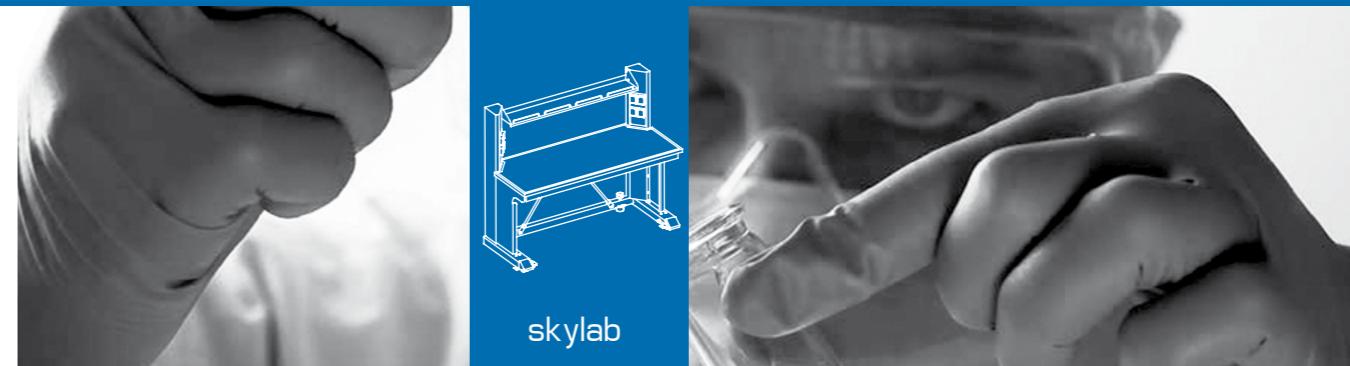
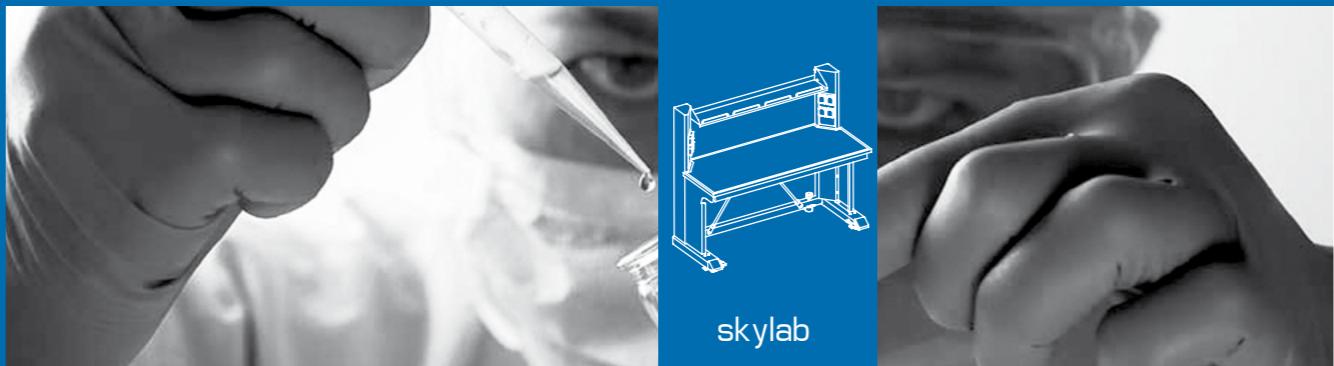
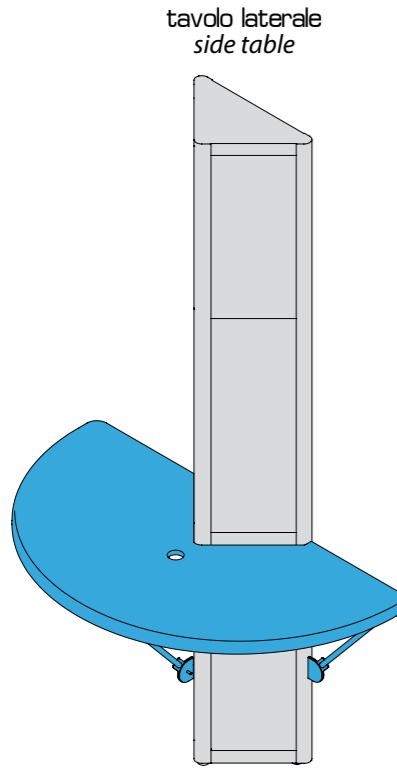
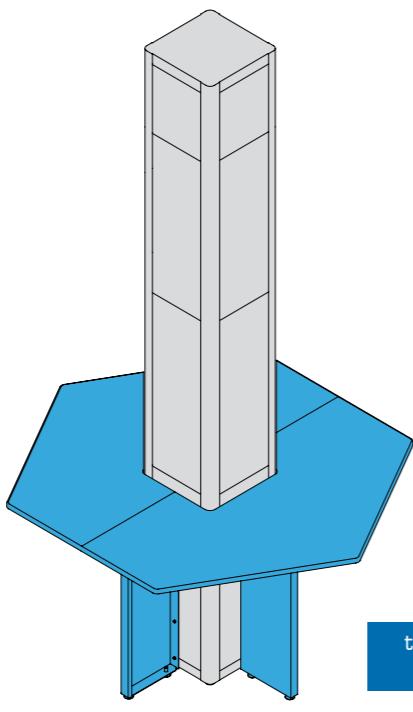
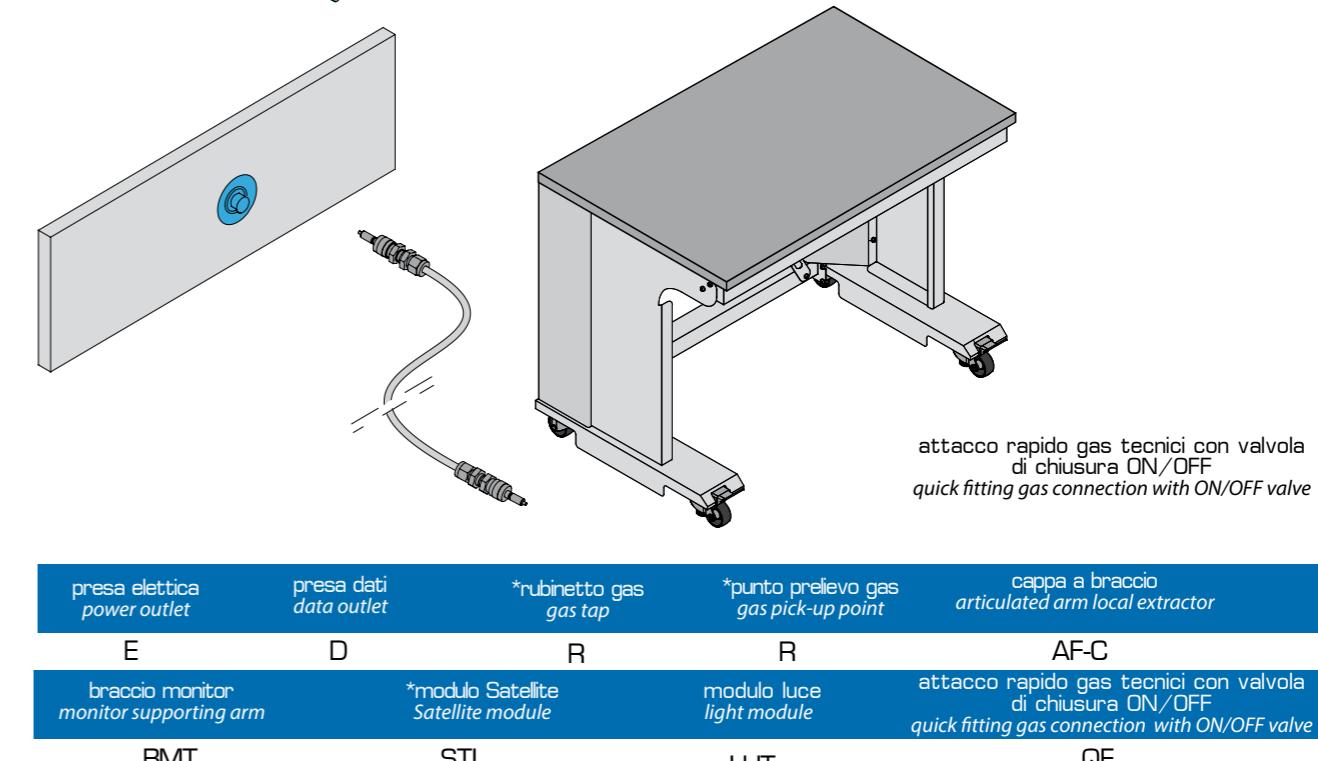
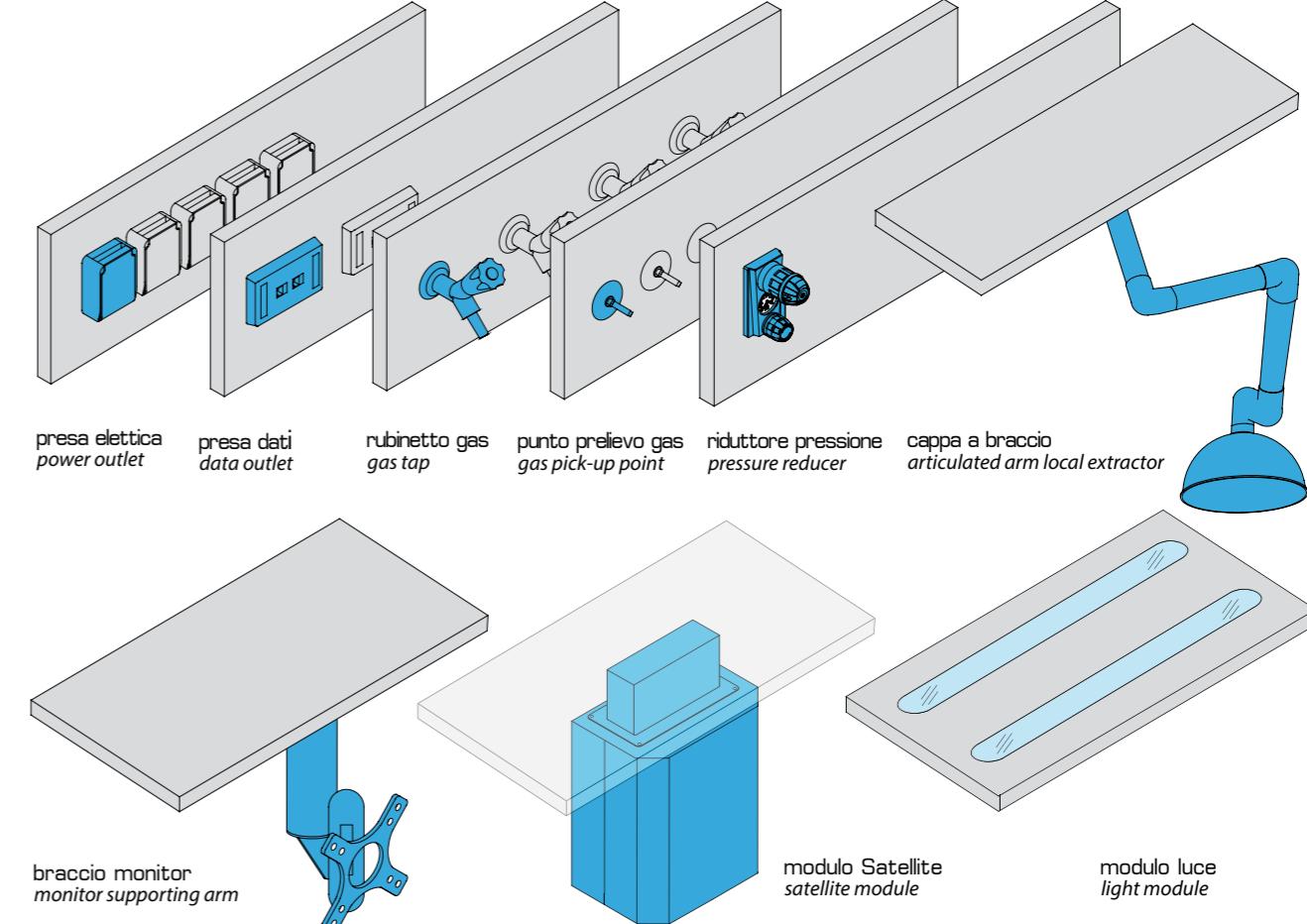
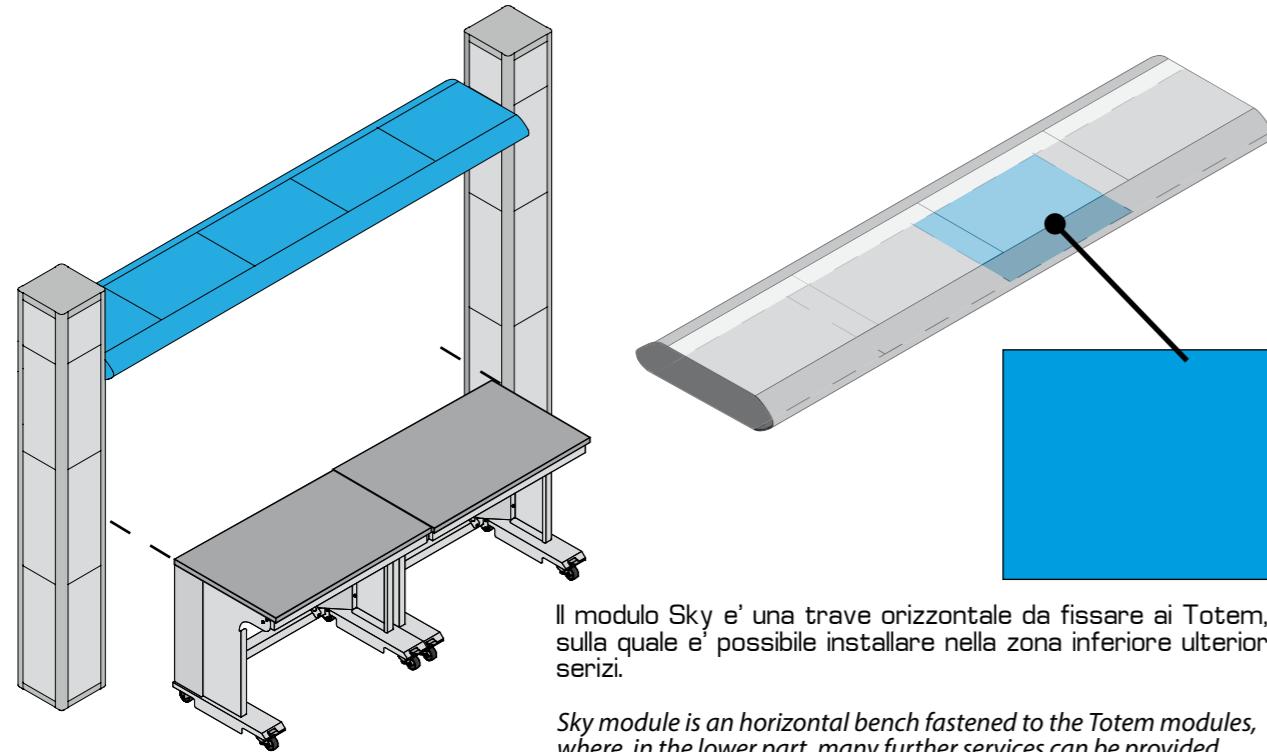
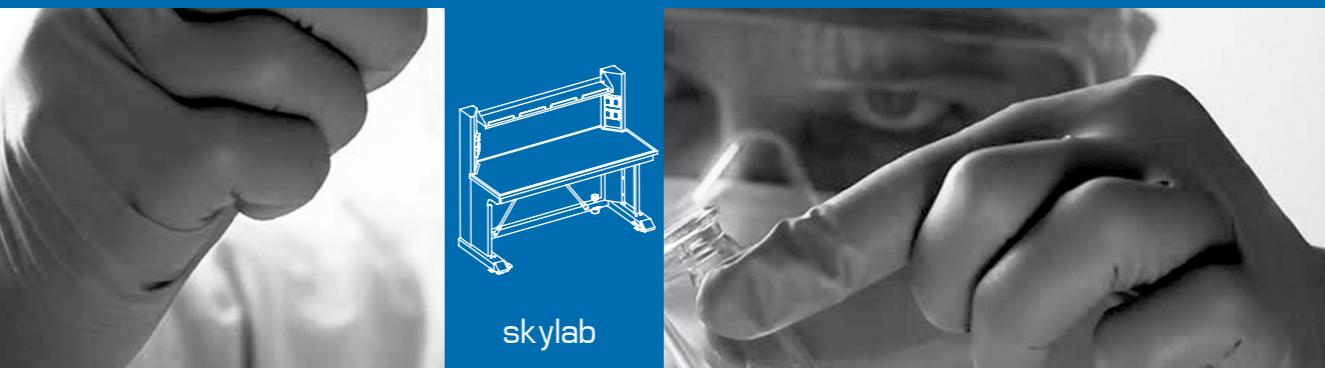
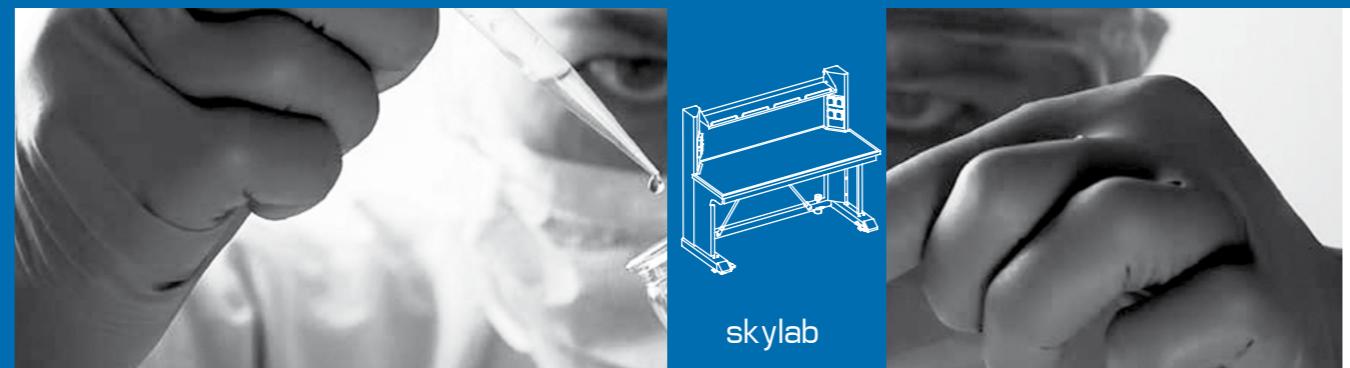
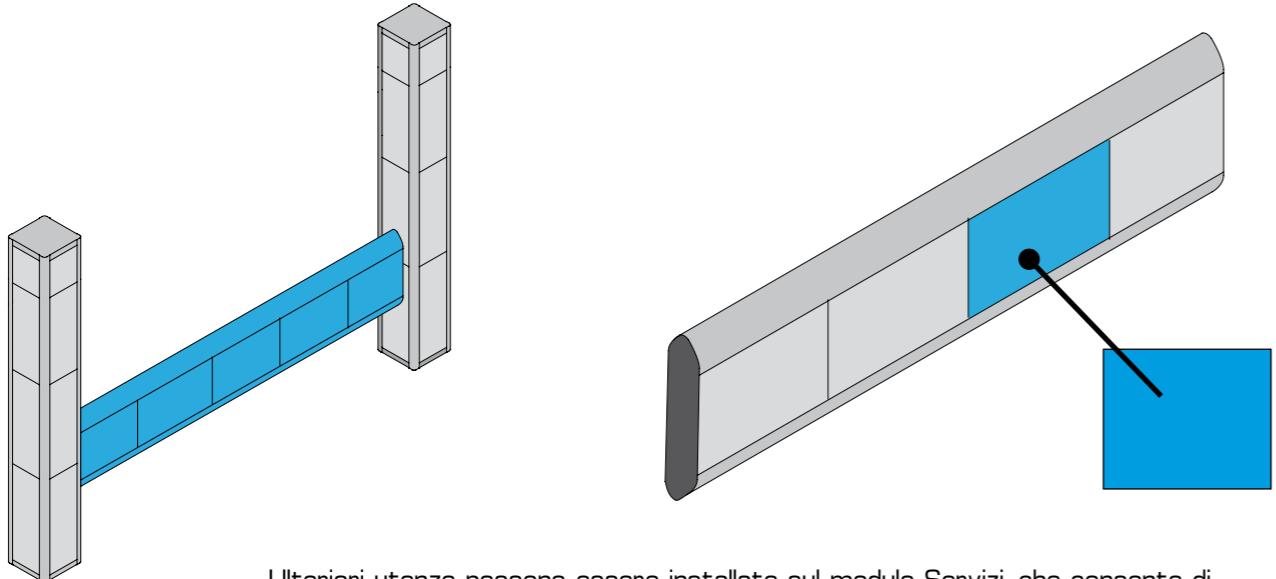
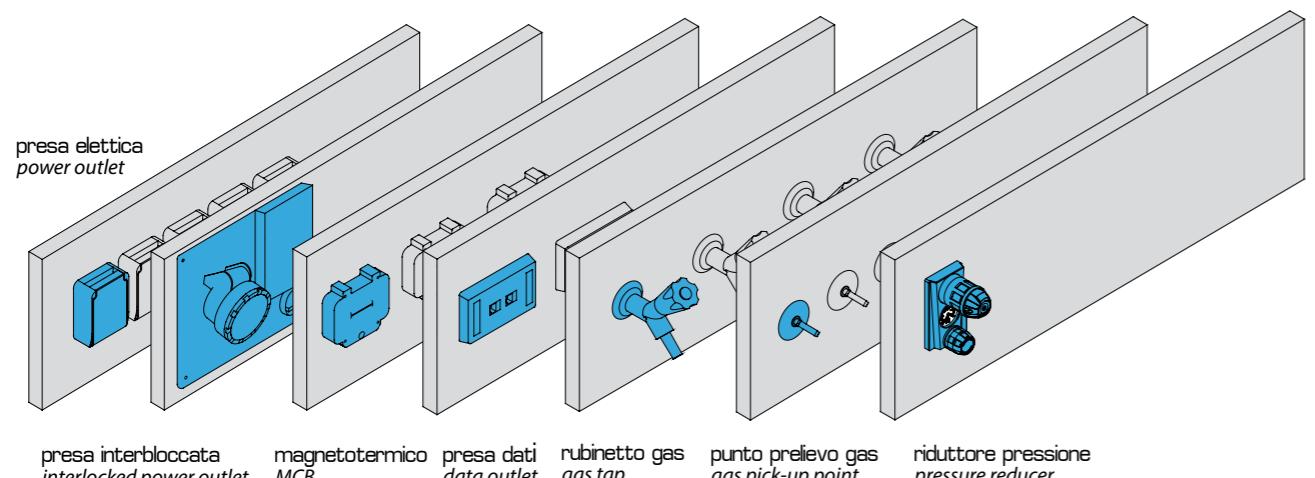
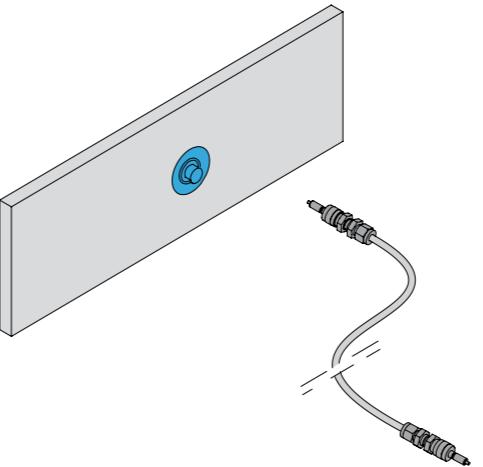
This module allows to obtain further services to the benches, and also grants the possibility of supplying electricity to trolley units located apart from totem modules.

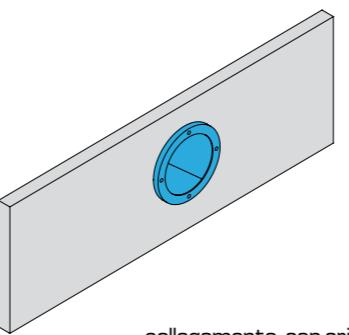
presa elettrica <i>power outlet</i>	presa interbloccata <i>interlocked power outlet</i>	magnetotermico <i>MCB</i>	presa dati <i>data outlet</i>	*rubinetto gas <i>gas tap</i>	*punto prelievo gas <i>gas pick-up point</i>
E	I	M	D	R	R

*riduttore pressione <i>pressure reducer</i>	attacco rapido gas tecnici con valvola di chiusura ON/OFF <i>quick fitting gas connection with ON/OFF valve</i>	collegamento asp. aria <i>connection for exhaust control panel</i>
PR	QF	EXH

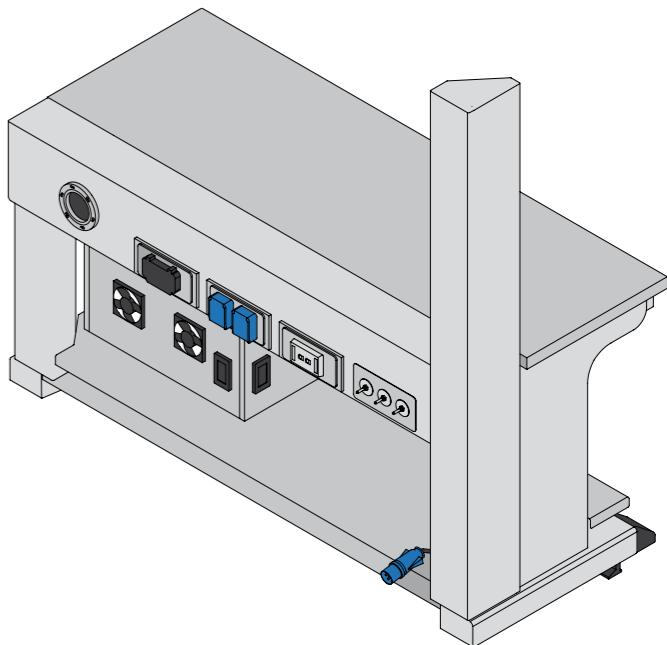
* codici relativi ai rubinetti gas a pag. 109-110
codes for gas taps at pag. 109-110

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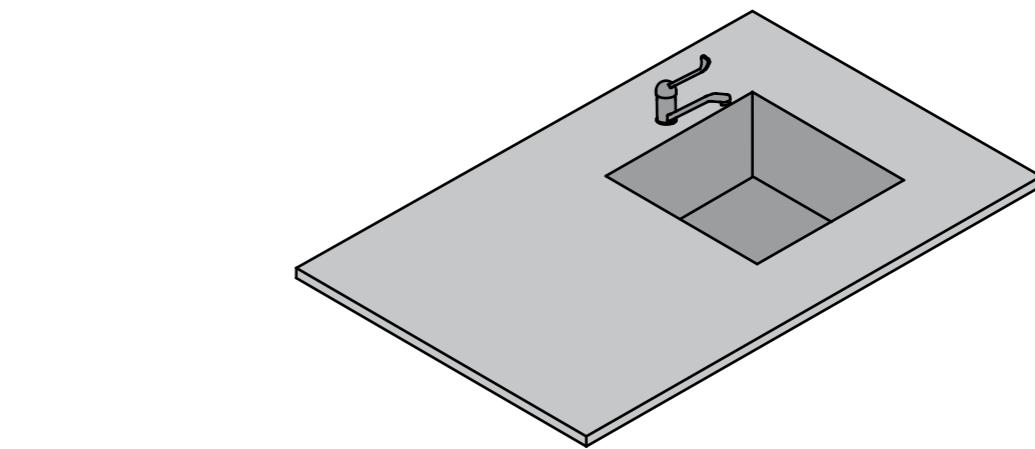
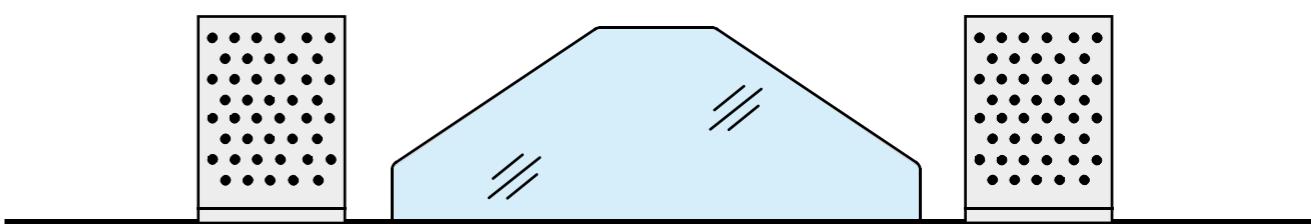
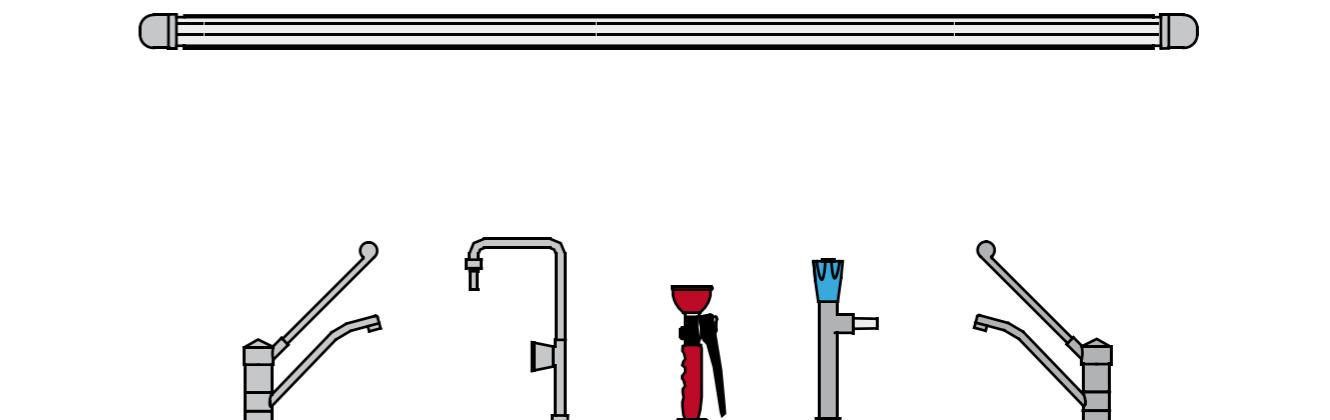
collegamento asp. aria
connection for exhaust

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*Thanks to the high flexibility and modularity of its production, **MOMOLINE** can equip benches with several accessories and items, in order to meet all customers's whishes.*

Accessories reported here, are the standard ones, but custom-made and specially designed solutions can be provided upon request, so that every customer can fulfill his own desires and expectations.

The possibility of providing tailor-made solutions applies to the type and number of installed accessories, as well as to the color of worktops and to the technical specifications of the services outlets.

MOMOLINE is always pleased to receive detailed and specific enquiries from customers, in order to provide the best solution for every occasion.

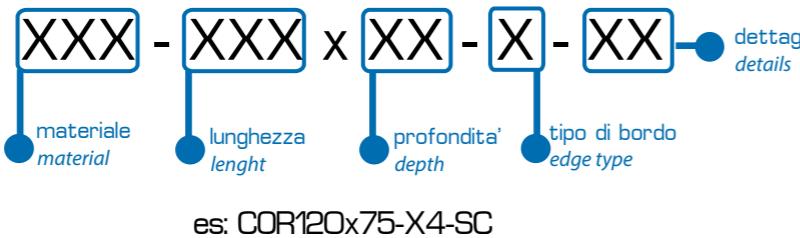
*The considerable experience of **MOMOLINE** in setting up laboratories for every field and application allows to support customers in the choice of the most suitable items, both from the aesthetic and technical point of view.*

applicazione / application

	COR	PP	HPL	LAM	WLAM	SW	EP	SP	SS	GR	ZL	GL
Utilizzo reagenti non aggressivi Employment of non-aggressive chemicals	●	●	●	●	●	●	●	●	●	●	●	●
Utilizzo reagenti molto aggressivi Employment of very aggressive chemicals	●	●	●	●	●	●	●	●	●	●	●	●
Utilizzo coloranti Employment of dyes	●	●	●	●	●	●	●	●	●	●	●	●
Settore petrochimico / Utilizzo solventi Oil&gas/Employment of solvents	●	●	●	●	●	●	●	●	●	●	●	●
Settore alimentare Food industry	●	●	●	●	●	●	●	●	●	●	●	●
Settore farmaceutico Pharmaceutical	●	●	●	●	●	●	●	●	●	●	●	●
Microbiologia Microbiology	●	●	●	●	●	●	●	●	●	●	●	●
Elevata igiene High hygiene standard	●	●	●	●	●	●	●	●	●	●	●	●
Appoggio strumentazione / Scrivania Equipment support / Desk	●	●	●	●	●	●	●	●	●	●	●	●

● idoneo
suitable● poco idoneo
not recommended● non idoneo
unsuitable

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materiale / worktop

COR PP HPL LAM WLAM SW EP SP SS GR ZL GL

Utilizzo reagenti non aggressivi
Employment of non-aggressive chemicals

● ● ● ● ● ● ● ● ● ● ● ● ● ●

AISI 316

Utilizzo reagenti molto aggressivi
Employment of very aggressive chemicals

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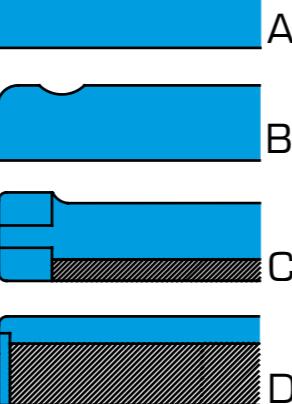
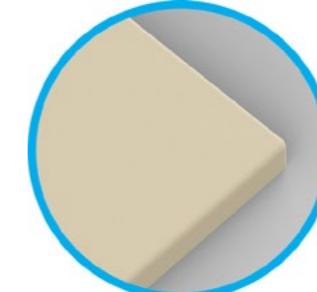
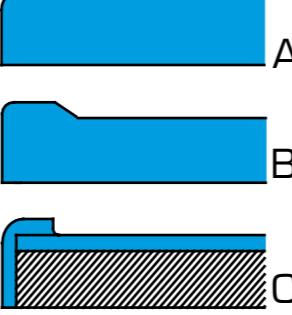
*Corian is a synthetic material, composed by 1/3 acrylic resin and by 2/3 of natural minerals. Corian has an incredibly uniform and homogeneous structure, does not delaminate, is not porous and is completely hygienic and stain-resistant. Corian also resists high temperature (up to 100°C) and most common chemicals. This material has excellent decontamination performances against viruses and bacteria, and its particular junction method allows to obtain virtually infinite surfaces with no evident junction grooves. For this reason, it is particularly suitable for microbiology and high-hygiene applications. **MOMOLINE** is part of Corian Quality Network, meaning all the products are officially recognized and guaranteed by DuPont, sole authorized manufacturer of Corian.*

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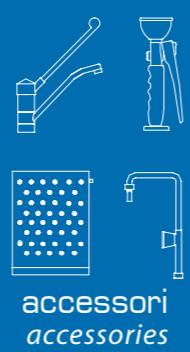
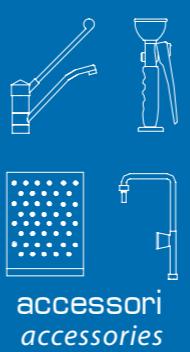
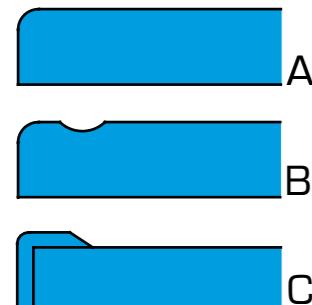
Polypropylene worktops are injection moulded and stand out for their extreme lightness. PP worktops provide very good resistance against acids, bases and alkalis.

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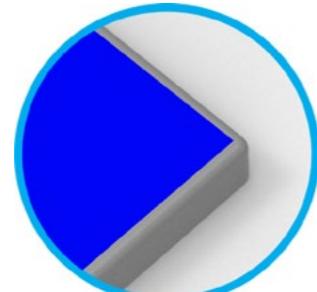
High Pressure Laminate worktops (such as Trespa) are composed by many cellulose fiber layers, soaked in phenolic resin and then heavily compressed through mechanical presses.

The surface of the panel is covered with cellulose fiber soaked in melaminic resin with wide choice of color. For aesthetic reasons the external surface can be specifically treated in order to obtain high resistance to chemicals, heat and scratches. Upon specific request, the phenolic resin can be additioned of special substances able to provide the worktop with flame retardant properties.

LAM

piano in LAMINATO PLASTICO/PLASTIC LAMINATE worktop

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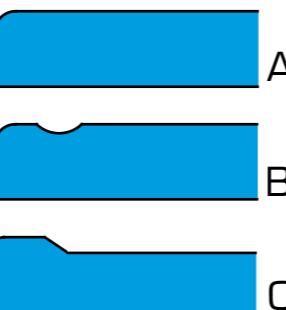
The core of this worktop is represented by a flame-retardant particle board panel, whose external faces are covered by flame-retardant, chemical resistant plastic laminate sheets.

The exposed edges are covered by ABS, for safety reasons and to provide further durability.

SW

piano in legno massello/solid wood worktop

SS

piano in acciaio INOX/stainless steel worktop

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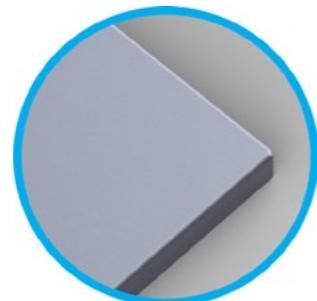
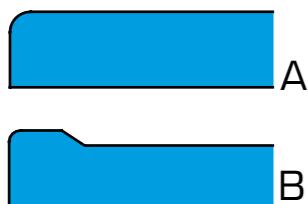
This worktop can be realized using different SS alloys (AISI 304 or AISI 316), according to the required application. SS worktops are realized using first choice steel sheets, cold bended, and can be provided with raised edge all around. The surface is completely seamless, perfectly shining and with no joints.

EP

piano in RESINA EPOSSIDICA/EPOXY RESIN worktop

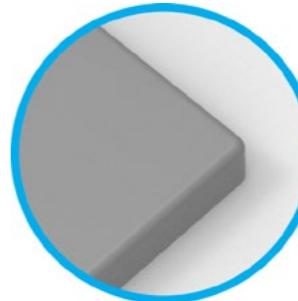
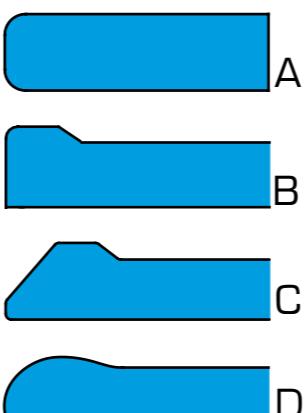
colori standard/standard colors

GR

piano in GRES MONOLITICO/CERAMIC worktop

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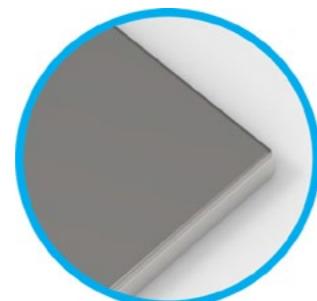
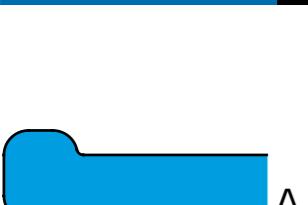
Epoxy resin worktops (such as Durcon) are obtained from molding and, then properly shaped using CNC (Computer Numerical Control) machines. These worktops are very easy to clean and decontaminate, are perfectly smooth and non-porous, very resistant to stains and to laboratory reagents.

Together with ceramic worktops, they represent one of the best choices when it comes to employment with very aggressive acids.

SP

piano in ACCIAIO PORCELLANATO/enamelled steel worktop

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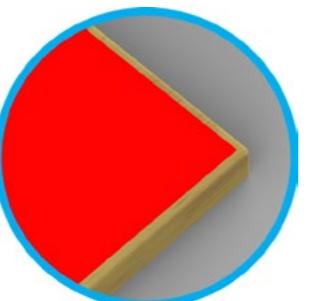
Enamelled steel is obtained by applying special ceramic enamel on a steel substratum. These worktops provide good resistance to most of common chemicals

ZL

piano in BETULLA LAMINATO/LAMINATED BIRCH worktop

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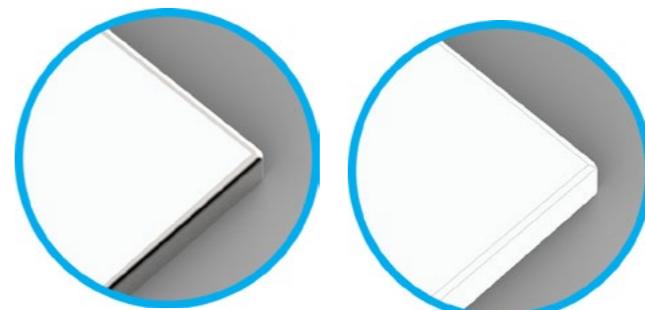
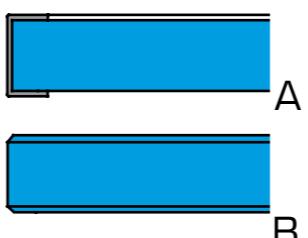
These worktops are made of bilaminated birch and they are ideal for desks.

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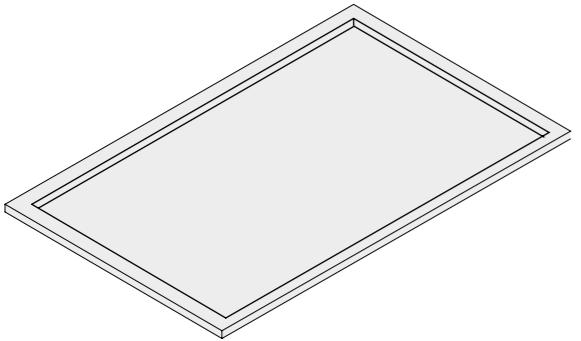
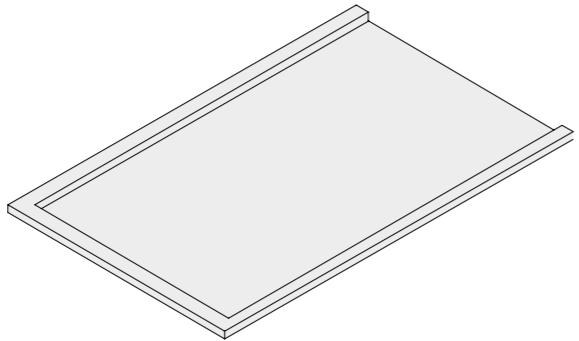
Worktop realized in tempered glass with optional AISI 304 stainless steel frame around its edges.

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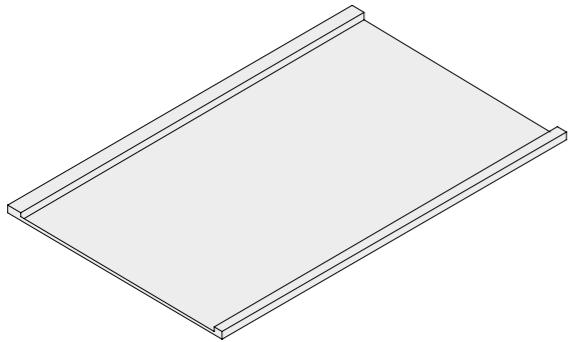
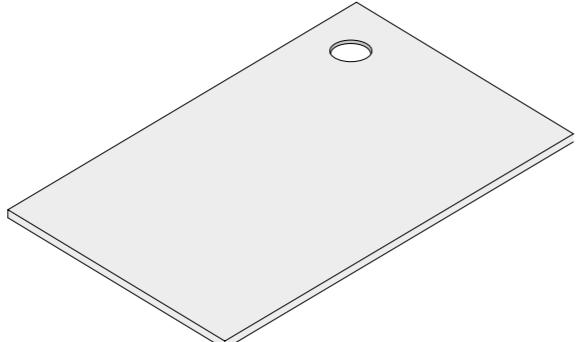
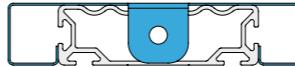
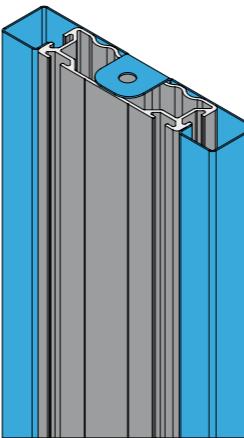
Pillar standard depth (100 mm) can be customized up to 150 mm.

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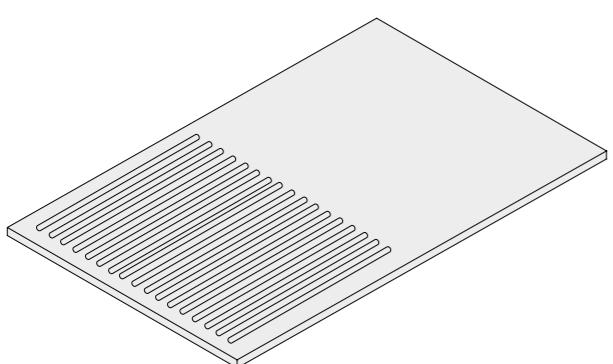
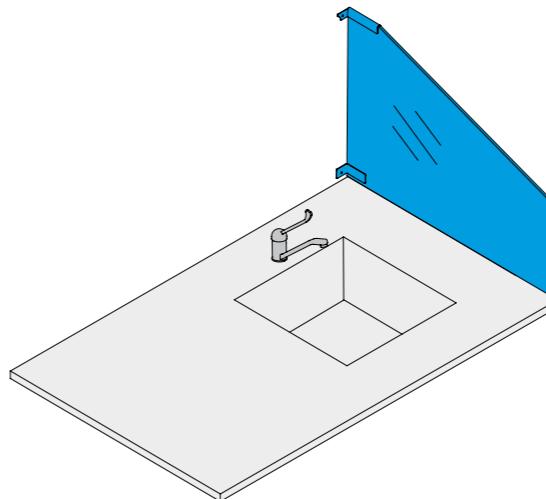
The lateral drain grooves are available for Corian, PP and ceramic worktops.

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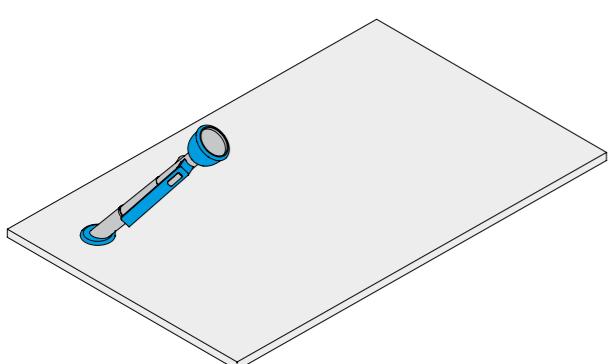
The lateral glass can act either as a splashback shield (if fitted near to a sink), or as a partition for benches and desks.

The panel can be transparent or opaque and can also be realized in anti-acid polycarbonate.

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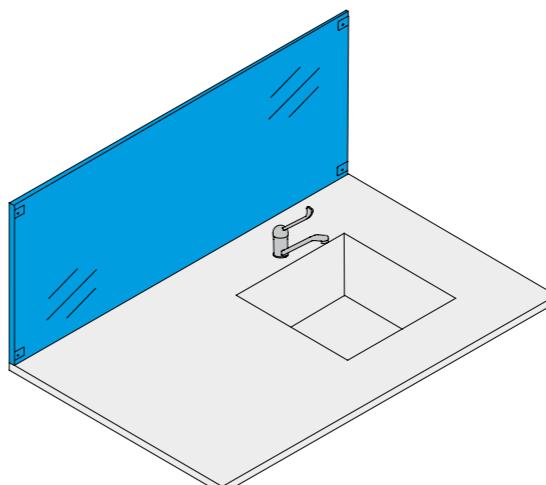
*Extractable eye-washer, fitted on the worktop.
The eye-washer gives first aid in case of accidental contact of the eyes with chemicals.
The water jet and the embedded regulator comply to DIN EN 246.*

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The rear glass can act either as a splashback shield (if fitted near to a sink), or as a partition for benches and desks.

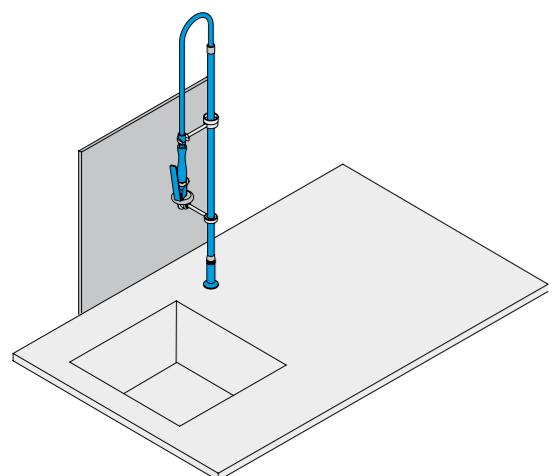
The panel can be transparent or opaque and can also be realized in anti-acid polycarbonate.

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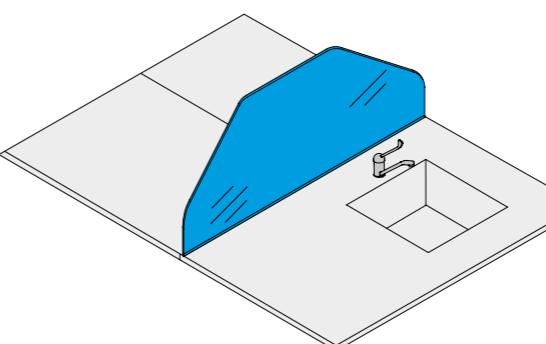
Entirely realized in chrome brass with stainless steel flexible hose and swivelling fitting.

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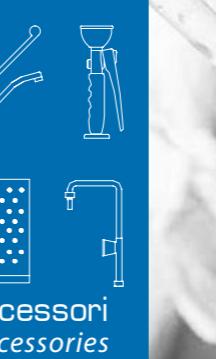
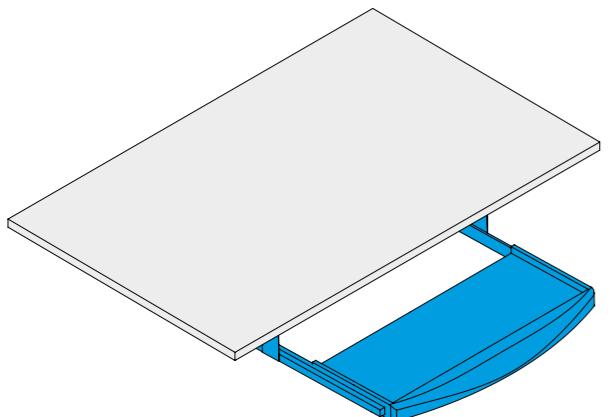
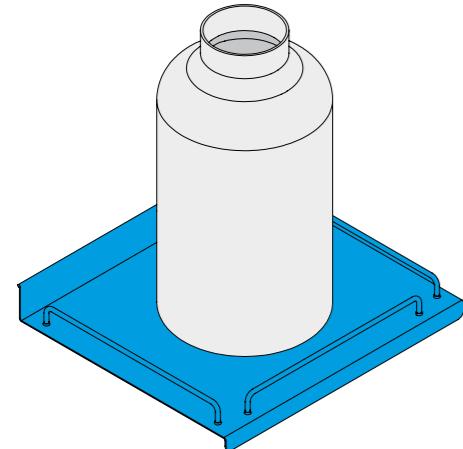
Pull-out keyboard tray fastened to the bench structure.

BT

portabottiglia / bottle holder

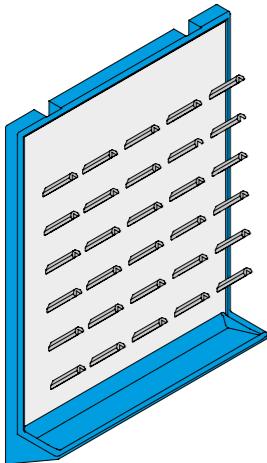
Portabottiglia in acciaio inox.

Stainless steel bottle holder.

CV

scolavetrerie / glassware drying rack

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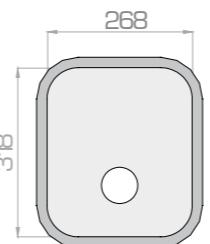
The drying rack for laboratory glassware is made of plastic material, and is fitted to the service spine. Plate is made up of 72 racks for standard size items, and 11 racks for smaller items. The rack is equipped with a drainage pipe. Also more than one rack can be installed on the service spine to obtain a complete wall unit.

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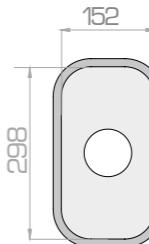
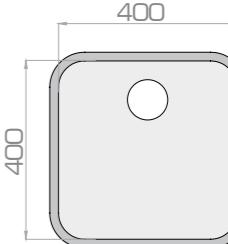
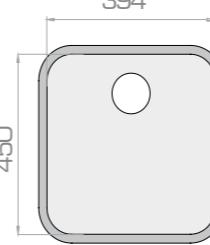
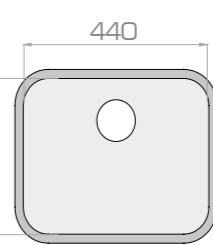
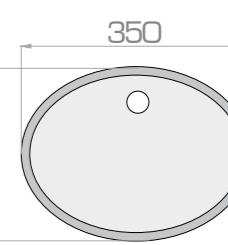
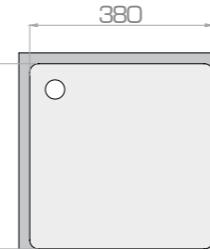
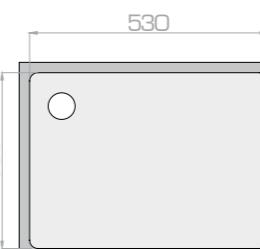
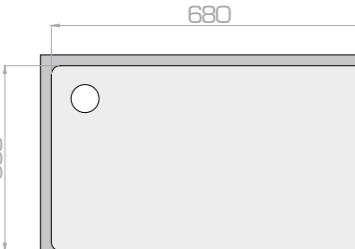
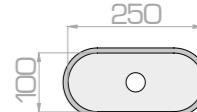
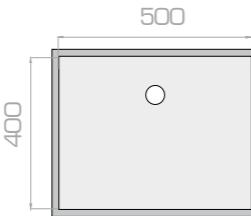
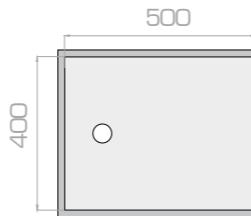
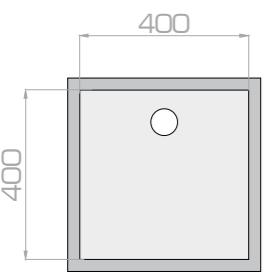
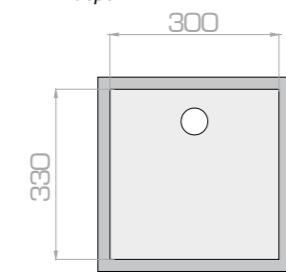
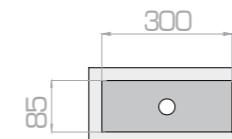
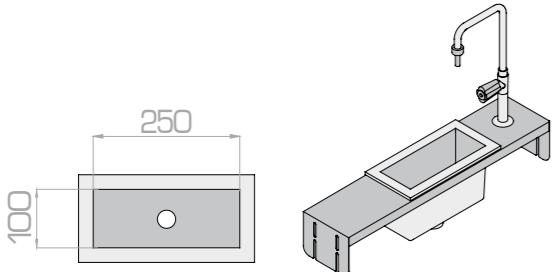
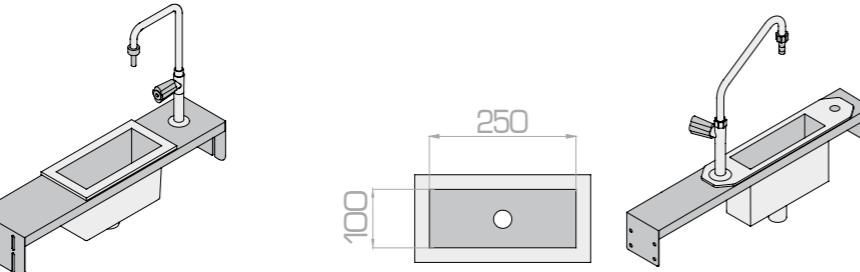
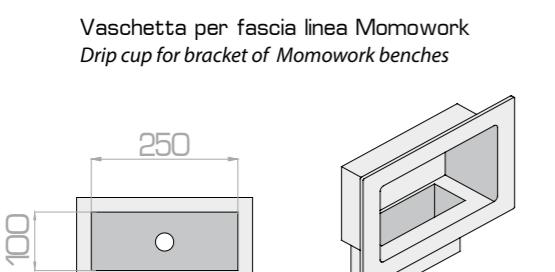
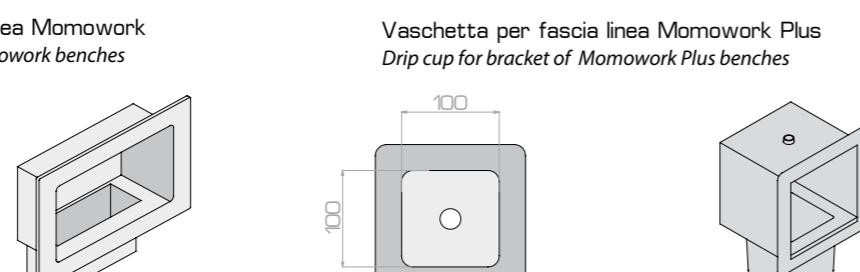
accessori
accessories

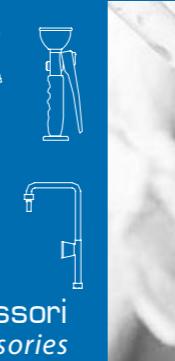
vasche in Polipropilene / Polypropylene sinks

Le vasche in Polipropilene sono accessori dei piani in Polipropilene, HPL, PVC e laminato. Polypropylene sinks are suitable for Polypropylene, HPL, PVC and laminate worktops.

codice / code
VSC-PP-1codice / code
VSC-PP-2codice / code
VSC-PP-3profondita' / depth
300 mmcodice / code
VSC-PP-4profondita' / depth
200 mmcodice / code
VSC-PP-5profondita' / depth
250 mmcodice / code
PS-MWcodice / code
PS-MWPprofondita' / depth
100 mmcodice / code
TW-MWcodice / code
TW-SKYprofondita' / depth
100 mm

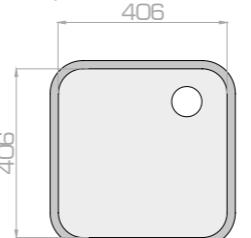
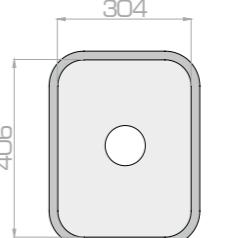
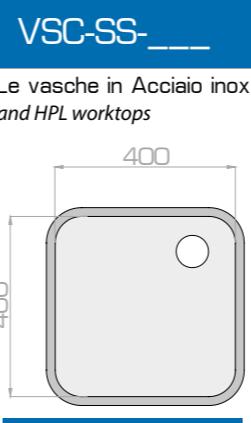
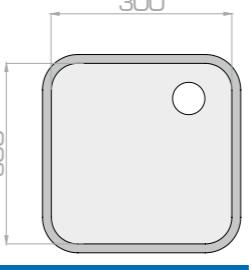
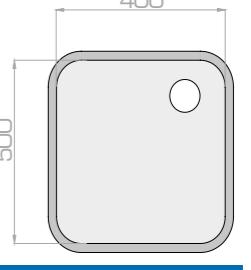
Vaschetta per fascia linea Momowork Plus
Drip cup for bracket of Momowork Plus benches

Vaschetta per Totem linea Skylab
Drip cup for Totem of Skylab benches

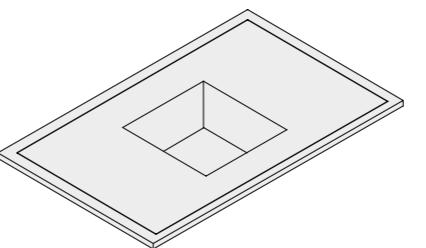
accessori
accessories

vasche in Resina Epossidica / Epoxy Resin sinks

Le vasche in Resina Epossidica sono accessori dei piani in resina epossidica. EpoxyResin sinks are suitable for Epoxy Resin worktops.

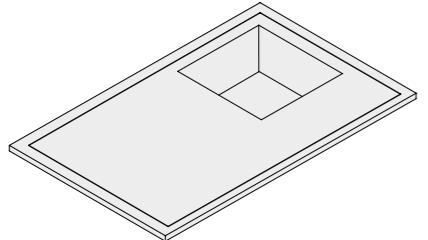
codice / code
VSC-DC-01profondita' / depth
210 mmcodice / code
VSC-DC-02profondita' / depth
222 mmcodice / code
VSC-SS-1profondita' / depth
250 mmcodice / code
VSC-SS-2profondita' / depth
180 mmcodice / code
VSC-SS-3profondita' / depth
300 mm

PP-VC

piano in Polipropilene con vasca integrata
Polypropylene worktop with embedded sink

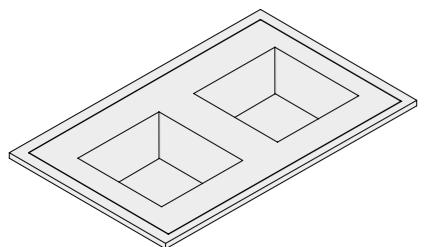
codice code	lunghezza piano worktop length	profondita' piano worktop depth	vasca l x w x p sink l x w x d
PP60X75-X4-VC	600	750	400x500x200
PP120X75-X4-VC-1	1200	750	300x300x200
PP120X75-X4-VC-2	1200	750	400x400x300
PP120X75-X4-VC-3	1200	750	400x500x300
PP150X75-X4-VC-1	1500	750	300x300x200
PP150X75-X4-VC-2	1500	750	400x400x300
PP150X75-X4-VC-3	1500	750	400x500x300
PP180X75-X4-VC-1	1800	750	300x300x200
PP180X75-X4-VC-2	1800	750	400x400x300
PP180X75-X4-VC-3	1800	750	400x500x300

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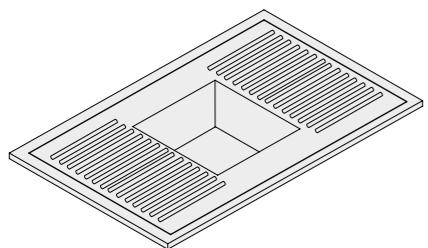
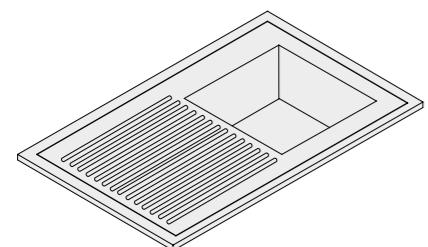
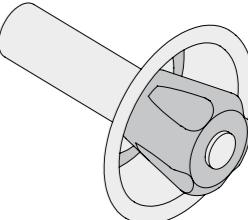
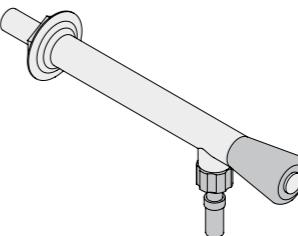
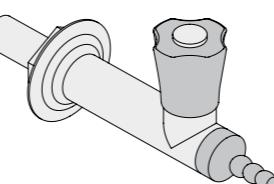
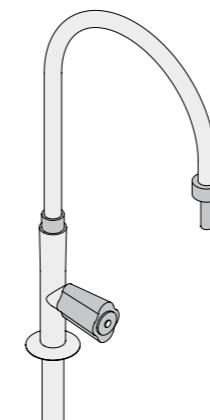
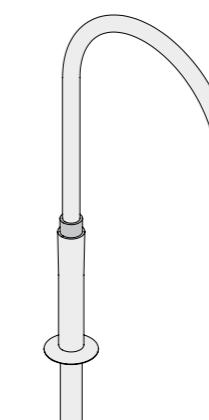
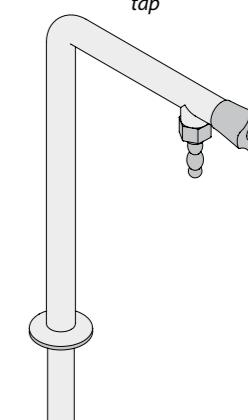
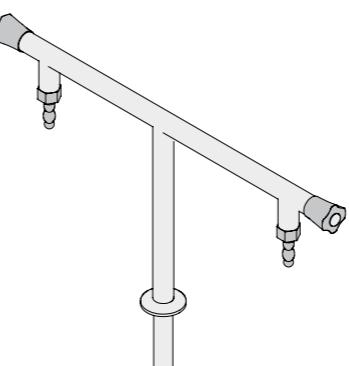
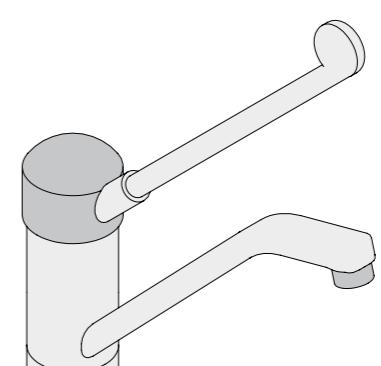
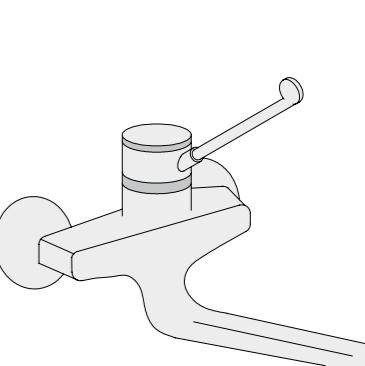
Q-____

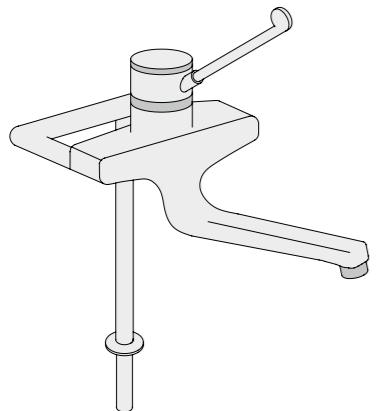
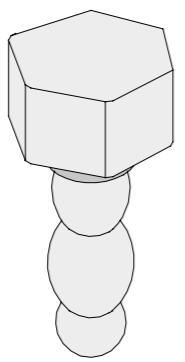
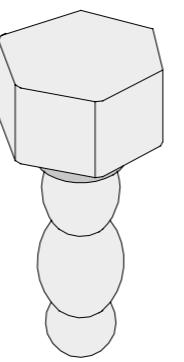
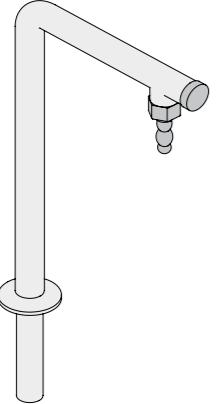
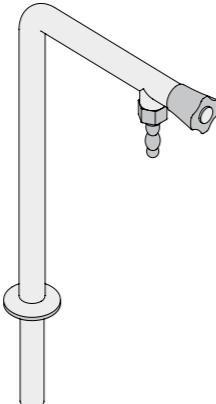
rubinetti e miscelatori acqua / water taps and mixers

codice / code
Q-01rubinetto con portagomma svitabile
straight tap with removable nozzlecodice / code
Q-02rubinetto 90° a parete
90° wall-mounted tapcodice / code
RC-01 / RC-02rubinetto di comando remoto acqua
/ demineralizzata
water / demineralized water remote control
tapcodice / code
Q-03rubinetto a collo di cigno, girevole
swivelling gooseneck spoutcodice / code
Q-04rubinetto a collo di cigno, girevole
swivelling gooseneck spoutcodice / code
Q-05rubinetto a "L"
"L" tapcodice / code
Q-06rubinetto a "T"
"T" tapcodice / code
Q-07miscelatore a leva clinica con areatore
mixer with long lever and aeratorcodice / code
Q-08miscelatore a parete, leva clinica e areatore
wall-mounted mixer, long lever and aerator

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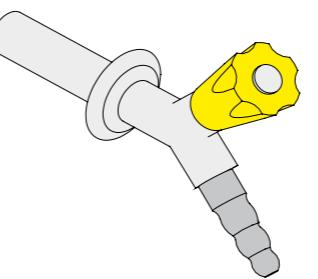
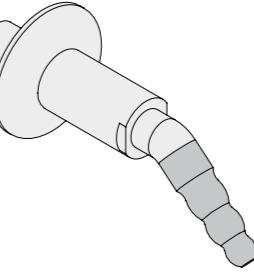
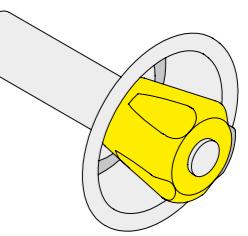
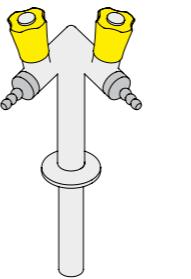
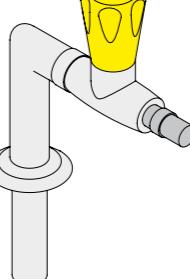
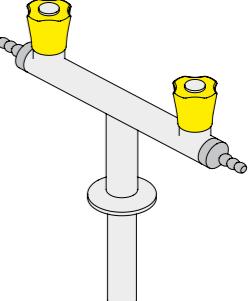
All taps are manufactured according to DIN 12918 norm.

The threaded demontable nozzle is manufactured according to DIN 12898 norm.

All handles are made of anti-acid polypropylene, and the colors comply to EN 13792 norm.

Maximum operating pressure is 7 bar. Temperature operating range 30°C to 150°C.

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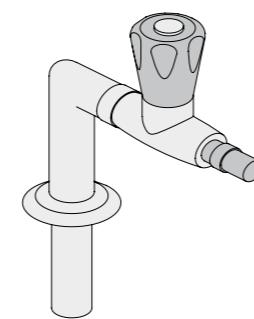
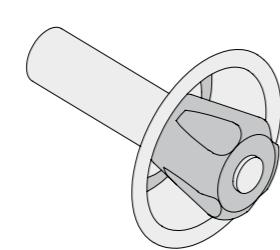
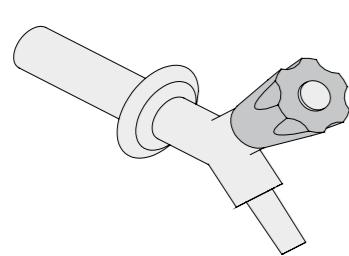
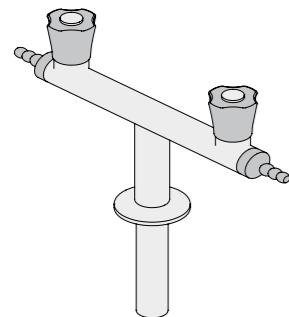
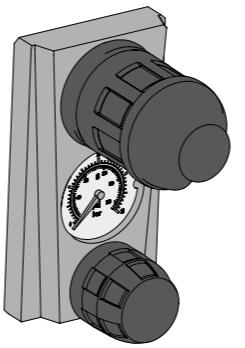
All taps are manufactured according to DIN 12918 norm.

The threaded demountable nozzle is manufactured according to DIN 12898 norm.

All handles are made of anti-acid polypropylene, and the colors comply to EN 13792 norm.
Maximum operating pressure is 7 bar. Temperature operating range 30°C to 150°C.

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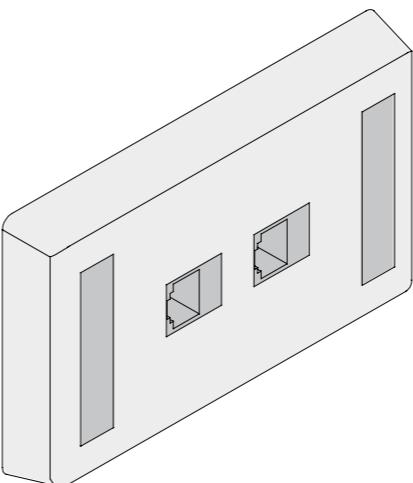
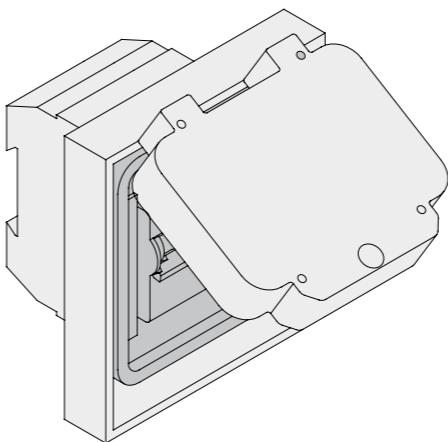
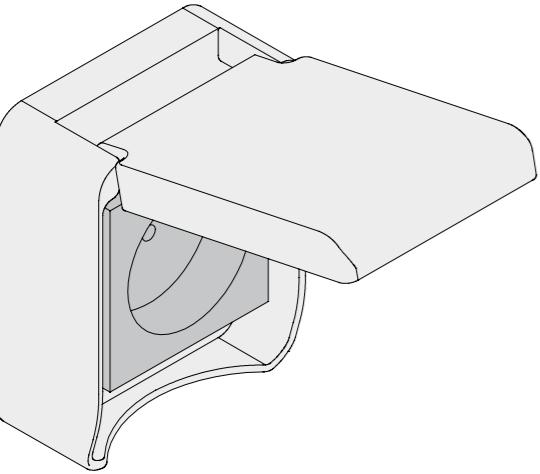
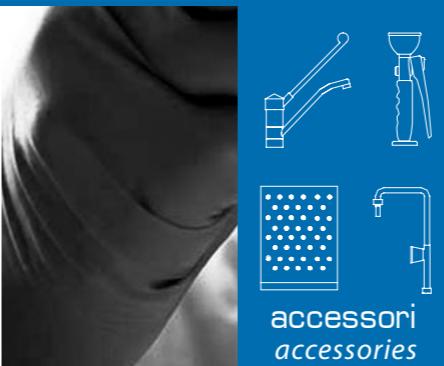
The provided MCB is certified by CE and IMQ for compliance to CEI 23-3(EN60898) and EN 60947-2 norms.

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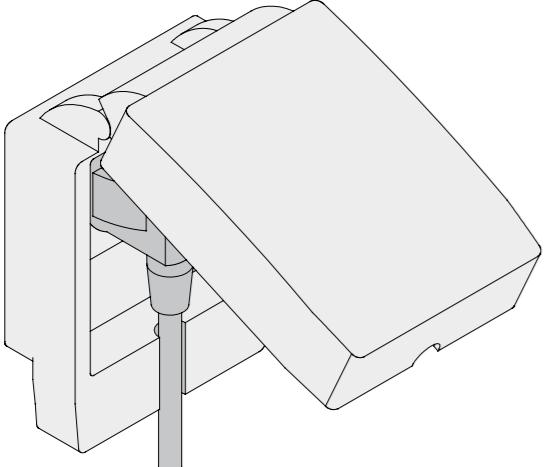
*IP65 Power outlet protection grade, according to IEC 60529 norm, IK09 protection grade against mechanical shocks, according to IEC 62262 norm; heat resistance 850°C, according to IEC 60695-2-1 norm.
10/16A, 250V.
Different socket standards available.*

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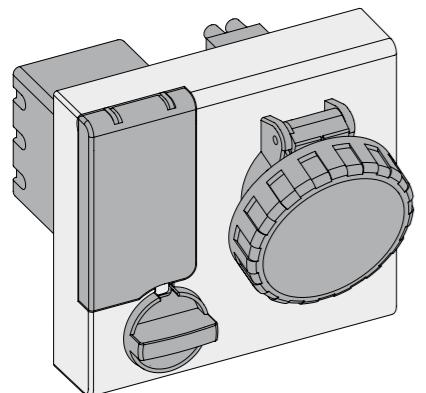
*IP66 Power outlet with fitted plug protection grade, according to IEC 60529 norm, IK08 protection grade against mechanical shocks, according to IEC 62262 norm; heat resistance 850°C, according to IEC 60695-2-1 norm.
10/16A, 250V.
Different socket standards available.*

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*IP67 interlocked power outlet, with embedded protection fuse and switch.
16 A, 50/60 Hz.
Sockets are marked IMQ and CE, and complies to EN 60309-1, EN 60309-2 norms.*

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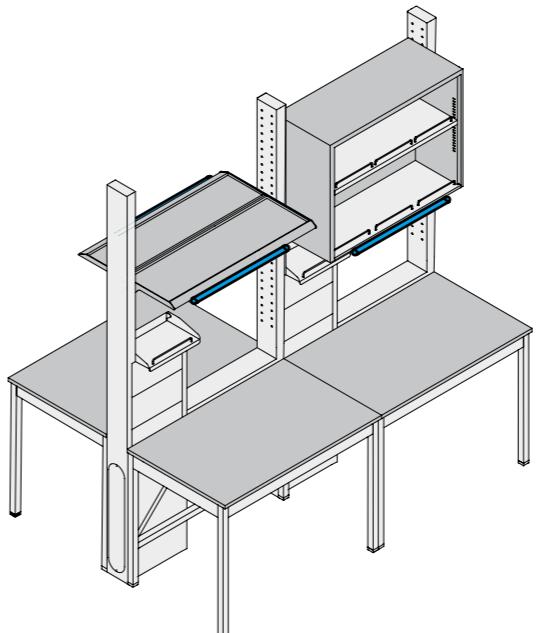
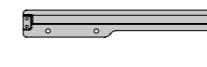
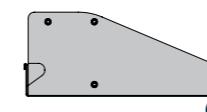
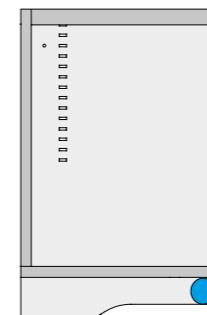
Lamps can be installed below the overhead cabinets, the storage shelves, and the reagent shelves, for all the benches lines.

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*Neon tube lamp 220-240V, 50/60Hz, IP20, G5/T5 fitting.
PVC case, polycarbonate covering, CE compliant.*

codice/code	per banco da/for bench length
NEON60	600
NEON80	800
NEON90	900
NEON100	1000
NEON120	1200
NEON150	1500
NEON180	1800

Lampada al neon 220V, 50Hz, IP67, 4000K. Omologazione VDE, conformita' CE.

*Neon tube lamp 220V, 50Hz, IP67, 4000K.
VDE omologation, CE compliant.*

codice/code	per banco da/for bench length
NEON60	600
NEON80	800
NEON90	900
NEON100	1000
NEON120	1200
NEON150	1500
NEON180	1800

Lampada a led 220V, 50Hz, IP65, 4500 K. Omologazione VDE, conformita' CE.

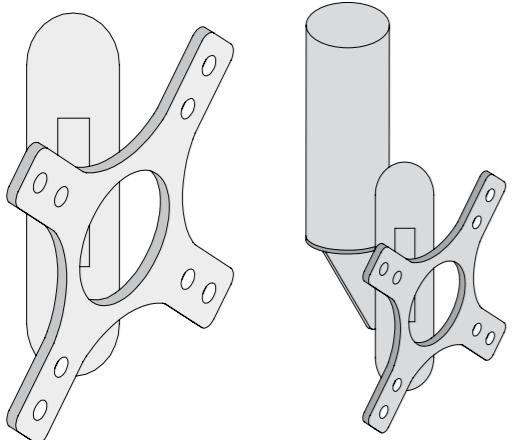
*Led lamp 220V, 50Hz, IP65, 4500 K.
VDE omologation, CE compliant.*

codice/code	per banco da/for bench length
LED60	600
LED80	800
LED90	900
LED100	1000
LED120	1200
LED150	1500
LED180	1800

BMT

braccio monitor / monitor supporting arm

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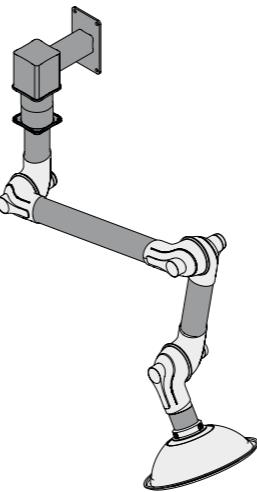
*The monitor supporting arm can be either suspended or wall-mounted.
It is suitable for 10" to 24" screens, having a maximum weight of 15Kg.
180° rotation, -5%+15° inclination.*

AF-C

cappa a braccio / articulated arm local extractor

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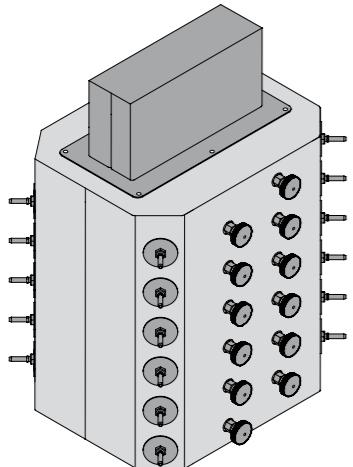
*The articulated arm local extractor is made up of anodized aluminium elements, with three joints.
The extraction bowl (350mm diameter) is made up of transparent polypropylene or aluminium.
The total length is 1500 mm. It can be wall mounted, ceiling mounted or fitted to the service spine. It can be equipped with extractor (M) and activated carbon filter box (P).*

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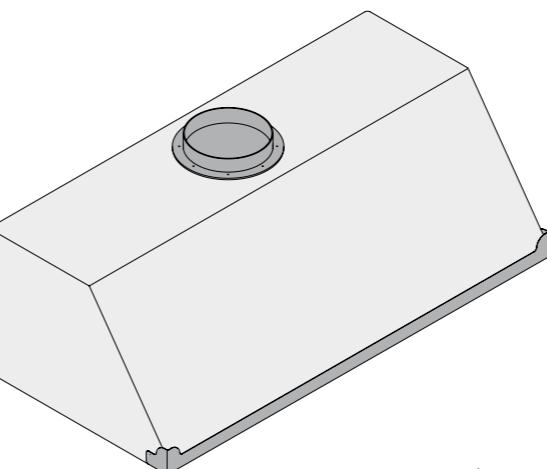
*Satellite module is designed to enhance and improve the performances of the gas distribution system.
Strongly recommended in case many gases are required, this module allows to centralize the second stage pressure reducer station, so that considerable savings can be achieved.*

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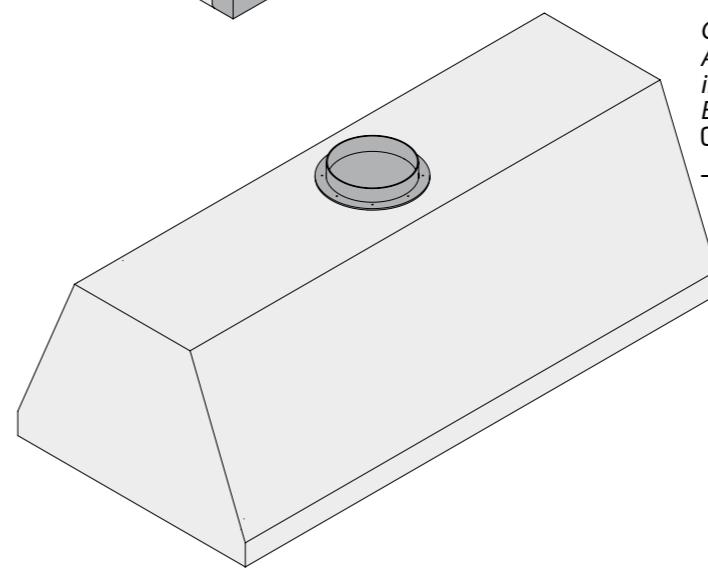
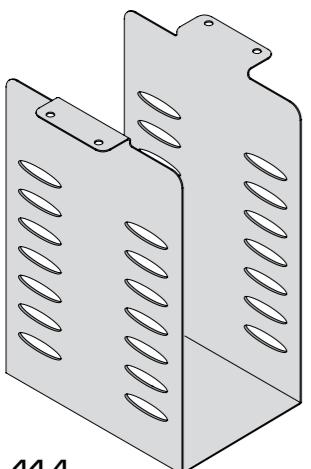
*Canopy hood for local extractor.
Available for fastening on bench pillars, on wall or hanging from ceiling.
Entirely realized in Anti-acid PVC thermo-welded. (code CH-__-PVC), or in AISI 304 Stainless Steel (code CH-__-SS).*

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Realized in epoxy-powder painted steel sheets.

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A fume cupboard (also called 'fume hood') is a very technologic machinery, employed in laboratory environments for safeguarding operators while handling potentially dangerous, toxic or hazardous substances.

Besides the simple protection from splashes, blasts or explosions, fume cupboards are intended to keep the noxious vapors away from the operator, in order to prevent their possible inhalation.

Fume cupboards are designed to generate a continuous forced ventilation flow through the frontal opening: forced-draught ventilation is produced and maintained by an electric fan, installed either onboard the fume cupboard or placed in a remote position, inside or outside the building (picture 1). The exhaust airflow is then expelled in the external environment.

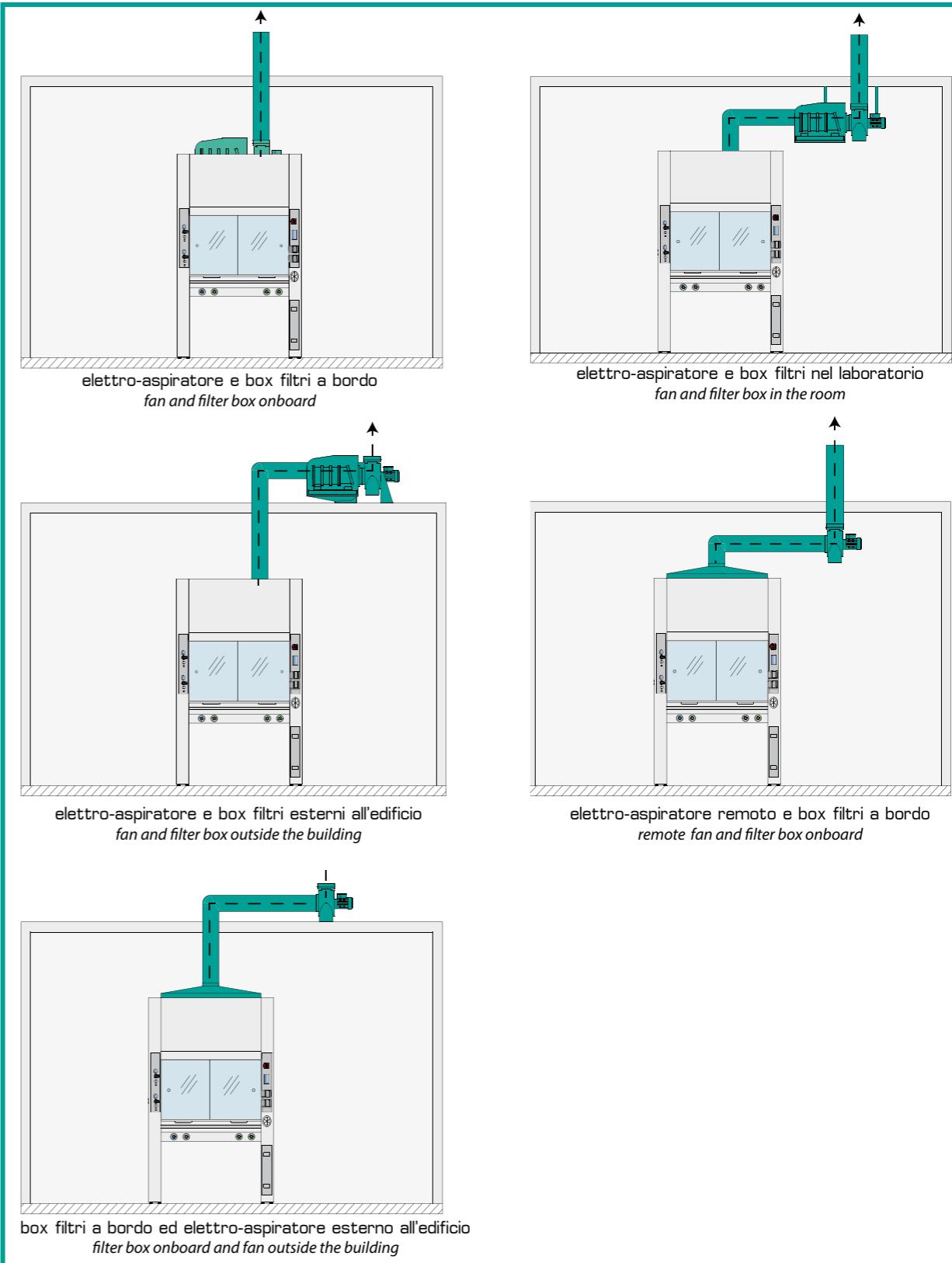
*The correct choice of the fan is strictly connected to the head loss along the exhaust duct, and is therefore a key issue for granting the proper working of a fume cupboard. Thanks to its long-time experience, technical staff from **MOMOLINE** is able to suggest the most adequate solution for every different application.*

Should the particular substances employed under the fume cupboard represent a risk for the environment, a filter box or a fume scrubber can be installed; different kinds of filters can be provided, according to the specific employed chemicals.

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Fume cupboards are generally divided into two main groups: ducted fume cupboards and recirculatory (or 'ductless') fume cupboards (picture 2).

Ducted fume cupboards are equipped with a dedicated exhaust duct, used to draw exhaust air to the outdoor environment.

Recirculatory fume cupboards do not require any exhaust duct, as exhaust air is fed back into the room after passing through some high-efficiency filters. In this case the fume cupboard is more compact and its installation is significantly easier, despite being not suggested in case particularly concentrated substances have to be employed, as well as in case of intensive use.

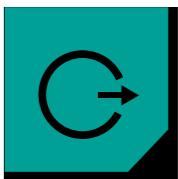
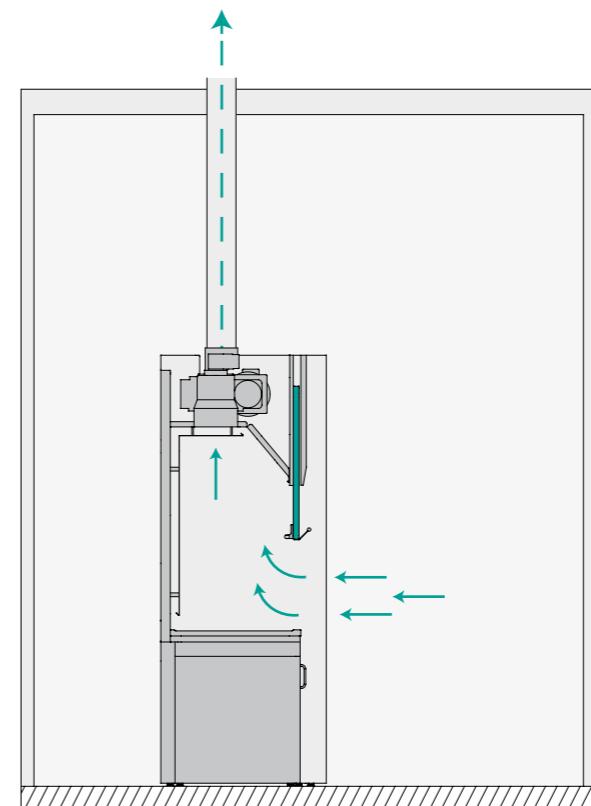
Performances and efficiency of a fume cupboard can be investigated, and therefore evaluated, relying the concepts of 'containment' and 'robustness' both defined in EN 14175 Norm : the first is defined as the capability of a fume cupboard to retain airborne contaminants within the work space, the latter as the capability to maintain containment during external front disturbances.

The efficiency of a fume cupboard is therefore defined as the capability of efficiently removing airborne contaminants from the work space.

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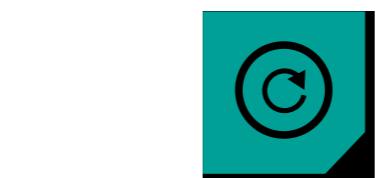
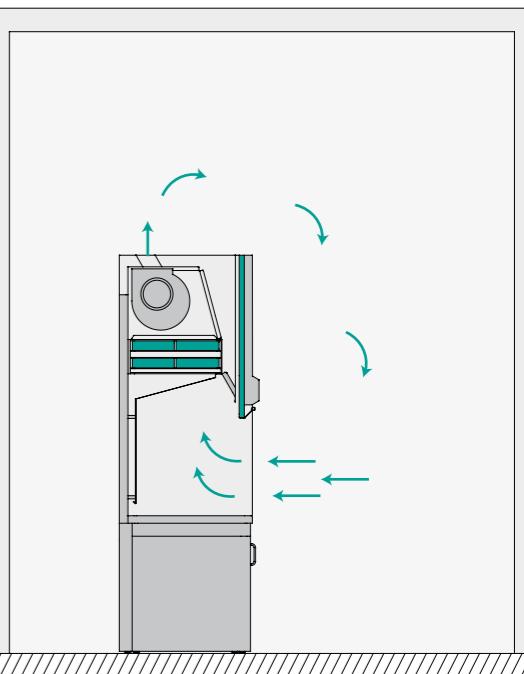
Actual performances of a fume cupboard strictly depend on the particular location where this is installed; for this reason, actual performances of fume cupboards shall be checked and certified not only after the installation, but even on a regular basis during their normal lifecycle, so that the operator's safety is always guaranteed.

*For this reason, after the installation of a fume cupboard, **MOMOLINE** staff can carry out a very strict testing procedure according EN 14175 Norm - Parts 3,4. After all the tests have been carried out, a Testing Certificate is issued, and all test reports are enclosed.*

Fume cupboards can be either from manual or automatic type. In the first case exhaust fan is simply controlled by an ON/OFF switch, and the actual airflow is not controlled by any electronic control.

On the other hand, automatic fume cupboards are fitted with an electronic controller continuously monitoring the airflow, so that their efficiency is constantly guaranteed.

Automatic fume cupboards can be optionally equipped with VAV (Variable Air Volume) system.

Furthermore, fume cupboards can be provided with specific alarms and active monitoring systems, able to inform the operator (in real time) of the actual efficiency and containment of the fume hood, in order to promptly intervene should the operator safety be endangered.

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The efficiency of a fume cupboard is strictly related to its 'face velocity', expressed in [m/s] of the air flow and defined as the velocity through its frontal opening, between the worktop and the sash.

This airflow is responsible for keeping noxious substances away from the operator and for preventing any accidental leak, in order to grant the robustness and the containment efficiency of the fume cupboard.

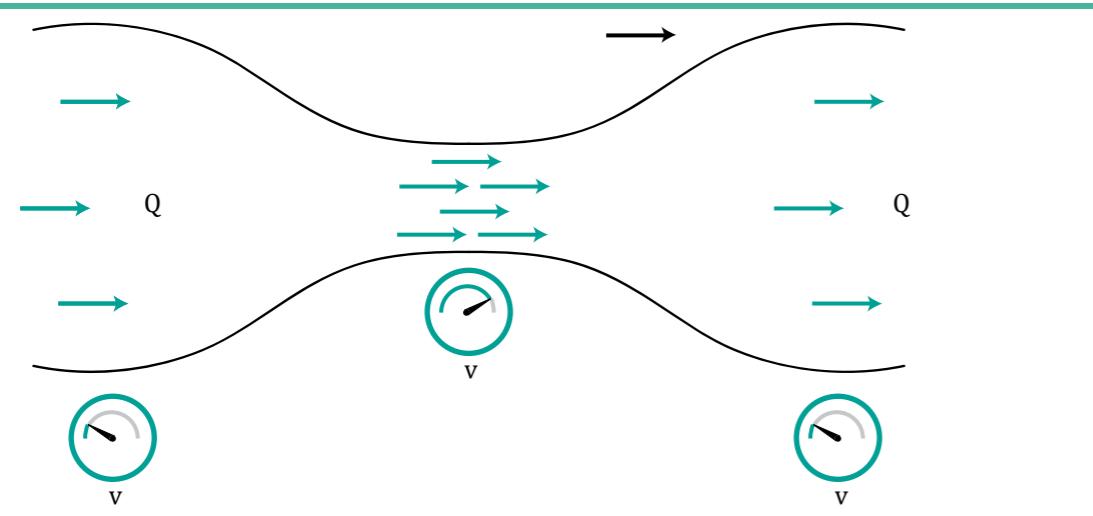
Many studies assume the face velocity of a fume cupboard as a reference parameter for determining its overall efficiency.

Major European and American Organizations agree that the face velocity should never be lower than 0.4 – 0.85 m/s, depending on the level of danger and toxicity of the specific substances employed.

Velocity (v) and total quantity (Q) of a fluid passing through a duct having cross section equal to S are related to equation A.

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The meaning of this relation can be easily understood with the help of the "Venturi Pipe" (picture 3).

If a certain quantity of fluid enters a duct, then this same quantity has to come out through the other end of the duct at any given time, provided that no other exit exists. For this reason, the Volumetric Flow rate Q is constant along the pipe.

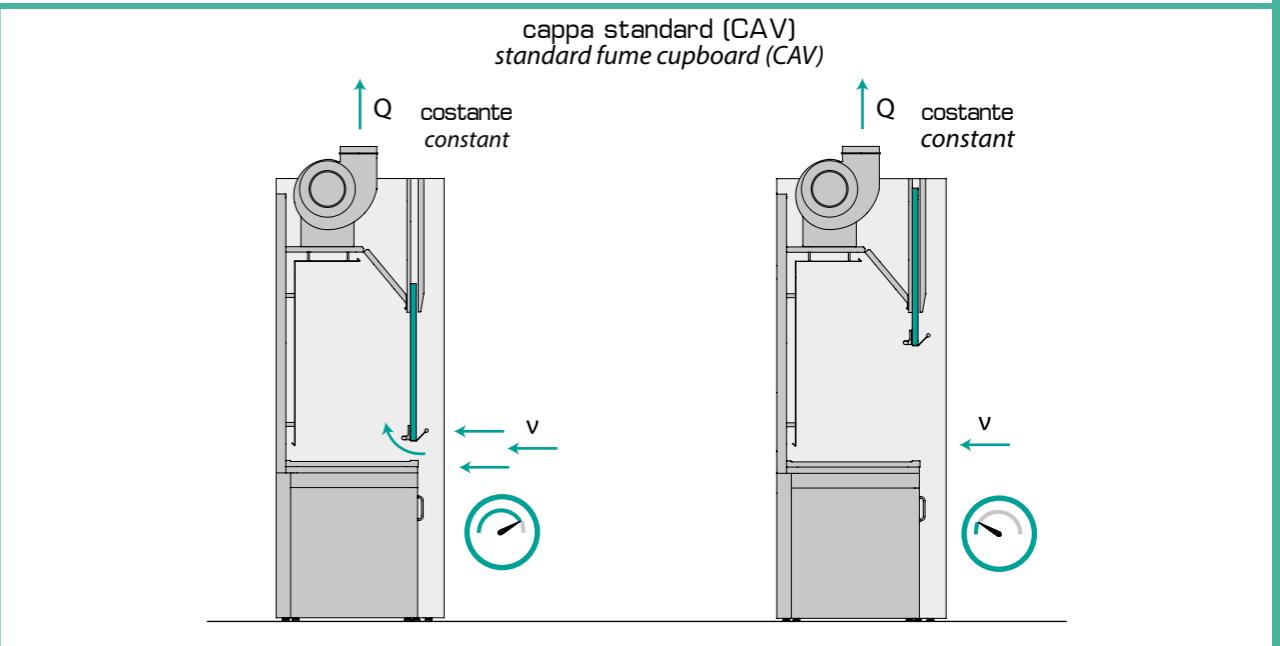
So, if the pipe narrows the fluid has to expedite, if the pipe widens it has to slow down, so that inlet and outlet flows always match.

This is exactly what happens in fume cupboards: when entering into the hood, the airflow is forced through the narrowing created between the worktop and the sash, like in Venturi Pipe.

In this case, the air flow rate Q is controlled by the rotational speed of the blower: if the blower rotates faster, then it is able to draw a larger air volume.

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Fans of conventional fume cupboards, without any airflow regulation, rotate at constant speed, and a Constant Airflow Volume is extracted (CAV system – picture 4).

In this case, when the sash is lifted up, then the S opening increases, and the inlet airflow slows down for the Venturi Principle. This may expose the operator to some risks, since the face velocity might decrease below the safety threshold.

The fan rotates at constant speed, so that the overall airflow Q is constant

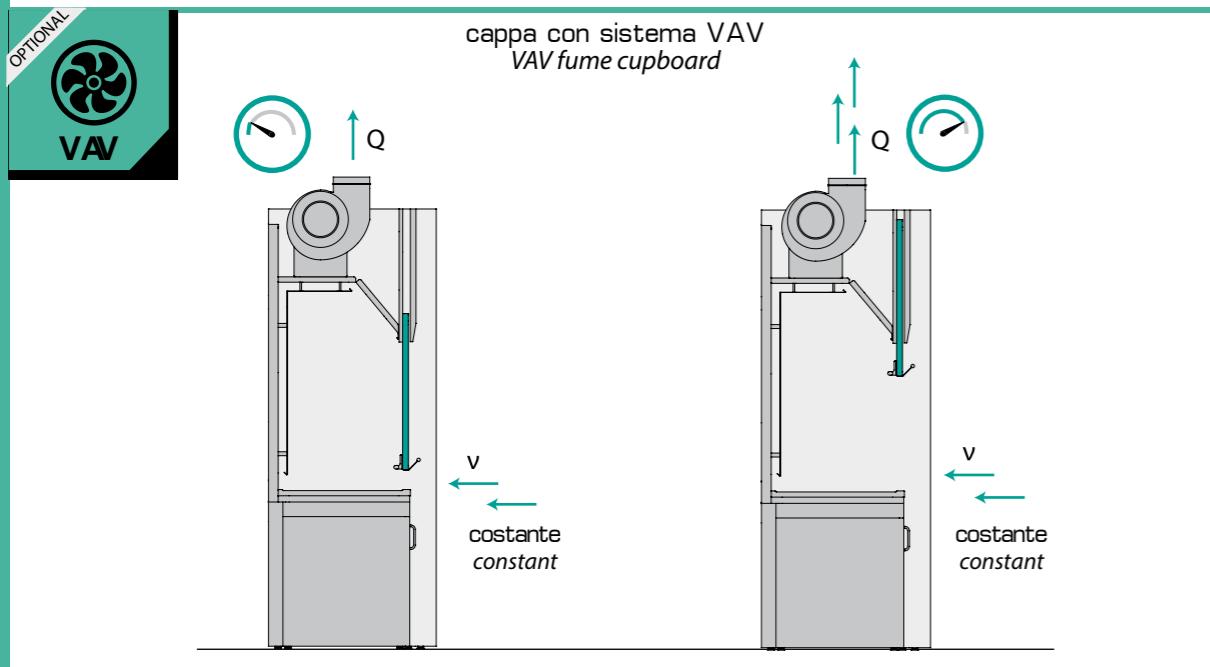
When the sash goes up, the S opening increases

To keep the airflow Q constant, the front air velocity v decreases

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VAV (Variable Airflow Volume) fume cupboards (picture 5), on the other hand, embed a control logic monitoring the airflow in real time: once the target face velocity has been set, this is kept constant with the sash opening, by changing the airflow Q.

The target value for the face velocity depends on the particular employed substances.

The most common regulation device for adjusting the Airflow makes use of an Inverter to change the rotational speed of the fan. If the sash goes up, the fan rotates faster to increase the airflow, so that velocity is kept constant and equal to the target value.

VAV system allows the operator to be always safeguarded, whatever is the sash position, since the fume cupboard adjusts its working conditions according to its movement.

The rotational speed of the fan can be adjusted in order to modify the inlet airflow Q

When the sash goes up, the S opening increases.

The blower rotates faster, so that the airflow Q increases and the air velocity v does not change.

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The same adjustment can also be obtained using an air damper installed in the exhaust duct, managing the airflow through the pipe.

The employment of VAV also grants considerable energy savings: VAV allows the fan to operate at high speed only when necessary (open sash) and makes it slows down if the sash is lowered.

In CAV fume cupboards, on the other hand, the fan must always rotate at high speed to grant the operator safety also in case of lifted sash, despite this is not necessary when the sash is closed.

Furthermore, when the sash is closed, an high airflow velocity can cause very harsh turbulences, making the work under the fume cupboard very difficult (little light elements might be sucked into the fume cupboard).

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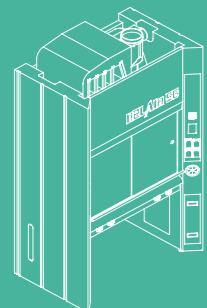
BELAIR56 is an automatic fume cupboard, standing as the last frontier of protection and safeguard for laboratory operators.

BELAIR56 has been designed and tested for many different applications, resulting able to grant full safety in every condition, even in case of heavy duty applications, when highly noxious chemicals are employed.

The aerodynamics of **BELAIR56** has been carefully designed to grant constant uniformity and efficacy of the extraction flows, so that containment performances are maximized.

The careful planning of aerodynamic flows has determined the implementation of really innovative technical solutions, as well as the application of cutting-edge technologies aimed to ensure top protection in every condition.

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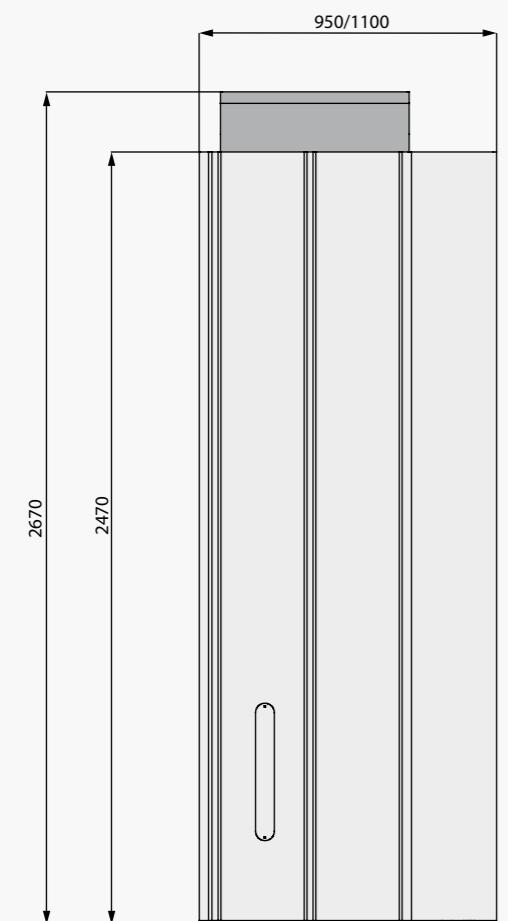
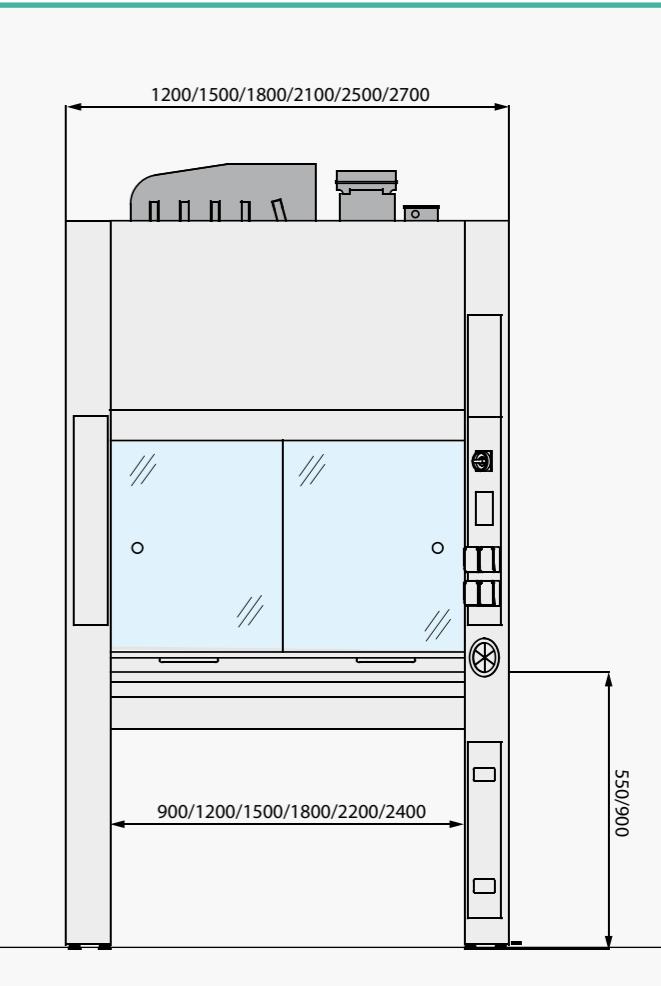
BELAIR56 has been certified for its compliance to EN 14175 Norm, and the certificate has been issued by Internationally recognized third bodies.

Besides the guarantee of safeguard and protection, **BELAIR56** fume cupboards empower every laboratory of a highly modern and impressive aspect, thanks to their soft and rounded lines and innovative design.

BELAIR56 fume cupboard is entirely realized using epoxy powder coated steel sheets, laser-cut and bended through cold processes.

*quote espresse in mm
dimensions expressed in mm

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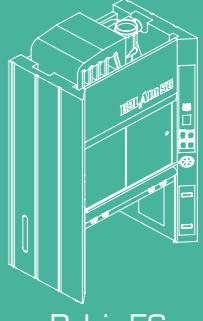
Steel sheets ensure a completely incombustible structure, so that **BELAIR56** is rated as 'Class 0' fire reaction. Furthermore, epoxy powder coating grants full protection from the most aggressive chemicals and from corrosion, making of **BELAIR56** a very exclusive product, suitable for installation in every kind of laboratory.

The exclusive employment of steel sheets makes **BELAIR56** a flagship product on the market, since common fume cupboards are often manufactured with wooden or plastic parts, offering much lower performances when it comes to durability and resistance to chemicals or fire attacks.

Upon specific request, and in case of necessity, **BELAIR56** fume cupboard can be entirely realized using Stainless Steel sheets: this ensures superior hygiene and decontamination performances.

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The position of services and utilities points has been carefully studied on the base of ergonomics principles, with the aim of making the use of the cupboard as pleasant as possible for the operator.

*Many different configurations are available for **BELAIR56** fume cupboard, so that every exigency or special necessity of customer can be fulfilled.*

Together with the wide choice of materials for worktop, many different accessories and optional items are available: motorized sash, double face fume cupboards or employment of explosion proof components are only some examples, but full details about all available optional items can be found in this Catalogue.

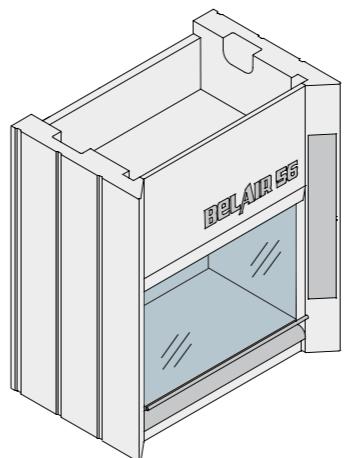
BELAIR56 fume cupboard is equipped with a highly technologic electronic control software, specifically designed for its management, and usable by the operator thanks to a very intuitive touch screen interface. The control software embeds a wide range of alarms and detecting probes, aimed to continuously monitor the working conditions of the fume cupboard, and ready to promptly intervene in case of necessity, so that operators are always safeguarded.

In addition to the standard software, Belair56 model can also embed some advanced airflow control features, designed to implement active safety and energy saving logics, such as the VAV, CLIMA and the TWIN SET systems.

Detailed specifications regarding all available accessories can be found in this Paragraph.

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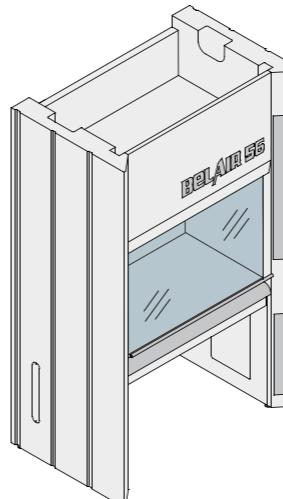
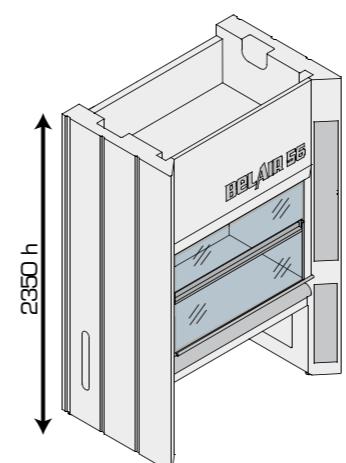
profondita' totale
total depth

950/1100 mm

codice code	larghezza piano worktop width	larghezza esterna external width
K90-S-BELAIR56	900	1200
K120-S-BELAIR56	1200	1500
K150-S-BELAIR56	1500	1800
K180-S-BELAIR56	1800	2100
K220-S-BELAIR56	2200	2500
K240-S-BELAIR56	2400	2700

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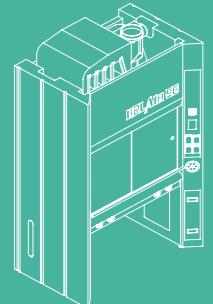
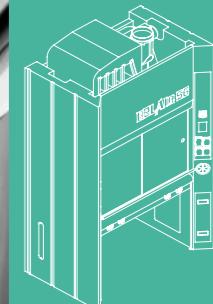
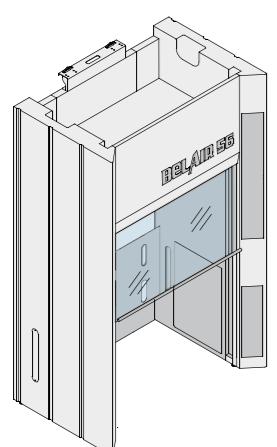
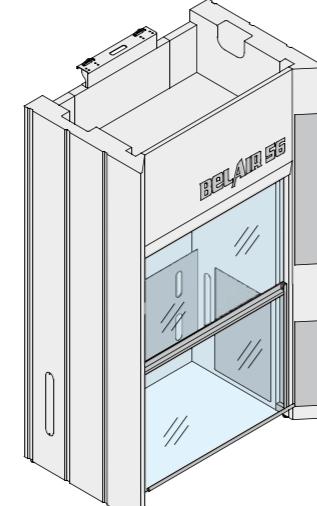
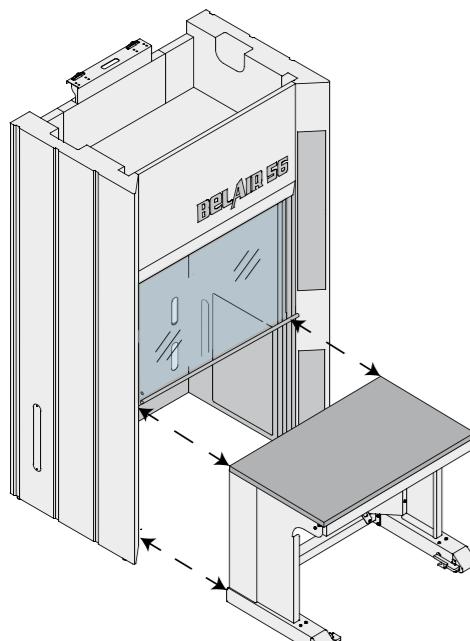
Walk-in version is not equipped with the conventional worktop to allow the insertion under the fume cupboard of tall equipment and trolley benches. Walk-in fume cupboards are often coupled to Shylab line trolley benches (code: BCAR), described in chapter 1 "lab benches".

cappa bifronte/double-sided fume cupboard

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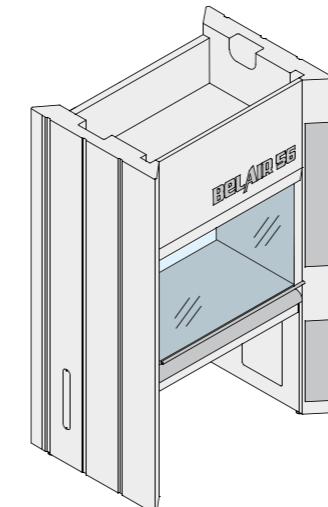
Double-sided fume cupboards can be used by two different operators at the same time. This fume cupboard is often located between different areas of the same laboratory, and is used to transfer dangerous samples.

K__-SB-BELAIR56

struttura bifronte base/double-sided basic structure

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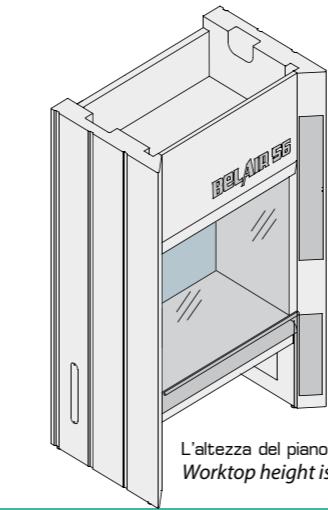
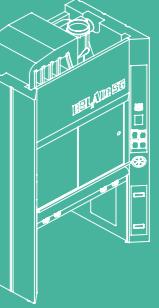
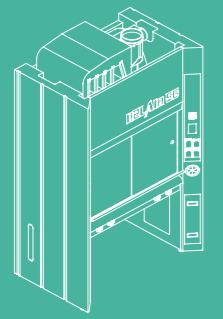
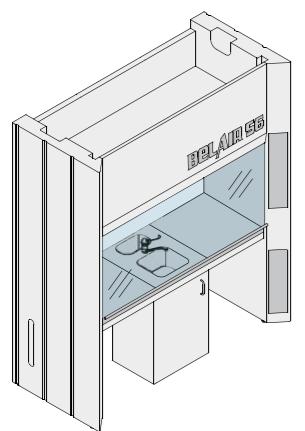
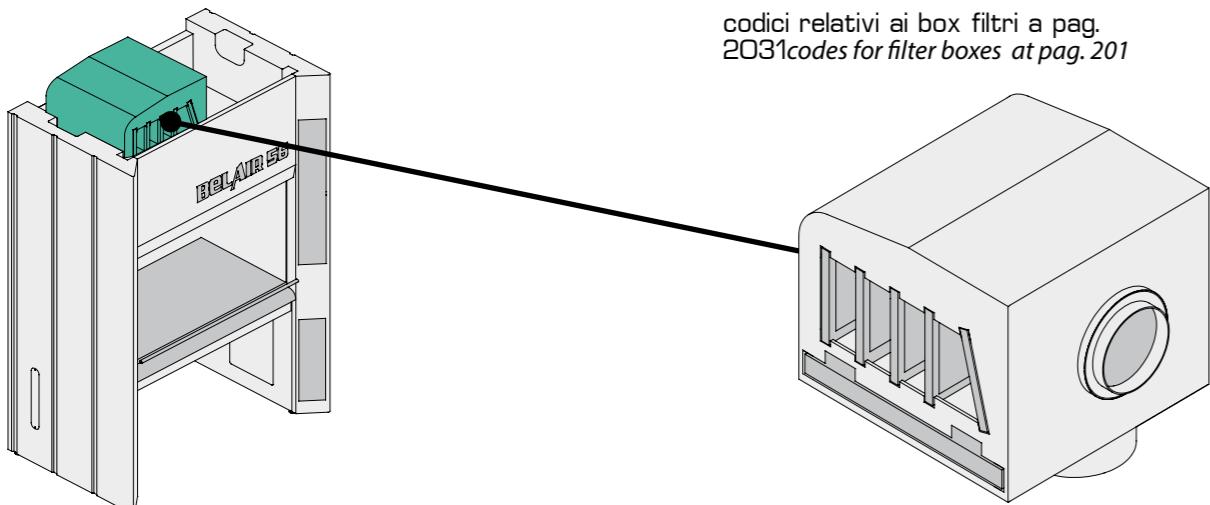
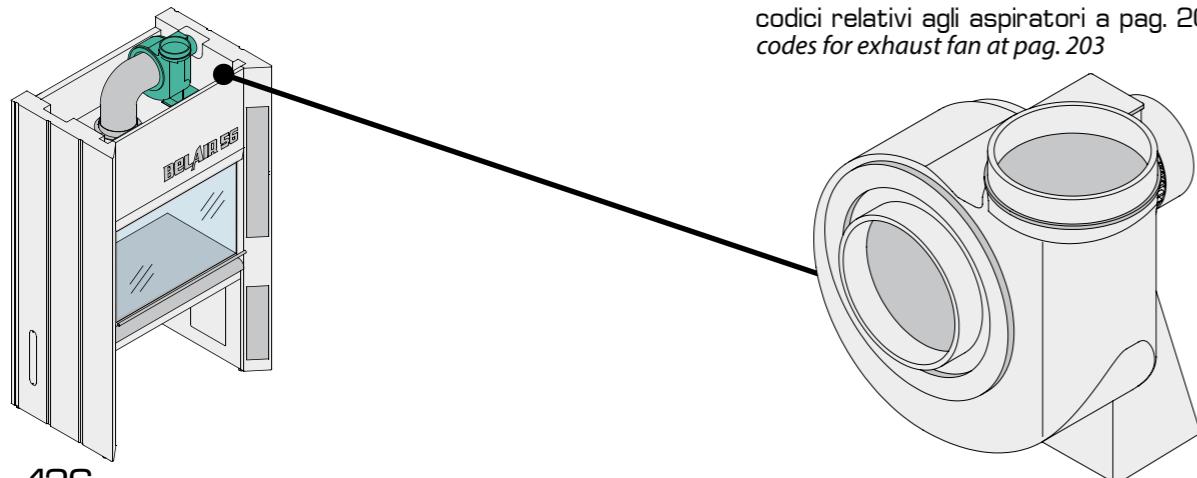
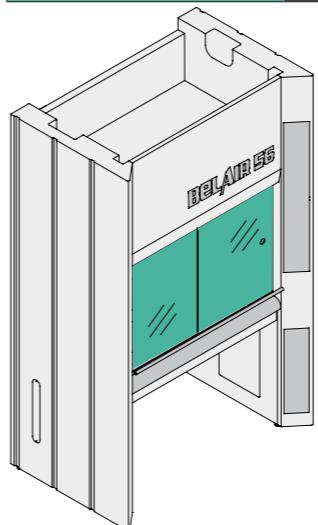
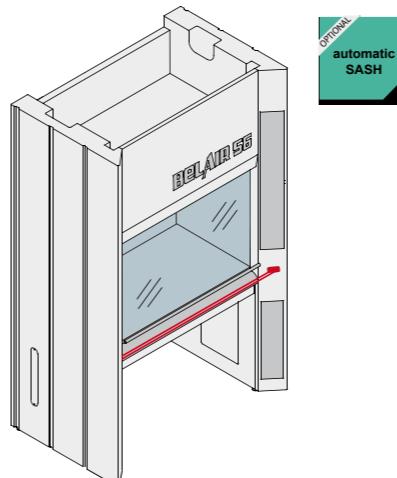
BELAIR56 is always equipped with a vertically sliding frontal safety glass, framed into a counter-balanced sash. Upon request, it is possible to provide a vertically sliding frame incorporating up to 4 glass panels, horizontally sliding.

Frontal sash is always equipped with a safety device granting its lock in case of failure of one or both its suspending chords, so that it is prevented from falling.

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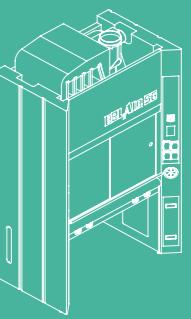
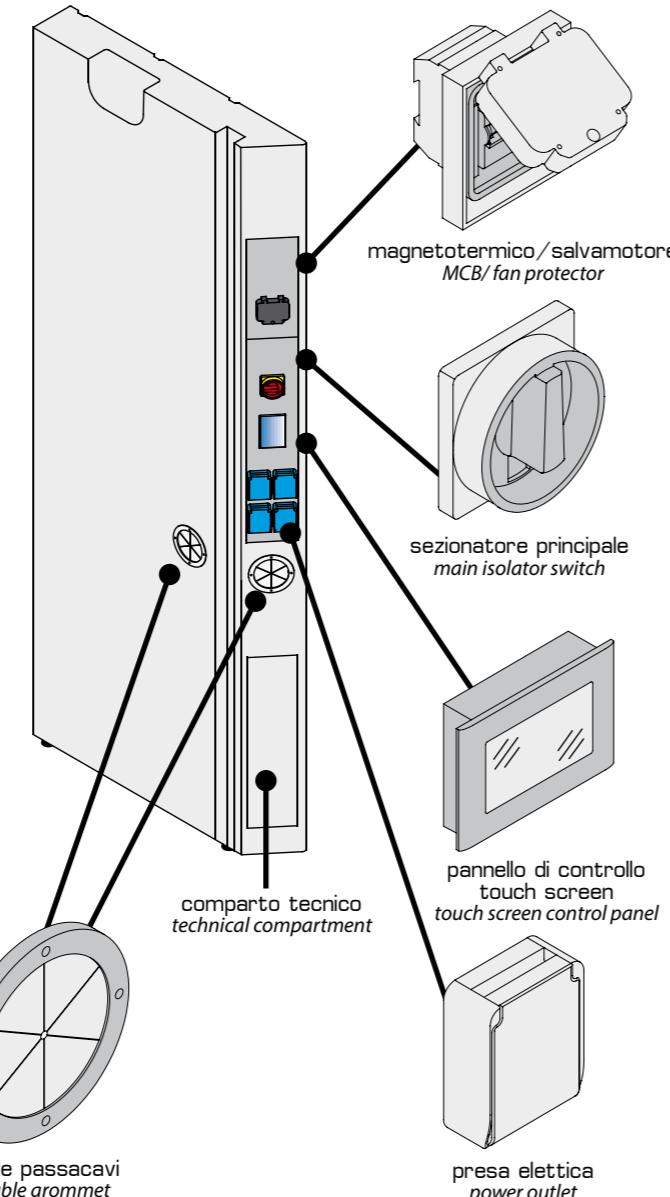
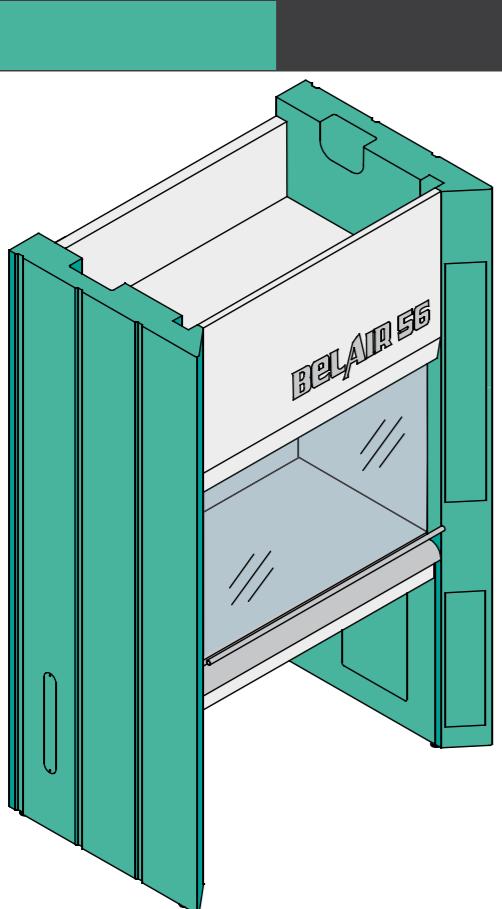
An electric engine can be provided to operate the movement of the frontal sash. The sash can be either operated manually (through dedicated buttons on the touch-screen panel) or automatically, through the operator's presence detection system. In case the operator's hands or any object is placed under the sash, the sensor automatically blocks the sash and raises it again. Even in case of motorized sash, the operator can always operate it manually.

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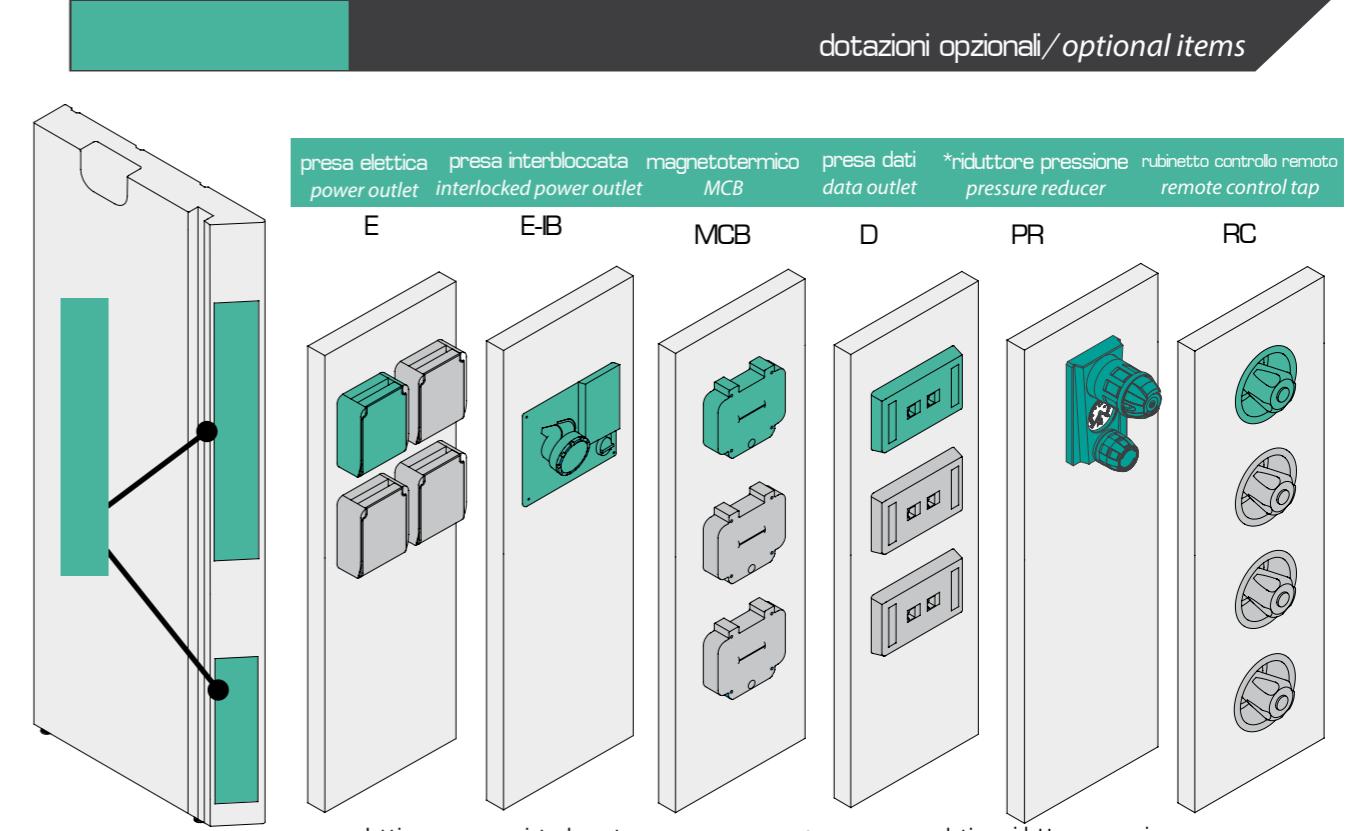
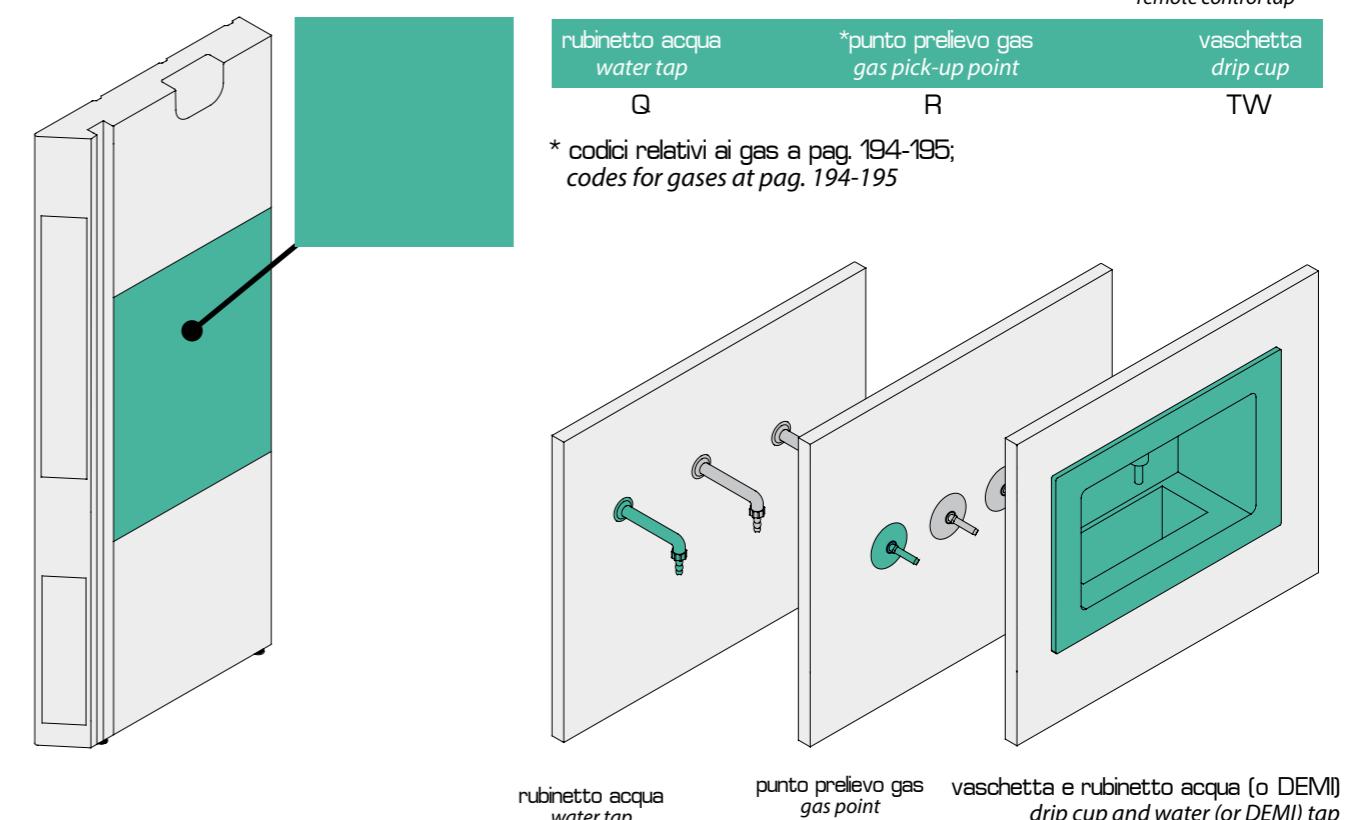
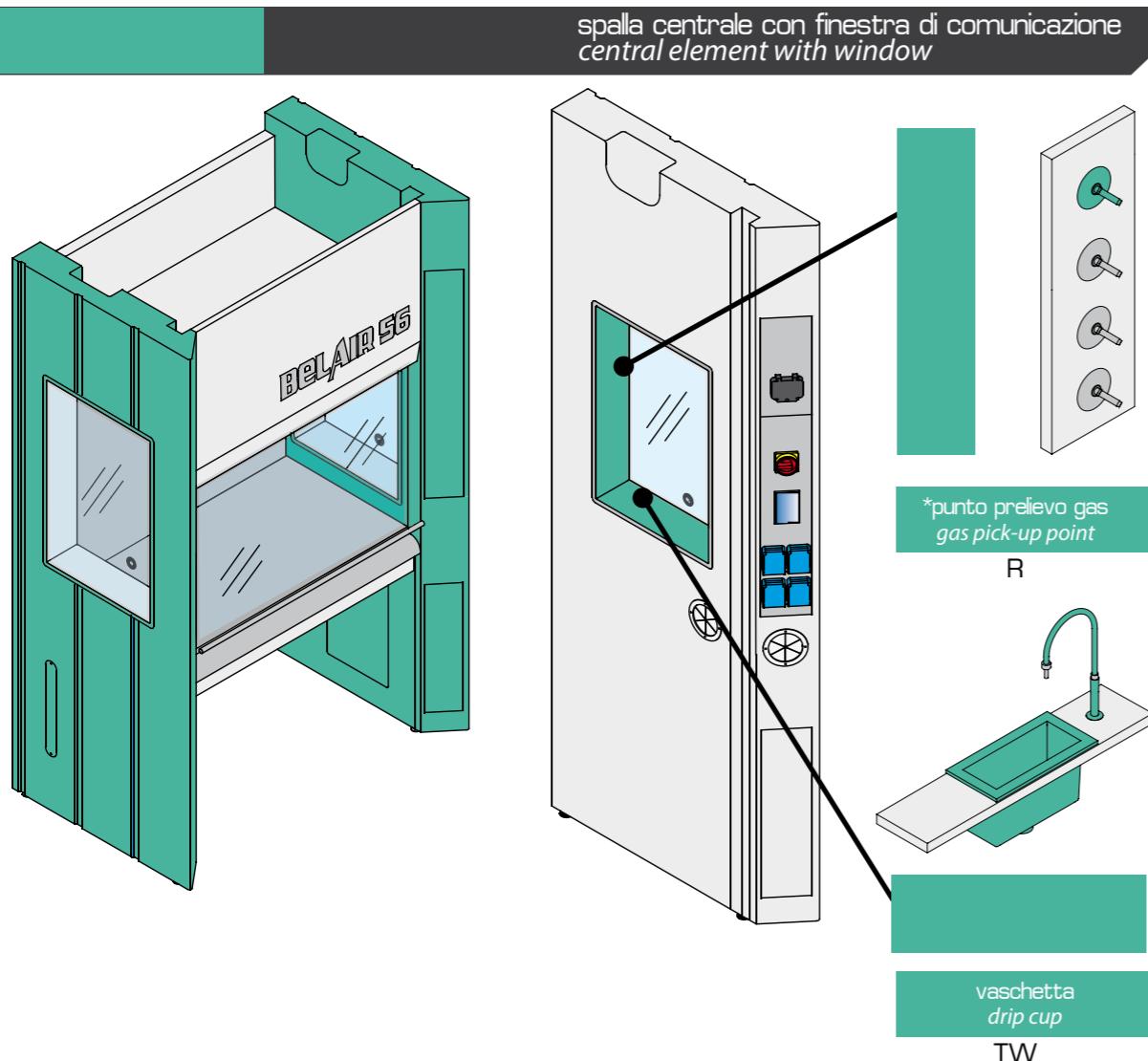
Side jambs of **BELAIR56** fume cupboard can be equipped with many accessories, as better amplified in the following page. In standard configuration the right jamb is equipped with electric services and the left jamb with fluid services. Upon request also left jamb can fit some electric utilities. Standard provisions for right jamb are: main isolator switch, MCB, 4 power outlets and a cable grommet, to be used for supplying power to all instruments placed on the worktop, with no risk of interfering with the frontal sash movement or with the extraction performances. Right jamb also incorporates a technical compartment to house the main control board.

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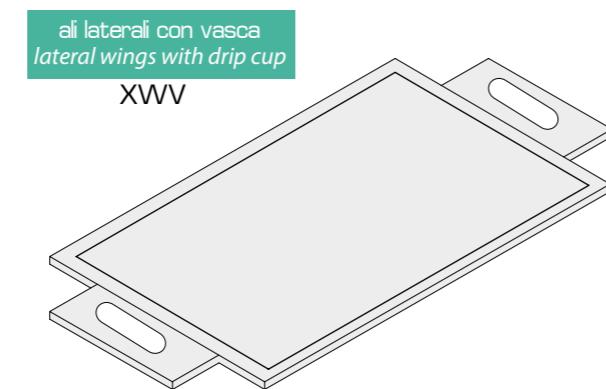
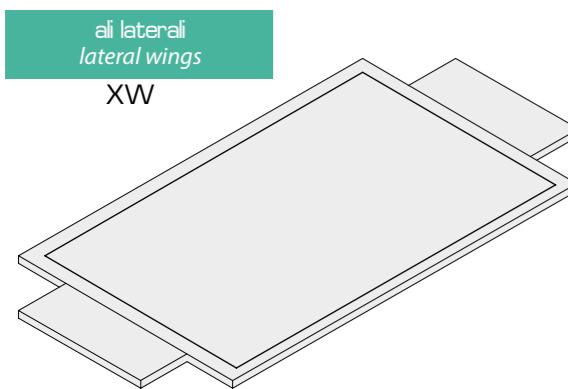
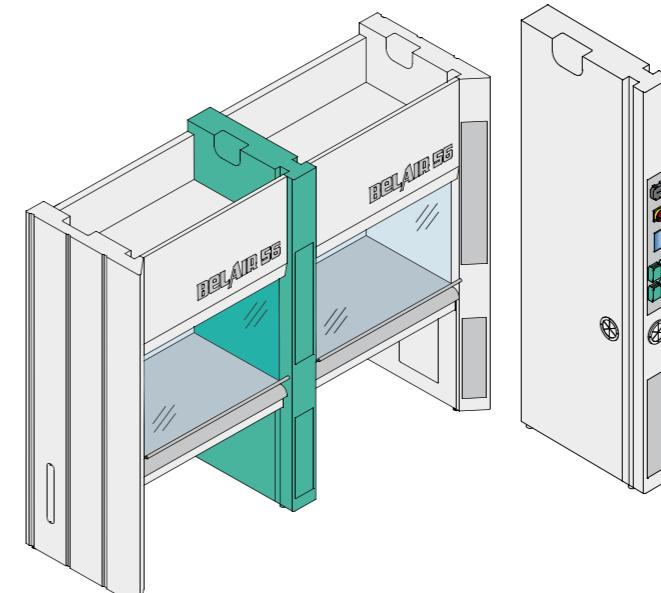
This version of **BELAIR56** can be provided with a worktop nested into the side elements of the fume cupboard, so that its width is equal to the external width of the fume hood.
The worktop can house a lateral drip cup.

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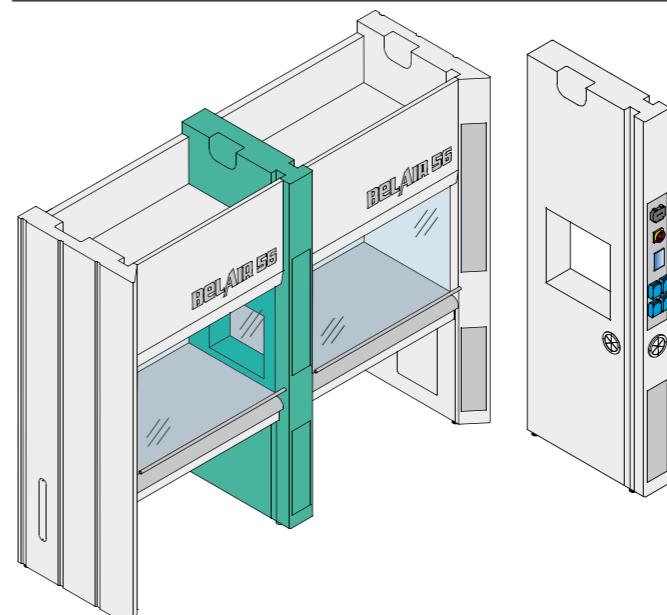
In case two different fume cupboards with inter-communicating working area are needed, a middle common jamb is provided. This solution is preferred either for passing samples between the fume cupboards, or for saving space. In this case the touh-screen control panel can be either common or separated. Standard provisions for the central jamb are: main isolator switch, MCB, touch-screen control panel, 4 power outlet and a cable grommet. Middle jamb can house the same accessories described in previous pages.

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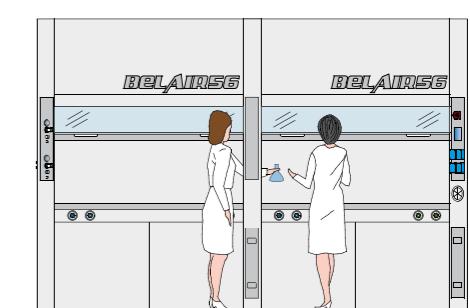
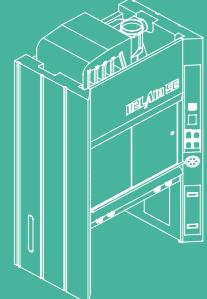
The middle jamb can be equipped with a glass window. In this case, it is not possible to install the dripcup (code TW).

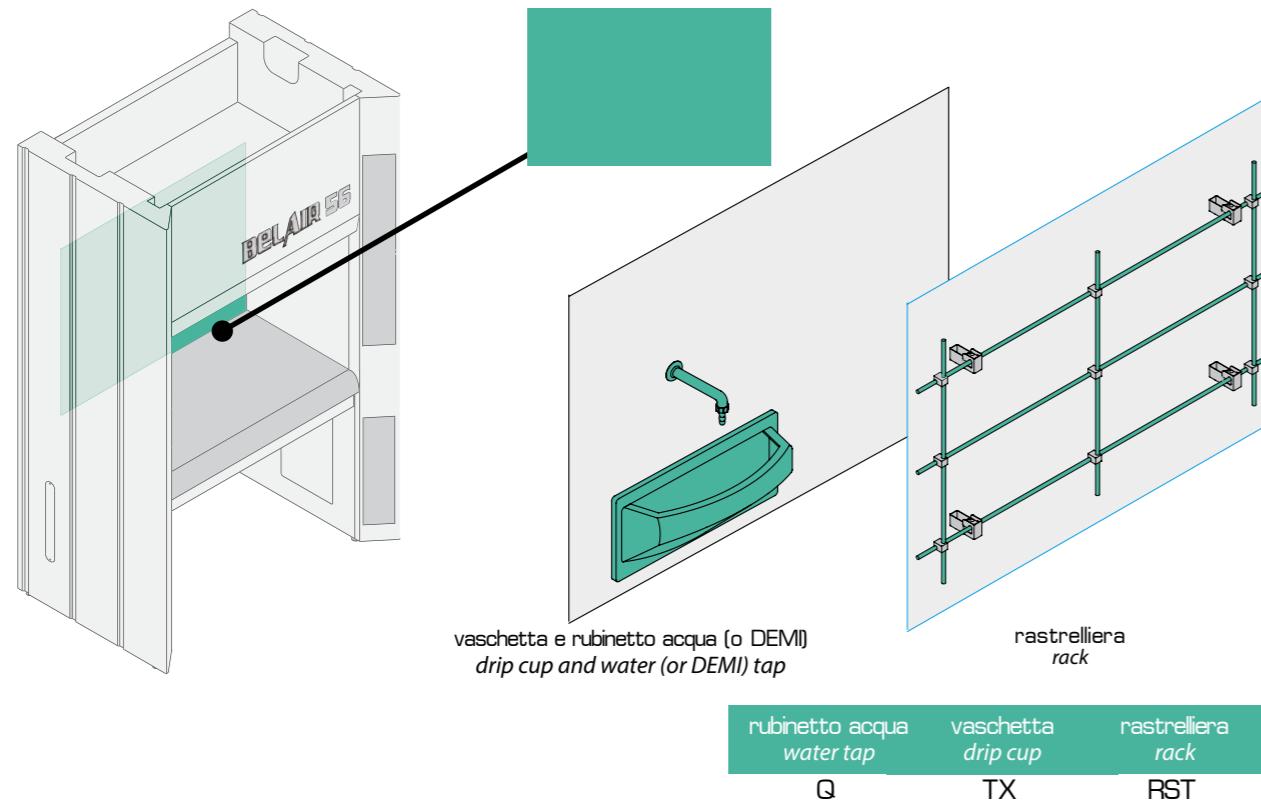
Belair 56

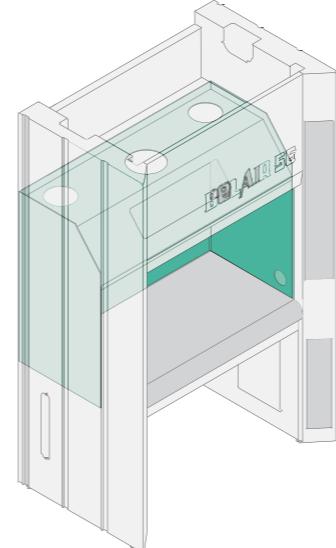
Belair 56

pannello posteriore/rear panel

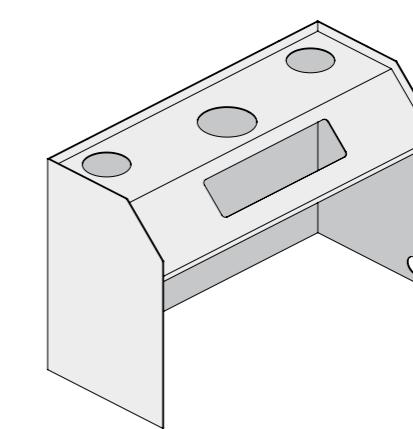
PLN-BELAIR56

rivestimento per alti aggressivi
internal covering for aggressive chemicals

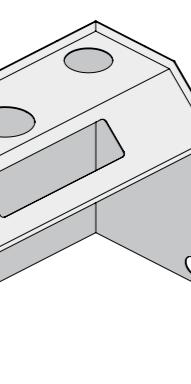
PLN-PP

rivestimento in PP
internal covering in PP

PLN-PVC

rivestimento in PVC
internal covering in PVC

PLN-GR

rivestimento in gres
internal covering in ceramic

PLN-HPL

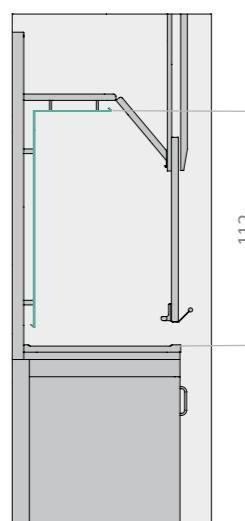
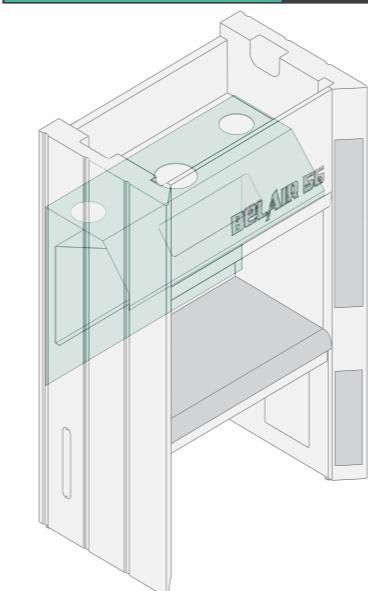
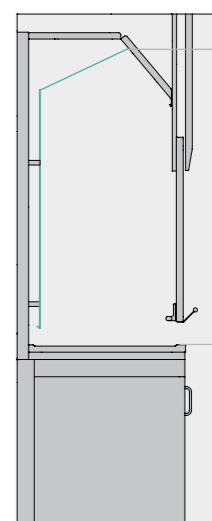
rivestimento in HPL
internal covering in HPL

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In case the fume cupboard has to be employed for manipulation of very aggressive chemicals (such as highly concentrated acids or acids at high temperature), a special internal covering can be provided. The internal covering panels can be made of anti-acid PVC, PP, epoxy resin or of ceramic, and protect the structure of the fume cupboard from corrosion.

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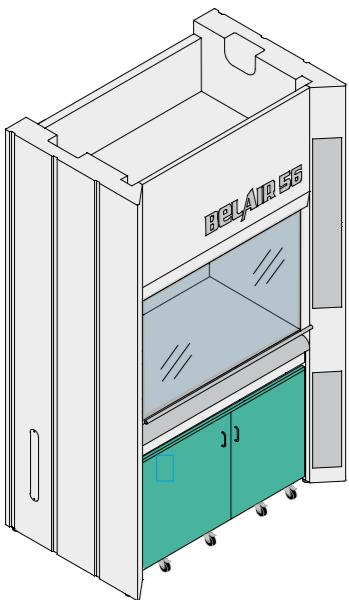
Upon specific request, the inner plenum of the fume cupboard can be provided with a greater height in order to offer further space and enough room for using tall instruments. In this case the fume cupboard can not be fitted with blower and filter box on board but these can only be located in remote position.

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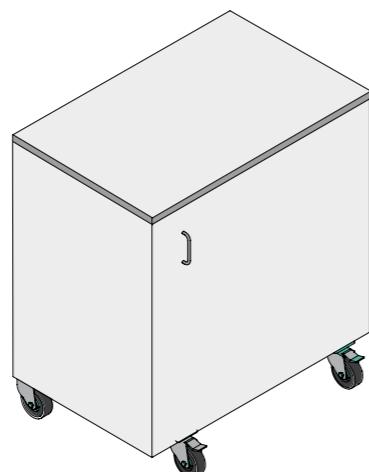
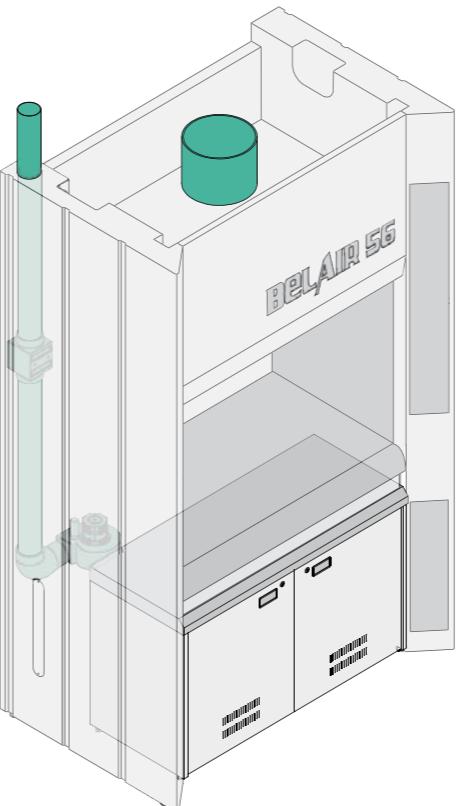
A wide range of storage units can be installed under the fume cupboard, for the user's convenience.

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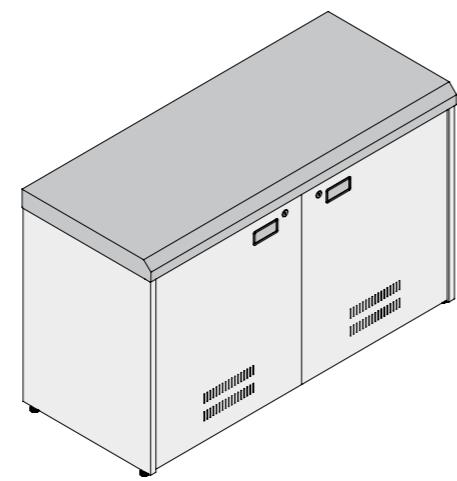
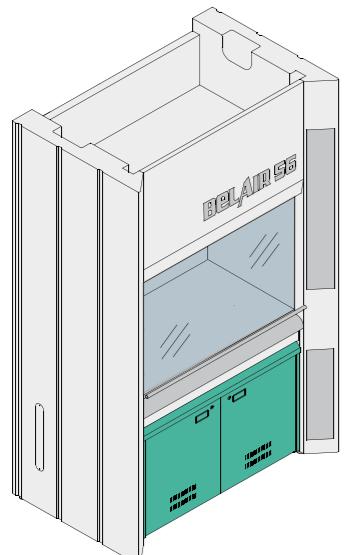
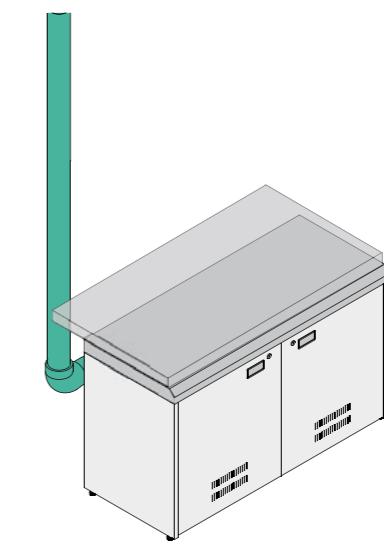
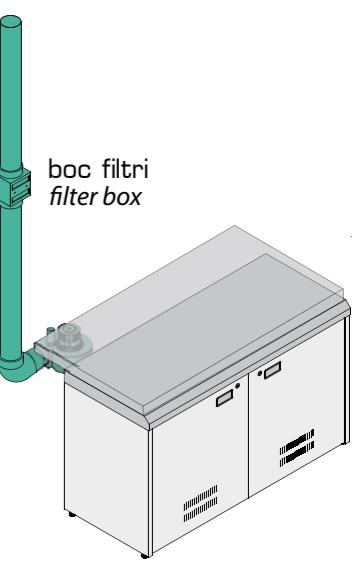
In case of necessity a safety cabinet for chemicals storage can be installed under the fume cupboard. The safety cabinet can also be equipped with extraction system upon request. For details and specifications of safety cabinets please refer to chapter 4 "safety cabinets".

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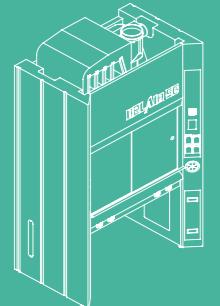
If equipped with VAV (Variable Airflow Volume), fume cupboards embed a controller monitoring their face velocity in real time: once the face velocity has been set, VAV system keeps it constant with the variation of sash position. VAV system can be obtained either using an inverter (cod. VAV-I-BELAIR56) or using an air dumper (cod. VAV-V-BELAIR56). More details at pag. 123

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Upon specific request, the fume cupboard can be equipped with an active temperature monitoring system, with alarms. Both the temperature of the working area and of the under-worktop units can be monitored. The latter is particularly important when safety cabinets or waste tanks for flammable liquids are positioned under the worktop.

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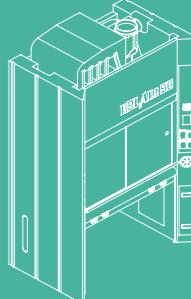
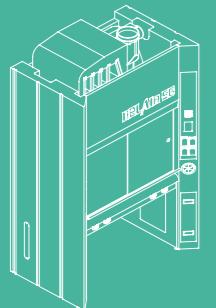
Upon specific request, the fume cupboard can be equipped with an automatic system for fire detection and suppression. The development of a fire inside the working area of the fume cupboard determines the sudden activation of the system (through a pressure rise in the probe duct) and the subsequent release of extinguishing agent through specific nozzles. The extinguishing agent is stored in a dedicated cylinder located outside of the fumehood and can be chosen according to the specific need (standard CO₂).

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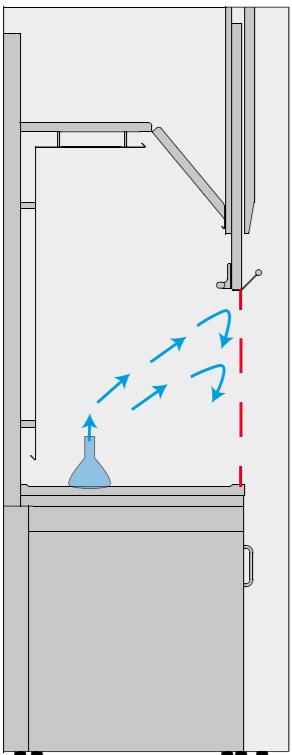
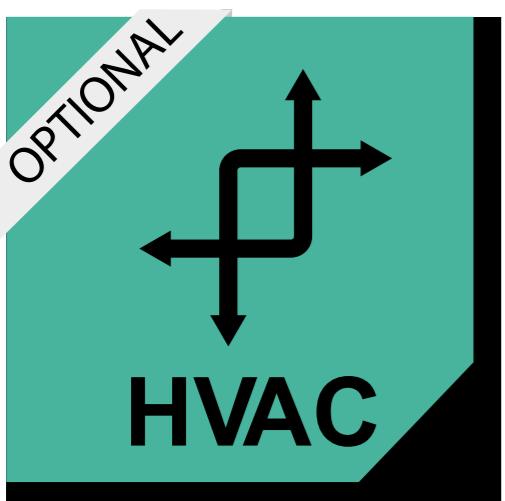
TWIN SET is an exclusive device able to integrate energy savings to operator's safeguard.

If TWIN SET system is installed, in case of necessity the operator can decrease the face velocity, in order to save energy and to reduce the total airflow extracted from the room.

In this case, to keep the containment performances unvaried, one or more additional blowers automatically activate, so that an air barrier is created.

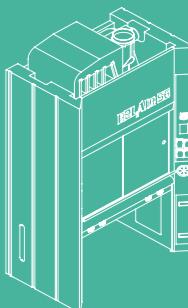
This barrier allows to completely shield the working environment under the fume cupboard, so that airborne contaminants are retained inside.

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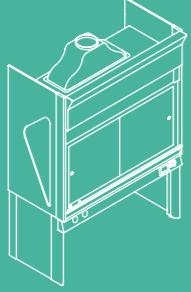
AK is an automatic chemical fume cupboard designed and tested to ensure full protection to lab operators in every situation, even the most challenging and advanced.

Thanks to its design, **AK** finds a place in every modern and innovative laboratory environment adding value and efficiency.

AK is characterized by a light appearance, thanks to a careful location of the utilities below the worktop, allowing a full employment of the available space, optimizing the efficiency and the overall usability of the fume cupboard.

Both the lateral jambs of **AK** are realized using anti-acid HPL sheets, providing at the same time stiffness and lightness to its structure, making it different from the conventional profiles available on market for chemical fume cupboards.

In case of fume hood for aggressive acids, its structure is entirely realized in PVC or PP.

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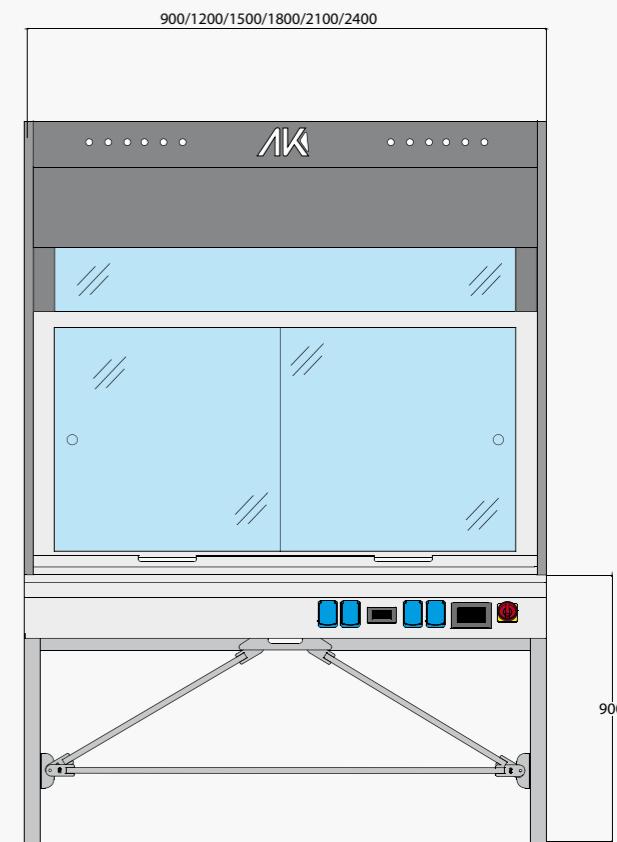
quote espresse in mm
dimensions expressed in mm

The lateral windows, standard fitting for all models, are realized in safety glass and contribute to make the structure of **AK** very light and sustainable, providing a smooth and futuristic design, together with an optimal illumination of the internal working area.

AK fume cupboard has been certified by a third party for its compliance to the Norm EN 14175:2003, and the certificate has been issued by Internationally recognized third authorities, so to ensure the efficiency and safety Standards are always met.

In order to obtain the EN 14175 Certification, **AK** has been tested in the most adverse and challenging conditions to grant full safety of the operators in all applications, proving superior features of containment and robustness of containment.

Together with the wide choice of materials for worktop, many different accessories and optional items are available: motorized sash, many automation optionals or employment of explosion proof components are only some examples.

AK fume cupboard is equipped as a standard with a highly technologic electronic control software, specifically designed for its management, and usable by the operator through a very intuitive touch screen interface. The control software embeds a wide range of alarms and detecting probes, aimed to continuously monitor the working conditions of the fume cupboard, and ready to promptly intervene in case of necessity, so that operators are always safeguarded.

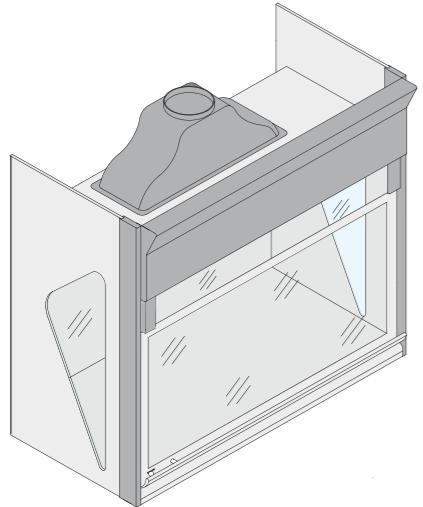
In addition to the standard software, fume cupboard can also embed some advanced airflow control features, designed to implement active safety and energy saving logics, such as the VAV, CLIMA and the TWIN SET systems.

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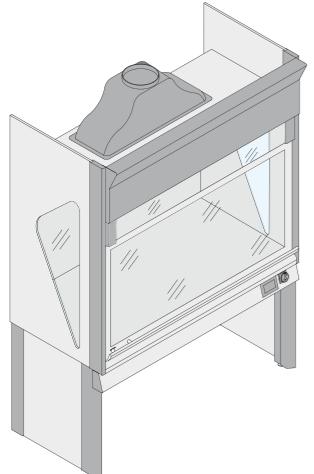
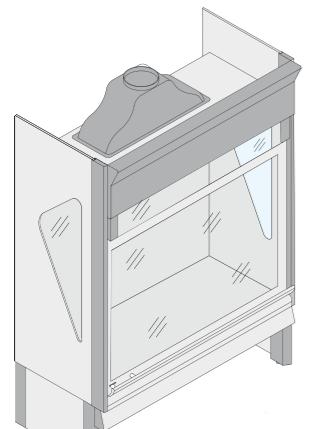
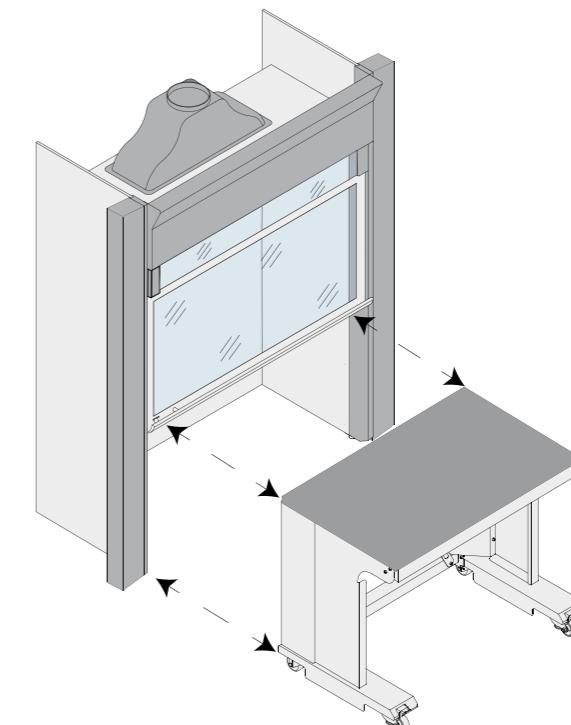
Equipments of **AK** walk-in can be positioned on side columns. Details at pag.162

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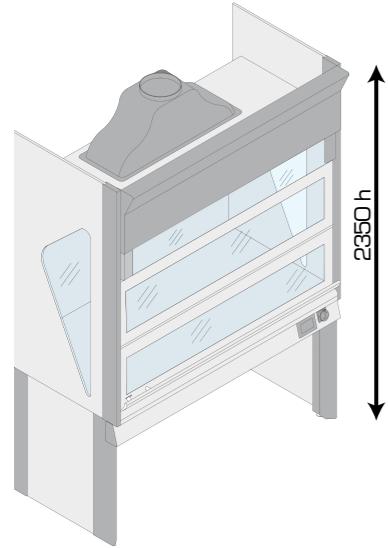
Walk-in version is not equipped with the conventional worktop to allow the insertion under the fume cupboard of tall equipment and trolley benches. Walk-in fume cupboards are often coupled to **Skylab** line trolley benches (code: BCAR), described in lab benches technical catalogue.

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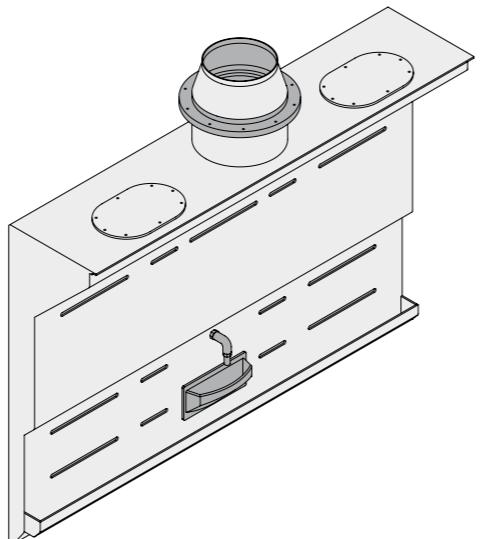
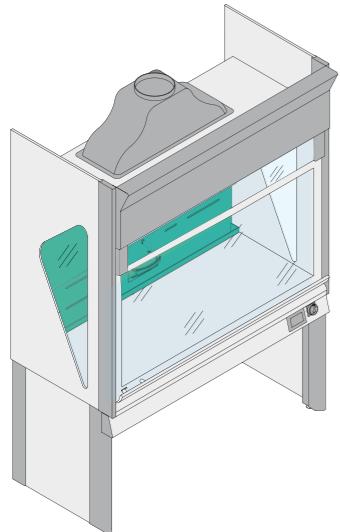
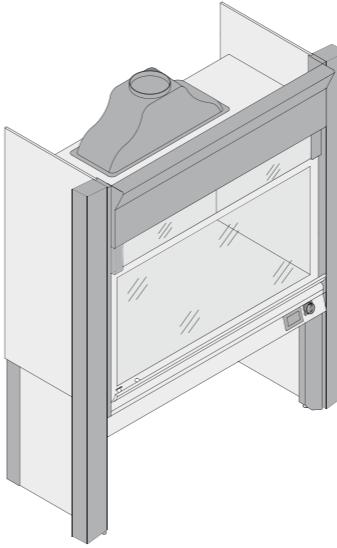
Upon request it is possible to equip fume cupboards with a fume washing system embedded in the plenum, in the back of the hood.

This system consist in a device spraying water in counter-current to the exhaust air, which is sent through some swirlers to increase the interface surface. Water can be collected into a tank and recirculated or directly properly disposed.

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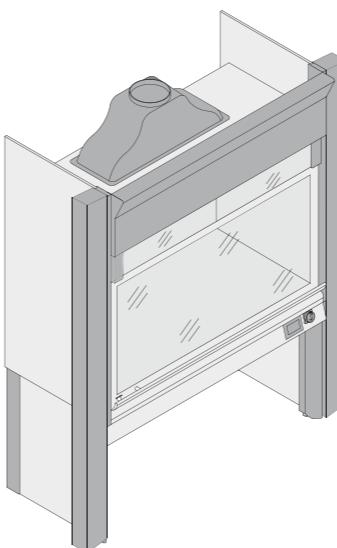
The upper part of this hood is chemical-resistant and entirely realized in PVC. It is suitable for application where highly acidic chemicals are needed. Sash sliding mechanism and his anti-fall safety device are positioned outside the working chamber in order to be protected from acid attacks.

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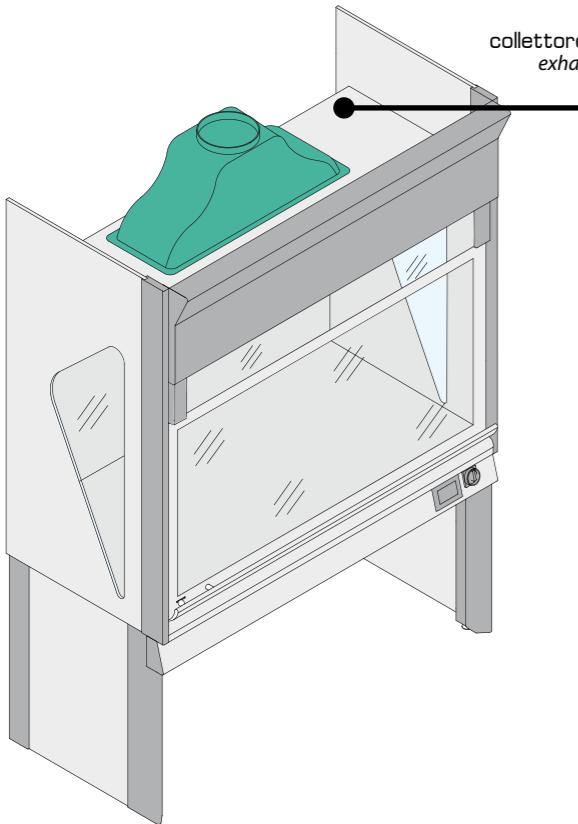
The structure for use of highly acidic and thermal load differs from the structure for use of highly acidic for the presence of a ceramic lining inside the chamber.

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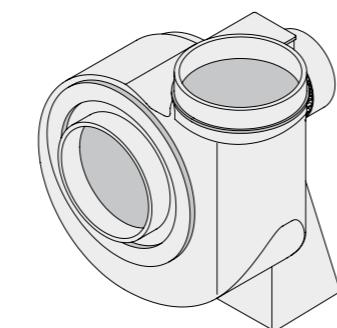
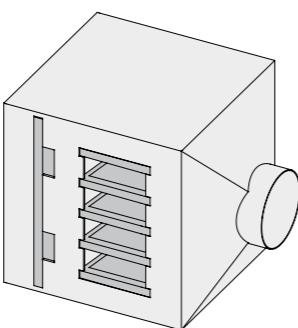
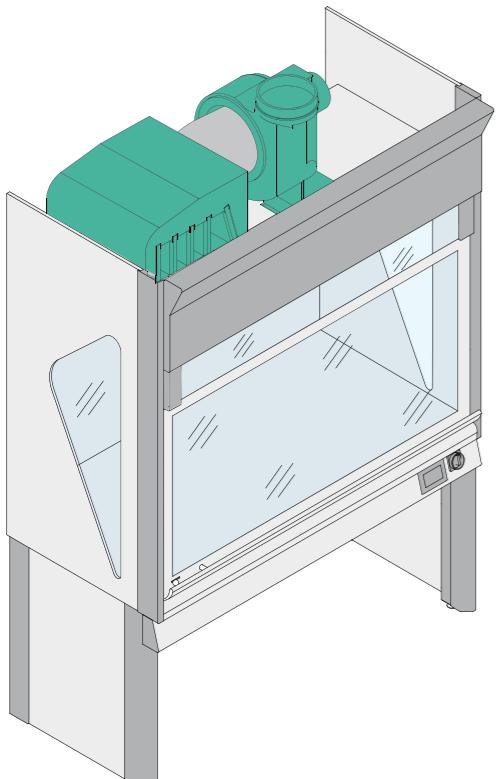
A-K is equipped, as a standard, with an exhaust collector maximising its containment performances.

If a filter box or an exhaust fan are provided, these are installed either on ceiling or on the roof.

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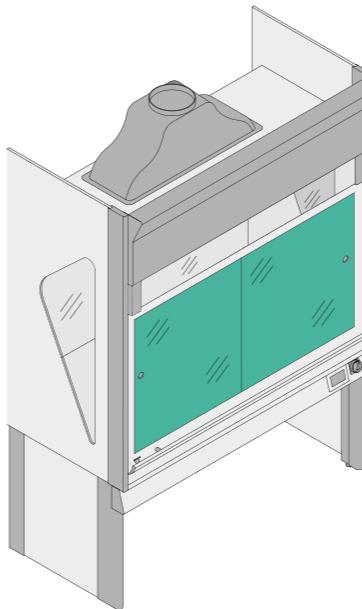
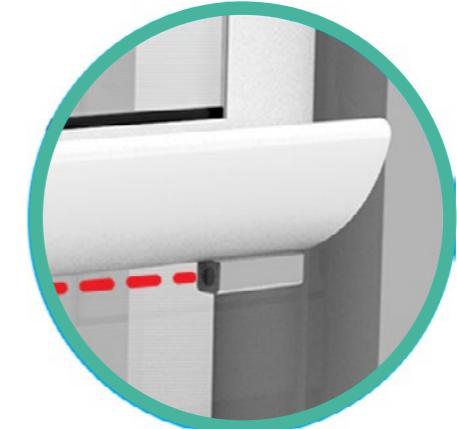
A-K is always equipped with a vertically sliding frontal safety glass, framed into a counter-balanced sash.

Upon request, it is possible to provide a vertically sliding frame incorporating up to 4 glass panels, horizontally sliding.

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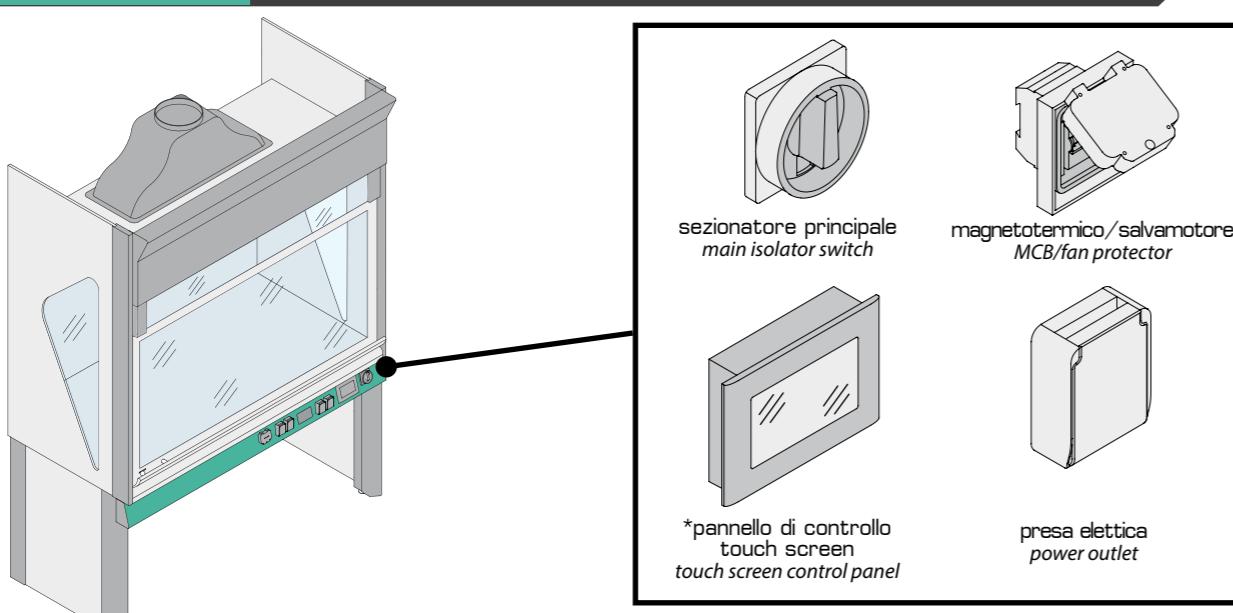
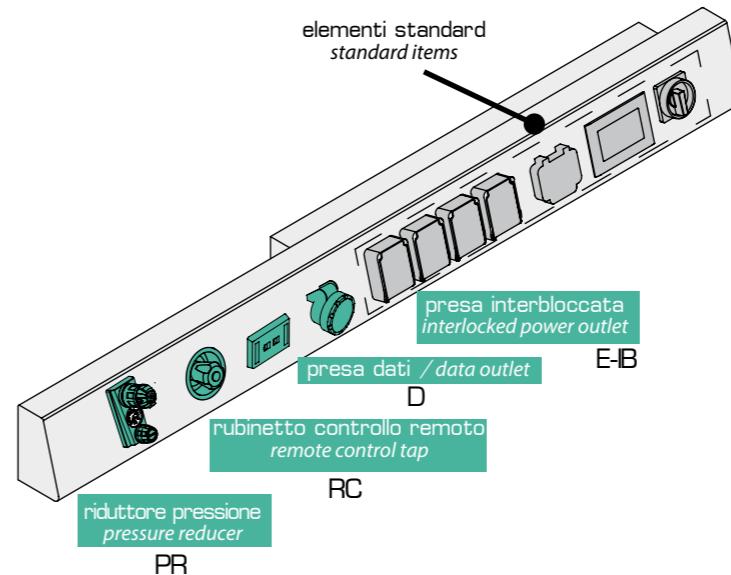
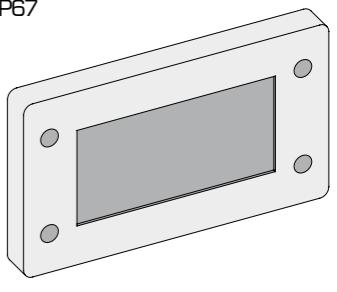
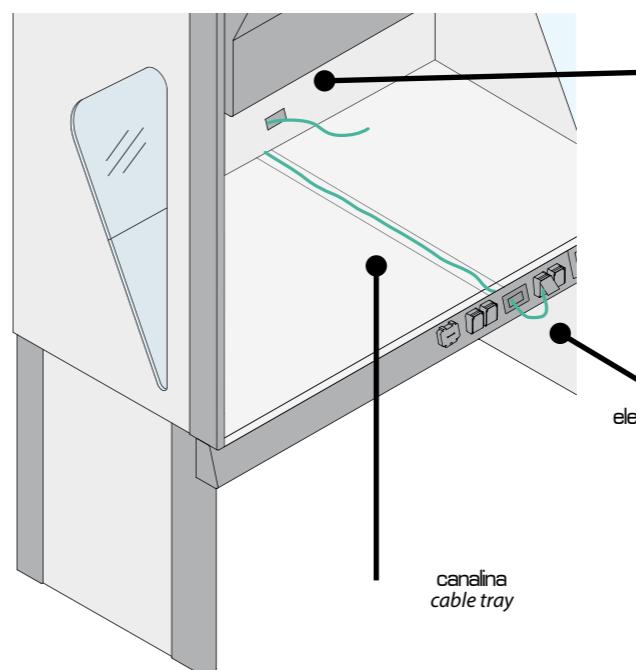
The frontal panel of AK fume cupboard can be equipped with many accessories, as shown above.
In standard configuration, basic provisions are: master control switch for turning ON and OFF the fume cupboard, MCB, 4 power outlets and a cable grommet, to be used for supplying power to all instruments placed on the worktop, with no risk of interfering with the frontal sash movement or with the extraction performances. Right side element also incorporates a technical compartment under the worktop area.

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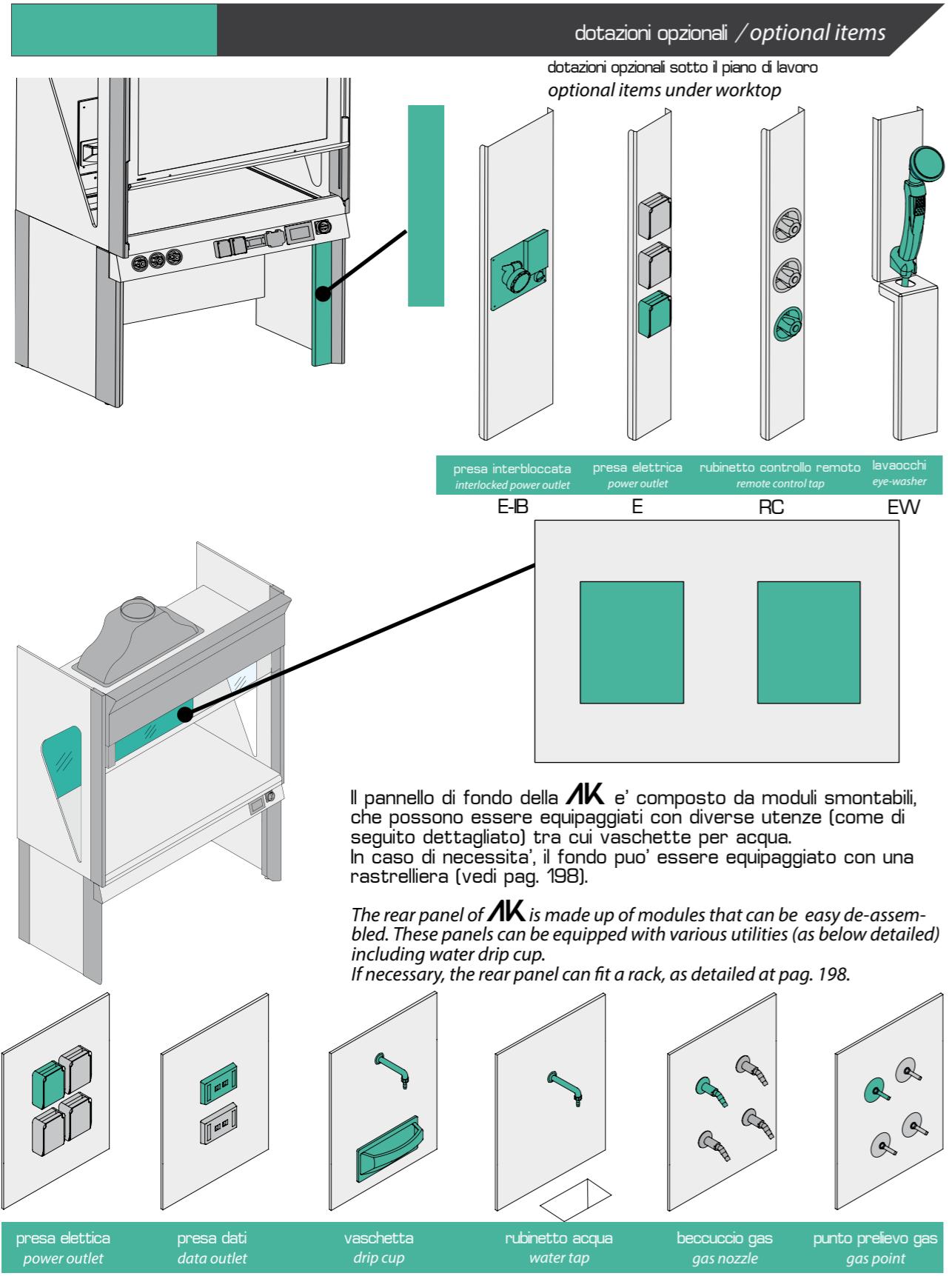
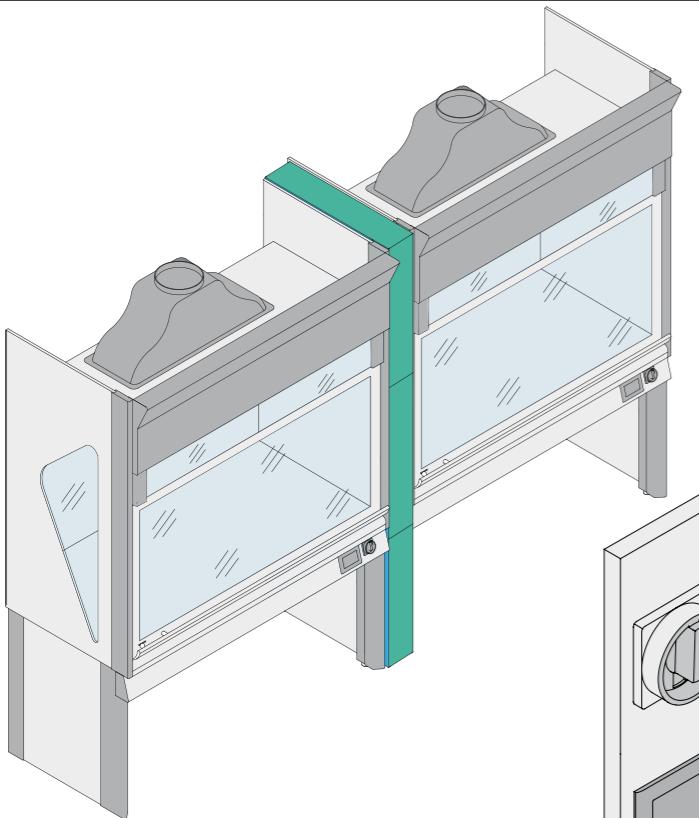
The rear panel of **AK** is made up of modules that can be easy de-assembled. These panels can be equipped with various utilities (as below detailed) including water drip cup.
If necessary, the rear panel can fit a rack, as detailed at pag. 198.

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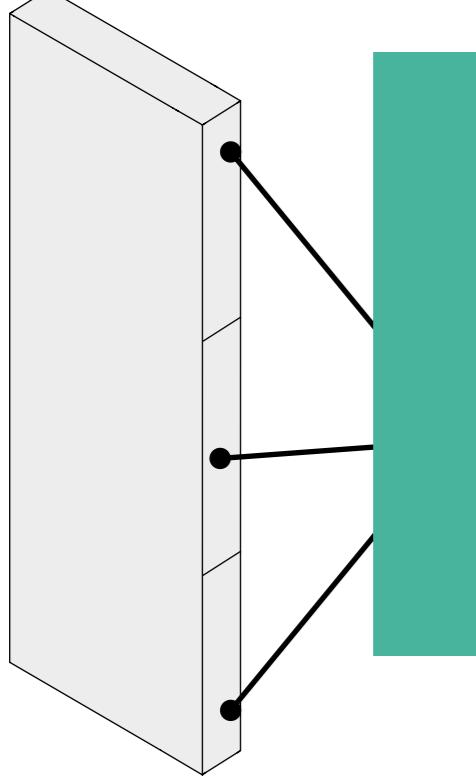
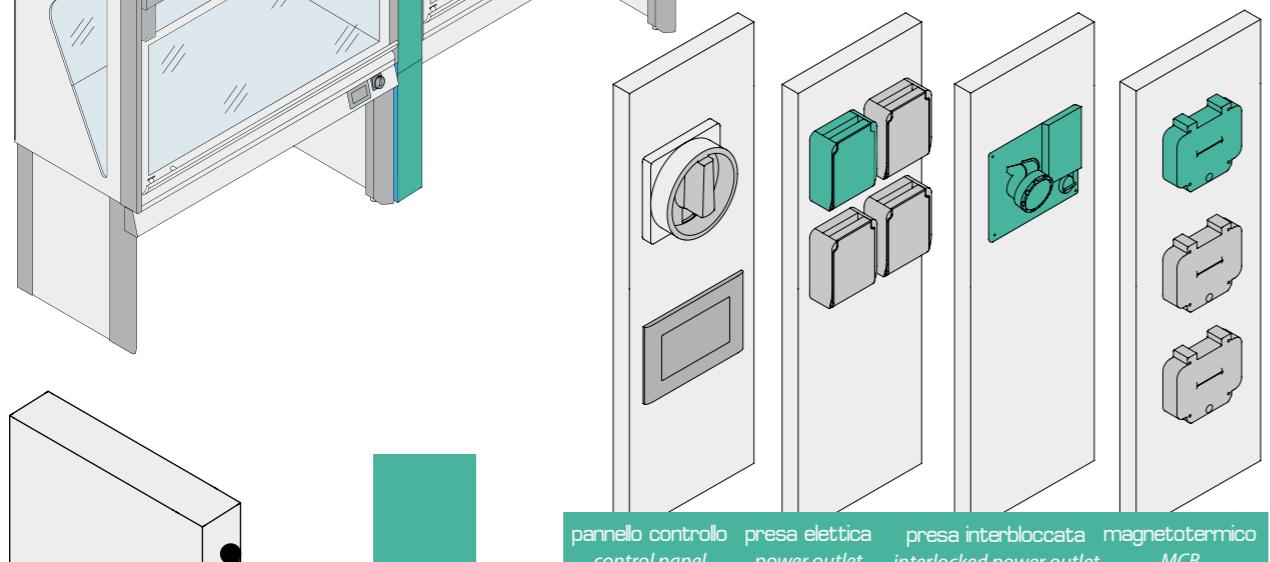
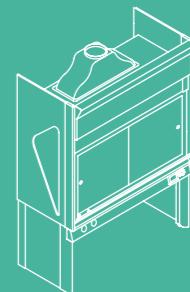
In case two different fume cupboard, with inter-communicating working area are needed, a central supporting element is provided. Middle element can incorporate up to 3 customizable panels.

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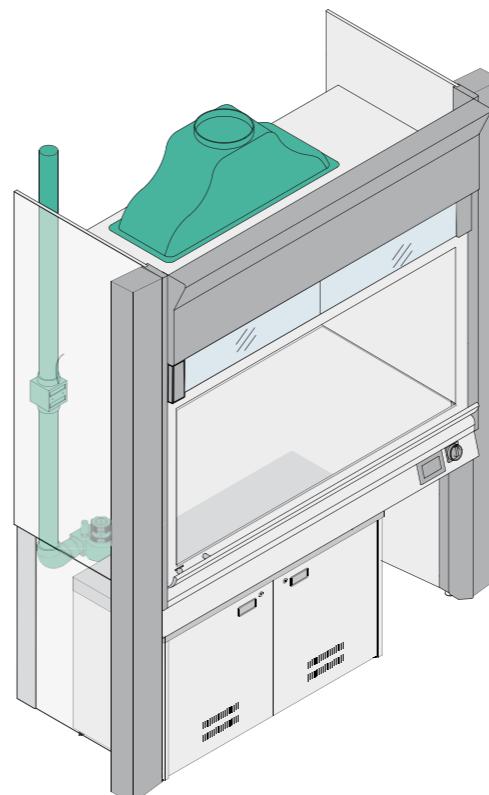
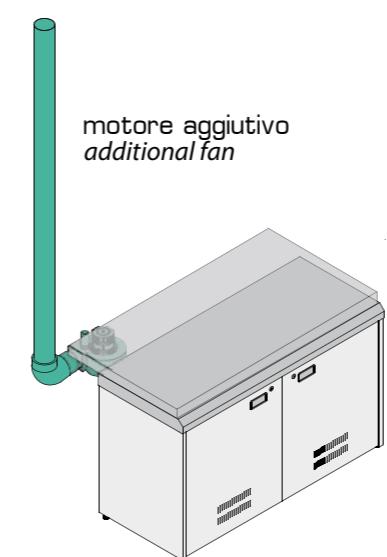
All safety cabinets installed under the worksurface can be connected to the exhaust system. Exhaust can be either common to the fume cupboard or independent. Upon specific request, also a dedicated filter embedded in the structure of the fume cupboard can be provided for safety cabinets.

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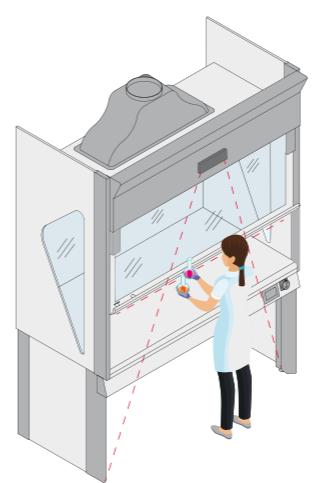
If equipped with VAV (Variable Airflow Volume) fume cupboards embed a controller monitoring their face velocity in real time: once the face velocity has been set, VAV system keeps it constant with the variation of sash position. VAV system can be obtained either using an inverter (cod. VAV-I-AK) or using an air dumper (cod. VAV-V-AK). More details at pag. 123

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Upon specific request, the fume cupboard can be equipped with an automatic system for fire detection and suppression. The development of a fire inside the working area of the fume cupboard determines the sudden activation of the system (through a pressure rise in the probe duct) and the subsequent release of extinguishing agent through specific nozzles. The extinguishing agent is stored in a dedicated cylinder located outside of the fumehood and can be chosen according to the specific need (standard CO₂)

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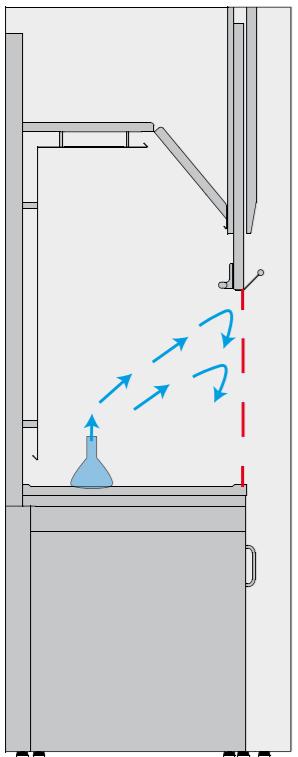
An electric engine can be provided to operate the movement of the frontal sash. The sash can be either operating manually (through dedicated buttons on the touch-screen panel) or automatically, through the operator presence detection system. In case the operator's hands or any object is placed under the sash, the sensor automatically blocks the sash and raises it again. Even in case of motorized sash, the operator can always operate it manually.

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ecoAIR is a manual fume cupboard, representing the optimal synthesis among efficiency, functionality and ease of use, at a reduced cost.

ecoAIRfume cupboard, though not equipped with a software for electronic control of extraction airflows, grants full safety to laboratory operators thanks to an aerodynamics planned and studied in details, granting superior containment.

ecoAIRfume cupboard is indeed designed and manufactured in full compliance to EN 14175:2003 Norm, as a guarantee of its efficiency

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ecoAIR fume cupboard is entirely realized using epoxy-powder coated steel sheets, laser-cut and bended through cold processes.

Steel sheets ensure a completely incombustible structure, so that **ecoAIR** is rated as 'Class O' fire reaction.

Furthermore, epoxy-powder coating grants full protection from the most aggressive chemicals and from corrosion, making **ecoAIR** absolutely suitable for installation in every kind of laboratory.

The exclusive employment of steel sheets makes also **ecoAIR** a flagship product on the market, since common fume cupboards are often manufactured with wooden or plastic parts, offering much lower performances when it comes to durability and resistance to chemicals or fire attacks.

Many different configurations are available for **ecoAIR** fume cupboard, so that every exigency or special necessity of customer can be fulfilled.

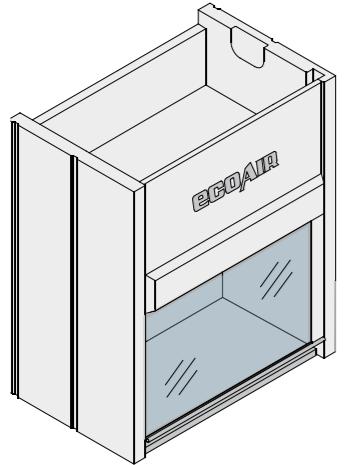
Together with the wide choice of materials for worktop, many different accessories and optional items are available, as better amplified in this Paragraph.

All services and utility points of the fume cupboard are placed in specific points, as a result of ergonomics studies, with the aim of making the use of the cupboard as pleasant as possible for the operator.

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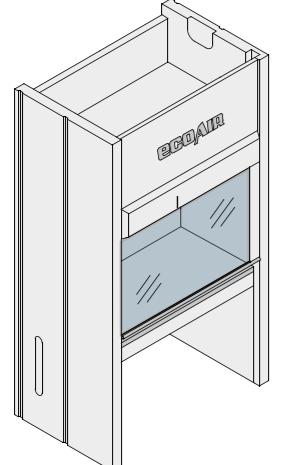
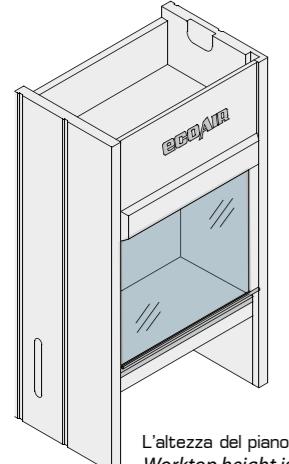
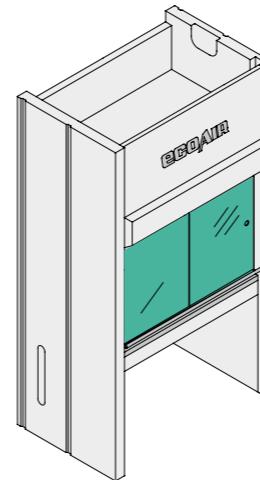
ECOAIR is always equipped with a vertically sliding frontal safety glass, framed into a counter-balanced sash. Upon request, it is possible to provide a vertically sliding frame incorporating up to 4 glass panels, horizontally sliding. Frontal sash is always equipped with a safety device granting its lock in case of failure of one or both its suspending chords, so that it is prevented from falling.

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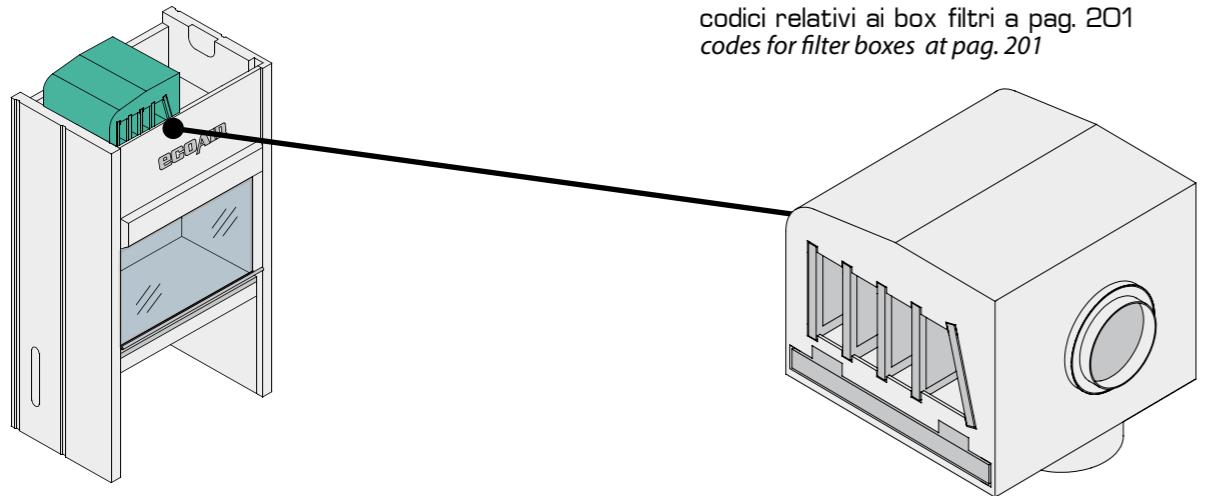
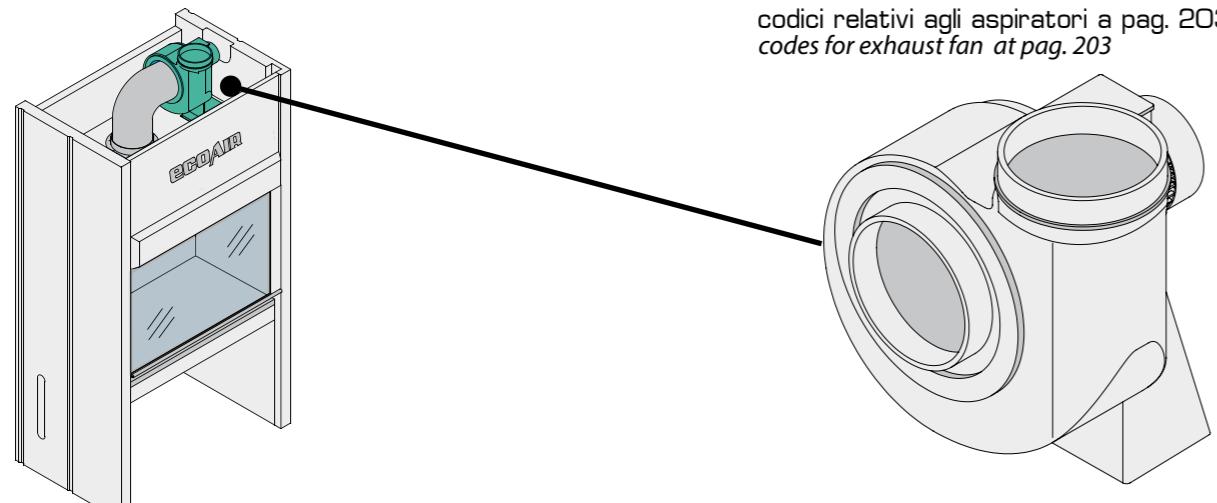
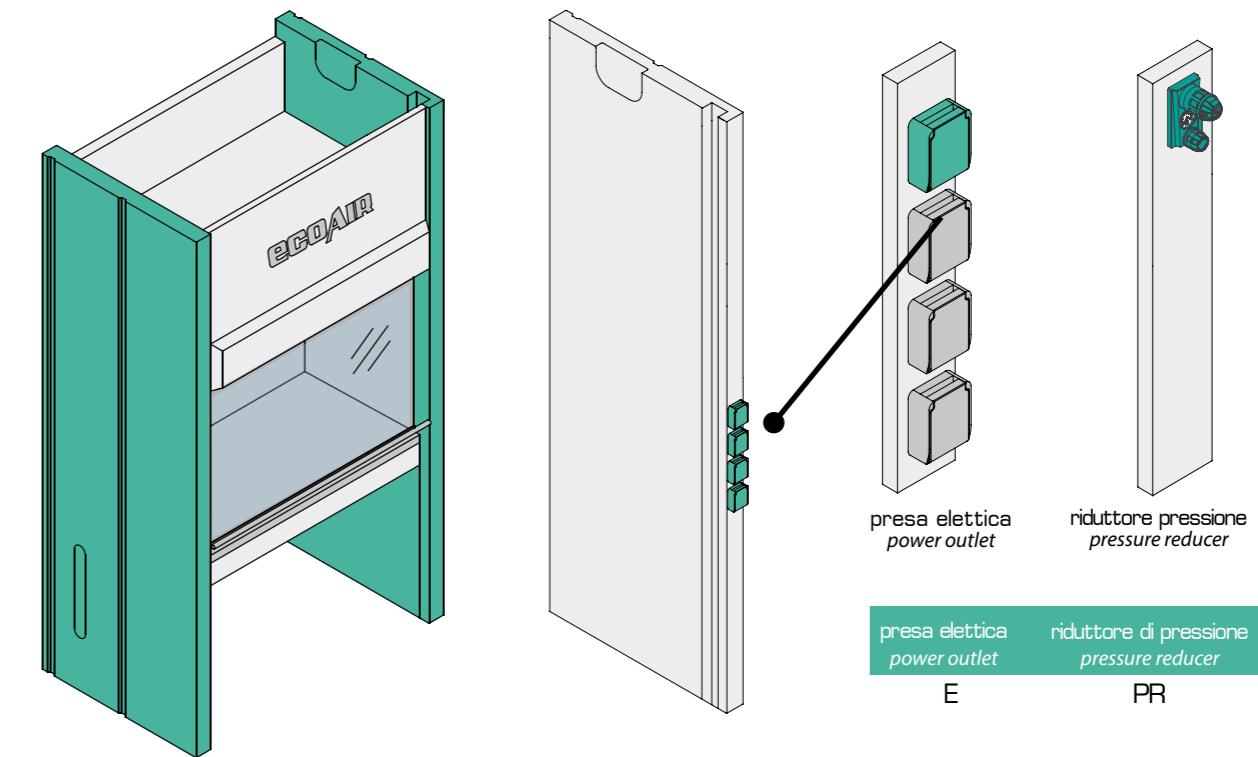
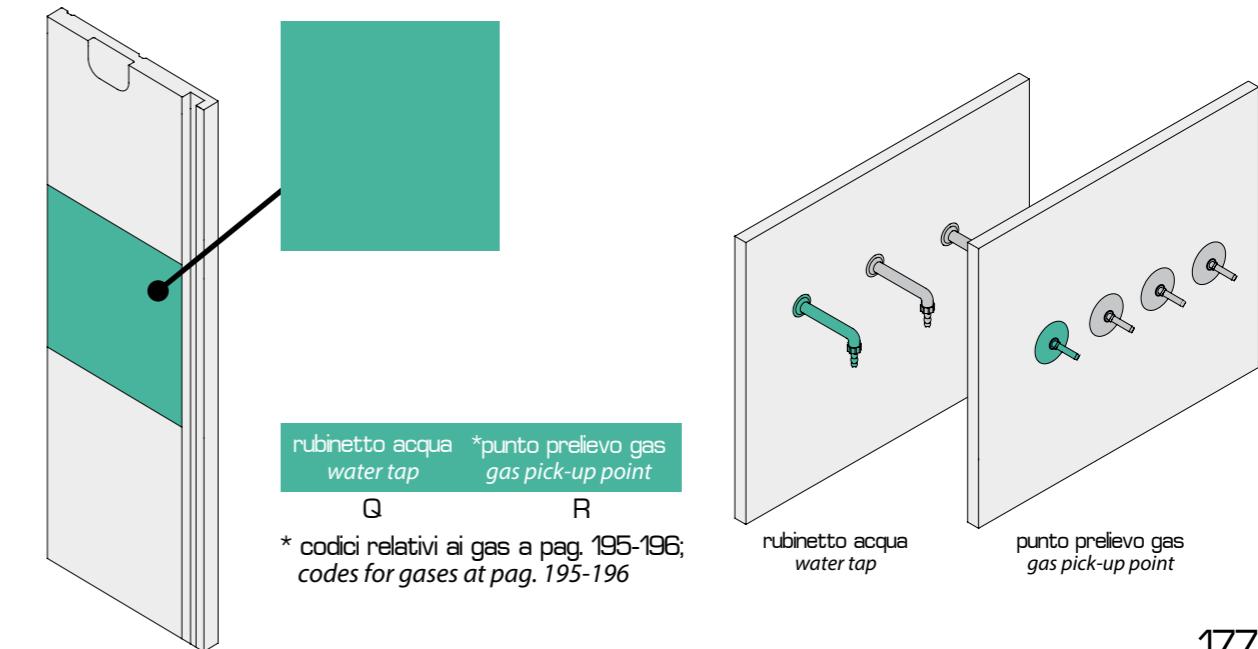
Side jams of **ECOAIR** fume cupboard can be equipped with many accessories. In standard configuration the right side is equipped with electric services and the left side with fluid services, but upon request also left side element can fit some electric services. Standard provisions for right side element are 4 power outlets.

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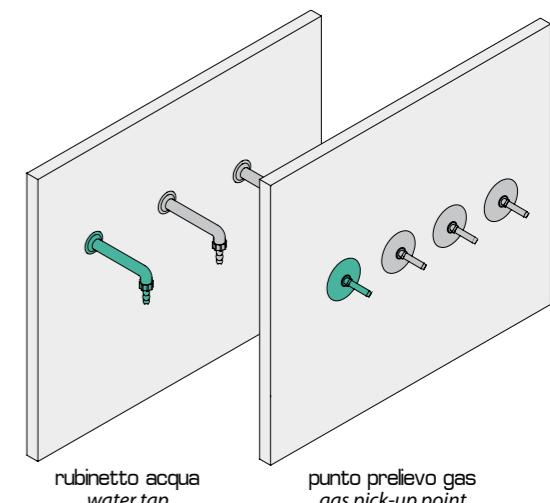
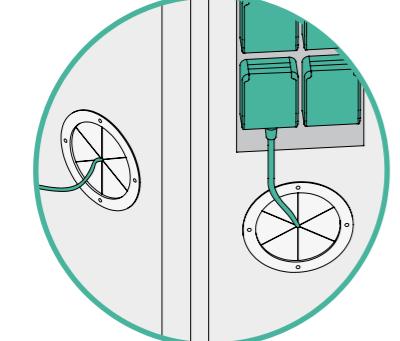
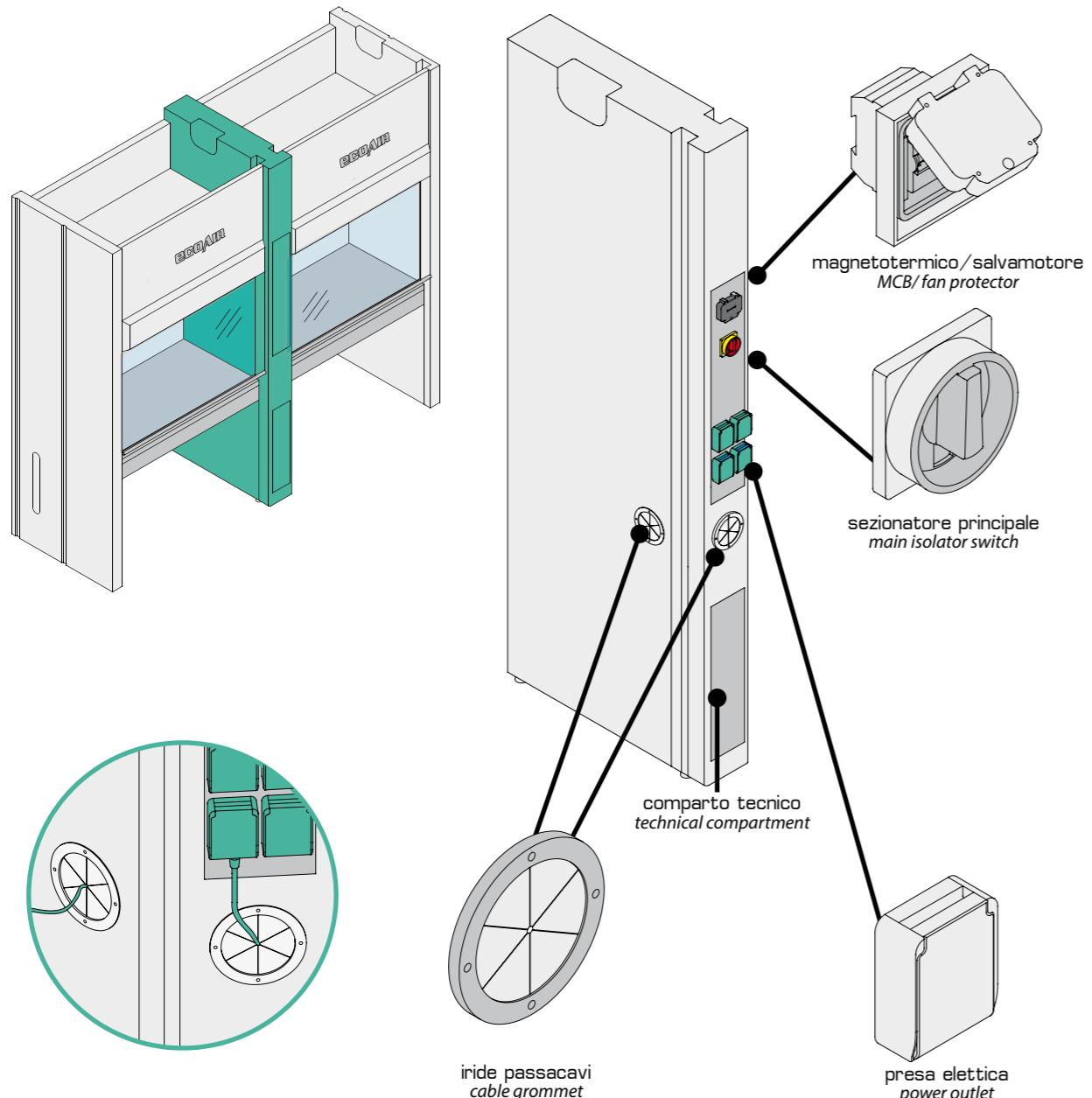
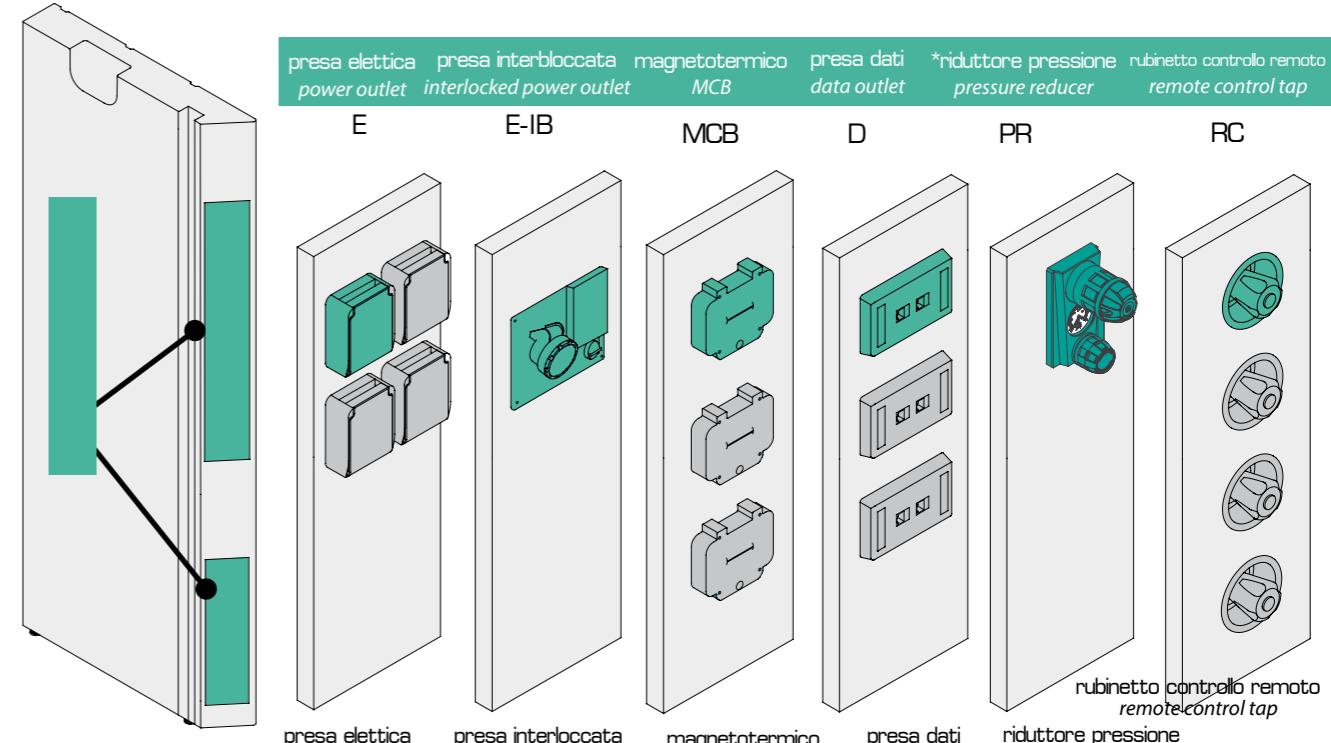
In case two different fume cupboards, with inter-communicating working area are needed, a common middle jamb is provided. This solution is preferred if samples have to be passed between the fume cupboards. Standard provisions for the central elements are: master control switch, MCB, 4 power outlet and a cable grommet. Furthermore the element can be provided with the accessories presented in the following page.

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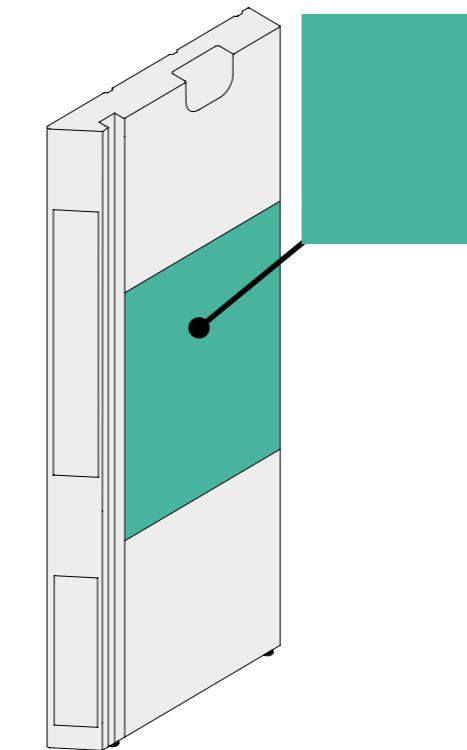
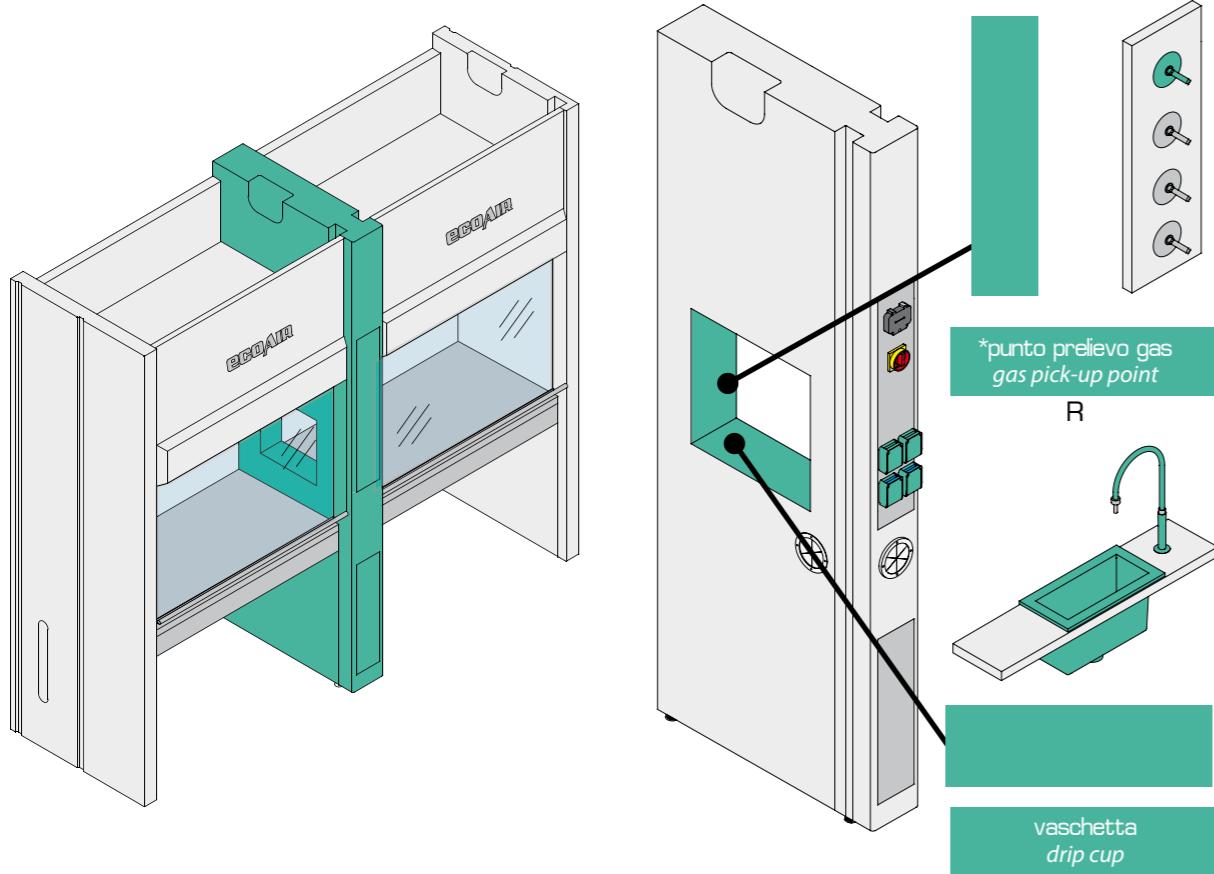
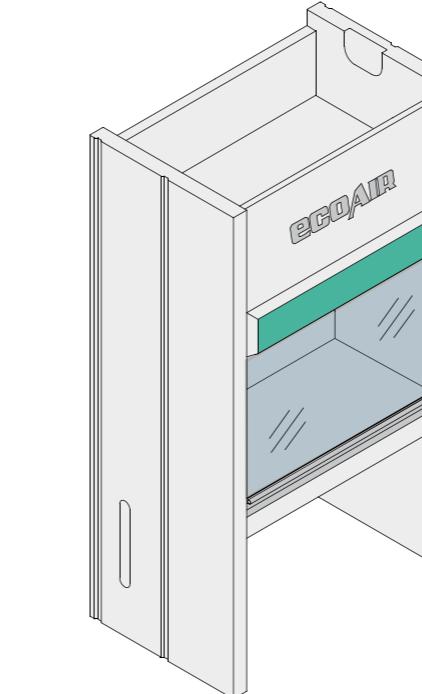
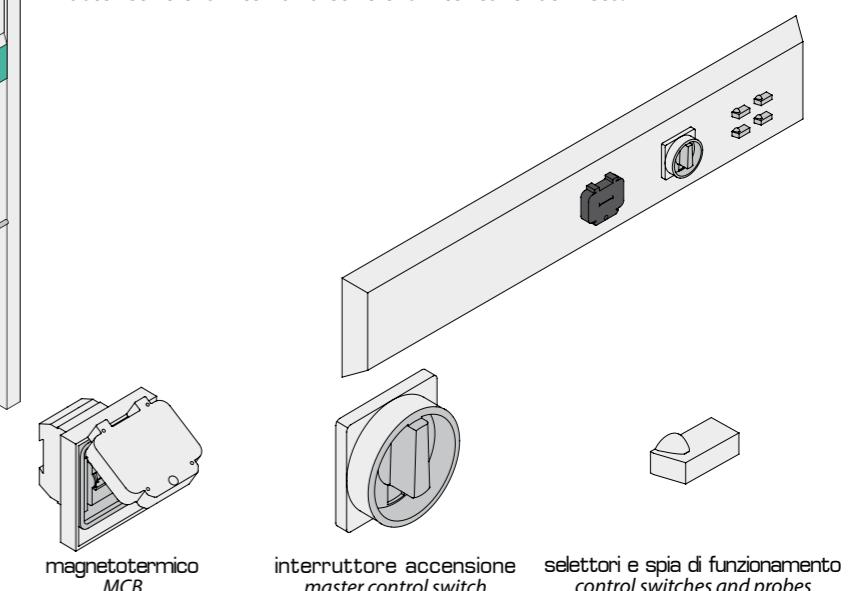
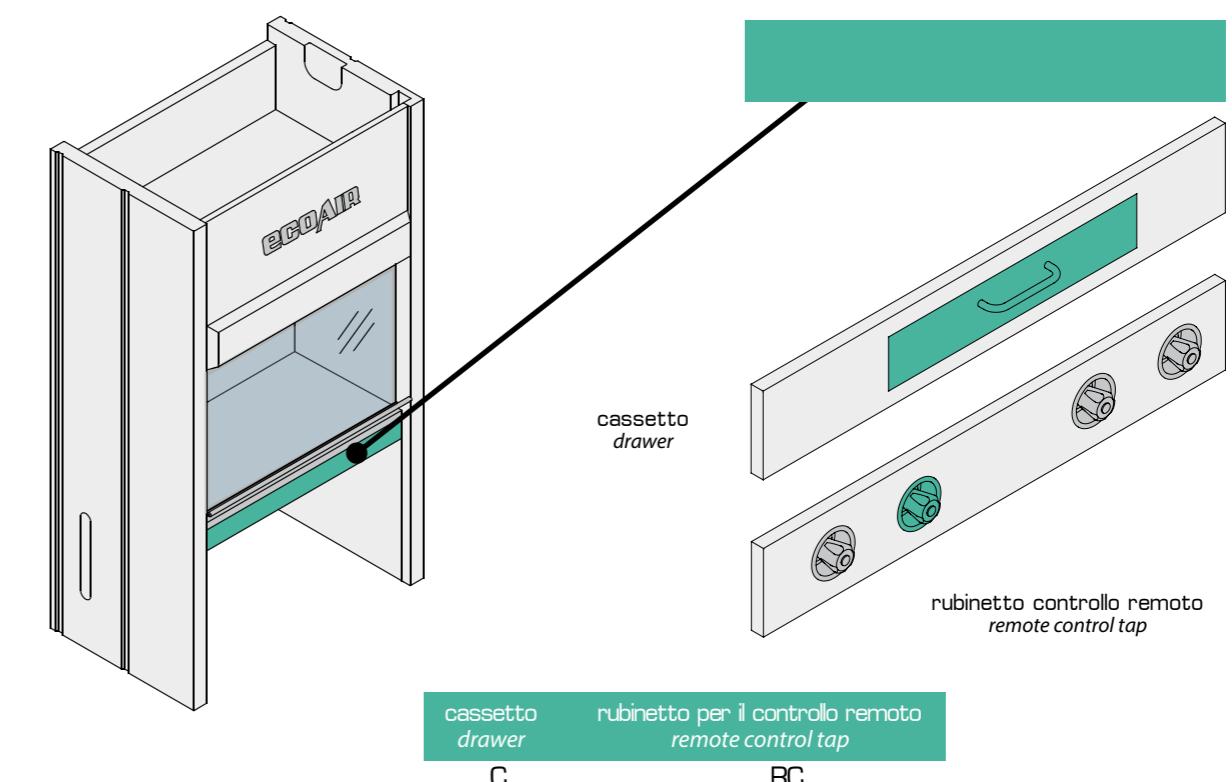
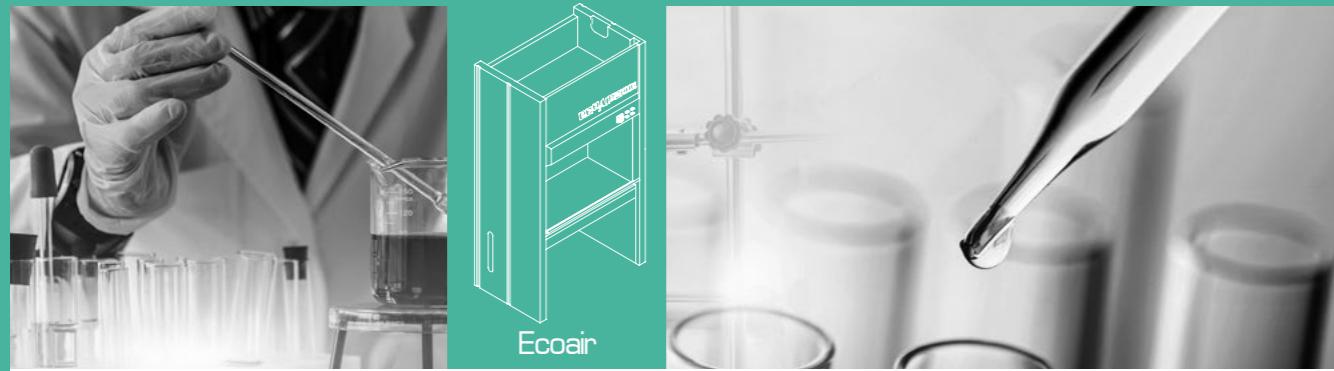
ECOAIR is provided with the following items on the top panel: MCB, master control switch and control switches for services.

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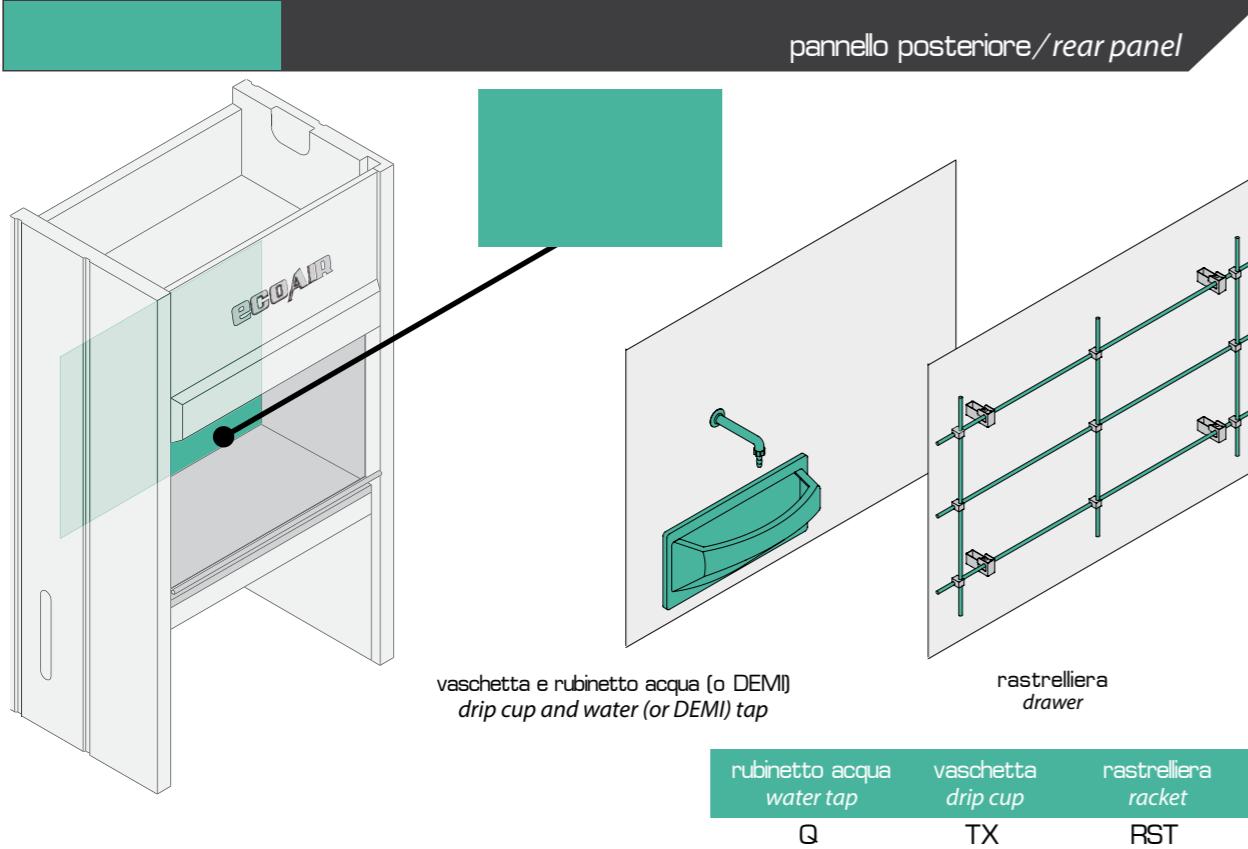
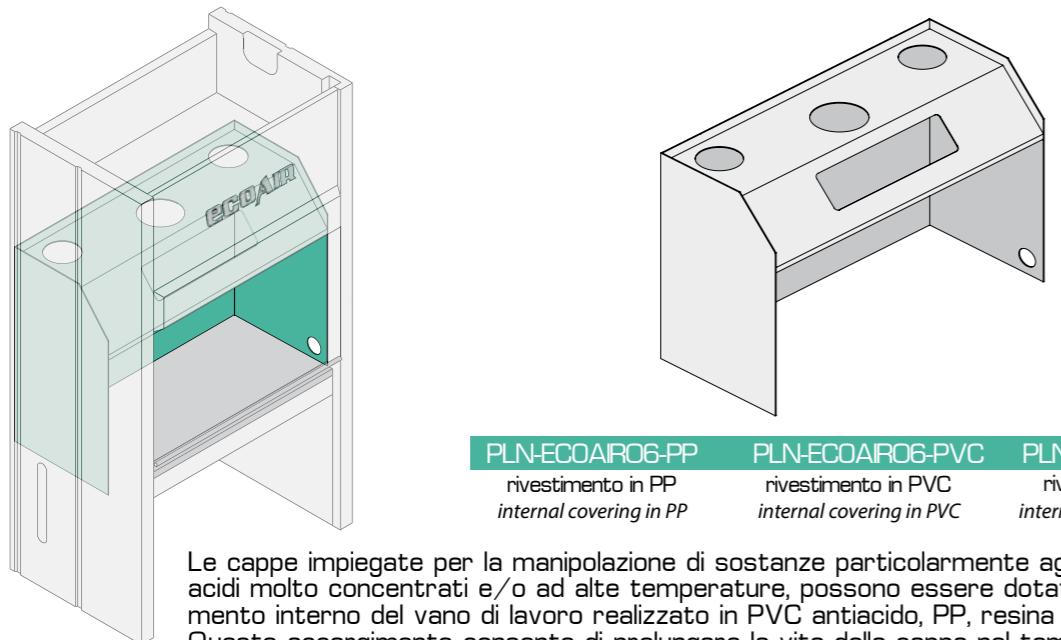
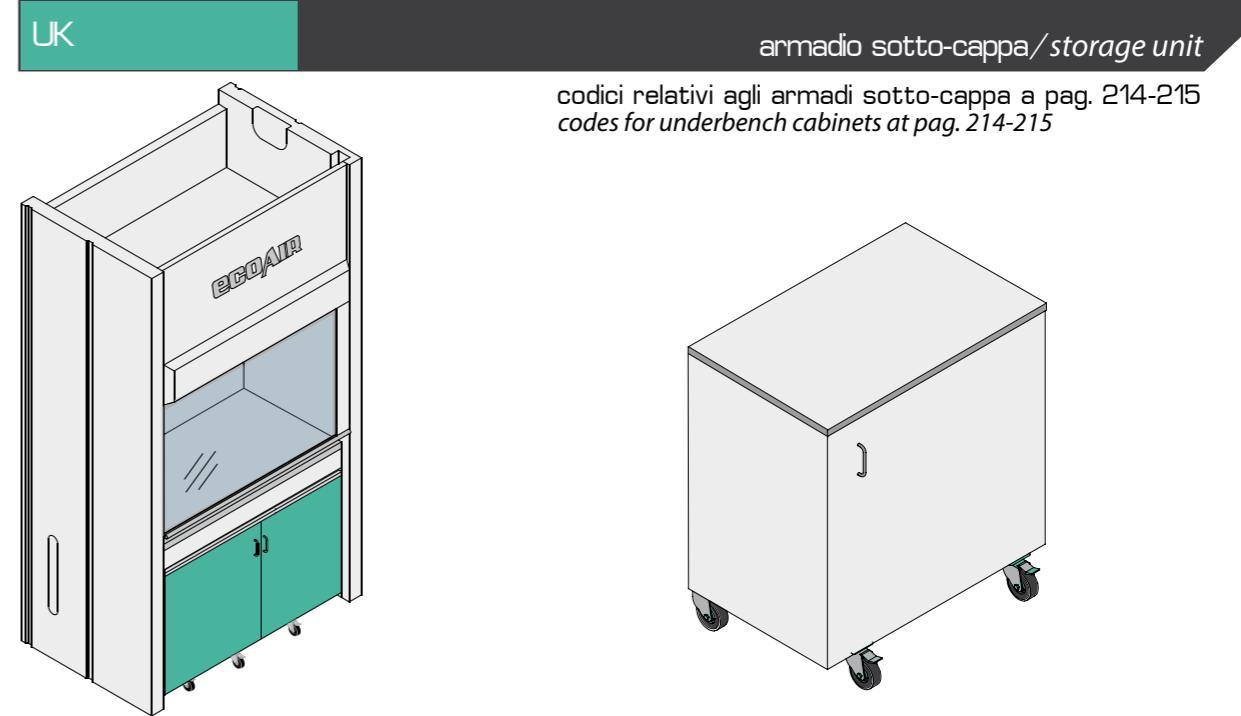
In case the fume cupboard has to be employed for manipulation of very aggressive chemicals (such as highly concentrated acids or acids at high temperature), a special internal covering can be provided. The internal covering panels can be made of anti-acid PVC, PP, epoxy resin, ceramic, and protect the structure of the fume cupboard from corrosion.

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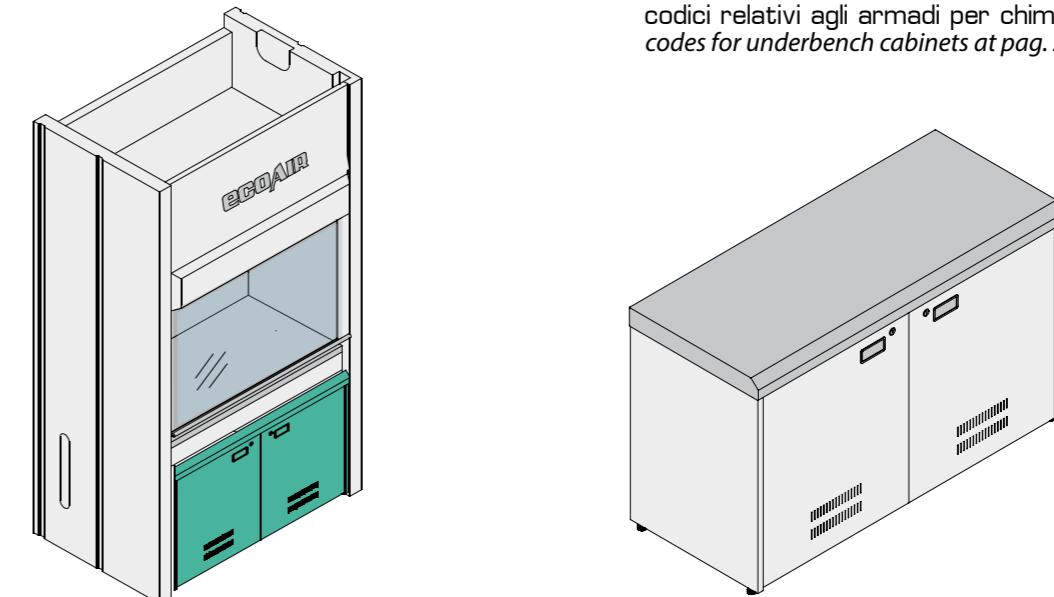
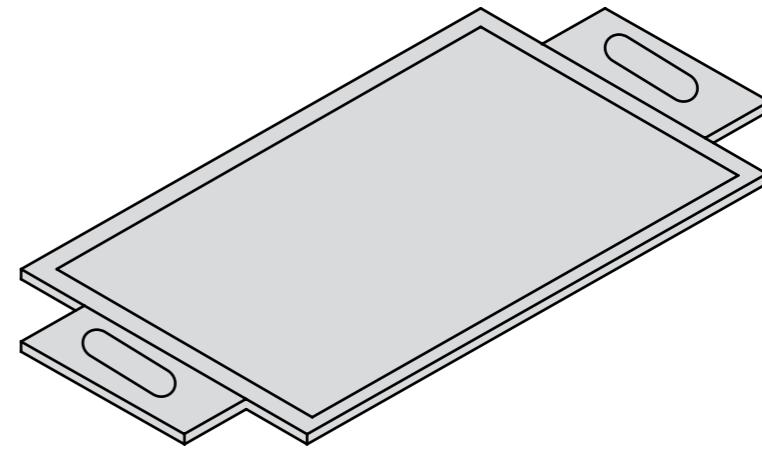
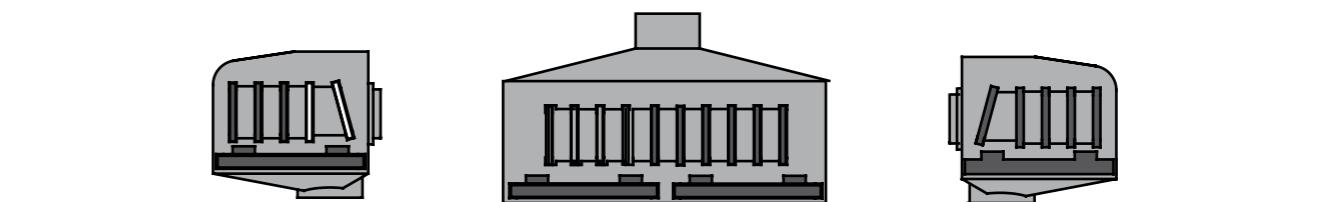
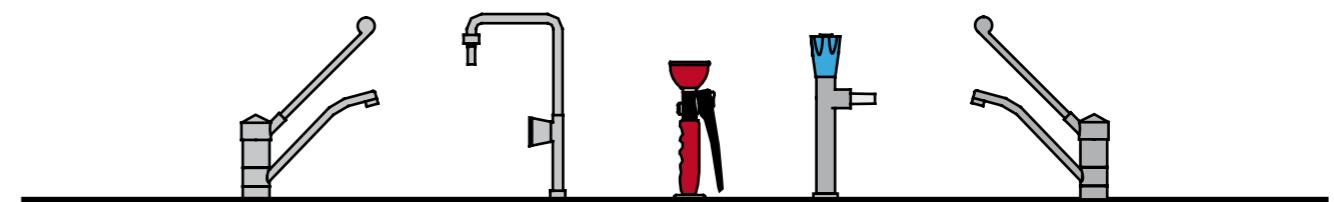
*The high flexibility and modularity of **MOMOLINE** production line allow to equip all fume cupboards with many different accessories and items, in order to meet all customer's wishes.*

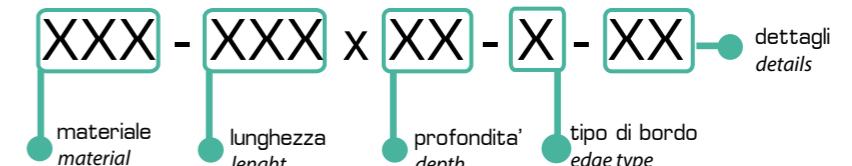
Accessories reported here are the standard ones, but custom-made and specially designed solutions can be provided upon specific request, so that every customer can fulfill its own desires and expectations from its new fume cupboard.

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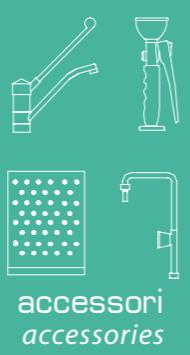
● idoneo
suitable ● poco idoneo
not recommended ● non idoneo
unsuitable

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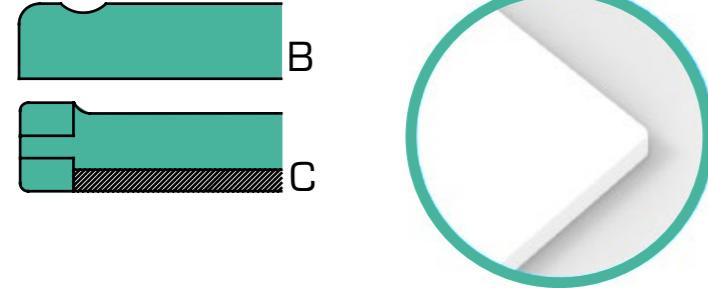
High Pressure Laminate worktops (such as Trespa) are composed by cellulose fiber layers, soaked in phenolic resin and then heavily compressed through mechanical presses. The surface of the panel is covered with cellulose fiber soaked in melaminic resin with wide choice of color. For aesthetic reasons the external surface can be specifically treated in order to obtain high resistance to chemicals, heat and scratches. Upon specific request, the phenolic resin can be additioned of special substances able to provide the worktop with flame retardant properties.

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Corian is a synthetic material, composed by 1/3 of acrylic resin and by 2/3 of natural minerals.

Corian has an incredibly uniform and homogeneous structure, does not delaminate, is not porous and is completely hygienic and stain-resistant. Corian also resists high temperature (up to 100°C) and most common chemicals.

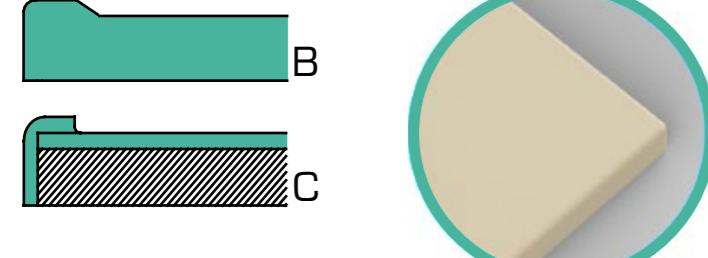
This material has excellent decontamination performances against viruses and bacteria, and its particular junction method allows to obtain virtually infinite surfaces, with no evident junction grooves. For this reason, it is particularly suitable for microbiology and high hygiene applications.

MOMOLINE is part of Corian Quality Network, meaning all the products are officially recognized and guaranteed by DuPont, sole authorized manufacturer of Corian.

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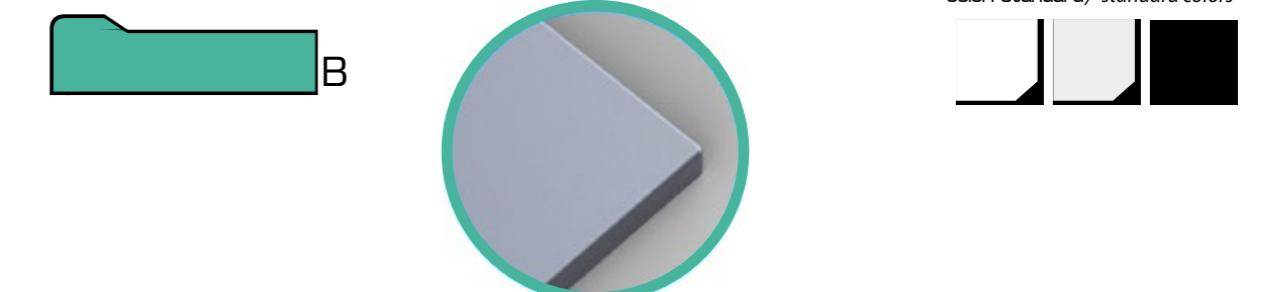
Polypropylene worktops are injection moulded and stand out for their extreme lightness.

PP worktops provide very good resistance against acids, bases and alkalis.

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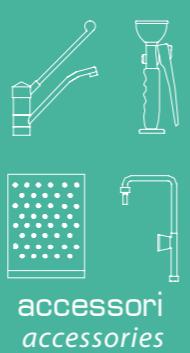
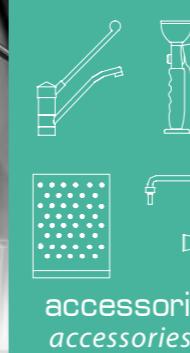
Epoxy resin worktops (such as Durcon) are obtained from molding and then properly shaped using CNC (Computer Numerical Control) machines.

These worktops are very easy to clean and decontaminate, are perfectly smooth and non-porous, very resistant to stains and to laboratory reagents. Together with ceramic worktops, they represent one of the best choices when it comes to employment with very aggressive acids.

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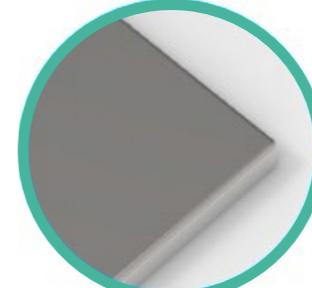
Enamelled steel is obtained by applying special ceramic enamel on a steel substratum. These worktops provide good resistance to most of common chemicals.

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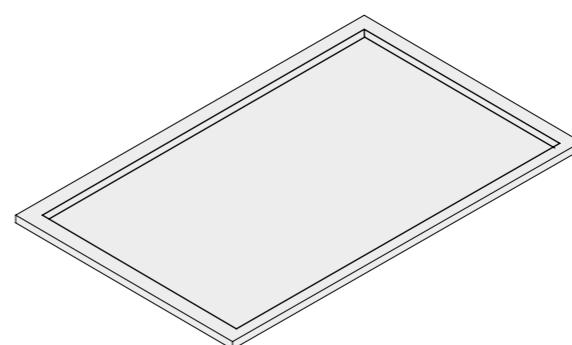
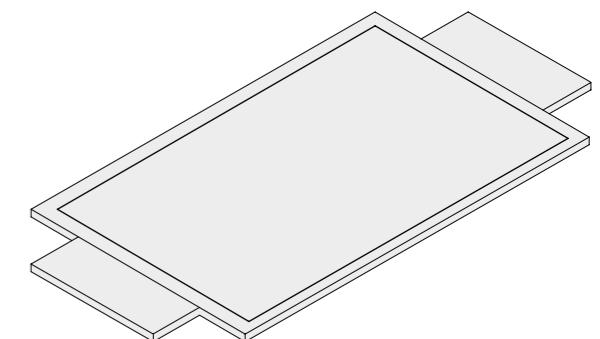
Ceramic worktops are manufactured through sintering of clay, feldspar and quartz, at a temperature higher than 1200°C. The worktops are perfectly seamless, so that an unique flat surface is obtained, with good resistance against mechanical shocks.

Ceramic worktops are recommended for all applications where a superior resistance to chemicals and heat is required.

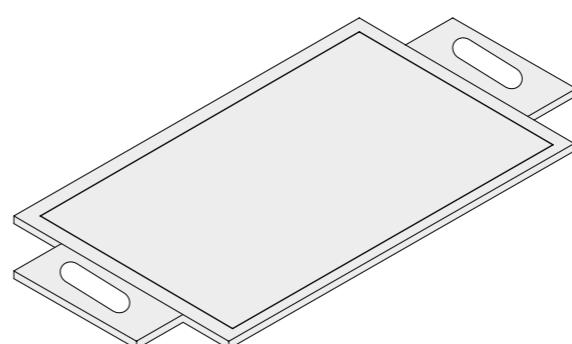
X4

bordo antidebordante 4 lati
4 sides rised edge

XW

ali laterali
lateral wings

XWV

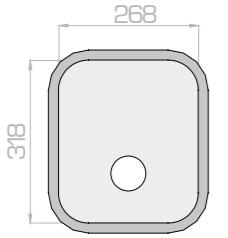
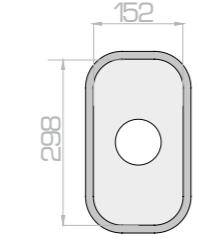
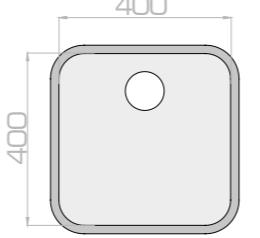
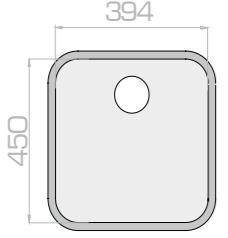
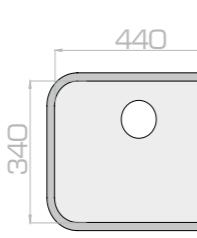
ali laterali con vasca
lateral wings with drip cup

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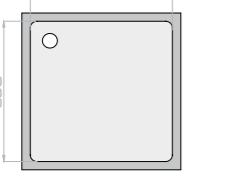
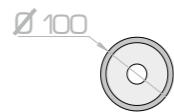
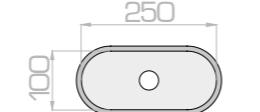
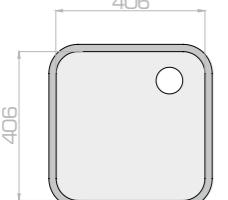
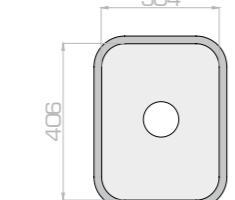
VSC-DC-____

vasche in Resina Epossidica/Epoxy Resin sinks

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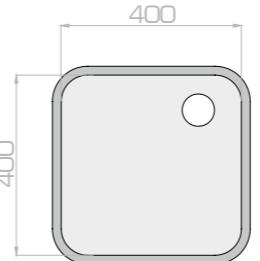
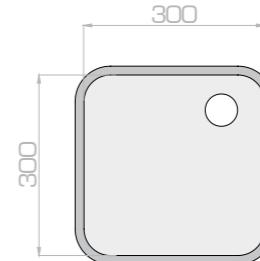
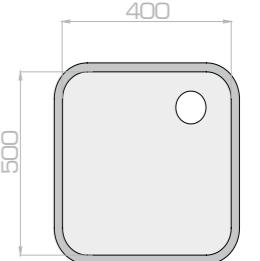
VSC-SS-____

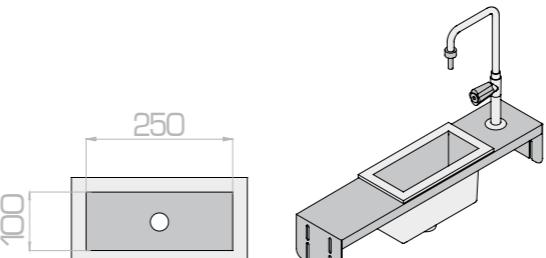
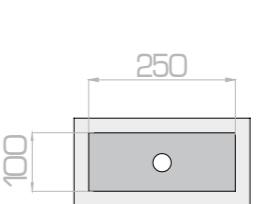
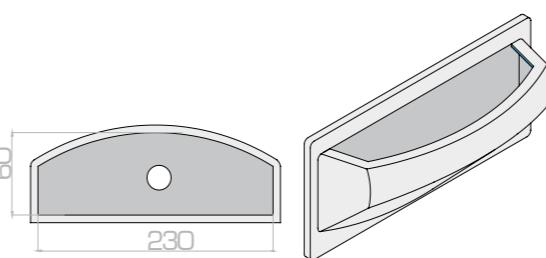
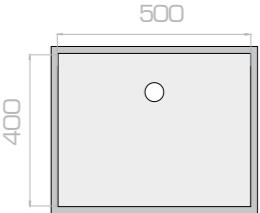
vasche in Acciaio inox/Stainless steel sinks

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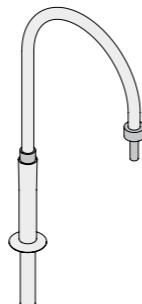
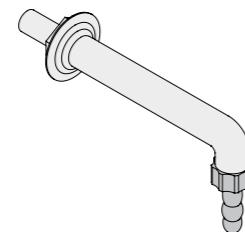
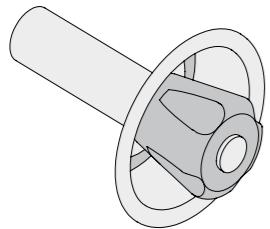
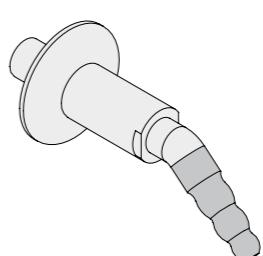
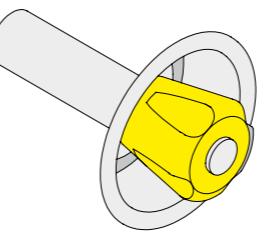
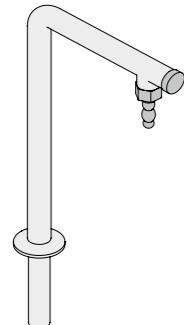
codice/code

Q-PP-2

beccuccio da piano in PP
PP worktop nozzle

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*All taps are manufactured according to DIN 12918 norm.**The threaded demountable nozzle is manufactured according to DIN 12898 norm.**All handles are made of anti-acid polypropylene, and the colors comply to EN 13792 norm.**Maximun operating pressure is 7 bar. Temperature operating range 30°C to 150°C.*

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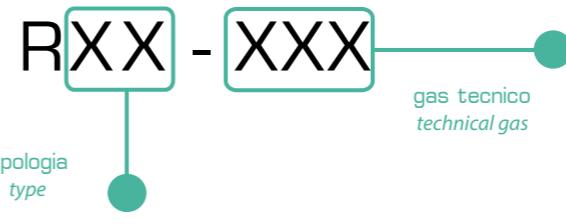
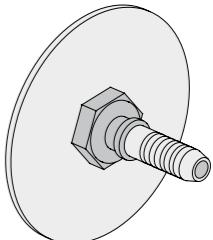
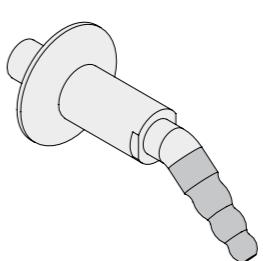
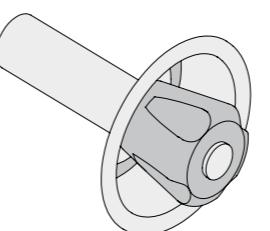
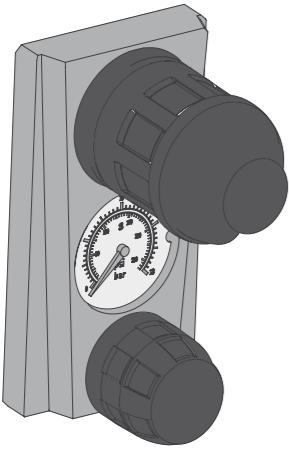
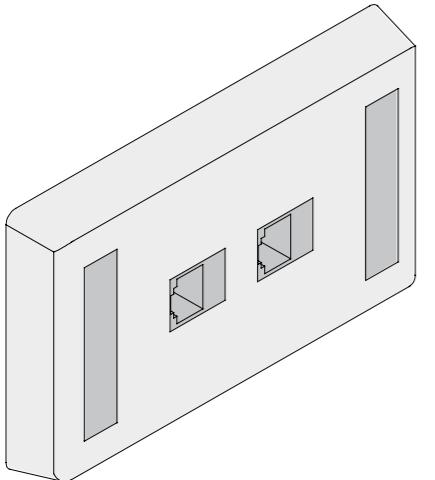
*Single stage pressure reducer for laboratory.
Flow pressure gauge compliant to EN 837-1 norm,
embedded needle type valves on outlet NY66.
Temperature operating range : -20/+60°C.*

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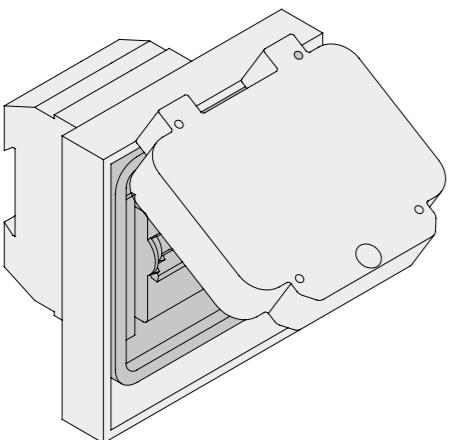
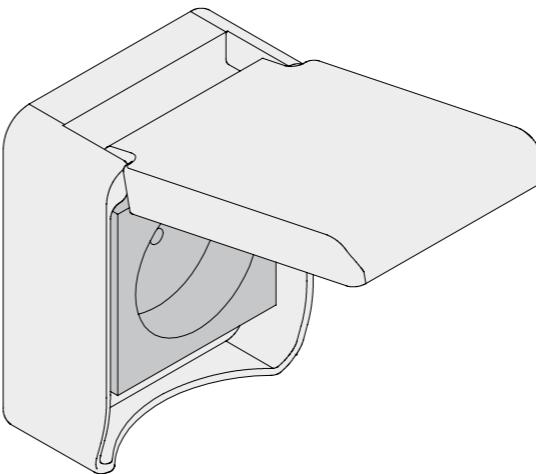
*The provided MCB is certified by CE and IMQ for compliance to CEI 23-3(EN60898) and EN 60947-2 norms.
Intervention curve C.
16A, 230V.
Interrupting capacity: 4,5kA a 230Vca (EN 60898).*

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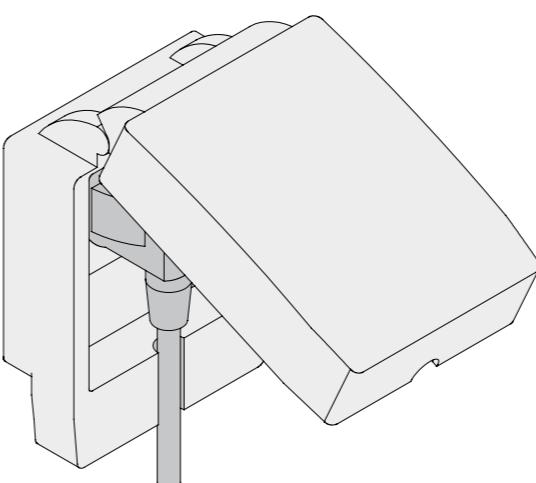
*IP65 Power outlet protection grade, according to IEC 60529 norm, IK 09 protection grade against mechanical shocks, according to IEC 62262 norm; heat resistance 850°C, according to IEC 60695-2-1 norm.
10/16A, 250V.
Different socket standards available.*

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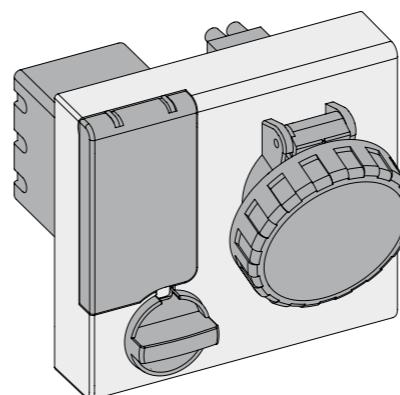
*IP66 Power outlet with fitted plug protection grade, according to IEC 60529 norm, IK 08 protection grade against mechanical shocks, according to IEC 62262 norm; heat resistance 850°C, according to IEC 60695-2-1 norm.
10/16A, 250V.
Different socket standards available.*

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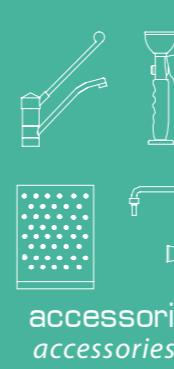
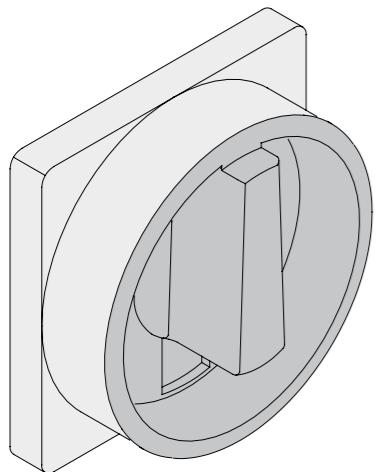
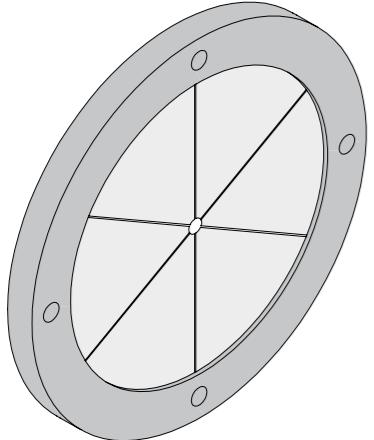
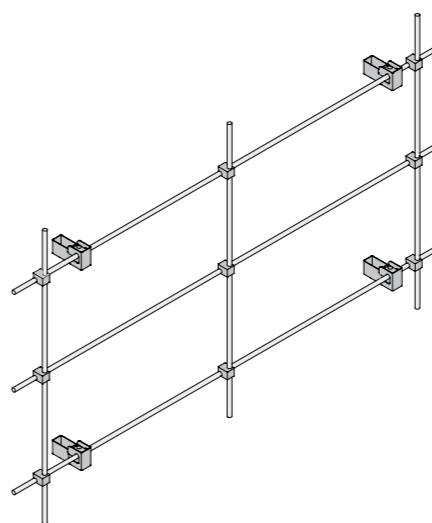
The cable grommet has a 100mm diameter, and can be used to connect all the equipments/instruments under the fume cupboard to a power supply, not interfering with the sash movement and with the airflow patterns. The cable grommet is entirely realized in soft anti-acid PVC, in order to grant a confortable and safe employmet, as well as to keep the containment performances unvaried.

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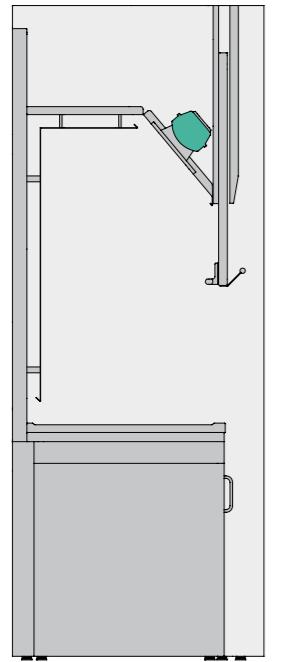
Stainless steel rack made up of properly shaped bars and fastened to the rear of the fume cupboard. The rack can be used for hanging/fastening objects or instruments, such as distillation equipment.

codice / code*	cappa da / cupboard length
RST-90	900
RST-120	1200
RST-150	1500
RST-180	1800
RST-220	2200
RST-240	2400

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Lamps are positioned outside the working area, so that the risk of triggering potentially explosive vapors is removed. The lamps are indeed separated from the working area by a dedicated wired safety glass.

Standard illumination makes use of a IP65 water-proof lamp with 2x18 W neon tubes having suitable power for granting approved enlightenment.

Upon specific request, other items can integrate or substitute the standard items:

- IP65 water-proof lamp with LED lamps;
- further adjustable light;
- UV lamp with timer.
- ex-proof lamp.

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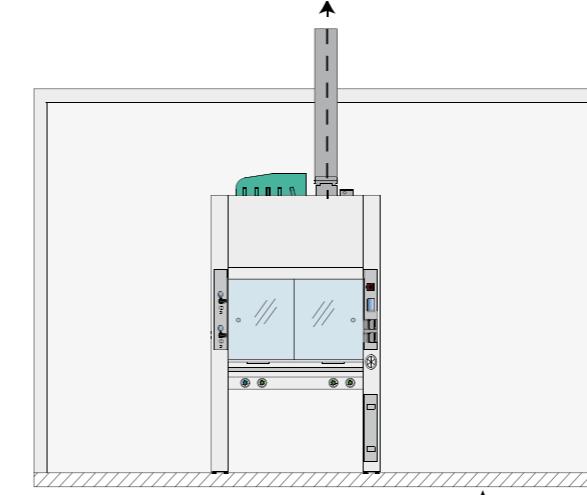
In CARBOCHEM model the lamps are positioned outside the working area, so that the risk of triggering potentially explosive vapors is removed. The lamps are indeed separated from the working area by a dedicated wired safety glass.

Standard illumination makes use of two led lamps 220V, 50Hz, IP65, 4500 K. VDE certification, CE compliant.

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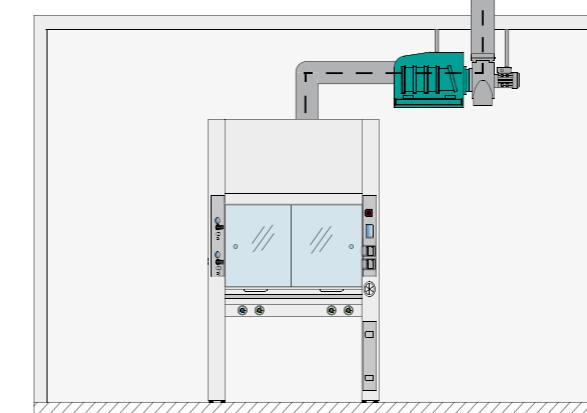
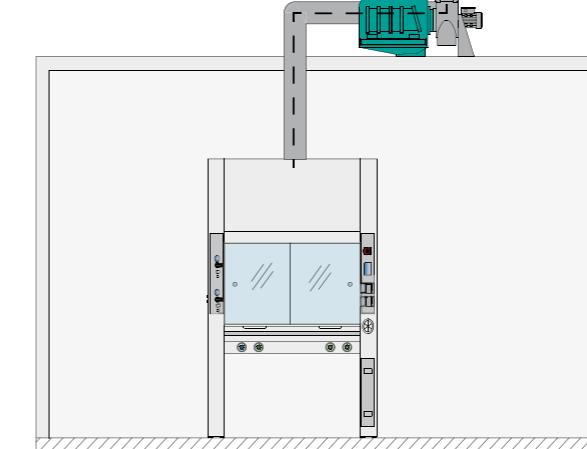
All hood models can be equipped, upon request, with filter boxes on board, wall/ceiling, or in a remote location outside the building.

The filter box is selected according to customer requirements and installation characteristics.

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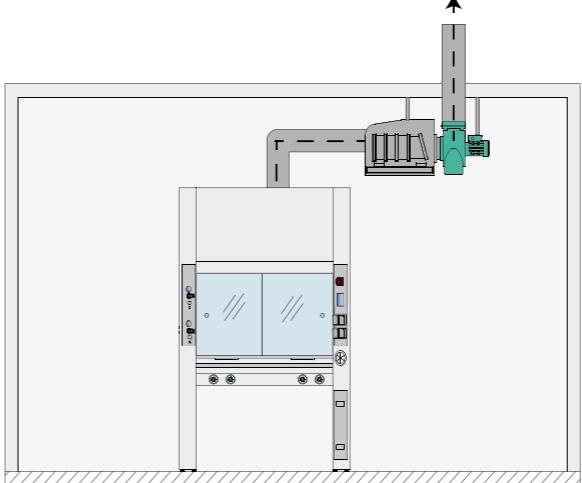
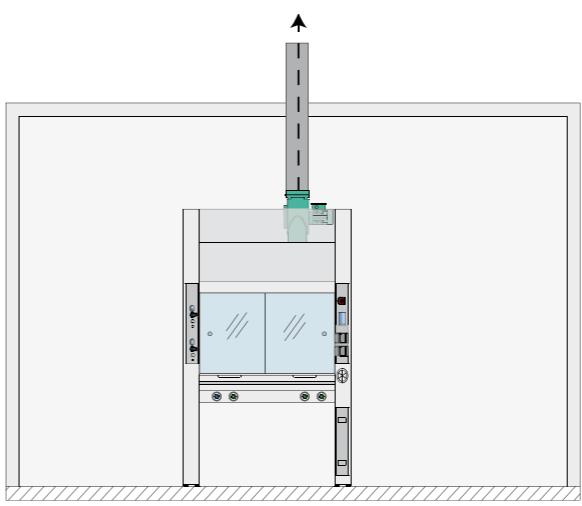
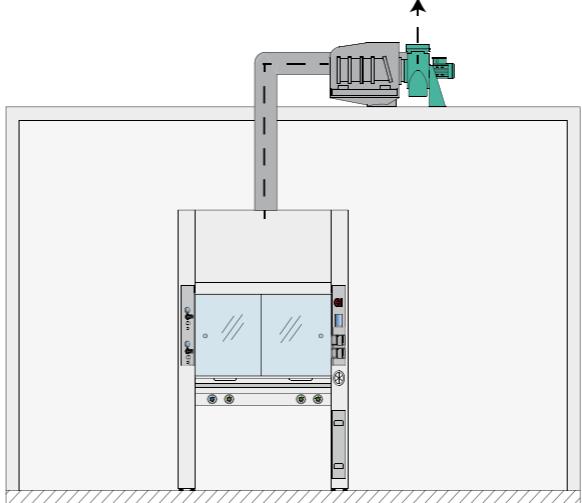
Exhaust fan is chosen according to the technical feature of the specific fume cupboard installed as well as on the base of the exhaust duct. The aspiration and containment parameters recommended by international norms are always guaranteed. ATEX (ex-proof) extraction system is available on request.

In case of blower installation outside the building, is available a kit inclusive of anchoring base, anti-weather covering, anti-birds net, IP65 electric breaker for maintenance operations (code ASP-EST).

ASP

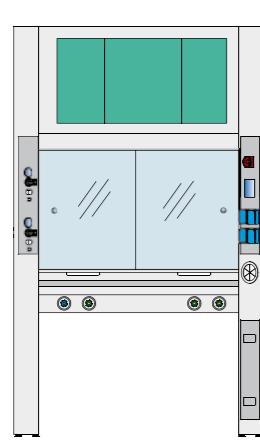
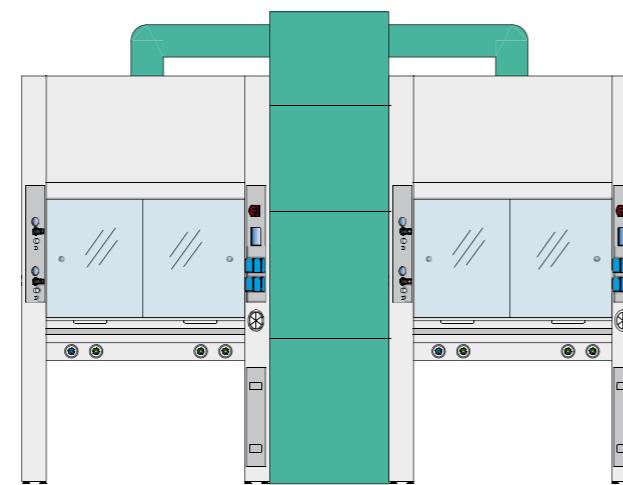
elettro-aspiratore / exhaust fan

codice / code
ASP-ONBmotore a bordo cappa
exhaust fan on the hoodcodice / code
ASP-WALLaspiratore a parete / soffitto
exhaust fan on the wall / ceilingcodice / code
ASP-ESTaspiratore all'esterno dell'edificio
exhaust fan outside the building

SCR

scrubber/scrubber

scrubber a bordo cappa
scrubber on the hoodscrubber a pavimento
floor scrubber

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The installation of a gas scrubber is recommended for use in laboratories where toxic, corrosive and odorous gases are released.

The laboratory gas scrubbers are used mainly for the absorption of strong acids. Inside the scrubber the contaminated exhaust air is brought into contact with an aqueous absorption solution.

To achieve a high efficiency it is important to ensure a large interfacial area between the air and the liquid. This is achieved by distributing the liquid over a packed bed.

All scrubbers are equipped with a level control. If the liquid falls below a certain level, fresh water is added automatically. The electrical conductivity of the washing solution is measured. If the adjustable set value is exceeded, a portion of the washing solution is discharged and replaced by fresh water.

Is also available with a pH measuring and control facility. By controlled addition of caustic soda or acid, the pH-value of the washing solution is kept at a constant value. This optimizes the performance of the scrubber.

Gas scrubber can be installed on the top of the fume cupboard, for dedicated employment, or on the floor, besides the fume cupboard.

Floor installation allows to increase the filtered airflow and is often employed for connecting more fume hoods to the same scrubber, optimizing the installation.

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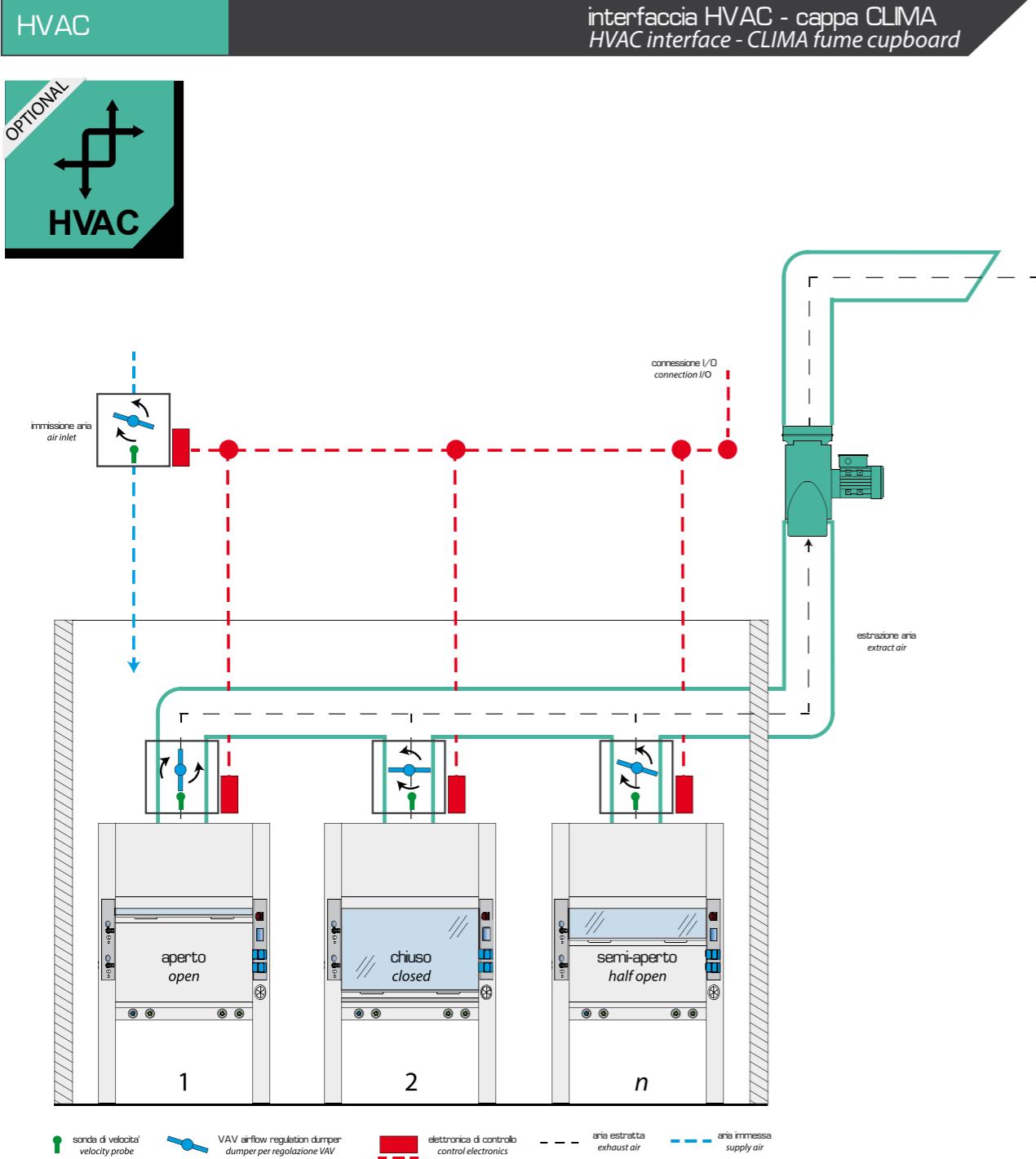
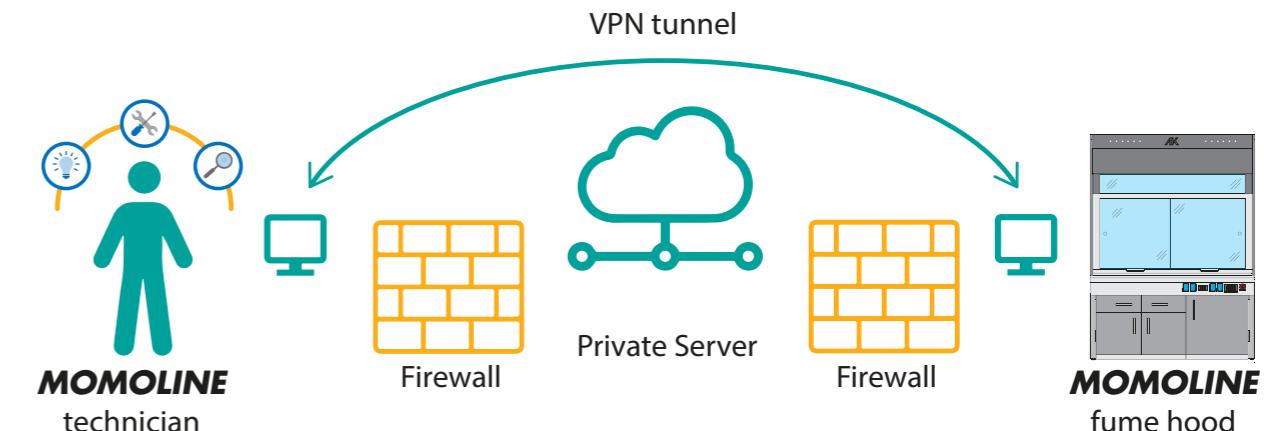
The control system of fume cupboard **BELAIRS6** can be connected to the air conditioning system (HVAC) of the building. The interface grants an optimal management of inlet and outlet airflows, ensuring at the same time the efficiency of the fume cupboards and an optimization of conditioned air volumes.

This solution can determine substantial energy savings.

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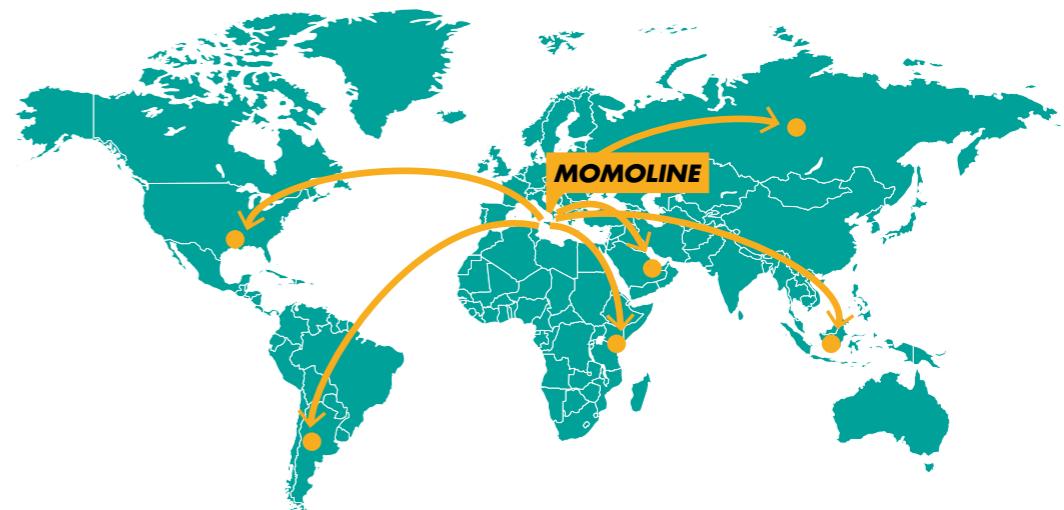
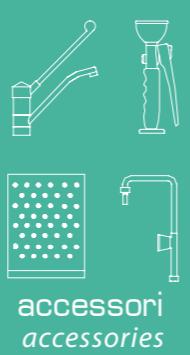
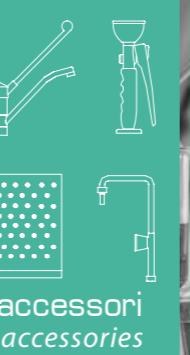
Fume hoods can be equipped with a remote assistance system to manage post-sales services such as technical assistance, software updating, maintenance, debug.

This solution enables **MOMOLINE** technicians to access the remote device (the single fume hood) connected to the network through a VPN based on proprietary technology having procedures with end-to-end sessions without any intermediate and optimized for industrial communication. Reliability and service continuity are granted thanks to a redundant and distributed server infrastructure.

The server infrastructure is globally distributed (Europe, United States, South America, Asia).

The system does not require additional hardware and allows technicians to operate on remote devices (after the authorization of the customer) eliminating the distance, reducing the need for on-site assistance, dramatically reducing postsale service costs.

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Within its facilities **MOMOLINE** has a modern test laboratory for fume cupboards realized according-
ly to EN14175-3:2019 and fully equipped to carry out all tests from standard EN and ASHRAE.

All models of fume cupboards sold by **MOMOLINE** have been tested (type test), and test results have been validated by Bureau Veritas, who has therefore issued third-party certification.

Upon request, specific tests on the single fume cupboard can be executed before the supply and test re-
port accordingly to EN14175 can be provided.

MOMOLINE is also certified by a third party according to EN14175-4 for the execution of in situ tests with the issuance of test reports in accordance with EN 14175-3:2019.

mobili-cabinets

indice • table of contents

mobili sottopiano / *underbench cabinets*

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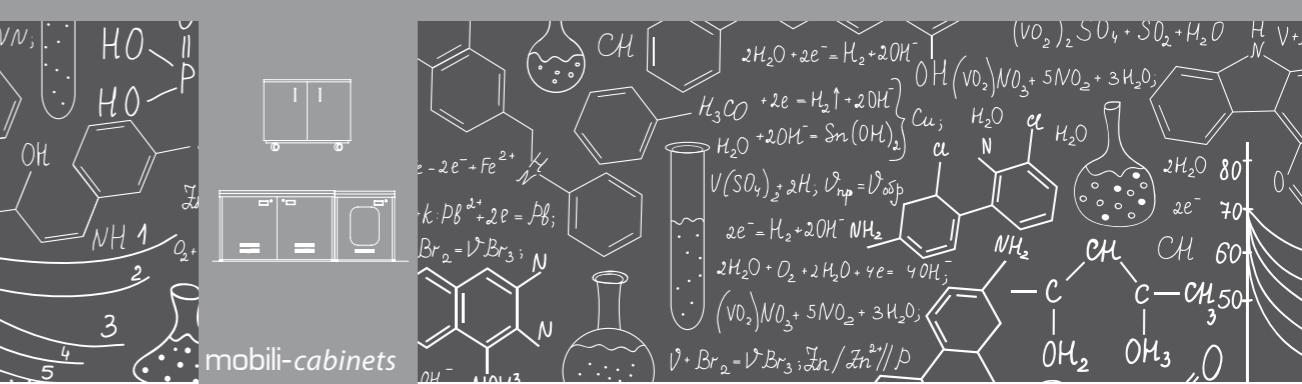
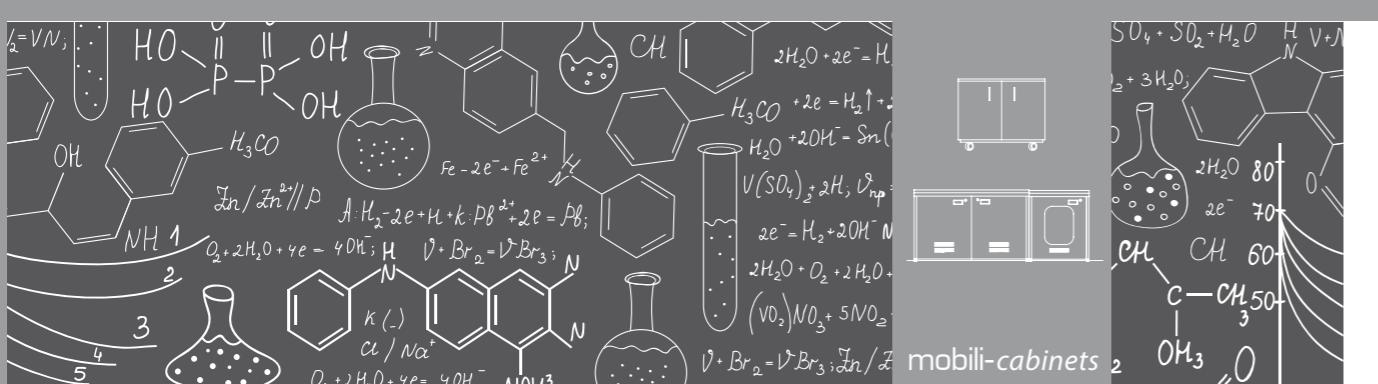
*Cabinets can be installed under all the lines of benches and hoods produced by **MOMOLINE**, or as stand-alone units.*

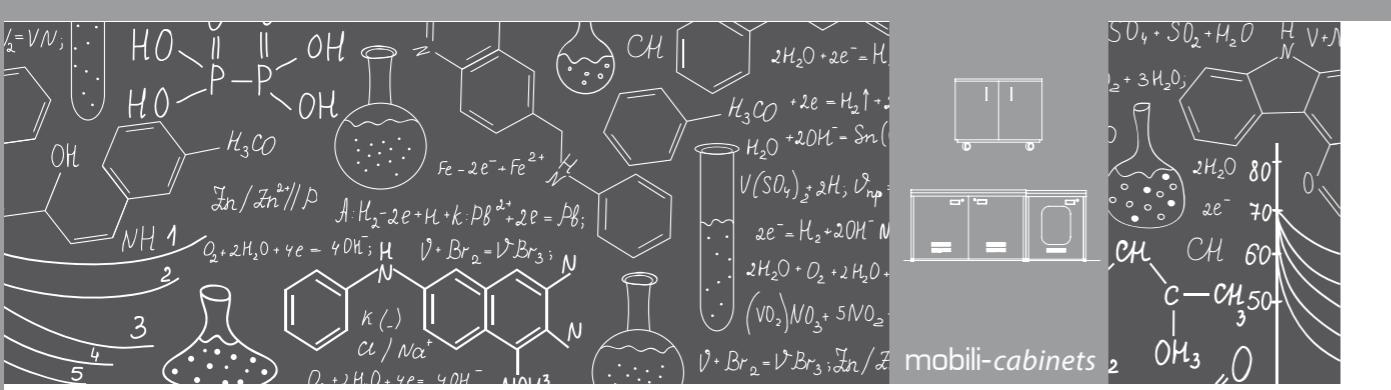
Cabinets can be realized in different versions and dimensions, with doors and/or drawers, and can be made either in wood or in metal, as better explained in this Paragraph.

Wooden cabinets are manufactured using bilaminated, chemical resistant and fire retardant panels (B-s1, d0 Fire Reaction Class, according to EN 13501-1:2009), obtained from 100% reclaimed wood, and marked E1 class for the low emission of formaldehyde.

The supporting base of wooden cabinets is realized using epoxy powder painted steel sheets, in order to grant further solidity and reliability to its structure and a better resistance against corrosion and fire.

Metal cabinets are manufactured using epoxy powder coated steel sheets, having a minimum thickness of 1 mm, bended and then coupled together to obtain a solid and reliable structure. Metal cabinets are classified as 'Class O' fire reaction, so that they are absolutely suitable for employment in every kind of laboratory, especially where flammable substances have to be handled. The epoxy powder painting process grants high resistance against corrosion, in compliance to all European Norms.

Upon specific request, it is possible to manufacture metal cabinets completely made of Stainless Steel, suitable for particular environments where a superior grade of hygiene is required. All the components fitted to the cabinets (such as its handles, hinges, rails, castors) are approved for laboratory use, and offer high resistance against corrosion. All cabinets can be equipped with lock and key, and cabinets with castors can be equipped with anti-capsize device, for safety reasons.

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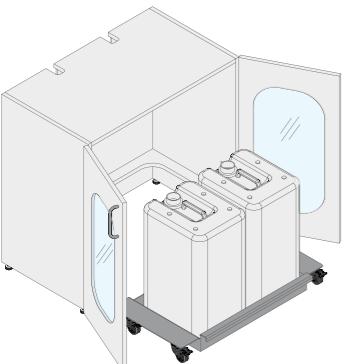
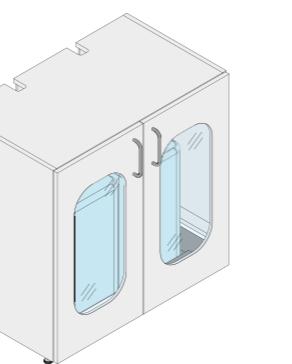
This special underbench unit is designed to allow the safe disposal and collection of chemical reagents or other waste liquids (such as oil).

Equipped with tank-holder trolley and one or more drainage collectors (funnels) above the worktop, it can also be equipped with overflow sensor/tank presence sensor (cod. SC-ALM).

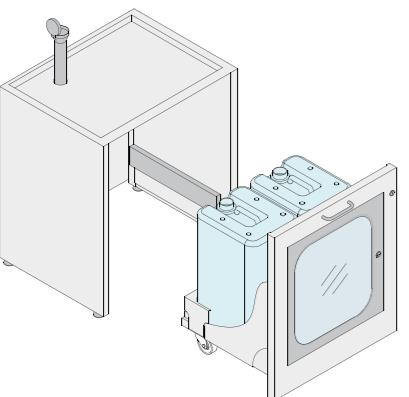
The unit can be realized in steel or HPL.

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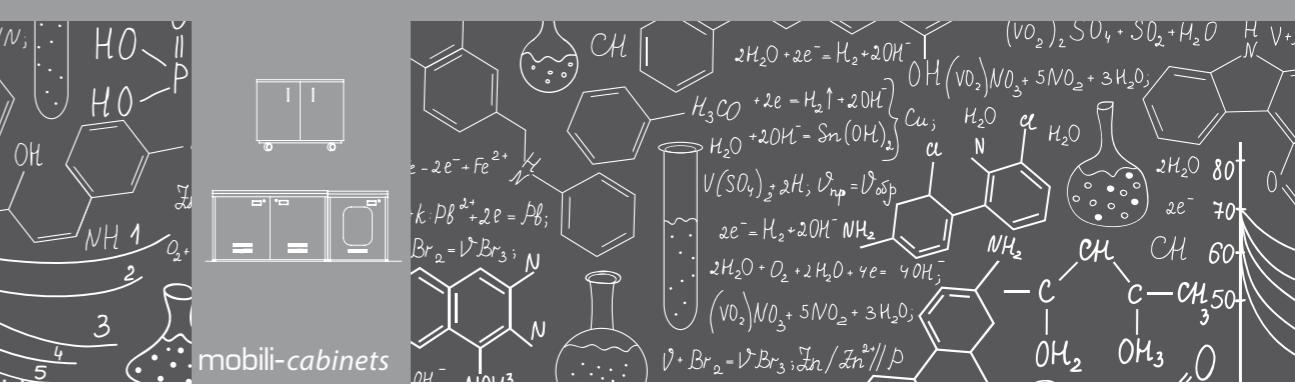
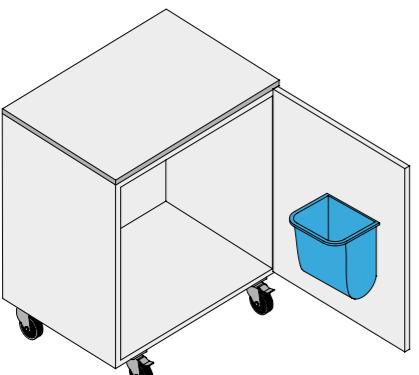
Plastic bin fastened to the door.

PG-___-WLL-

pensile a muro a giorno
open shelving overhead cabinet for wall fastening

profondita'
depth

400 mm

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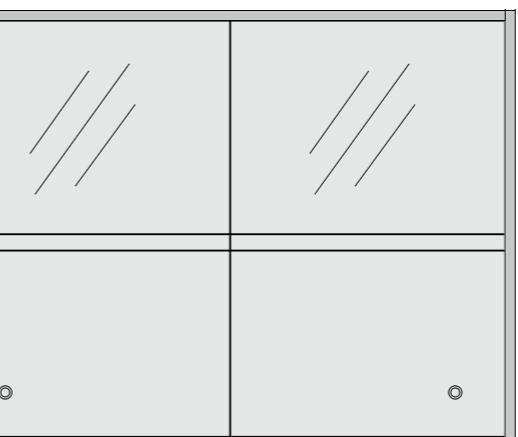
600/800 mm

codice/code per banco da/for bench length

PG60-WLL-X	600
PG80-WLL-X	800
PG90-WLL-X	900
PG100-WLL-X	1000
PG120-WLL-X	1200
PG150-WLL-X	1500
PG180-WLL-X	1800

PV-___-WLL-

pensile a muro con vetri scorrevoli
overhead cabinet with sliding glass doors for wall fastening

profondita'
depth

400 mm

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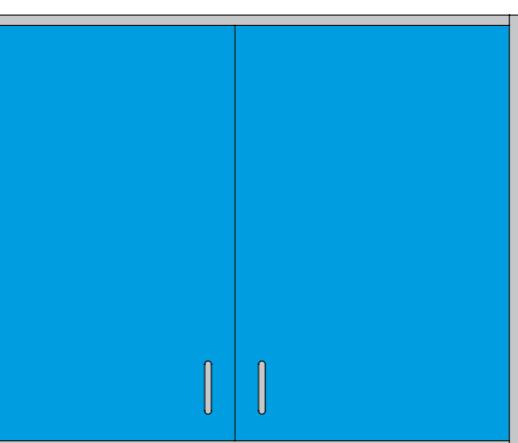
600/800 mm

codice/code per banco da/for bench length

PV60-WLL-X	600
PV80-WLL-X	800
PV90-WLL-X	900
PV100-WLL-X	1000
PV120-WLL-X	1200
PV150-WLL-X	1500
PV180-WLL-X	1800

PA-___-WLL-

pensile a muro con ante
overhead cabinet with doors for wall fastening

profondita'
depth

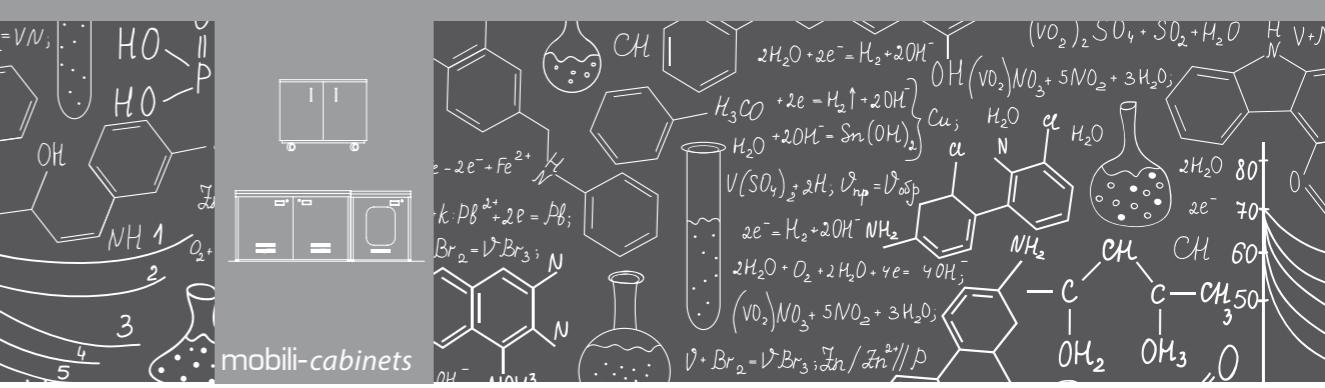
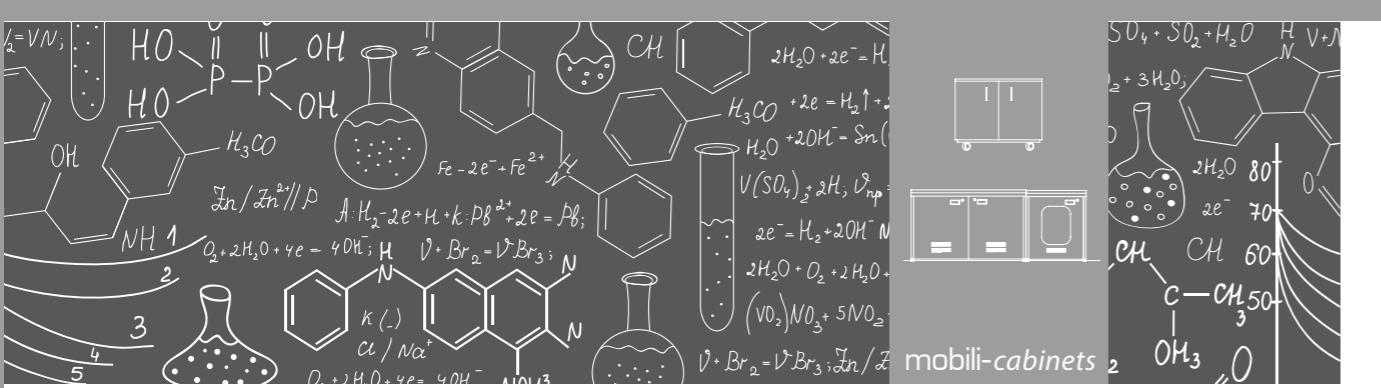
400 mm

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600/800 mm

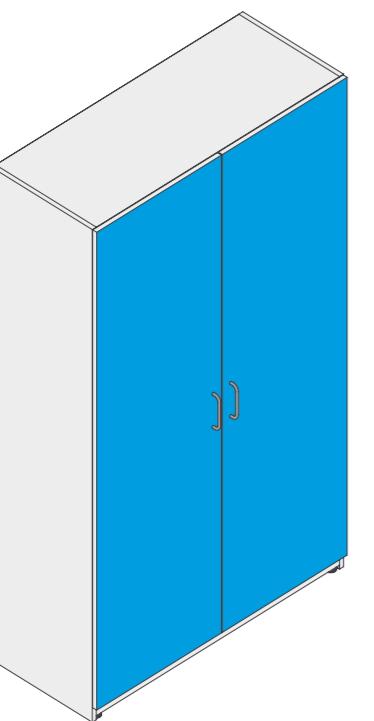
codice/code per banco da/for bench length

PA60-WLL-X	600
PA80-WLL-X	800
PA90-WLL-X	900
PA100-WLL-X	1000
PA120-WLL-X	1200
PA150-WLL-X	1500
PA180-WLL-X	1800

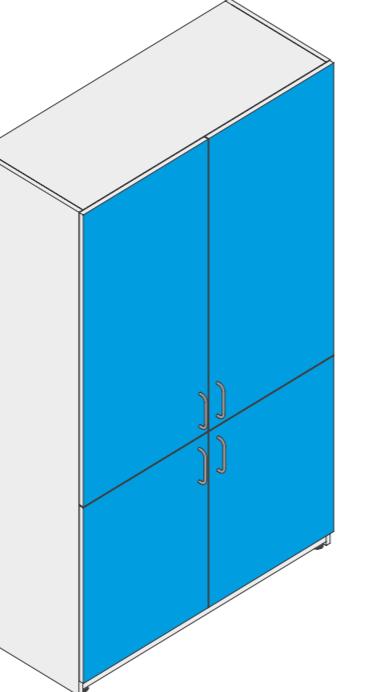
TALL-C

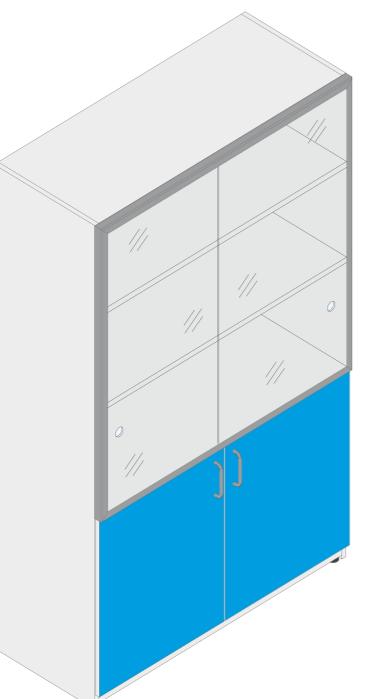
armadi alti/*tall cabinets*

codice code	lxpxh lxdxh
TALL-C-90-L/U	900X500X1900
TALL-C-120-L/U	1200X500X1900

codice
code lpxh
lxdhx
TALL-C-120-L/U-4A 1200X500X1900

codice <i>code</i>	lpxlh <i>lxdh</i>
TALL-C-90-V-L/U	900X500X1900
TALL-C-120-V-L/U	1200X500X1900

mobil
cabinets

3

OMO I

mobili
cabinets

MOMOLINE

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safety cabinets

indice • table of contents

armadi per reagenti chimici
cabinets for chemicals storage

222 - 225

armadi per reagenti infiammabili
cabinets for flammable storage

226 - 227

armadi per bombole gas
cabinets for gas cylinders storage

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Chemiguard is a wide line of safety cabinets designed for the storage of hazardous chemical substances in all laboratory environments. These cabinets are essential in all kind of laboratories in order to safely stock the necessary reagents, in full accordance to all law requirements.

Safety cabinets are available in many different sizes and configurations, and in different versions: structure in epoxy-powder coated steel sheets, in acid-proof Polypropylene sheets or in PVC. All Chemiguard cabinets are furthermore available in the following configurations: "BASE": Simple storage cabinets, with no filters nor air extraction system, but already set-up for connection to extraction systems; "ECO": Storage cabinets equipped with carbon filters, with no air extraction system, but already set-up for connection to extraction systems; "PLUS": Complete storage cabinet, equipped with carbon filters and independent air extraction system.

Main features:

- main structure with high chemical resistance insulating rubber feet allowing height adjustments
- internal sliding and height-adjustable shelves, tray-shaped to retain possible liquids' spillage
- bottom tray for additional safety liquids' spillage containment (available Stainless Steel upon request)
- front grid for fresh air intake and proper ventilation, with dust filter
- carefully designed and planned airflow patterns for best extraction efficiency
- possibility of separate storage for Acids and Bases (double door and internal partition wall), with possibility of separate filter boxes and separate extraction fans
- technical compartment for filter box and fan separate from storage area, and easily accessible for maintenance operations.
- switch to control fan operation, with MCB and thermal protection (only for metal cabinets)
- possibility of integration between Chemiguard safety cabinets for chemicals and a safety cabinet for flammables

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armadi per reagenti chimici / security cabinets for chemicals storage

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DÜPERTHAL as the German quality and technology leader, sets the standards in the field of system solutions for process-optimized storage of hazardous goods.

Type 90 safety cabinets meet the requirements of EU standards EN 14470 (Part 1 for the storage of flammable liquids, Part 2 for the storage of gas cylinders) and EN 16121 and 16122 (on the durability and longevity of container furniture) and therefore are also labelled with the GS mark. Furthermore, DÜPERTHAL safety cabinets have been awarded with the High Quality certificate by TÜV Süd for their durability, user-friendliness and upscale features.

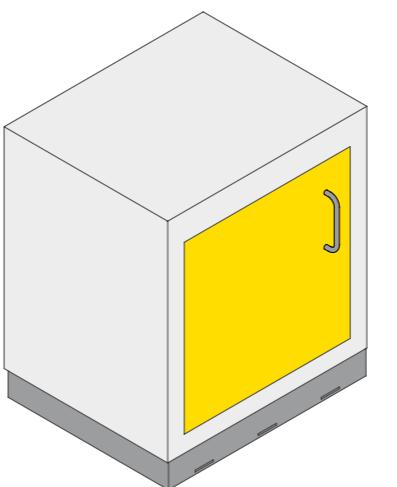
The cabinets are available in various sizes and equipment and are prepared factory-made for connection to an exhaust air system and the grounding. For maximum ease of use, the cylinder lock is located at handle height.

The following features are available for the storage of flammable hazardous materials:

- Hinged or folding doors
- Storage shelves or pull-out trays
- Single-handed operation or two separate door leaves
- Optional stainless steel interior

In case of fire, the doors are self-closing. All cabinets are equipped with a collection tray and the tall cabinets are also equipped as a standard with a base that can be underdriven.

codice code	lxpxh lxdhx
DUP-29-200667-030	594X612X2085
DUP-29-200967-030	894X500X1900
DUP-29-201267-030	1194X500X1900

codice code	lxpxh lxdhx
DUP-29-060657-001	601X493X631
DUP-29-060957-001	888X500X631
DUP-29-061157-001	1100X500X631
DUP-29-061457-001	1400X500X631

armadi per bombole gas / security cabinets for tanks storage

armadi per bombole gas da esterno
outdoor security cabinets for tanks storage

MOMOLINE

armadi per bombole gas da interno
indoor security cabinets for cylinders storage

GASG

armadi per bombole gas da esterno outdoor security cabinets for tanks storage

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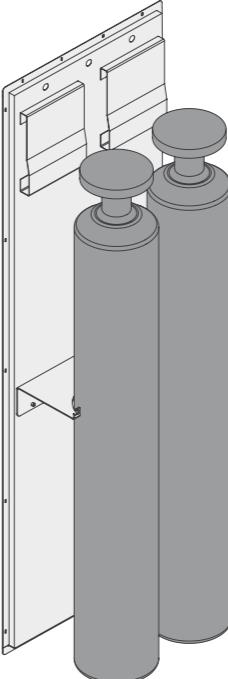
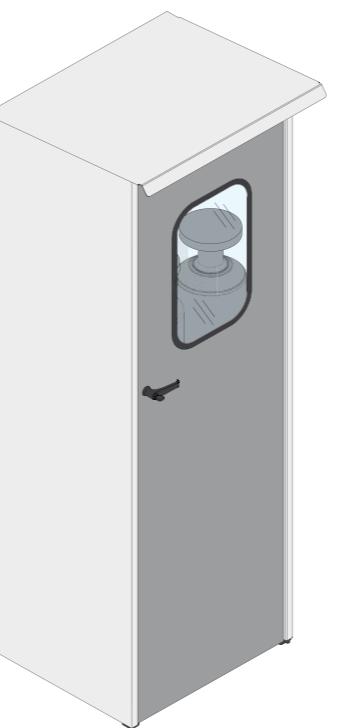
Gas Guard is a line of safety cabinets specifically designed for the storage of compressed gas cylinders. These cabinets are entirely realized using epoxy powder coated steel sheets, and can be specifically manufactured in AISI304/316 Stainless Steel upon request.

Main features:

- Adjustable leveling feet on the base
- Magnetic device granting full closure of the door
- Wing door with 110° opening hinges
- Closure handle with locker and key
- Housing slot for cylinders with fastening chain
- Pre-arranged fastening slot for manifold (1st stage pressure reducers)
- Extractable aluminum ramp allowing the loading of cylinders
- Air intake grids on the sides

Available Optional Items:

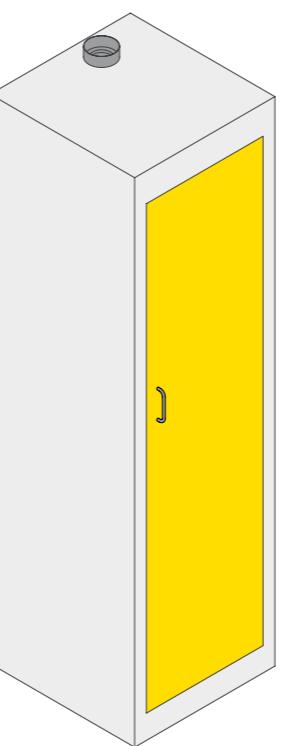
- Anti-weather covering (for outdoor installations)
- Transparent window on the door
- Manifold (1st stage pressure reducer)
- Flexible connection hose for cylinder

DUP

armadi per bombole gas da interno indoor security cabinets for tanks storage

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DÜPERTHAL Security cabinets for storage of pressurised gas cylinders include many practical features that facilitate the handling of the gas cylinders:

- Robust roll-in flap for easy insertion of the cylinders
- Individually adjustable door opening angle
- Standard adjustable mounting rails for the installation of fittings

These cabinets meet the requirements of EU standards EN 14470 Part 2 for the storage of gas cylinders.

	codice code	lxpxh lxpxh
1 bombola gas / 1 gas cylinders	DUP-79-200660-021	595X610X2070
2 bombole gas / 2 gas cylinders	DUP-79-200960-021	895X610X2070
3 bombole gas / 3 gas cylinders	DUP-79-201260-021	1195X610X2070
4 bombole gas / 4 gas cylinders	DUP-79-201460-021	1395X610X2070

Note

Notes

A blank 10x10 grid of black lines on a white background. The grid consists of 9 horizontal rows and 9 vertical columns, creating a total of 81 equal-sized squares. The lines are thin and black, with no additional markings or text.

Note

Notes

A blank grid consisting of 20 horizontal rows and 10 vertical columns, created by black lines on a white background. The grid is intended for a crossword puzzle, with empty squares available for letter placement.

graphic design by
Matilde Merciai